




making work easy

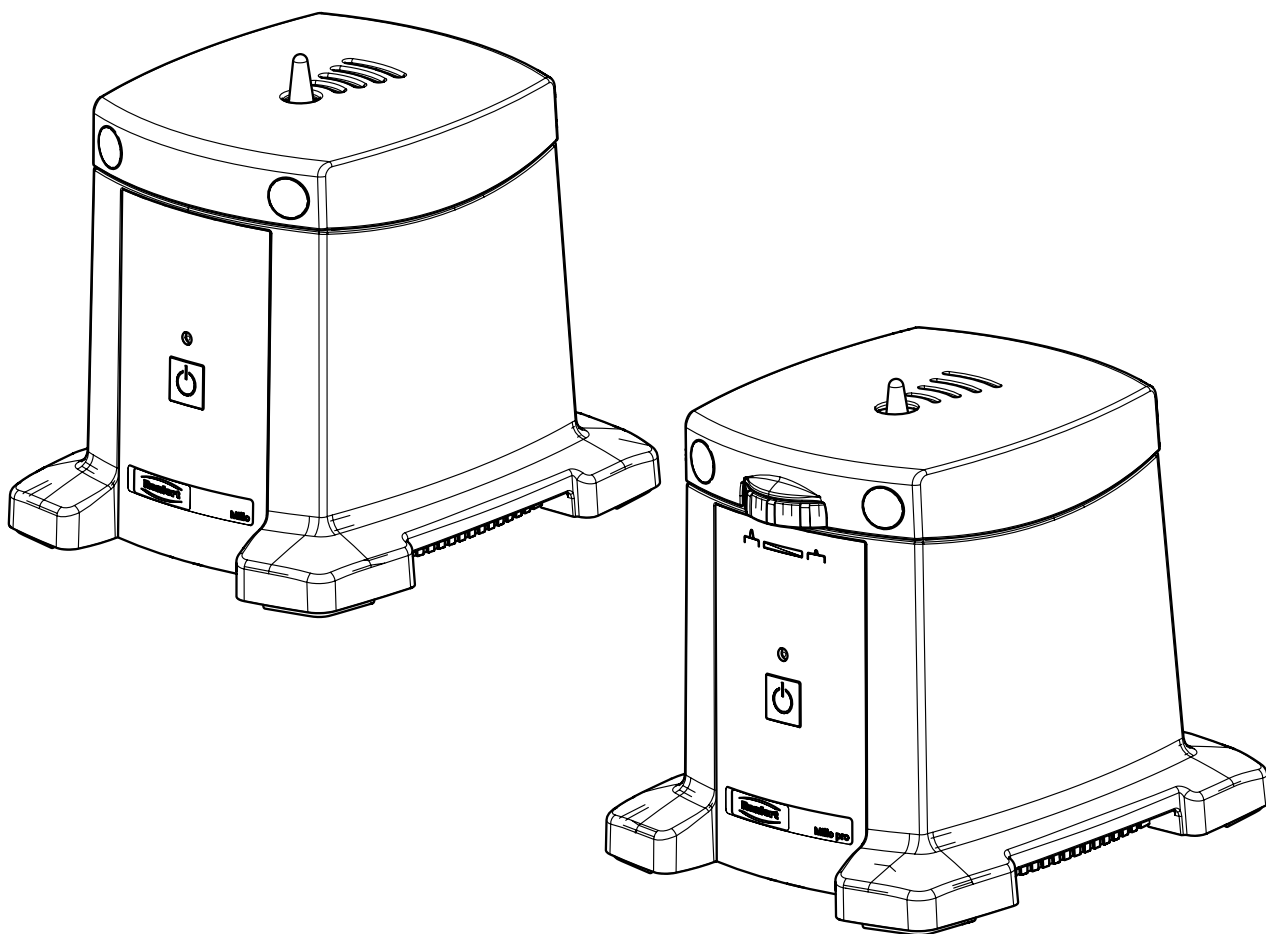

Millo / Millo pro

Service Instructions

Made in Germany

21-6849 05122025 / A

Contents

90003 3385

Spannbuchse • Clip socket • Douille de serrage • Boccola di serraggio
Casquillo de fijación • Зажимная втулка

90003 3428

Schutzkappe • Protection cap • Capuchon protecteur • Calotta protettiva
Caperuza protectora • Защитный колпачок

90003 3416 / 90013 4185

Sicherungsset • Set of fuses • Jeu de fusibles • Kit di fusibili • Juego de fusibles
Комплект предохранителей

90003 3430

Arbeitstisch Millo pro • Millo pro work table • Table de travail Millo pro
Piano di lavoro Millo pro • Mesa de trabajo Millo pro • Столик Millo pro

90003 3429

Arbeitstisch Millo • Millo work table • Table de travail Millo • Piano di lavoro Millo
Mesa de trabajo Millo • Столик Millo

90003 3552 / ..3 / ..5 / ..6 / ..7

Adapter Schuko • Schuko adapter • Adaptateur Schuko • Adattatore Schuko
Adaptador Schuko • Schuko Адаптер

90003 3425 / 90003 3426

Netzkabel • Power cable • Câble de réseau • Cavo di rete • Cable de red
Сетевой кабель

90003 3438 / 90003 3443

Leiterplatte • Circuit card • Carte imprimée • Circuito stampato • Placa de circuito
Электронная плата

90003 3409 / 90003 3410

Leiterplatte • Circuit card • Carte imprimée • Circuito stampato • Placa de circuito
Электронная плата

90012 1402 / 90012 1403

Motor • Engine • Moteur • Motore • Motor • Мотор

90002 1402 / 90002 1403

Motor • Engine • Moteur • Motore • Motor • Мотор

90003 3399 / 90003 3401

Motorkondensator • Engine condenser • Condensateur du moteur
Condensatore del motore • Condensador del motor • Конденсатор мотора

90003 3384

Gerätefüße • Rubber bases • Pieds de l'appareil • Piedini dell'apparecchio
Pies de l'aparato • Ножки прибора

90003 3386



Höhenverstellung kpl • Height adjustment complete • Régleur d'hauteur complet
Regolatore d'altezza completo • Ajusta de altura completo • Установка по высоте

90003 3405



Überstromschutzschalter • Overcurrent circuit breaker
Protecteur contre les surintensités • Protezione da sovratensioni
Protección de sobrecargas • Защитный выключатель при перегрузке тока

90003 3385


Spannbuchse • Clip socket • Douille de serrage • Boccola di serraggio
Casquillo de fijación • Зажимная втулка

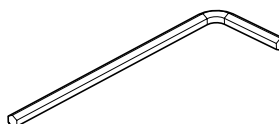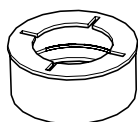


1 Millo / Millo pro

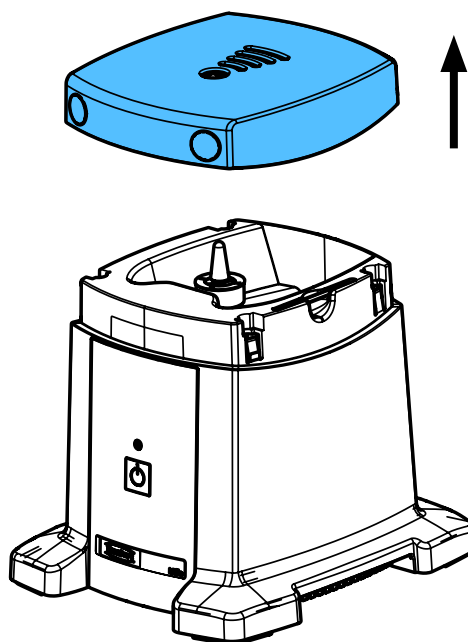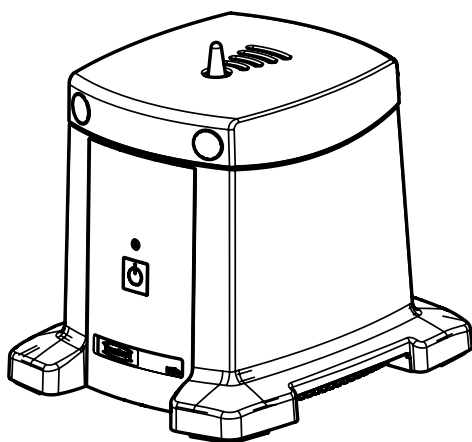


2

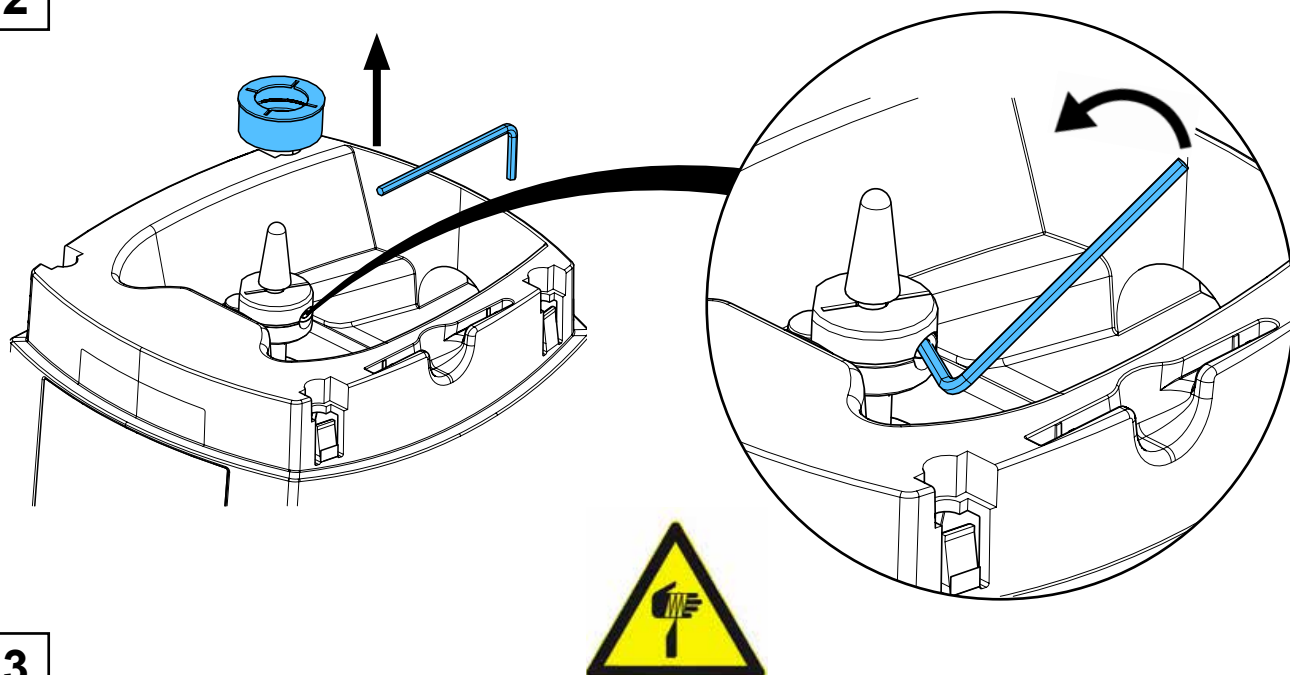




3

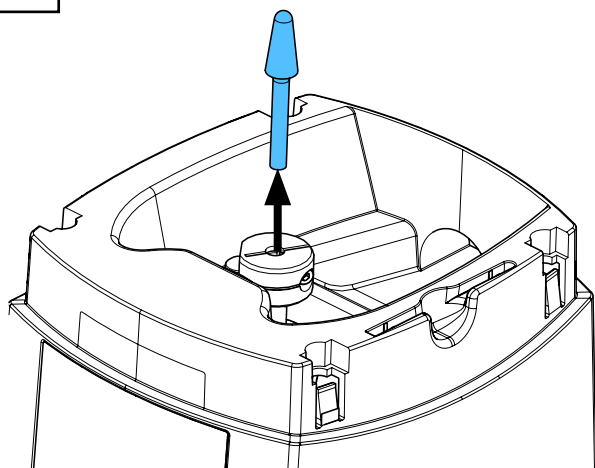



4

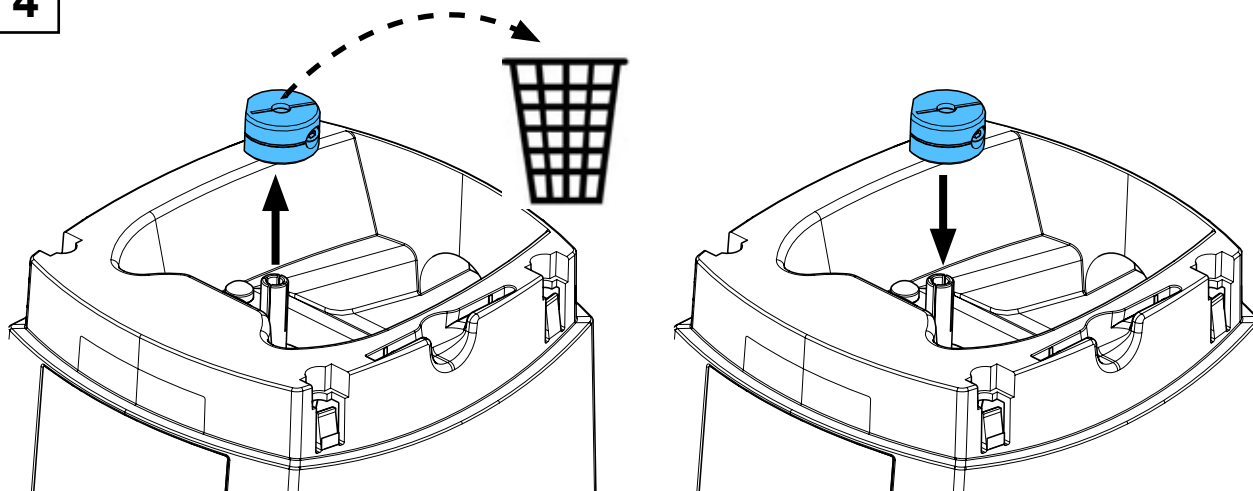


5

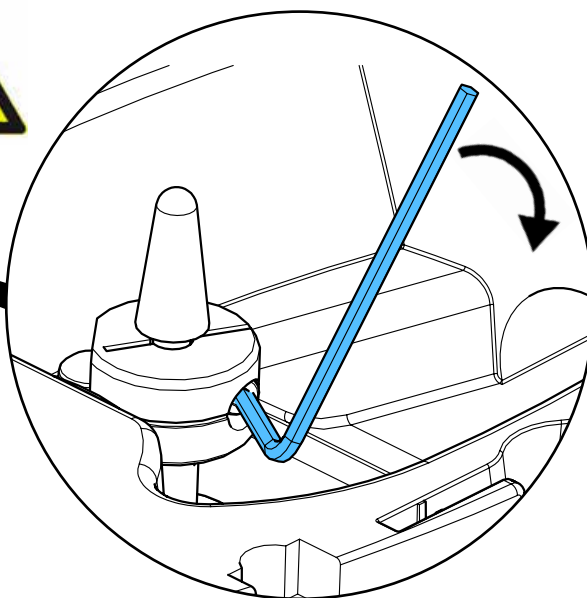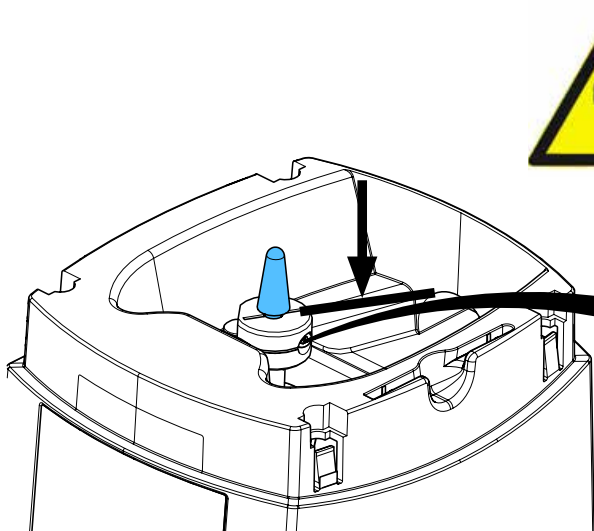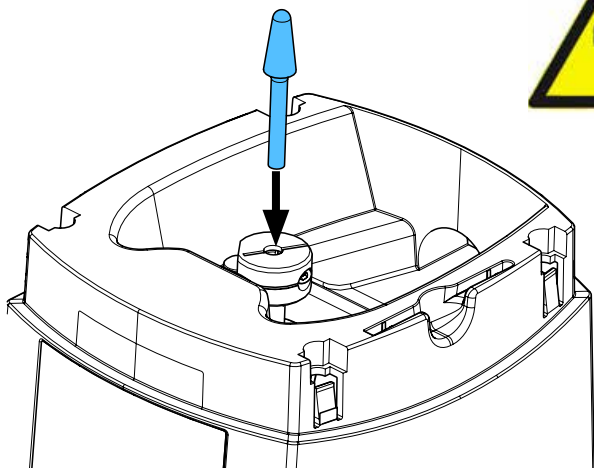


6

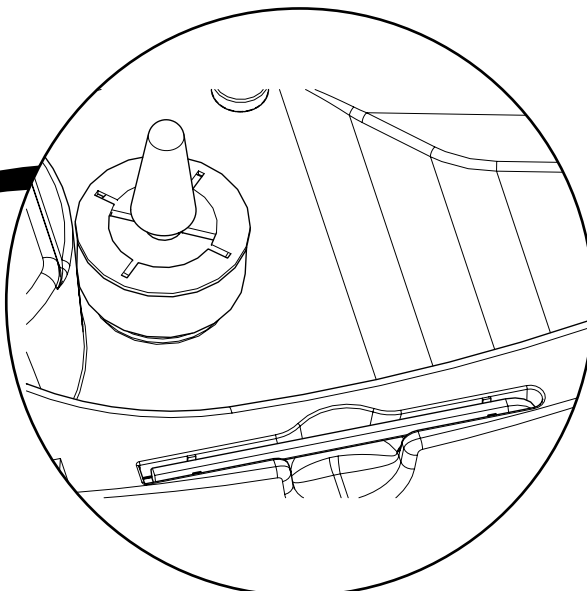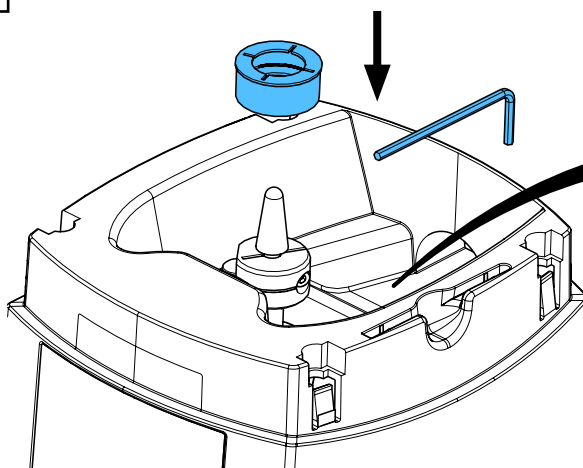

7

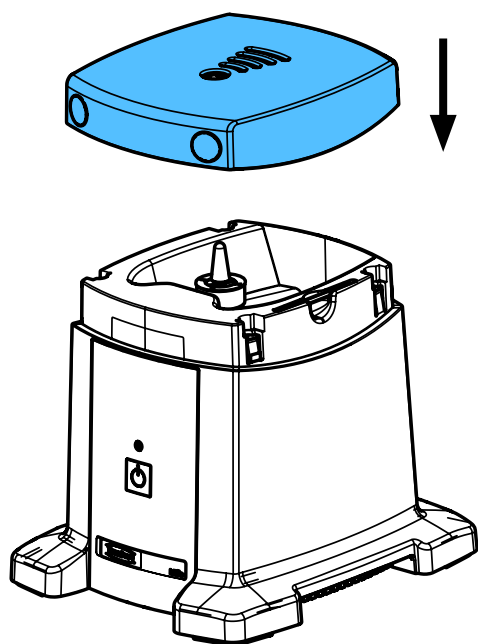


8

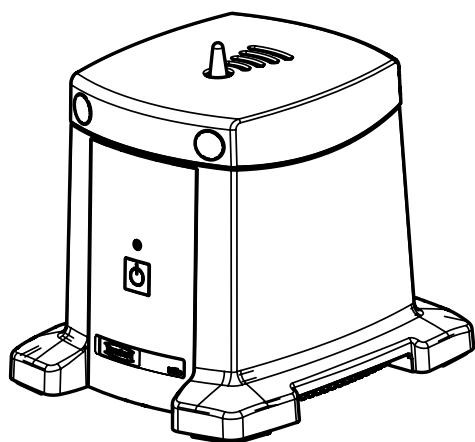




90003 3428

Schutzkappe • Protection cap • Capuchon protecteur • Calotta protettiva
Caperuza protectora • Защитный колпачок


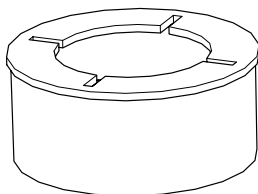


1


Millo / Millo pro

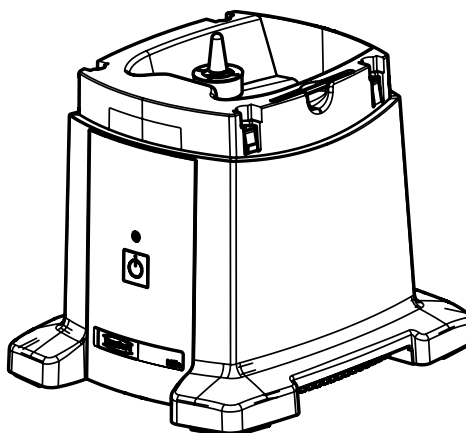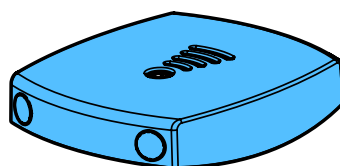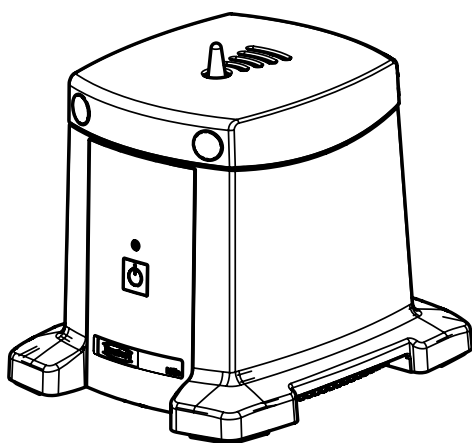



2

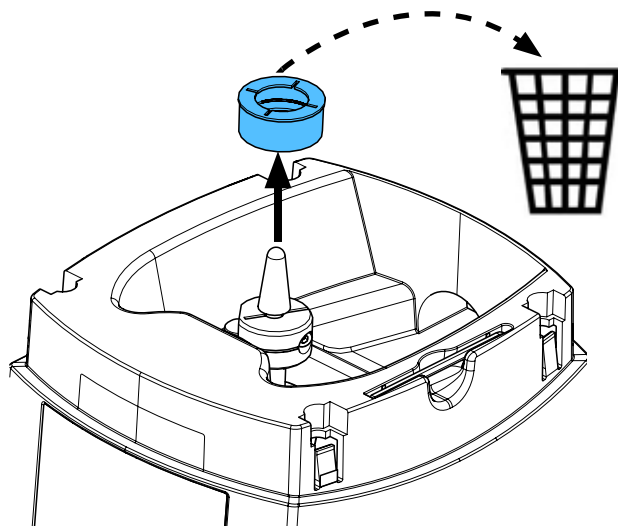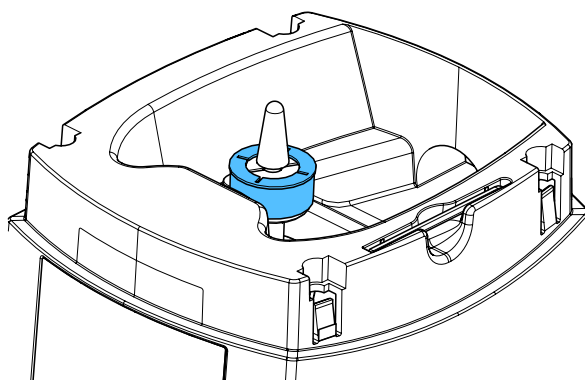

3


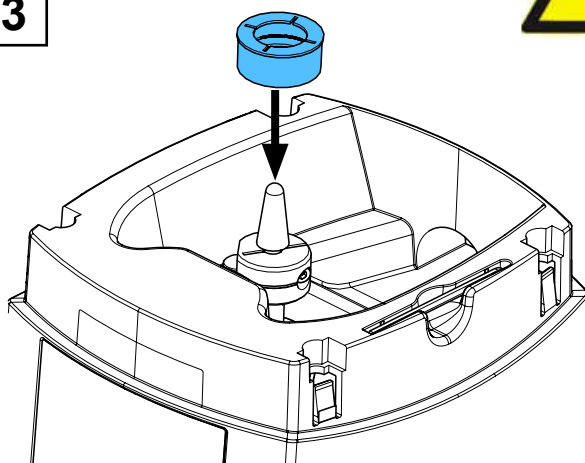


4

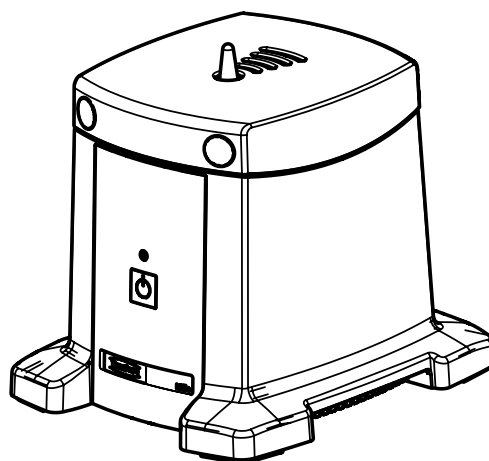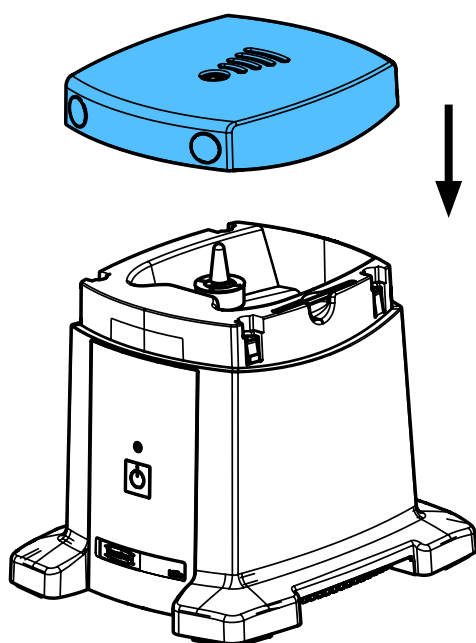


90003 3416 / 90013 4185



Sicherungsset • Set of fuses • Jeu de fusibles • Kit di fusibili • Juego de fusibles
Комплект предохранителей

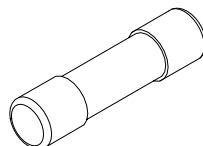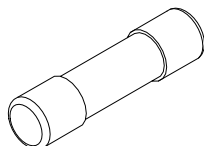


90003 3416 = 230V
90013 4185 = 120V

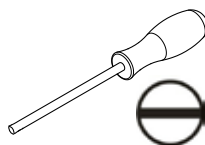



1

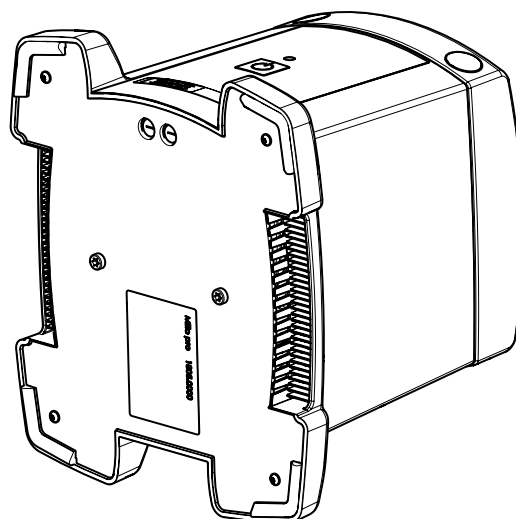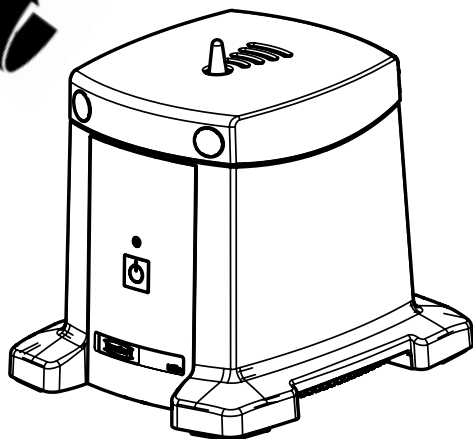



2

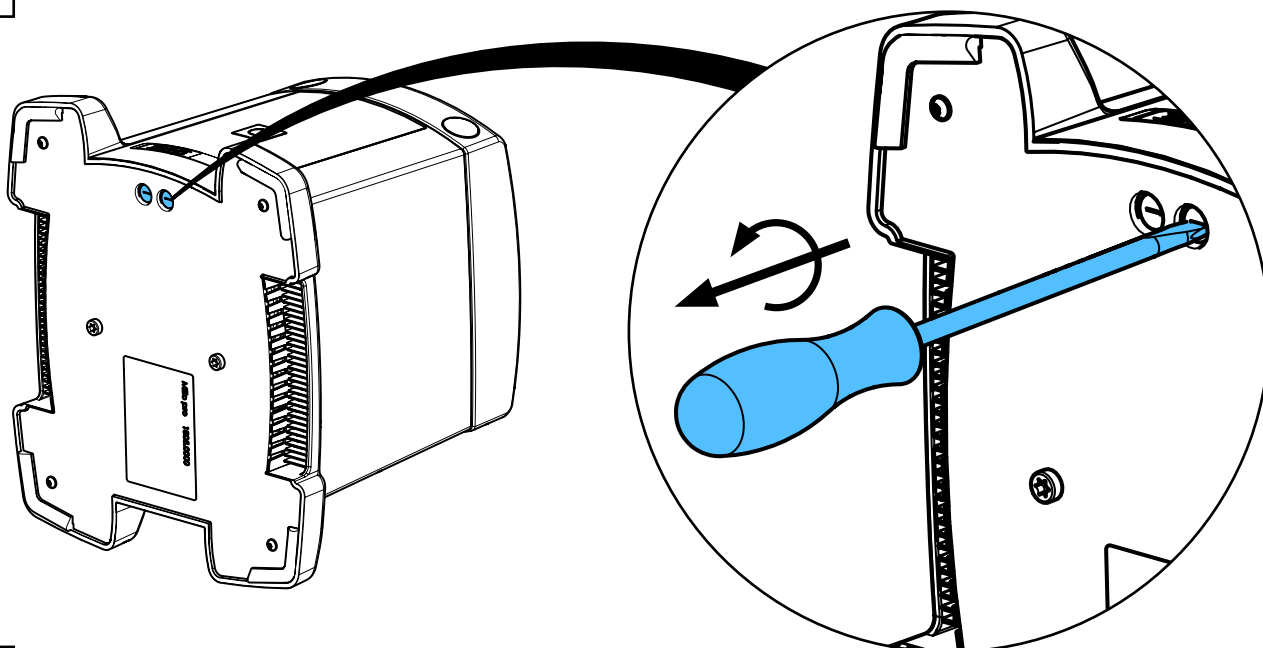

3

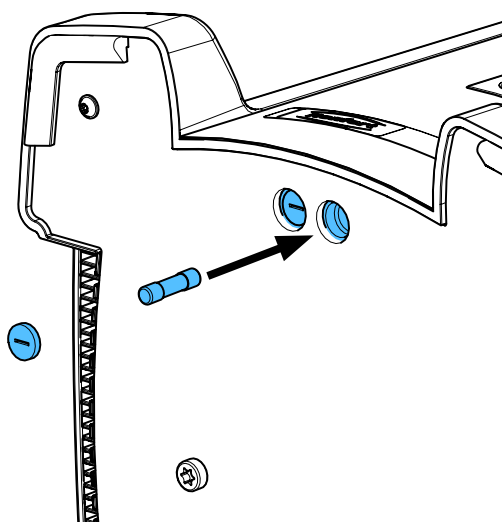
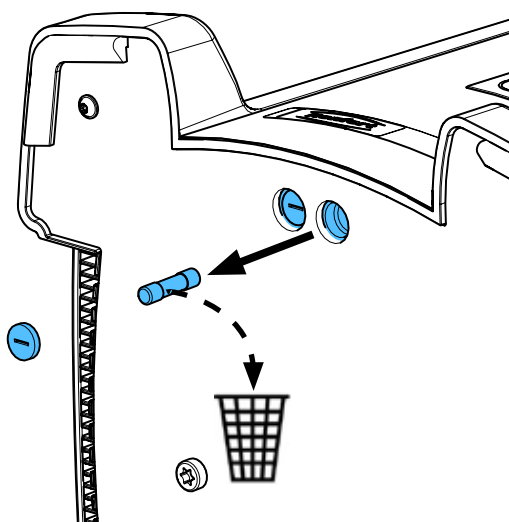


4

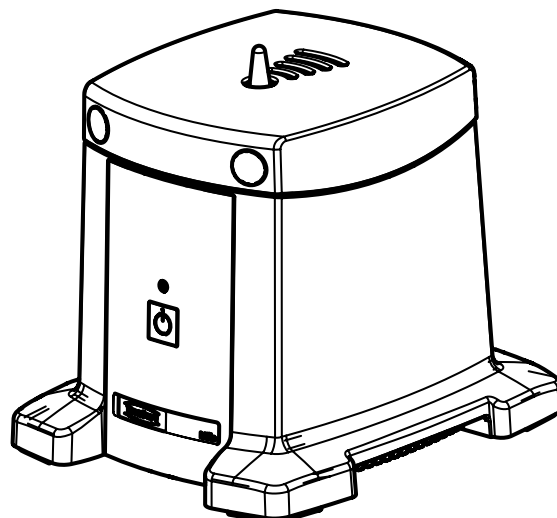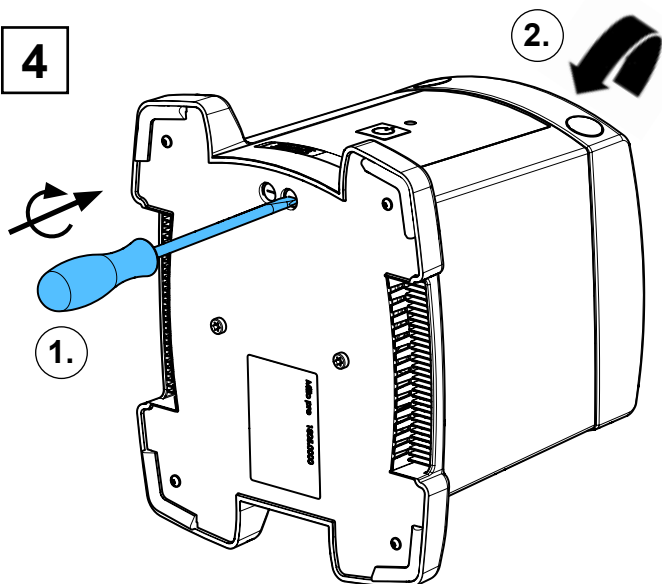


90003 3430


Arbeitstisch Millo pro • Millo pro work table • Table de travail Millo pro
Piano di lavoro Millo pro • Mesa de trabajo Millo pro • Столик Millo pro

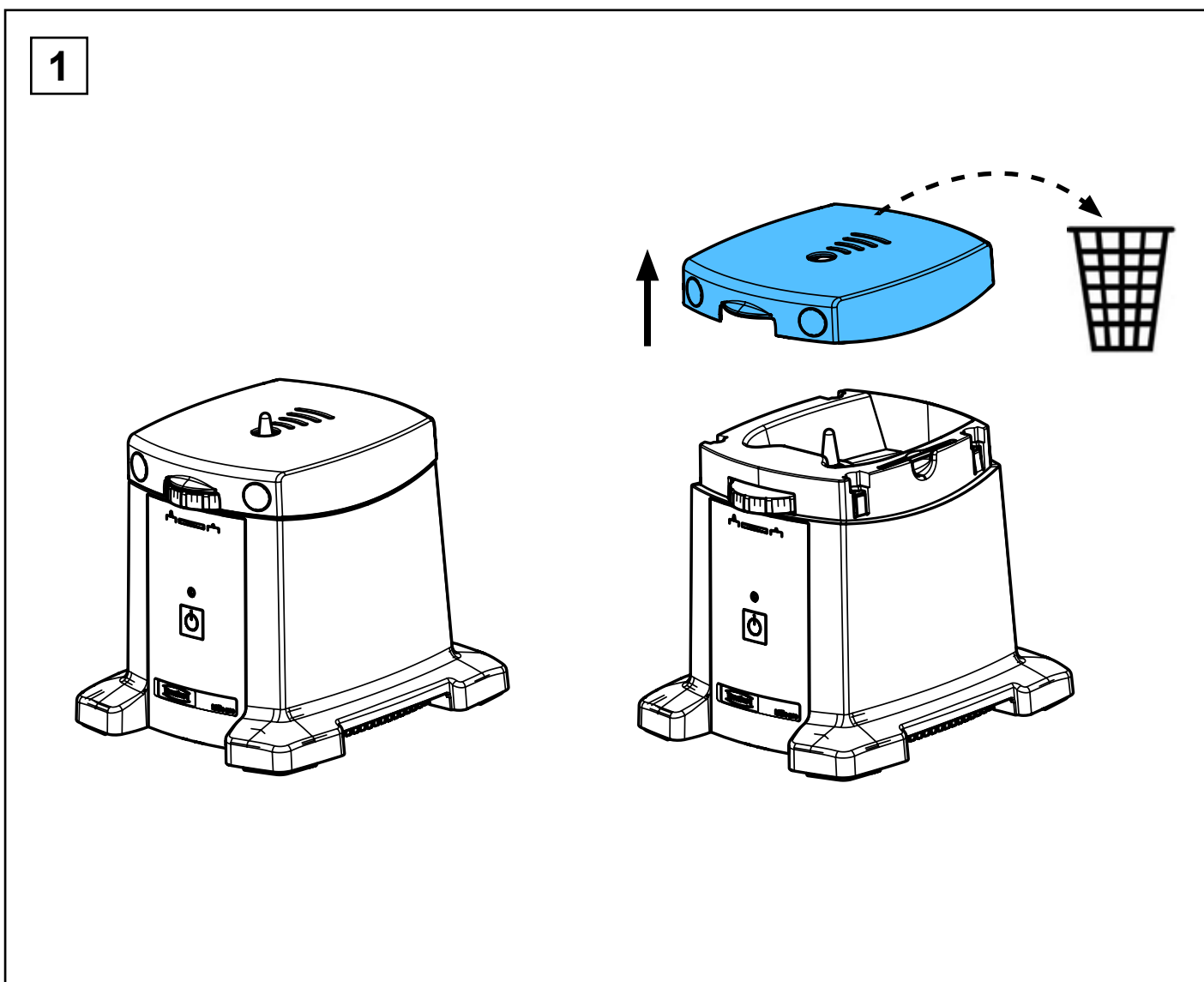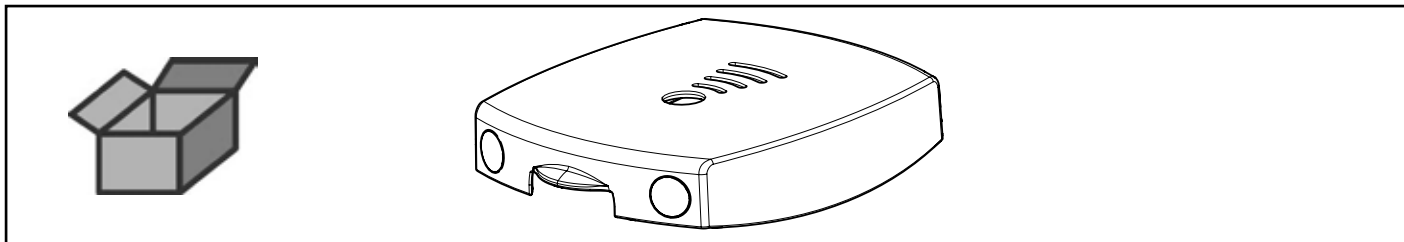

2


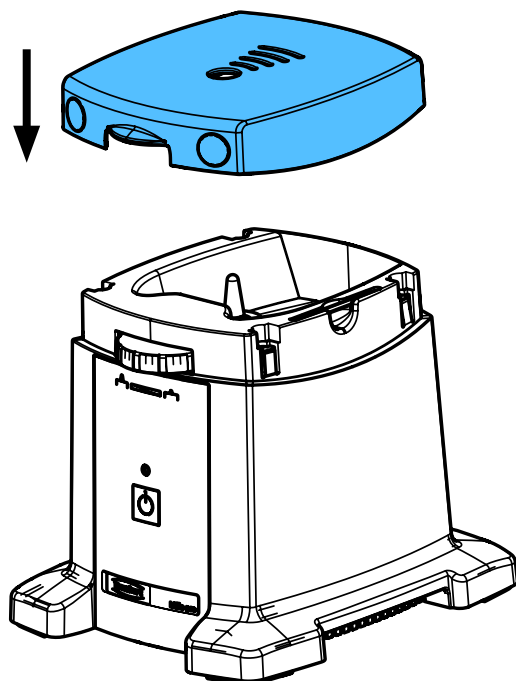


3

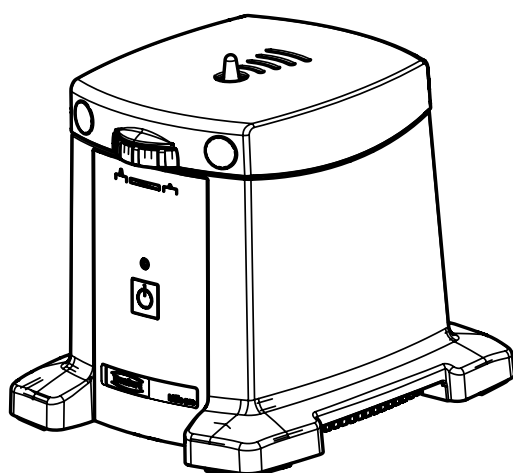


90003 3429


Arbeitstisch Millo • Millo work table • Table de travail Millo • Piano di lavoro Millo
Mesa de trabajo Millo • Столик Millo

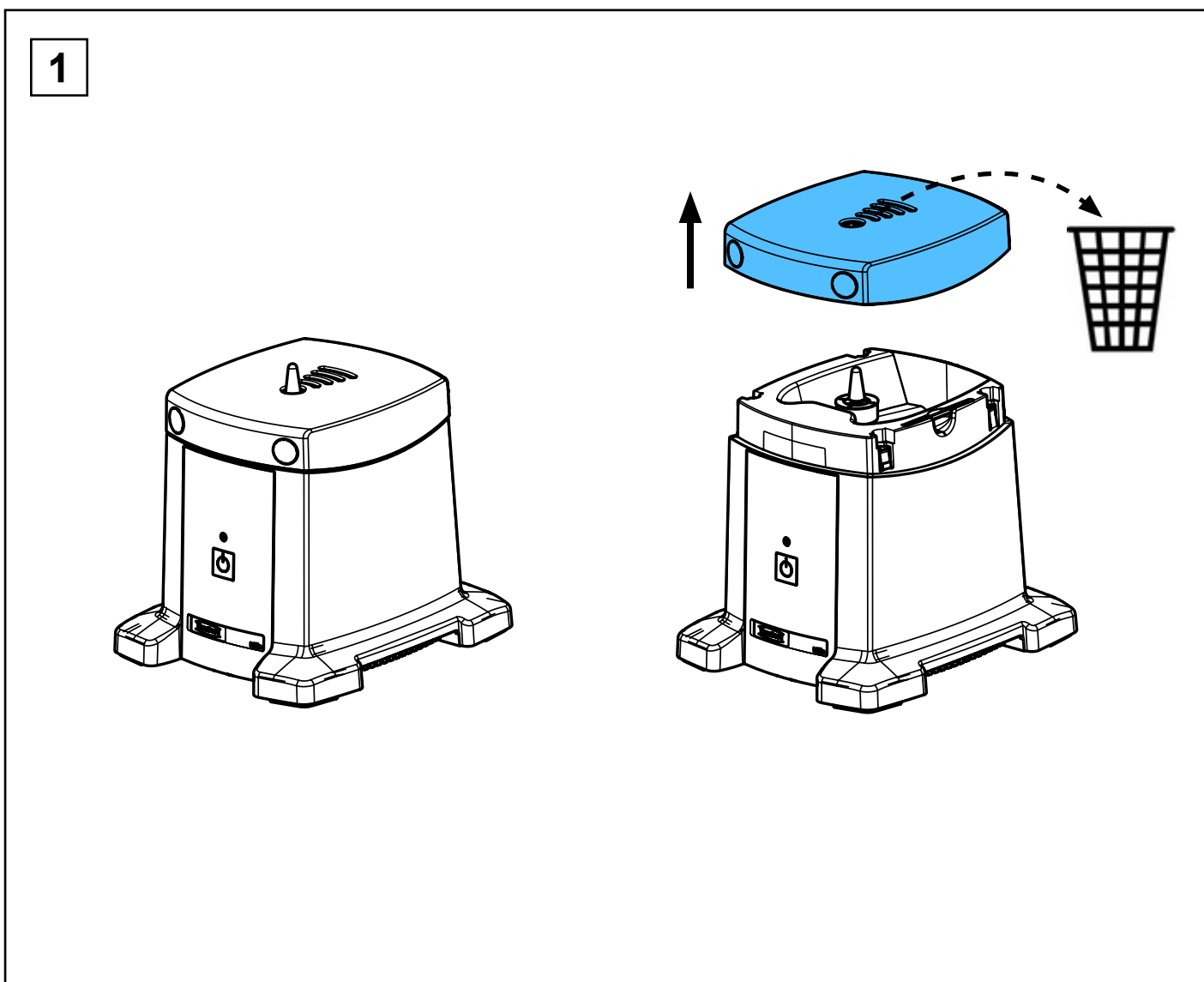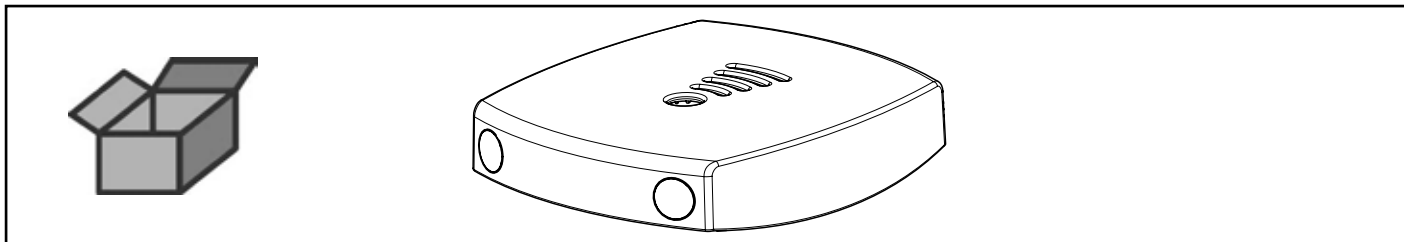

2

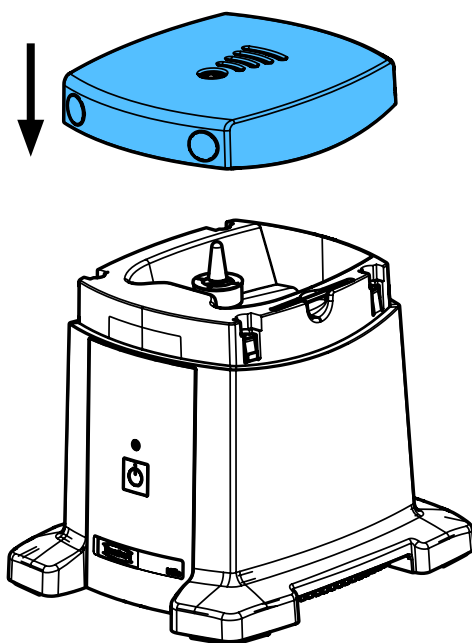


3

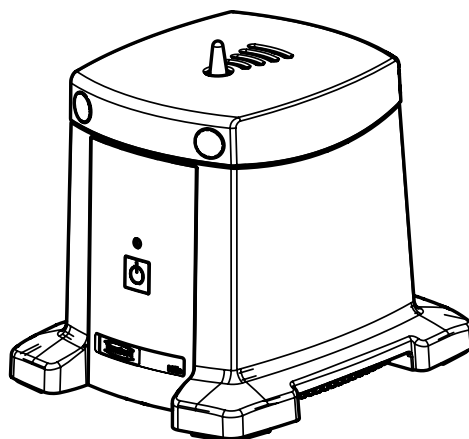


90003 3552 /..3 /..5 /..6 /..7

Adapter Schuko • Schuko adapter • Adaptateur Schuko • Adattatore Schuko
Adaptador Schuko • Schuko Адаптер

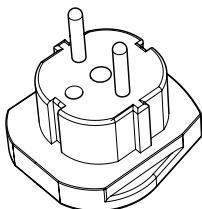


1

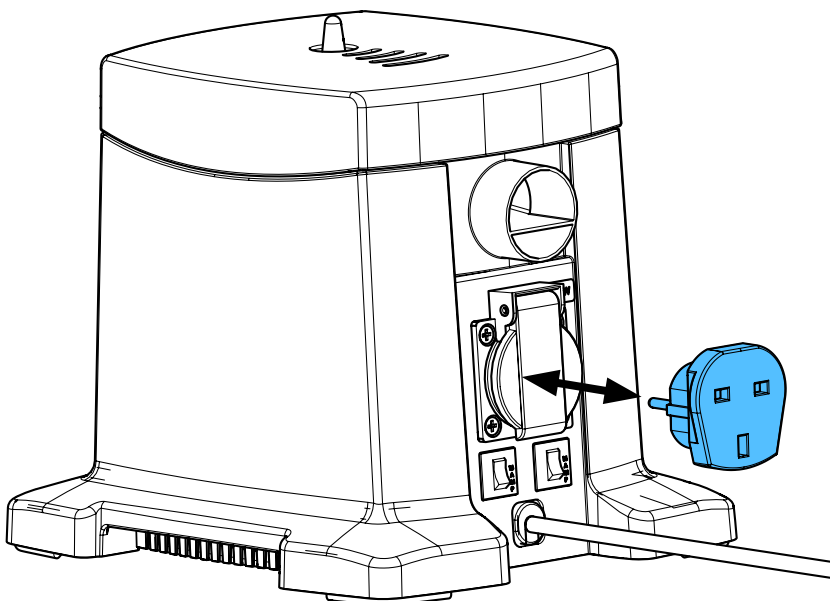


90003 3425 / 90003 3426

Netzkabel • Power cable • Câble de réseau • Cavo di rete • Cable de red
Сетевой кабель

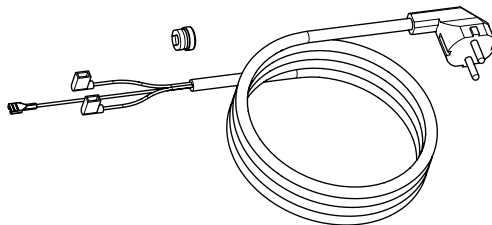



90003 3425 = 230 V

90003 3426 = 120 V

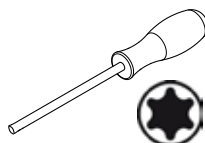


T10


1 Millo 1 - 9 / Millo pro 10 - 18

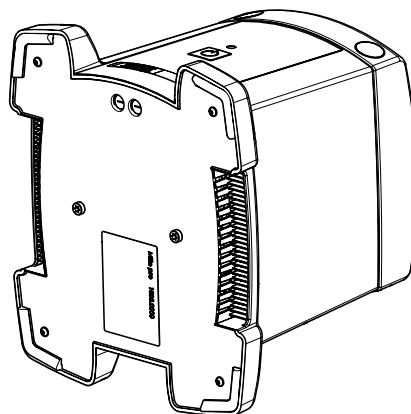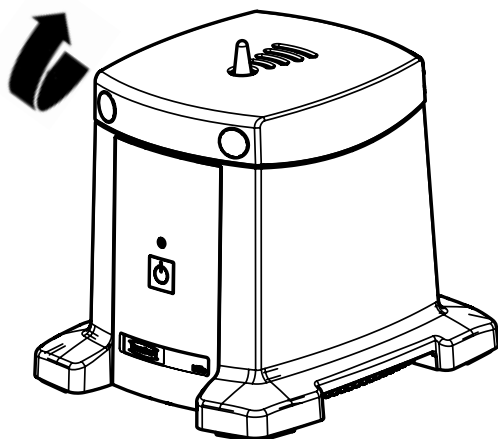


2

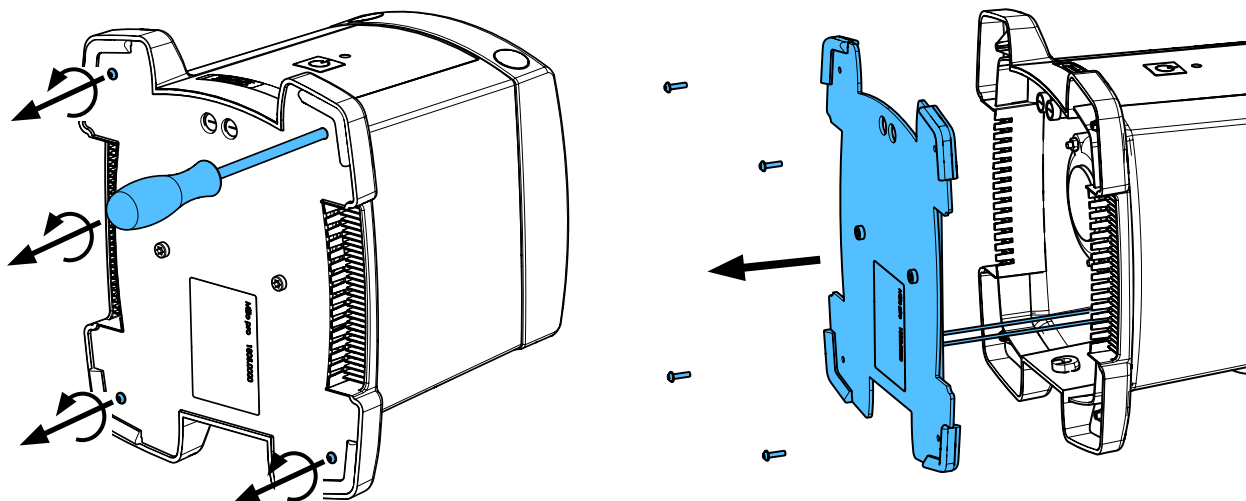



3

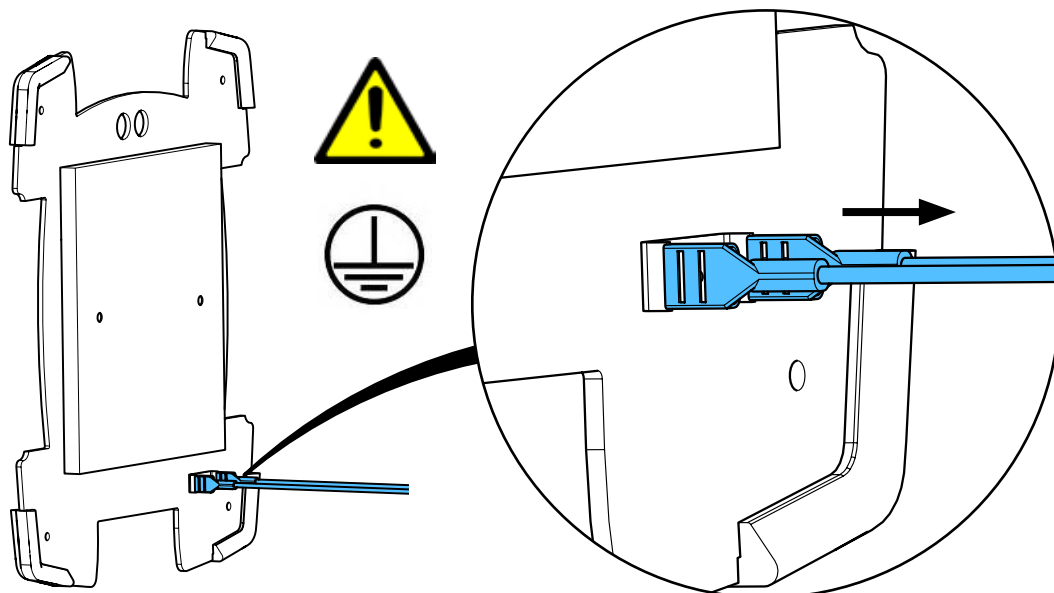



4

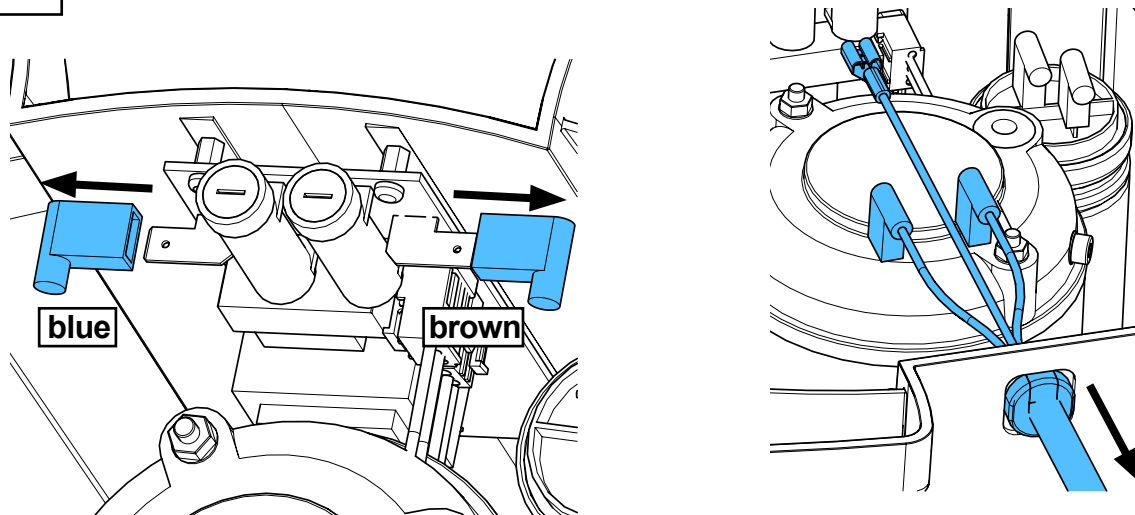



5

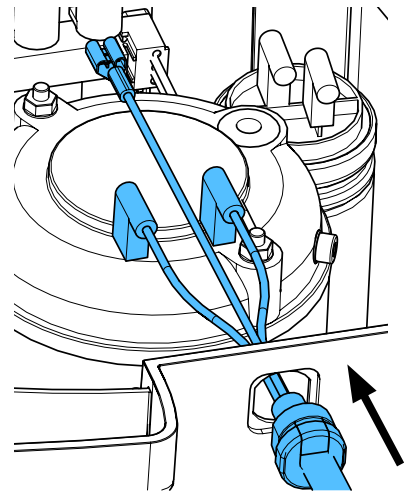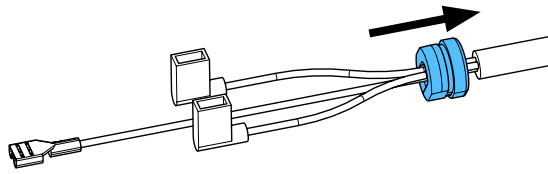



6

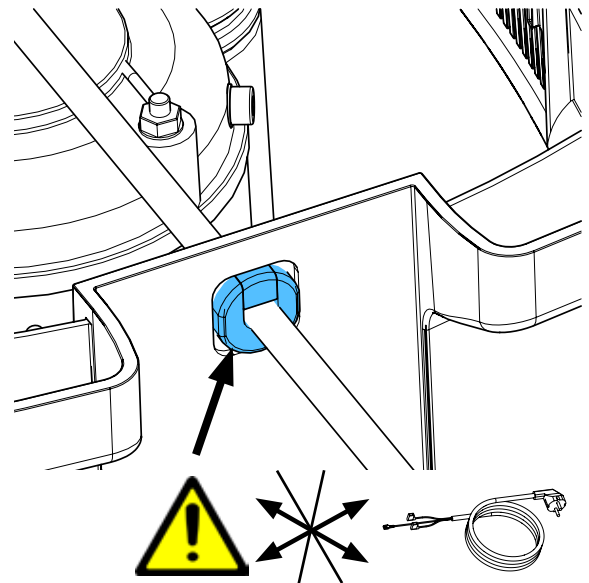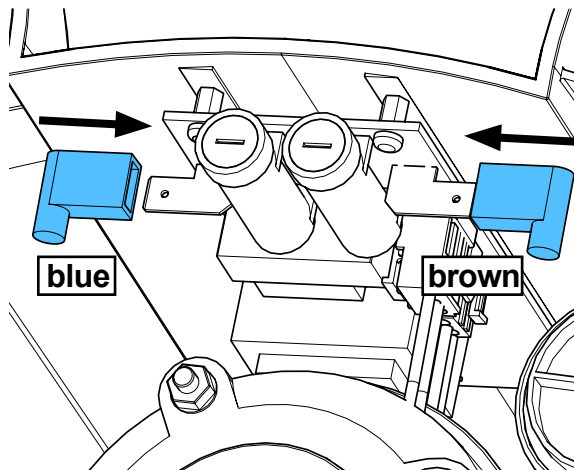


7

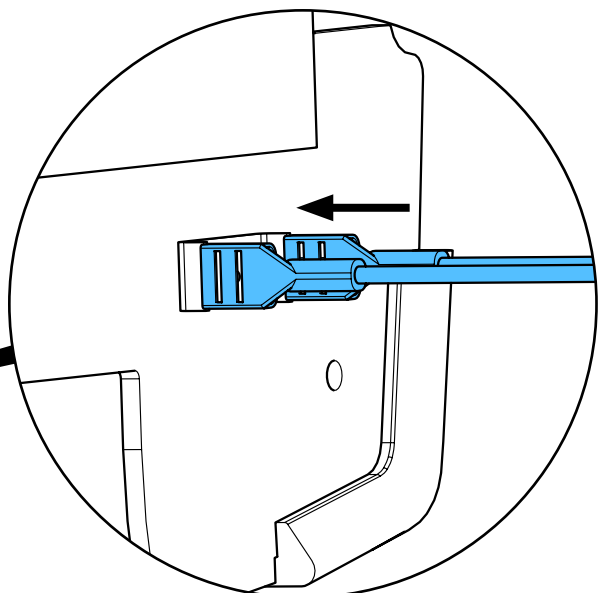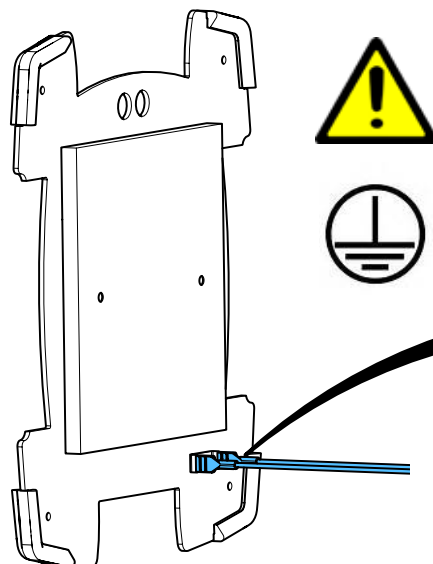

8

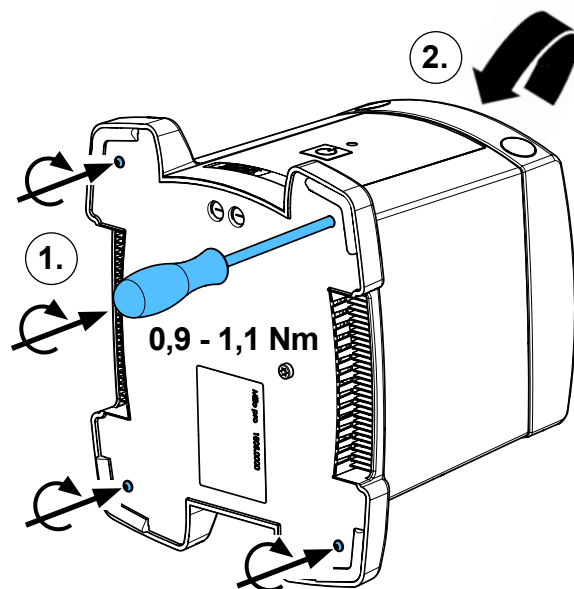
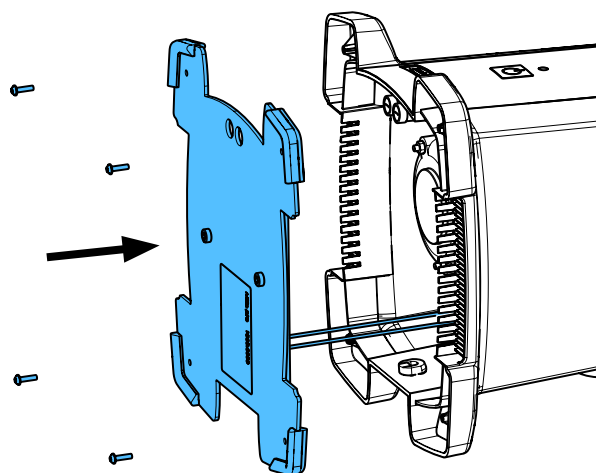


9


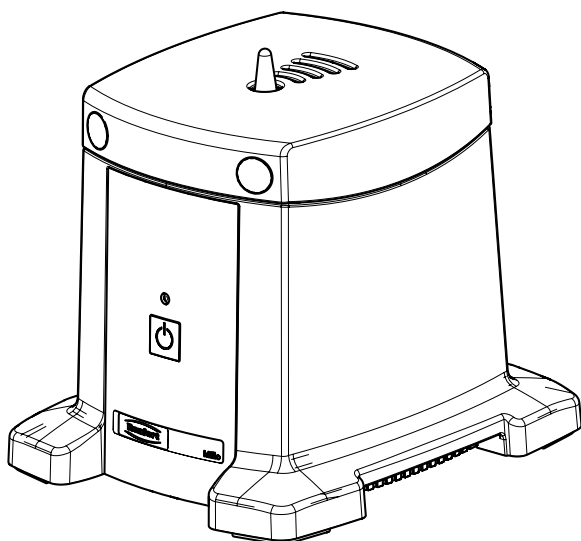


10



Millo pro 10 - 18

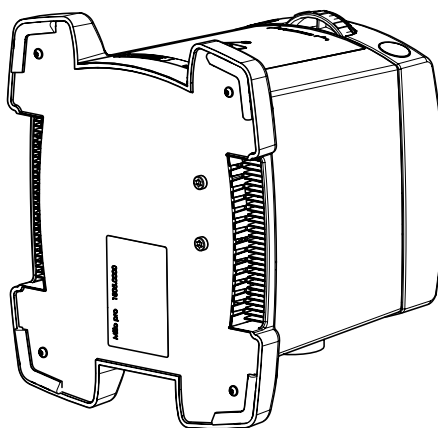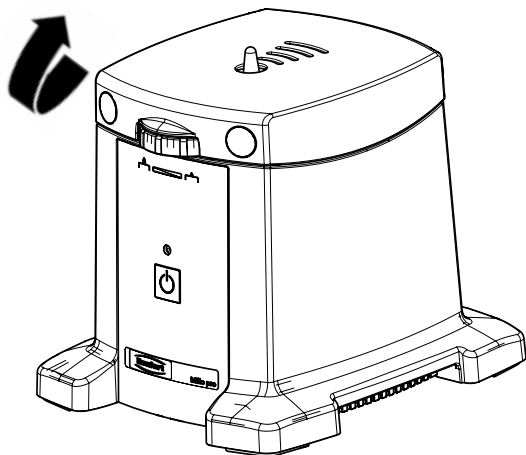



11

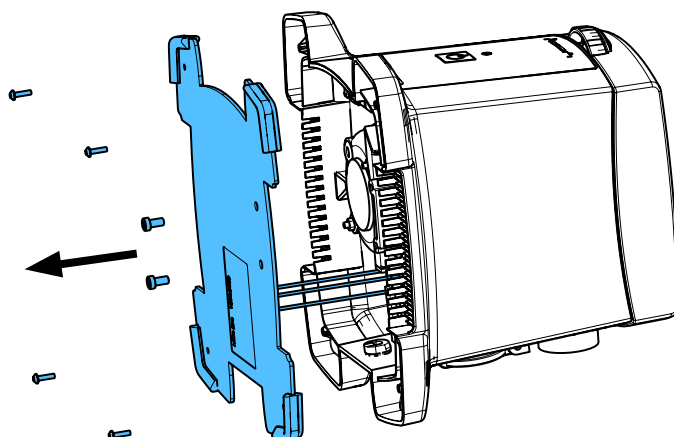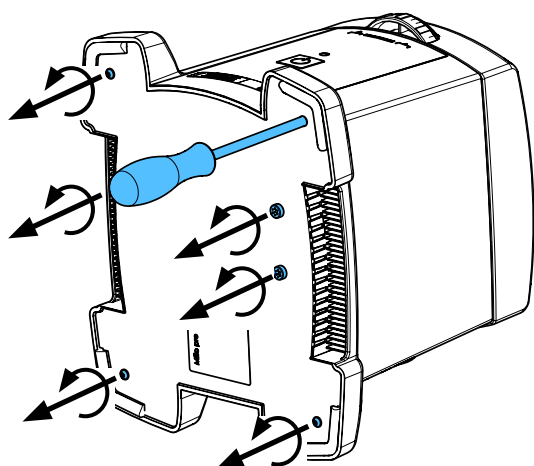



12

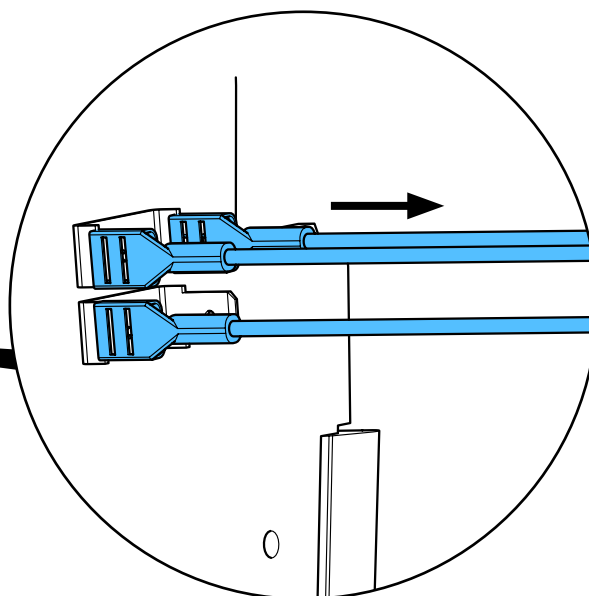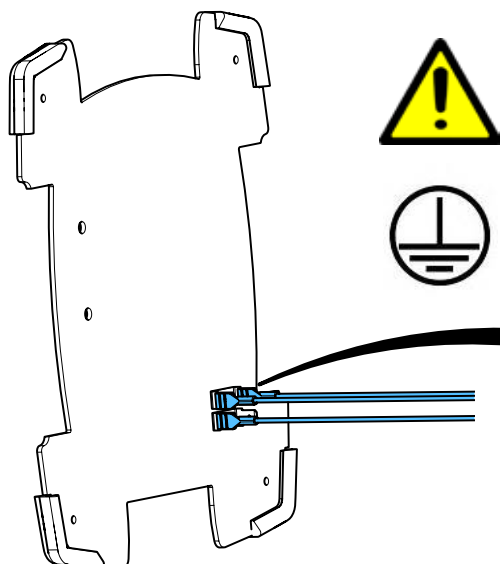




13

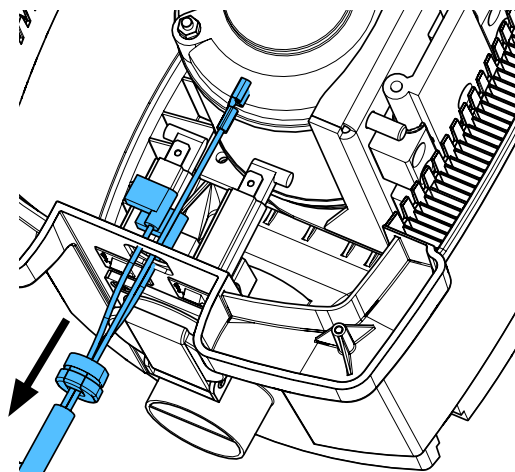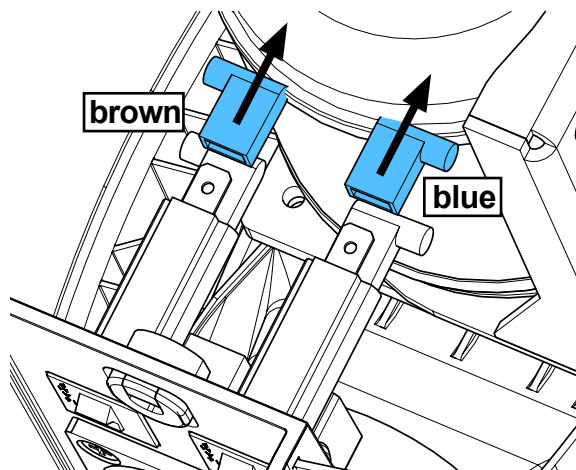


14

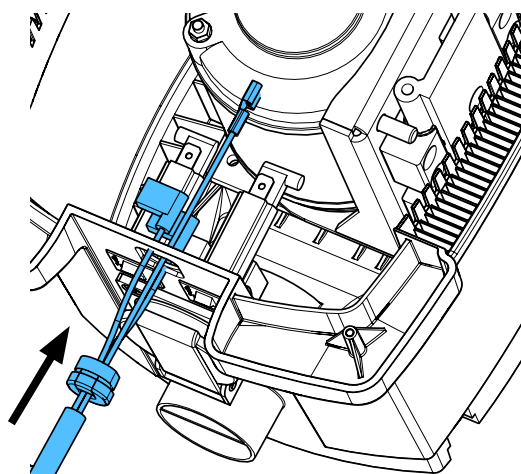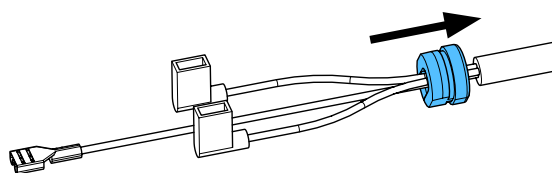



15

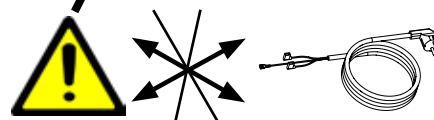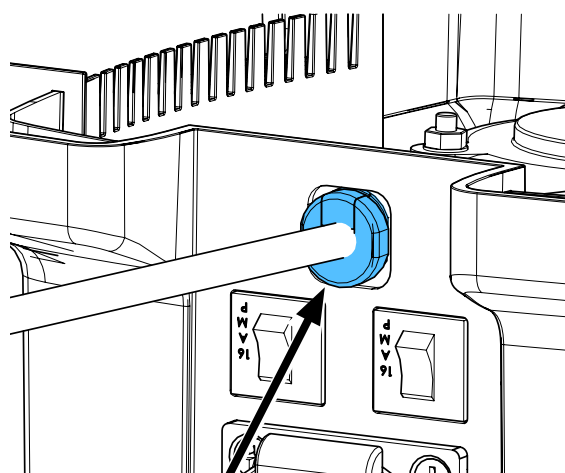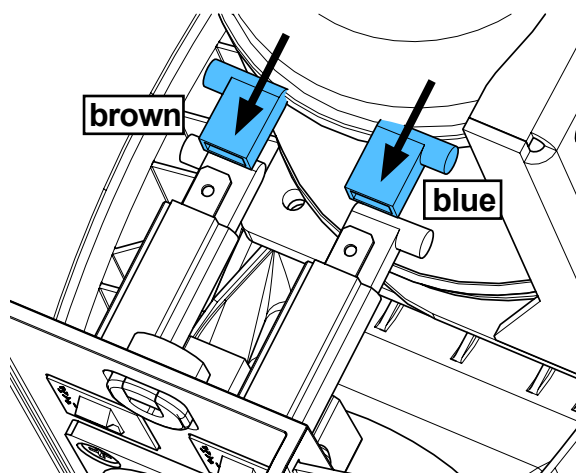


16

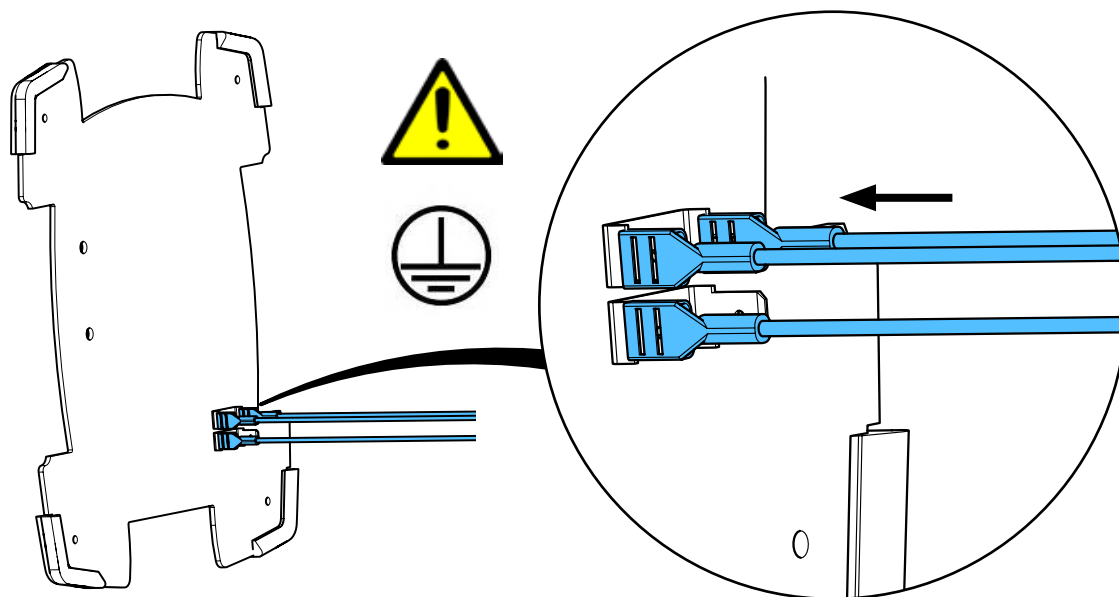




17

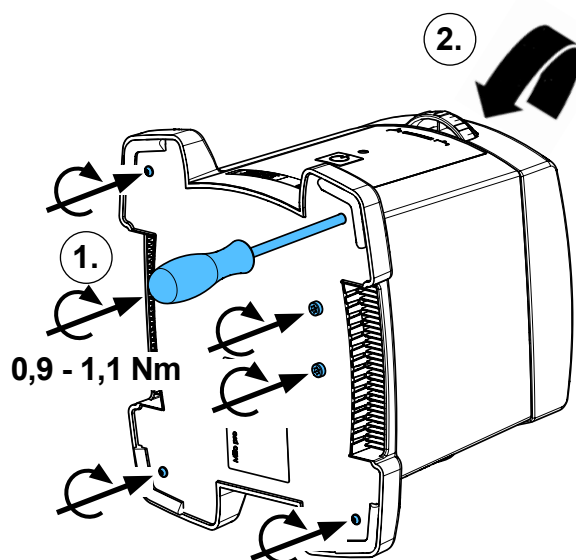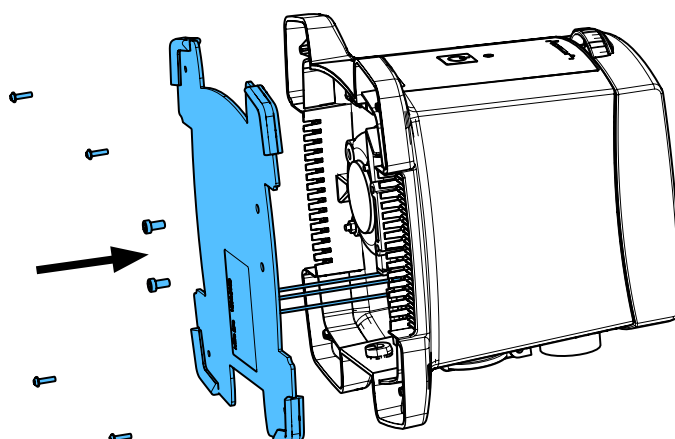

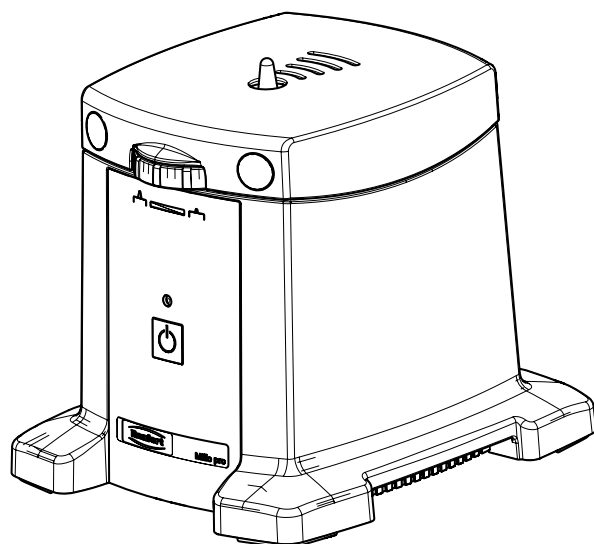


90003 3438 / 90003 3443


Leiterplatte • Circuit card • Carte imprimée • Circuito stampato • Placa de circuito
Электронная плата

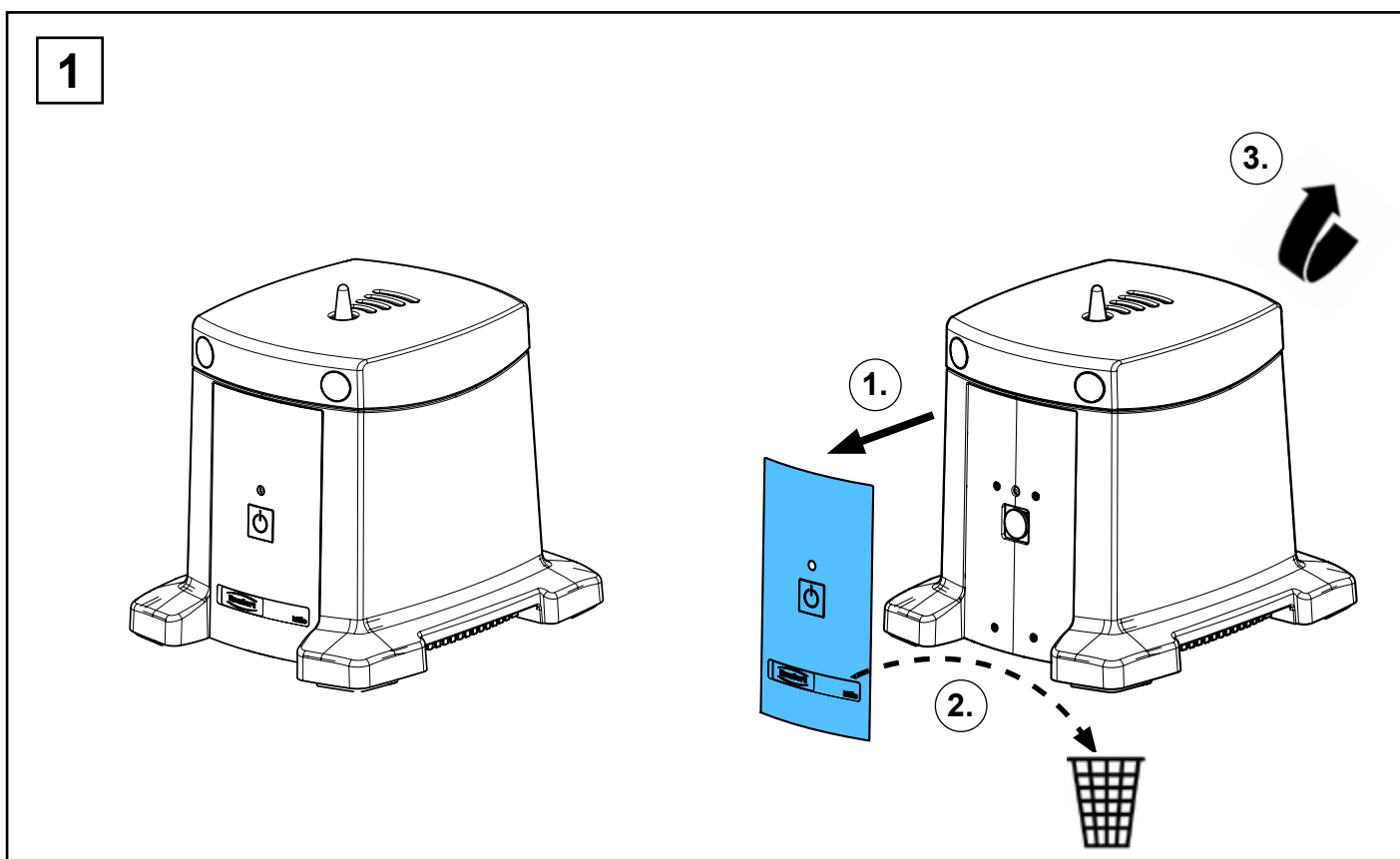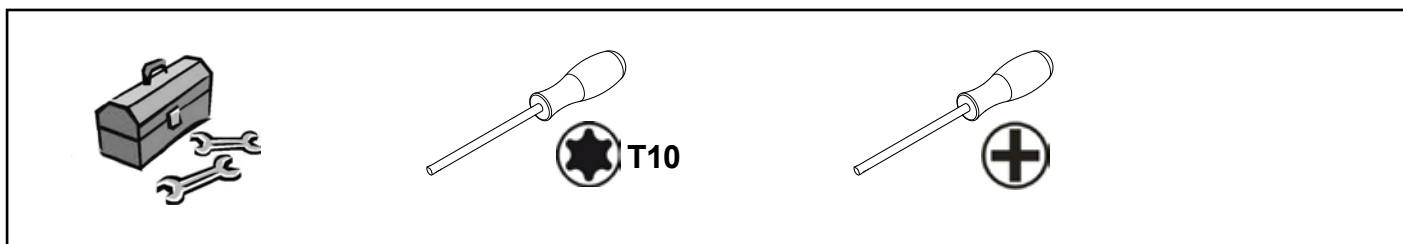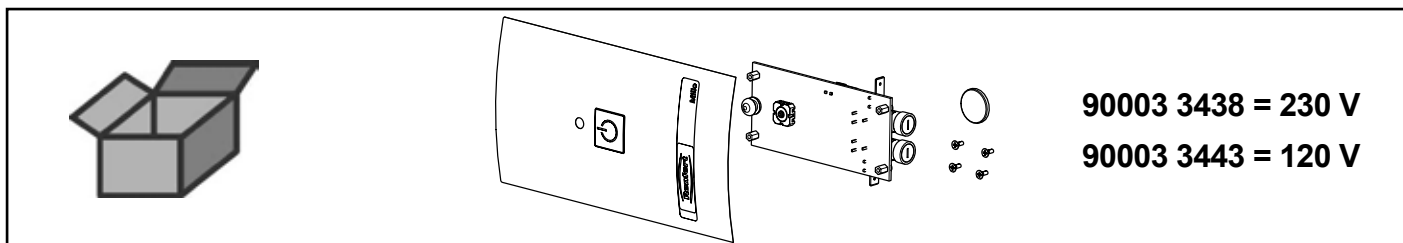


2

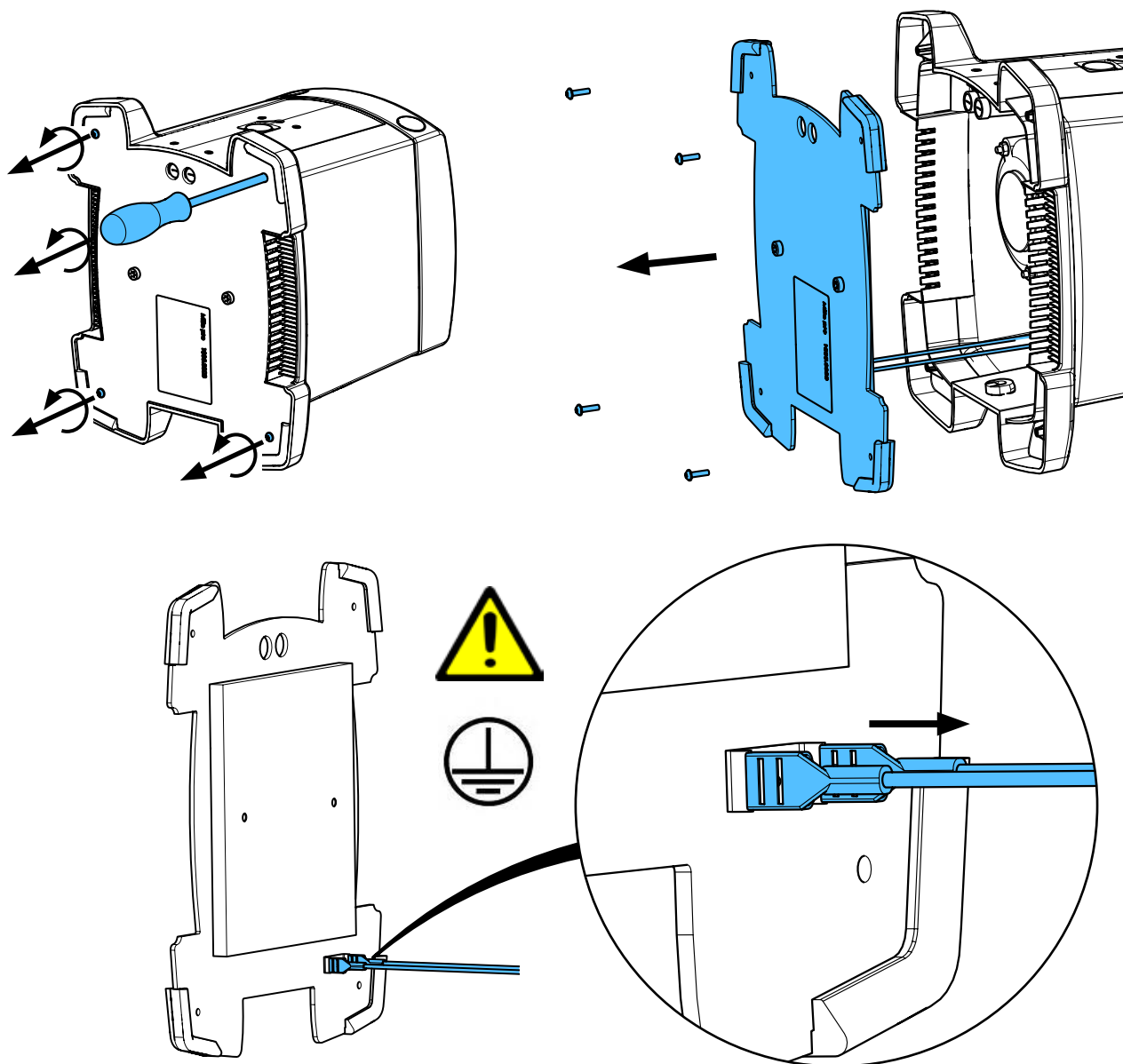


3

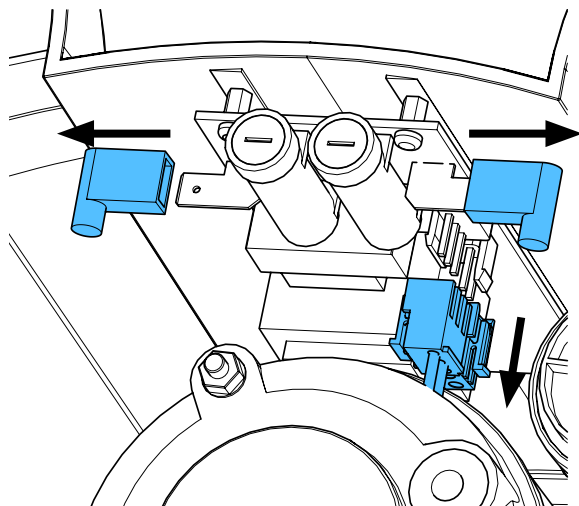


4

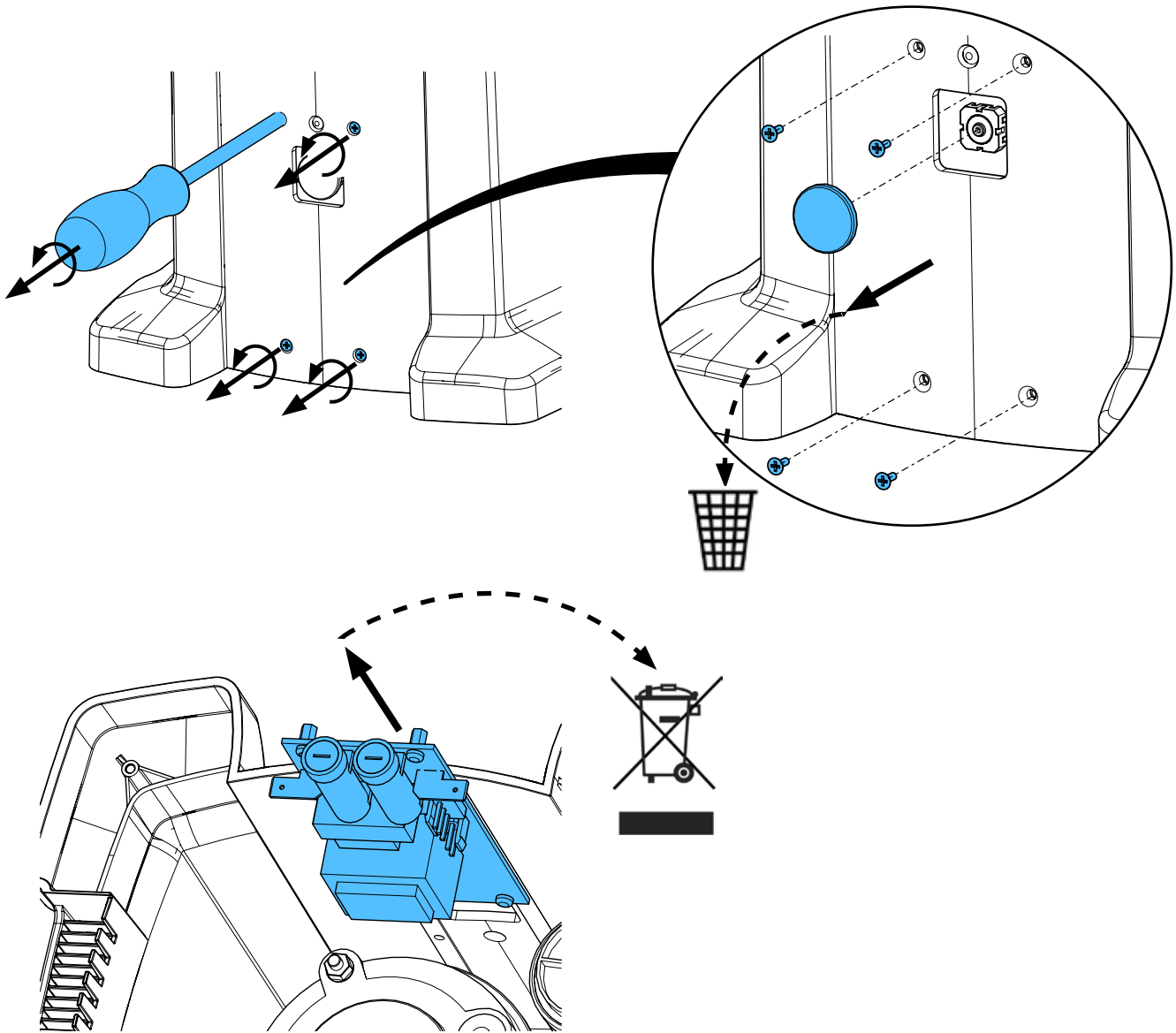




5


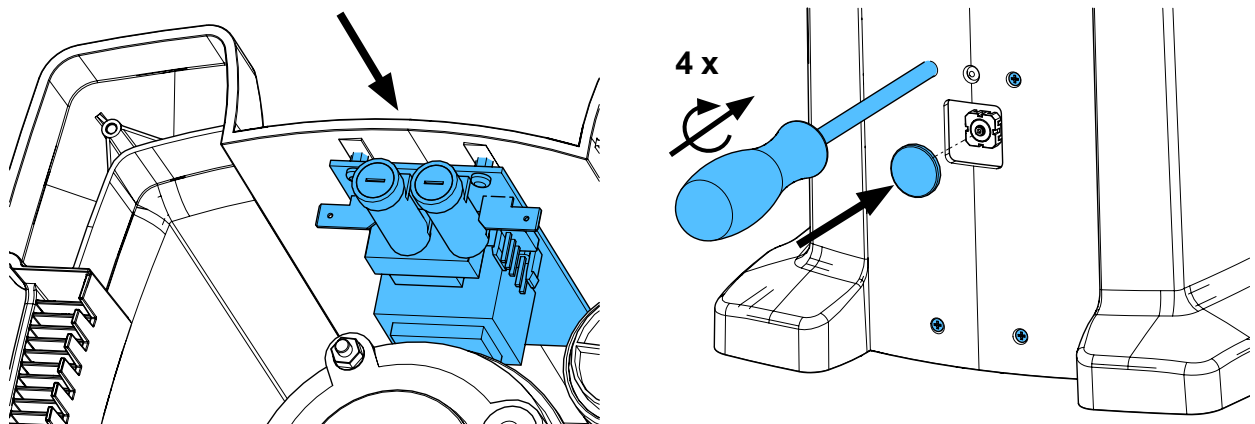



6

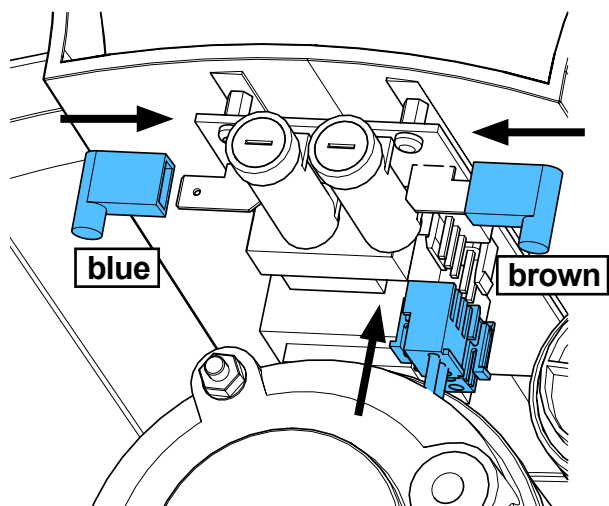

7

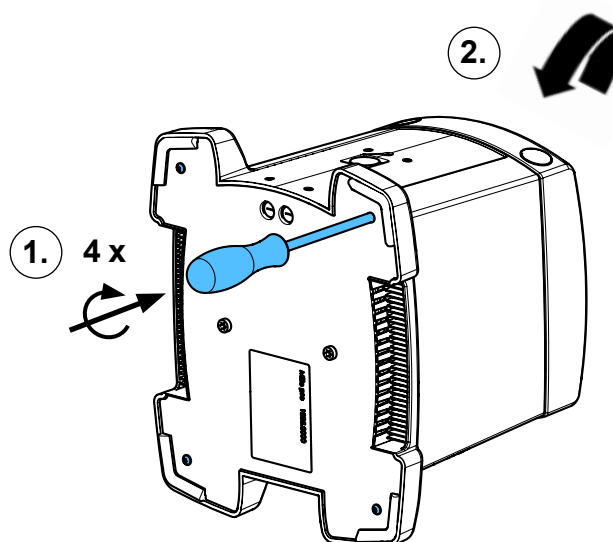
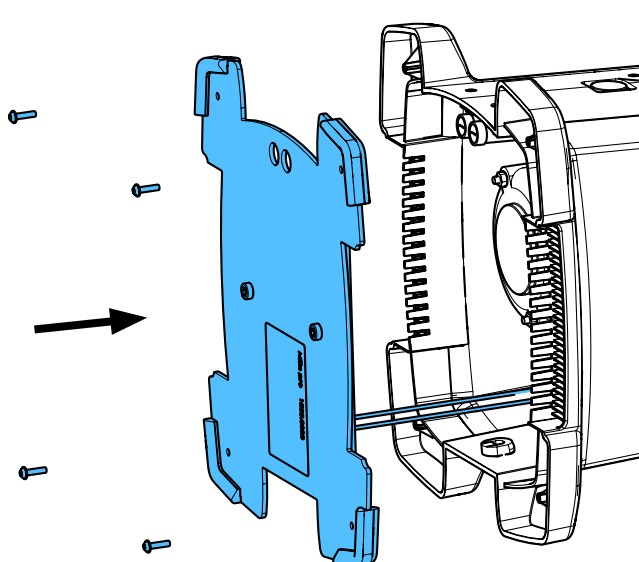
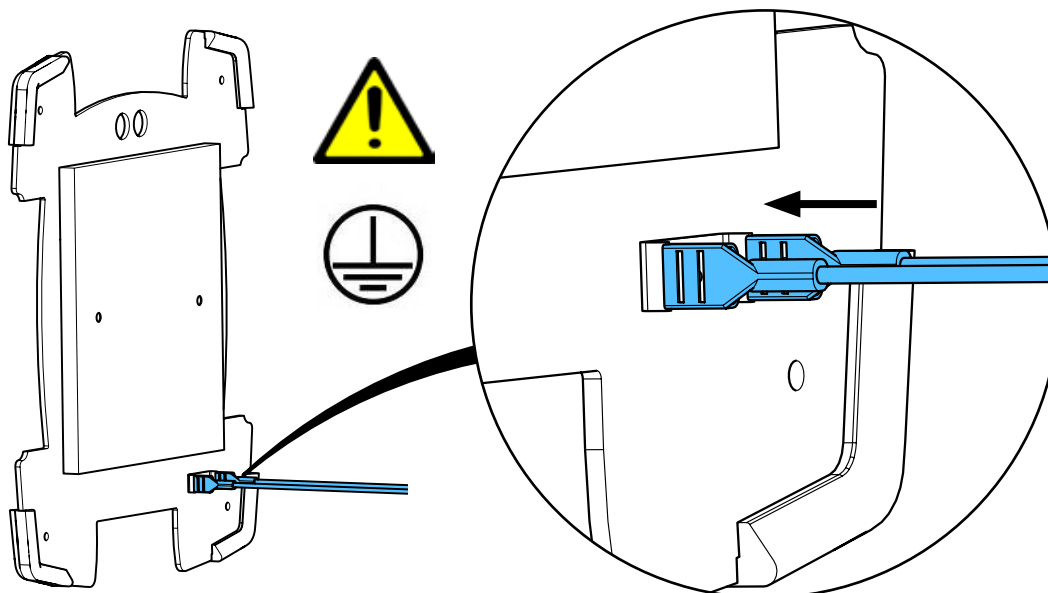


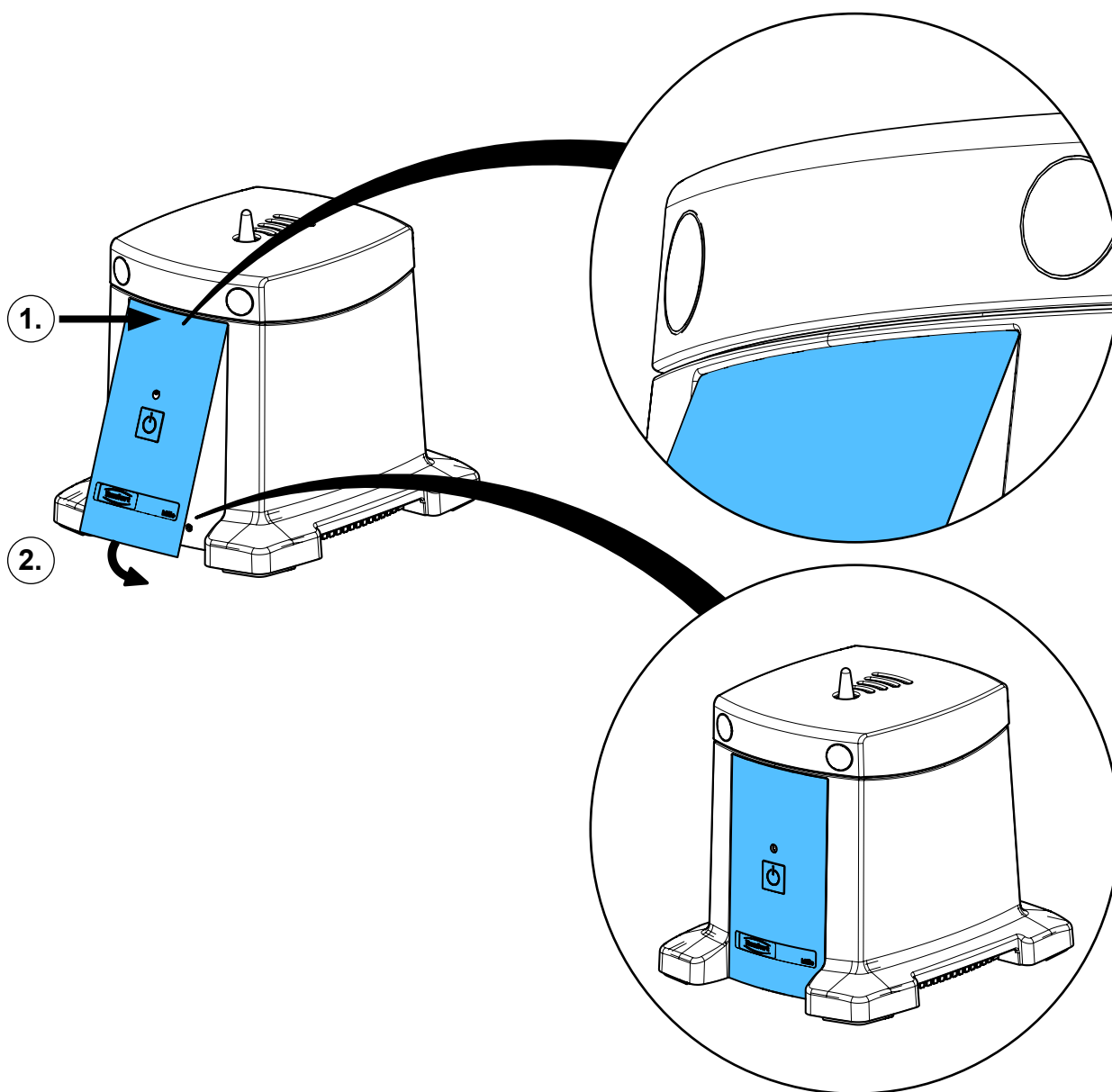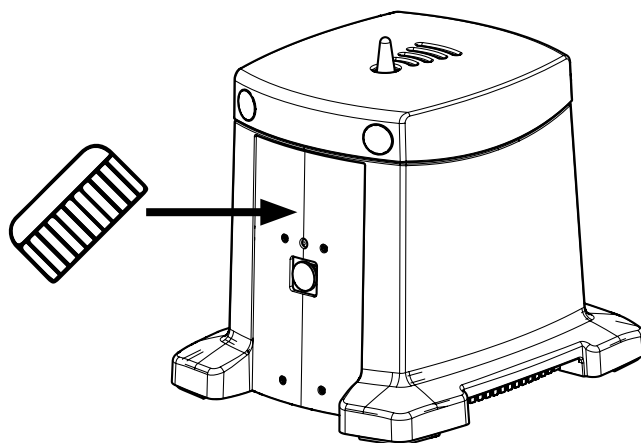




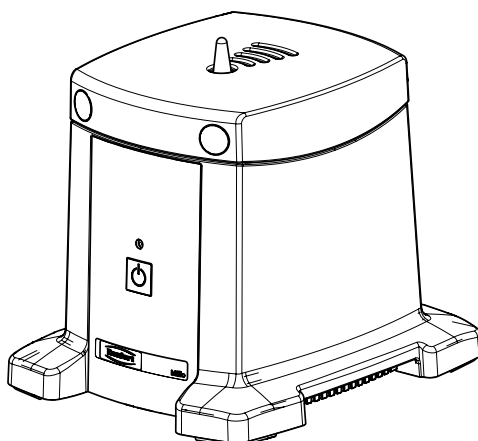


90003 3409 / 90003 3410


Leiterplatte • Circuit card • Carte imprimée • Circuito stampato • Placa de circuito
Электронная плата

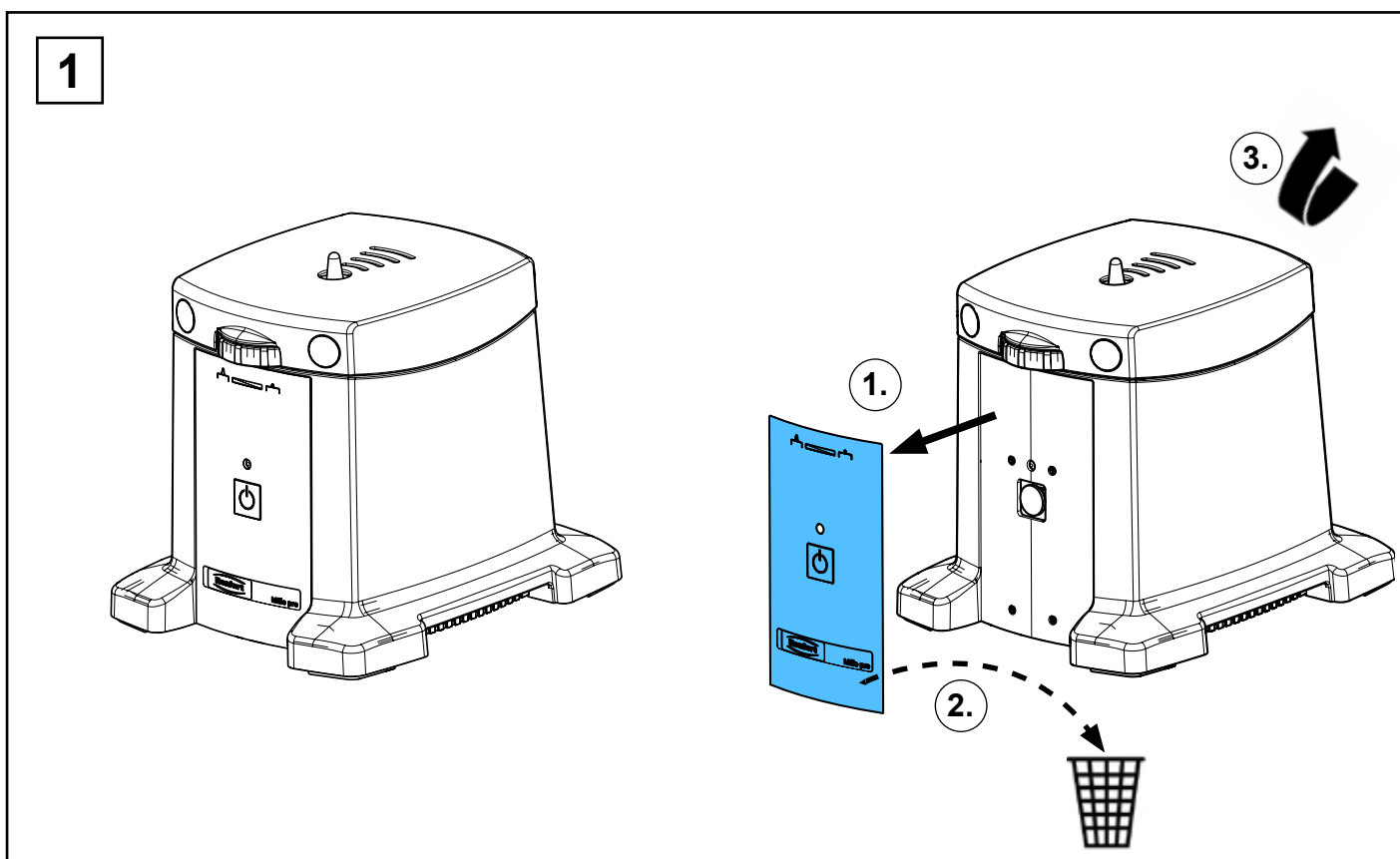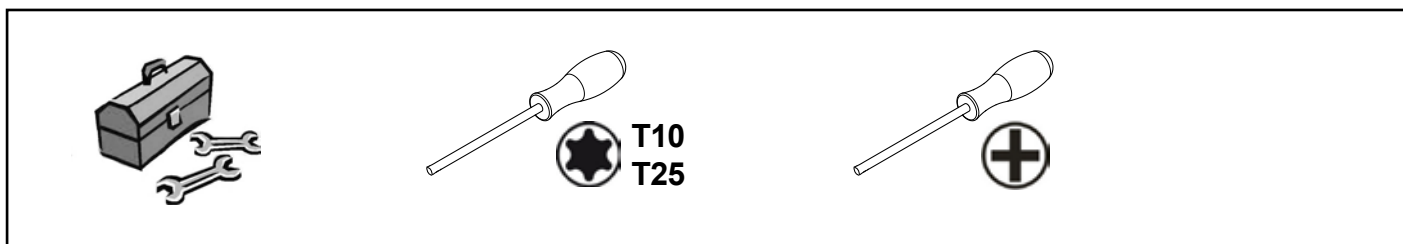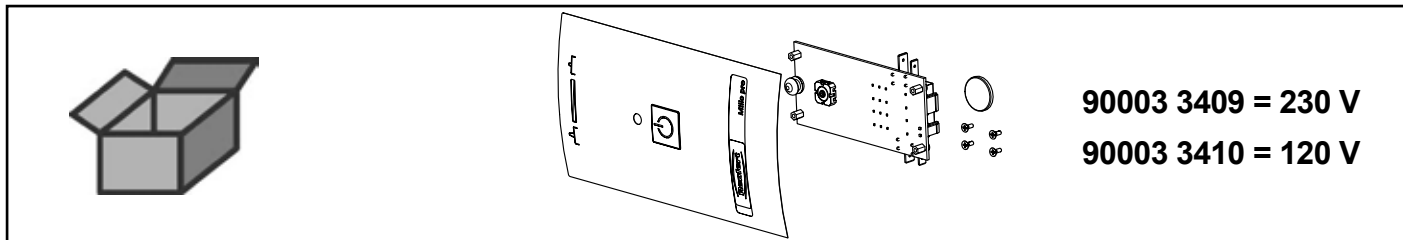


2

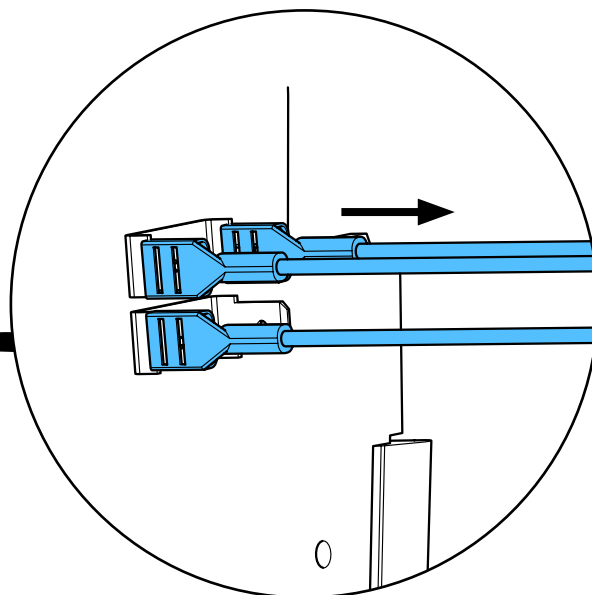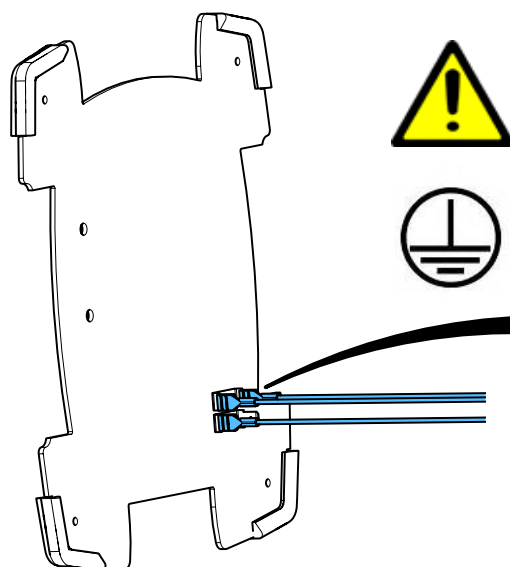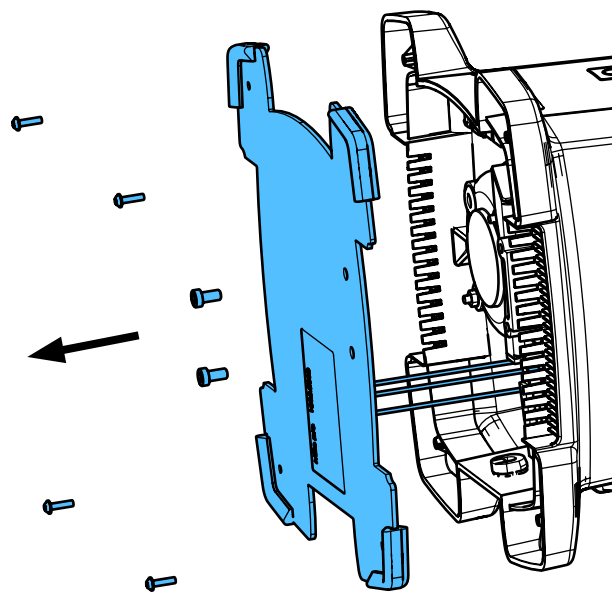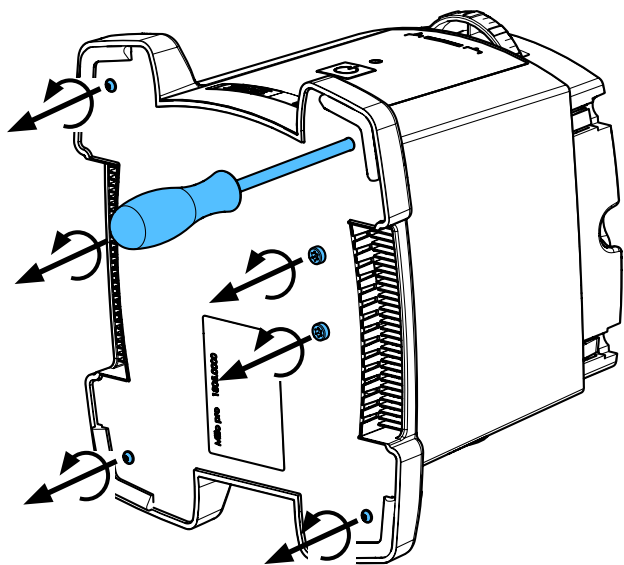


3

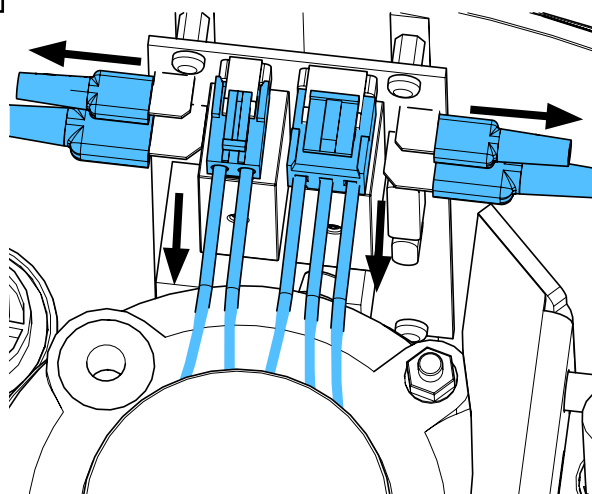


4

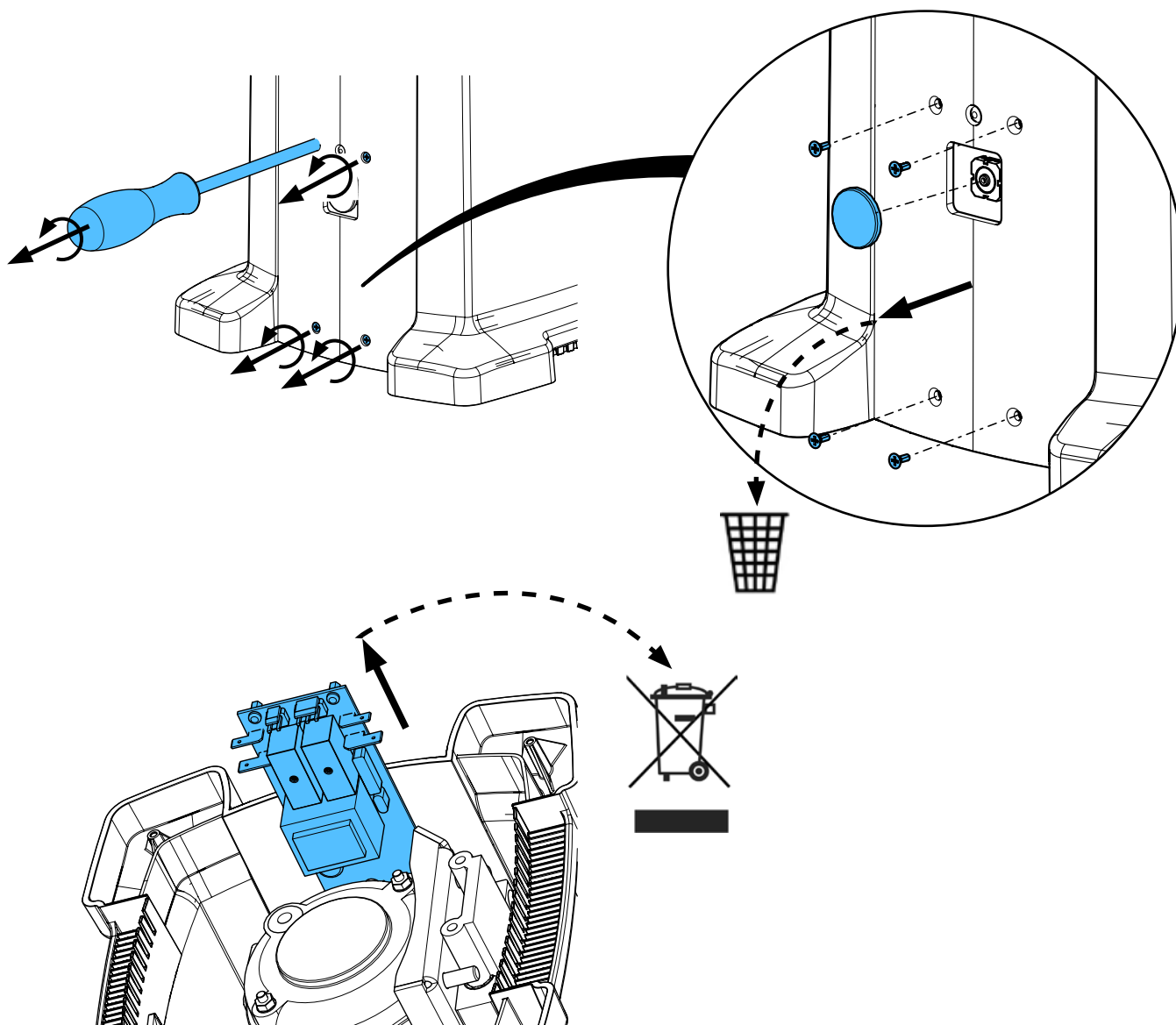



5

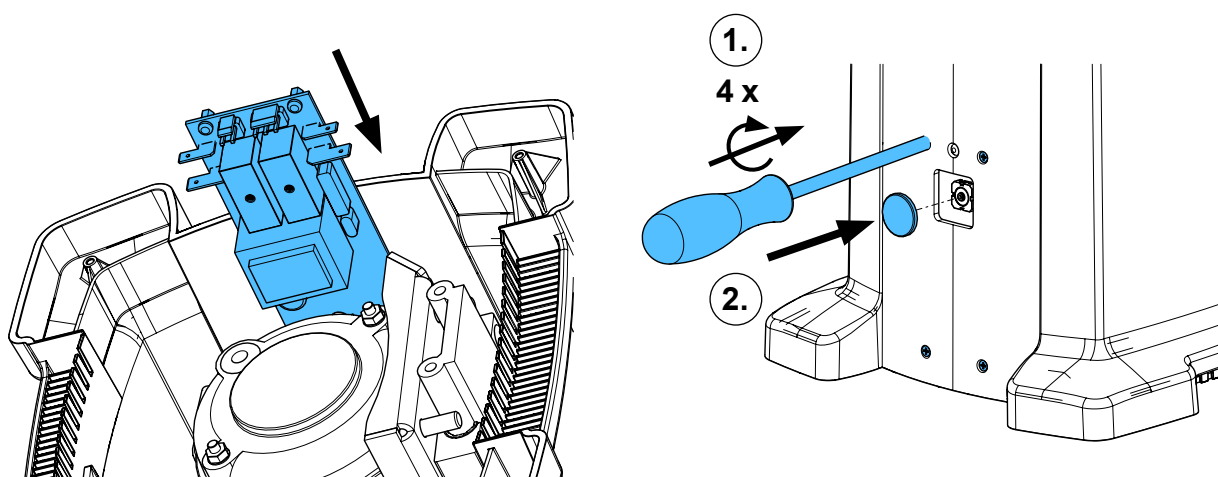



6

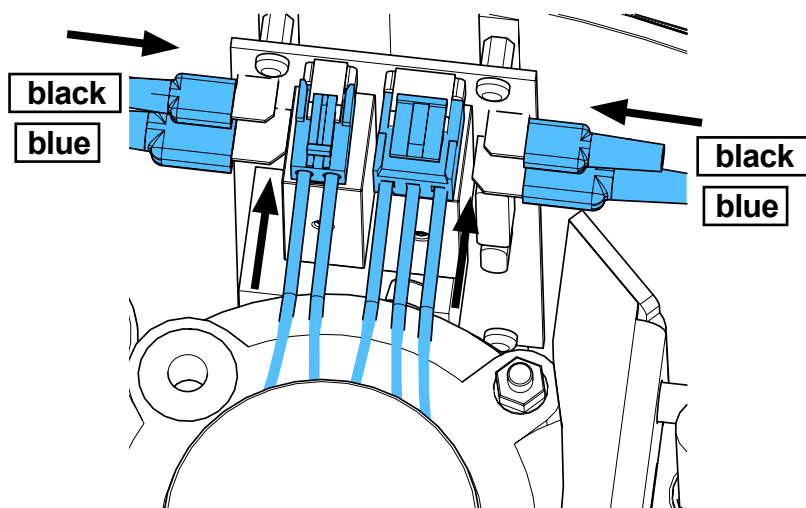

7

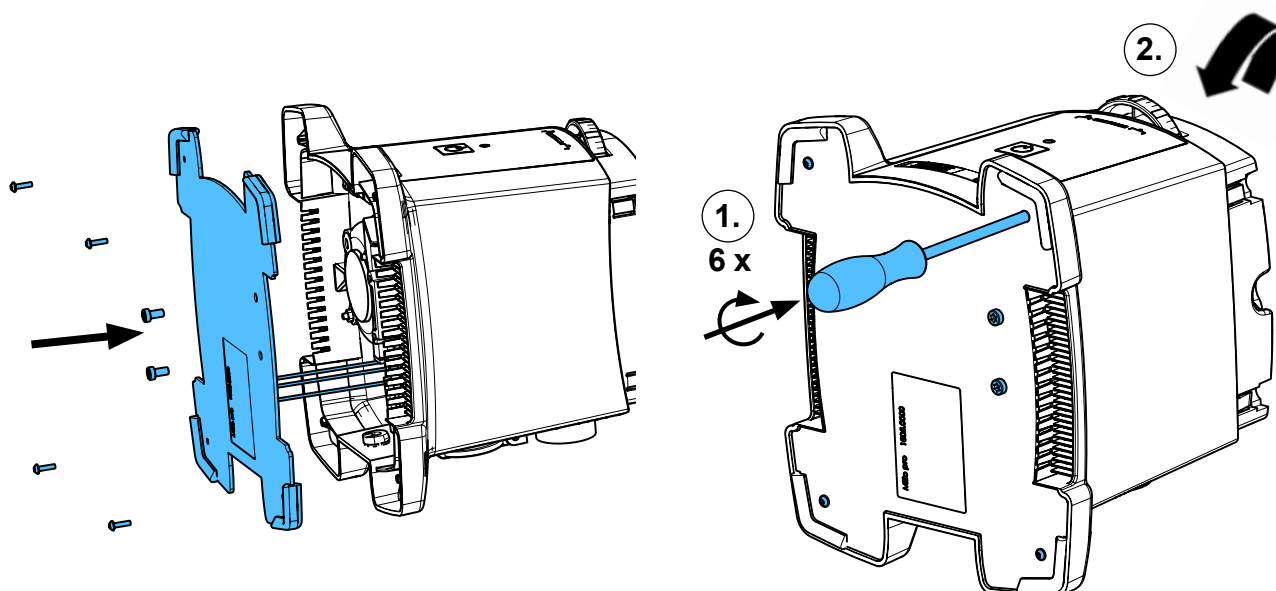
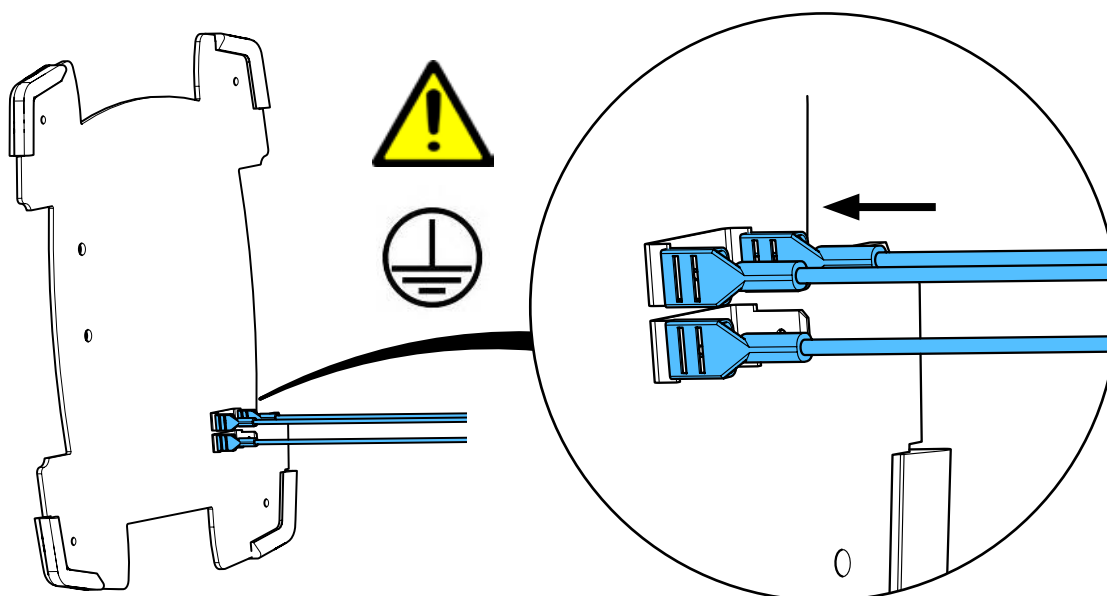


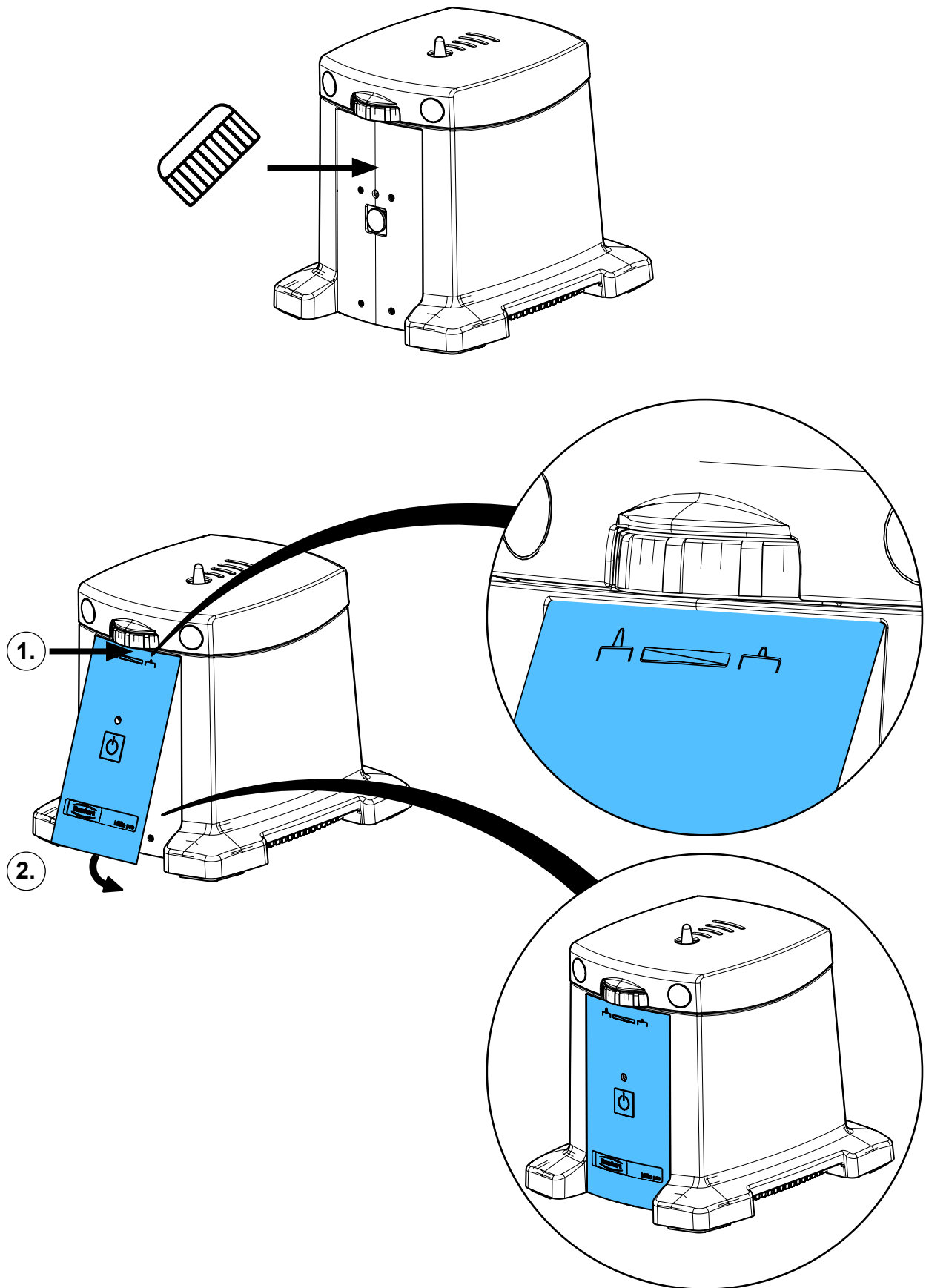

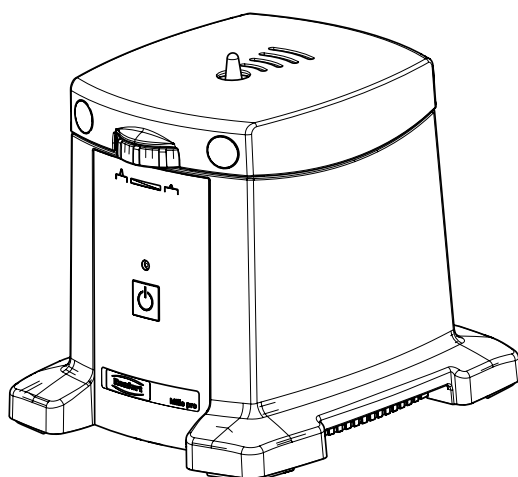


90012 1402 / 90012 1403


Motor • Engine • Moteur • Motore • Motor • Мотор

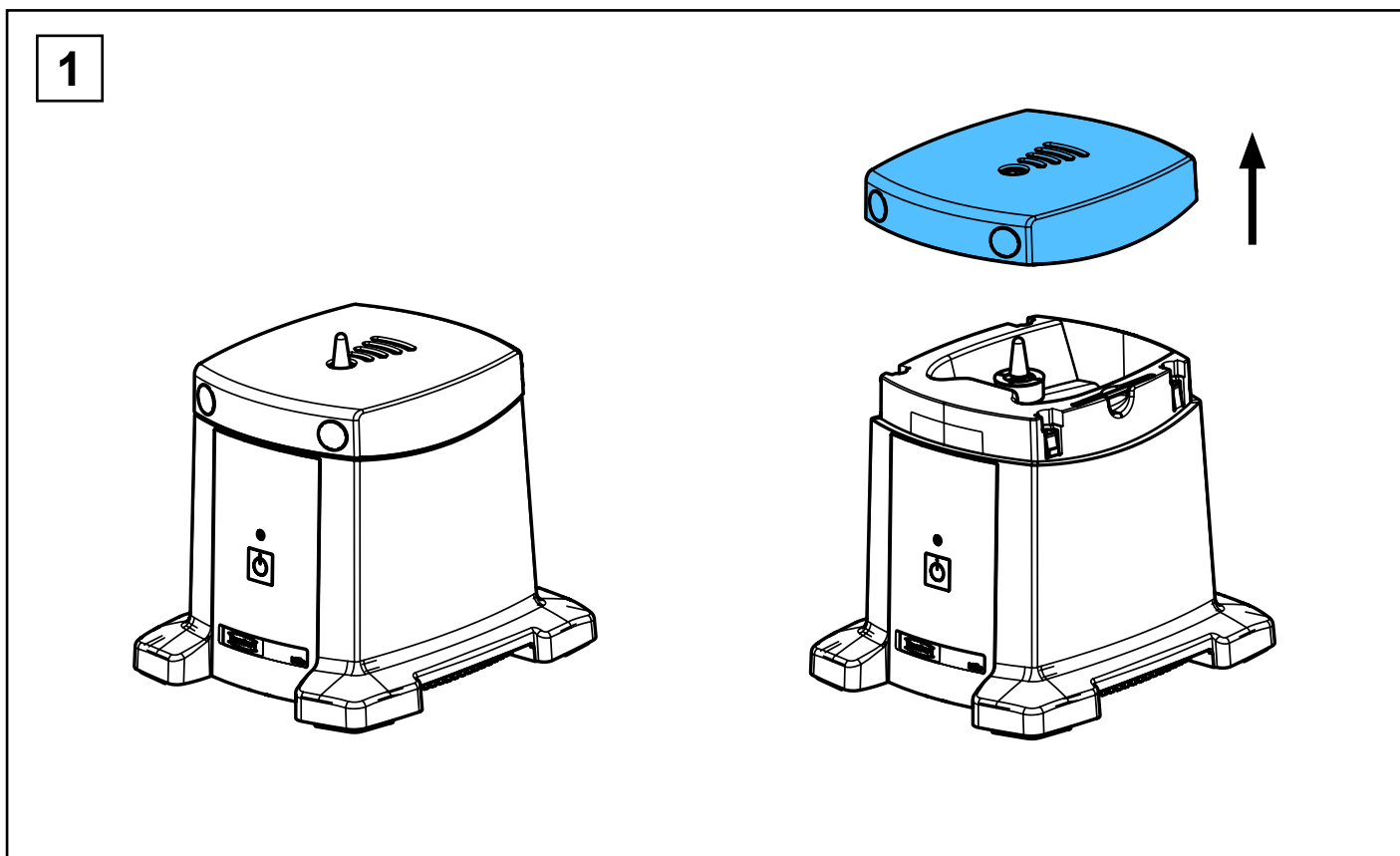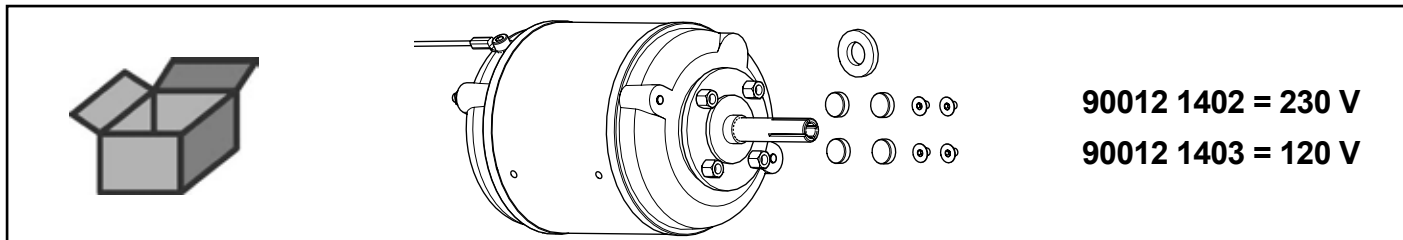


2

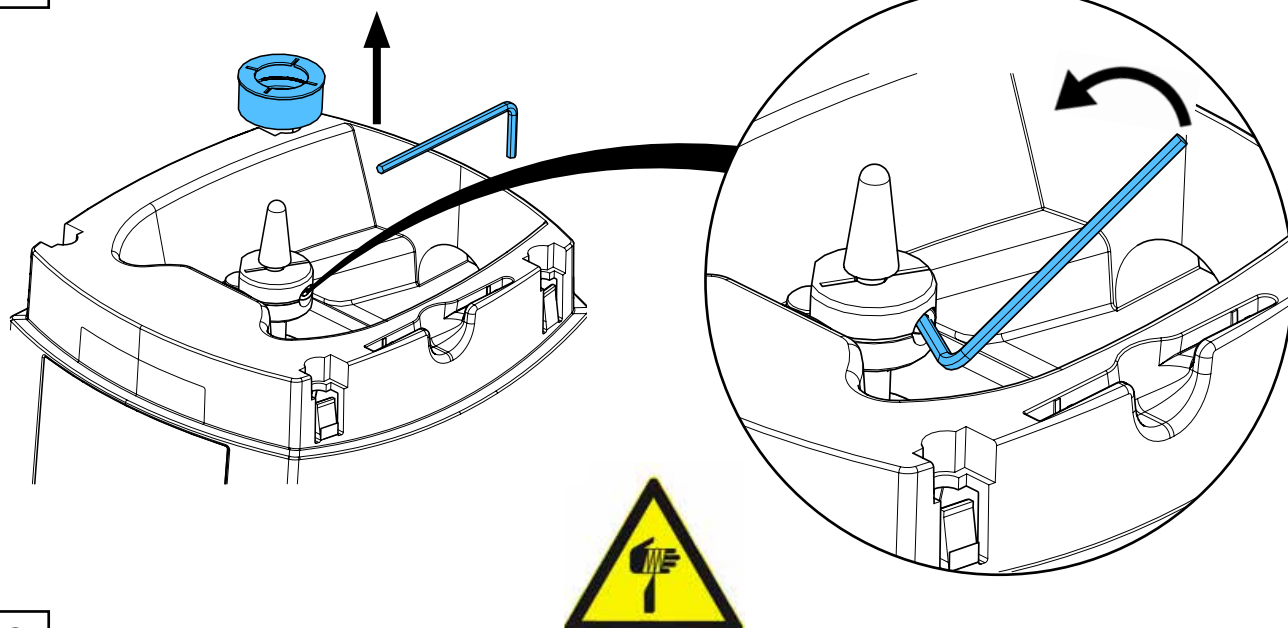



3

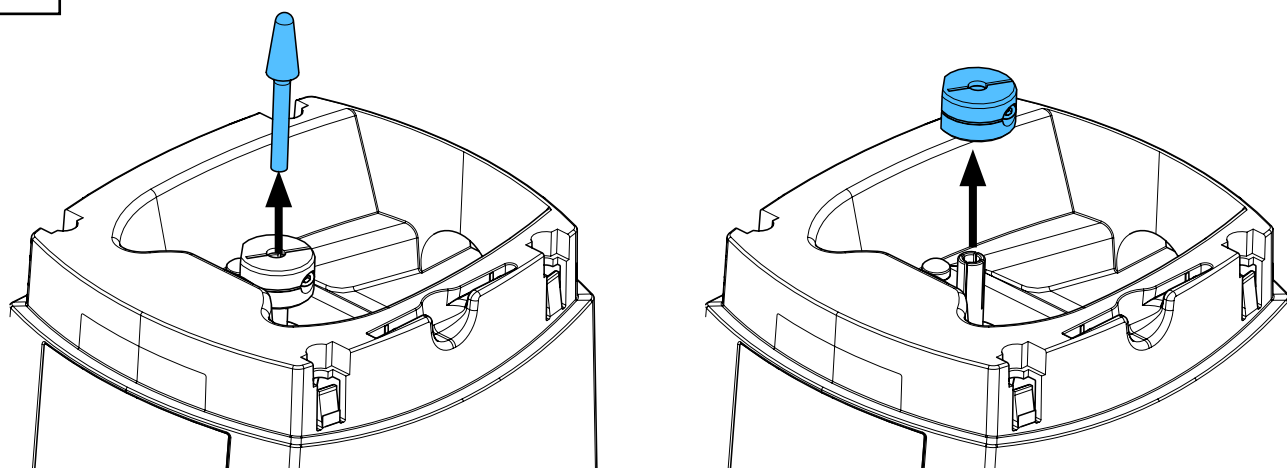



4

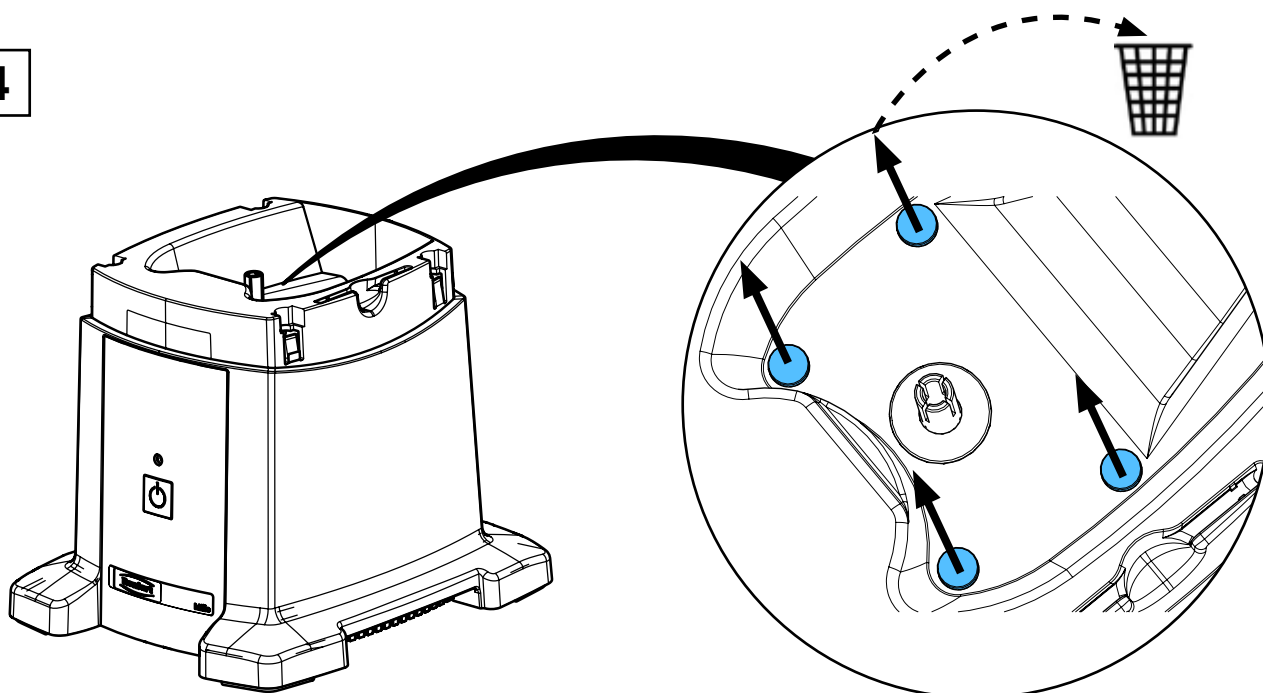



5

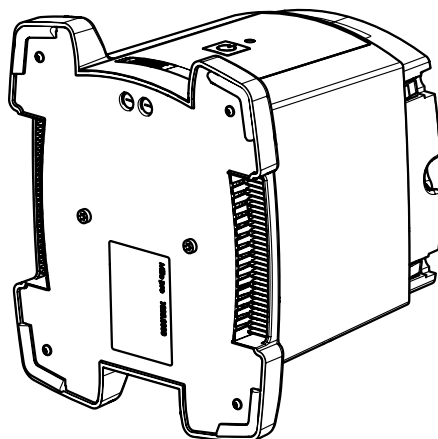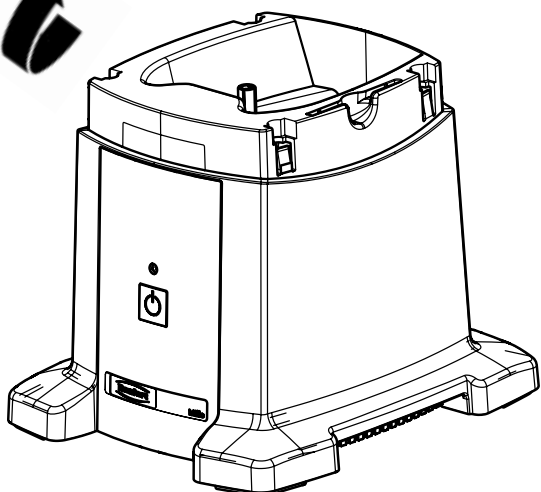


6

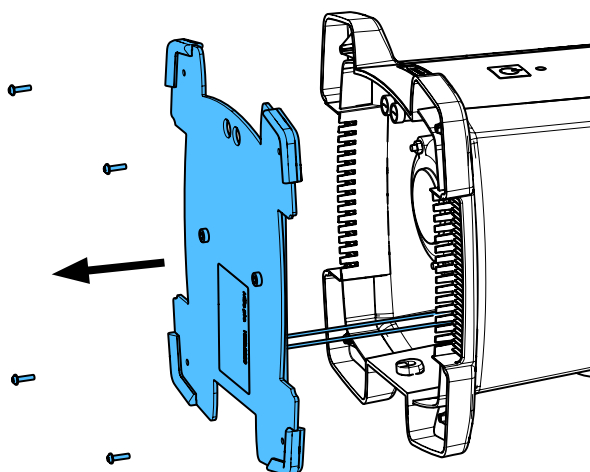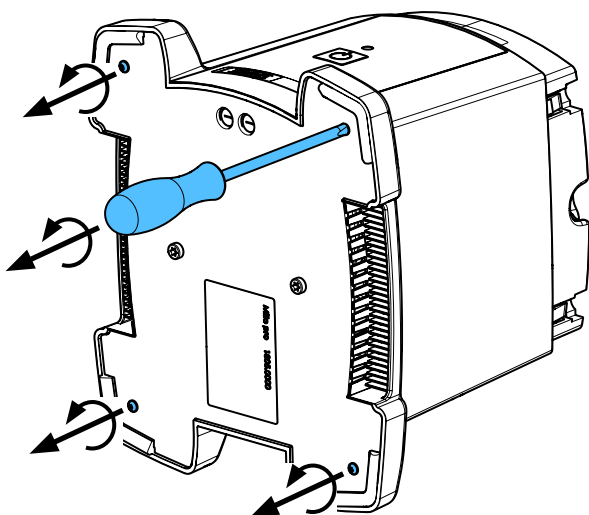


7

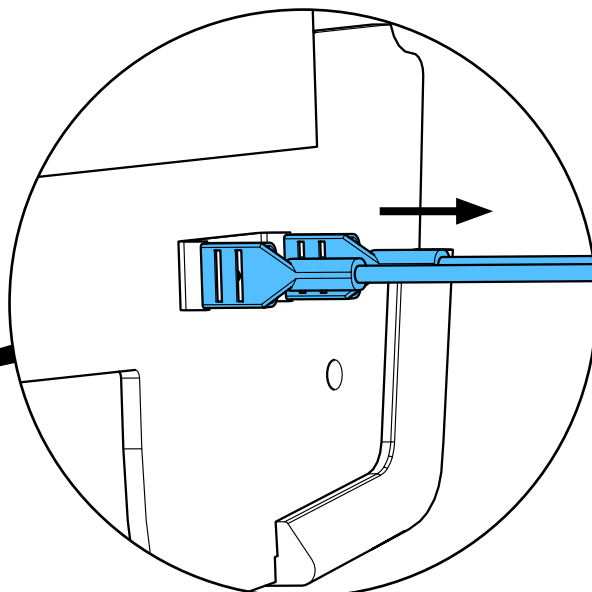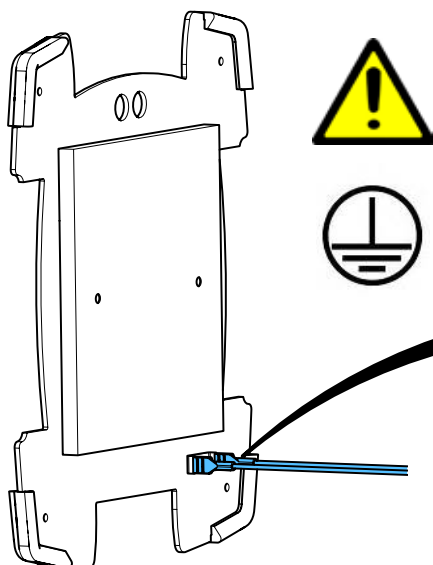


8

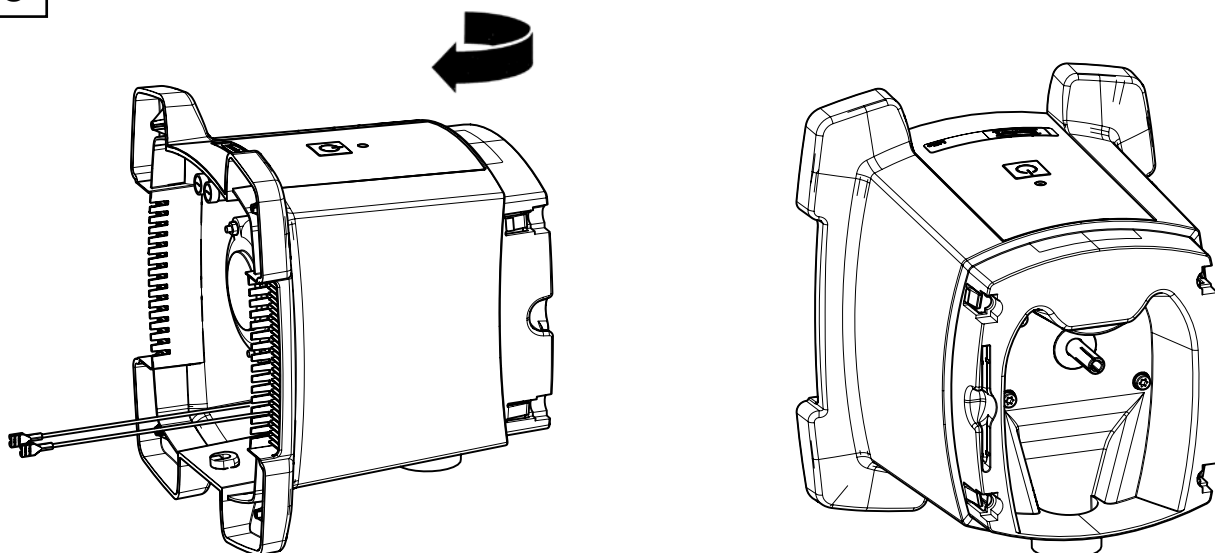

9

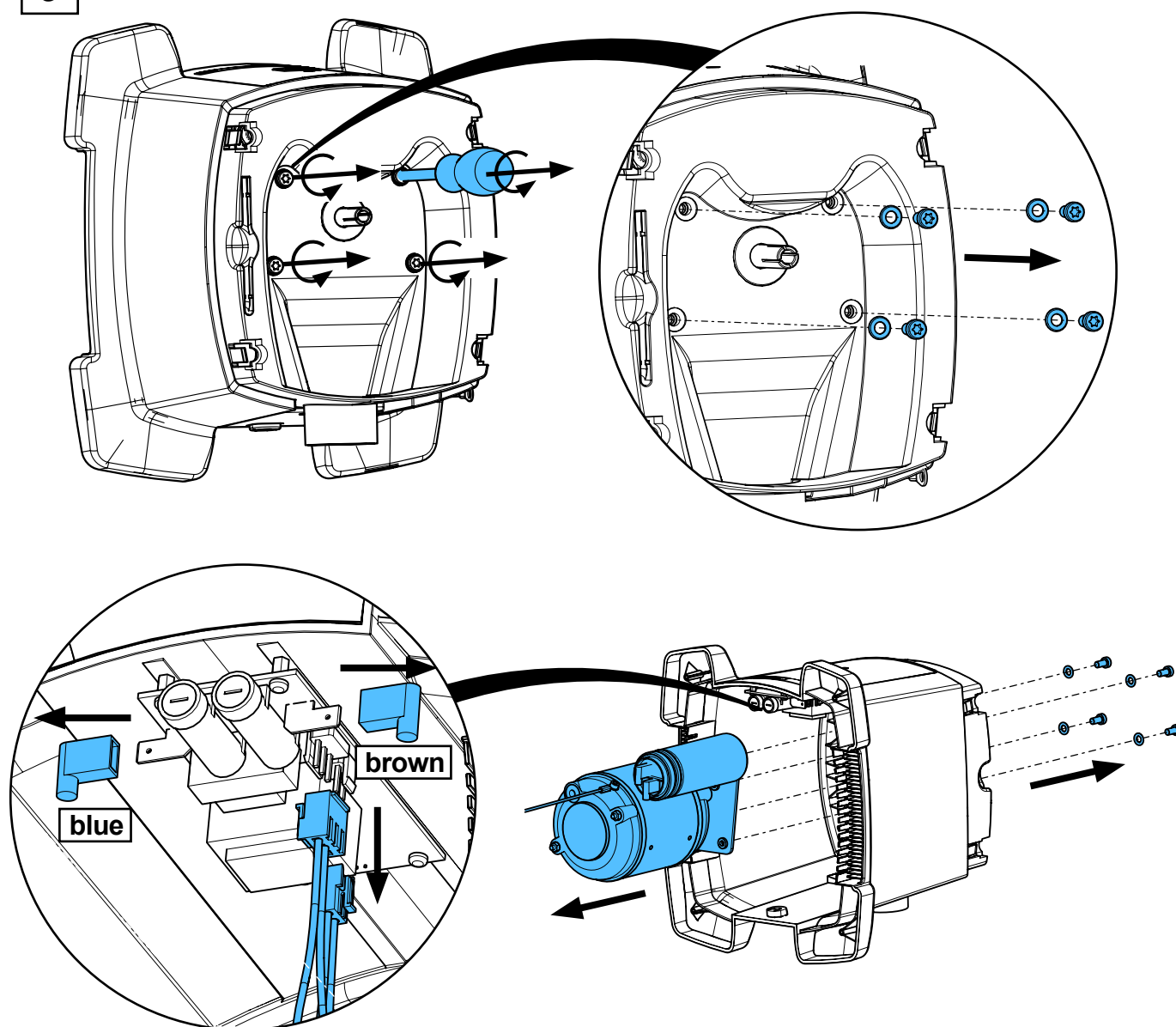







10

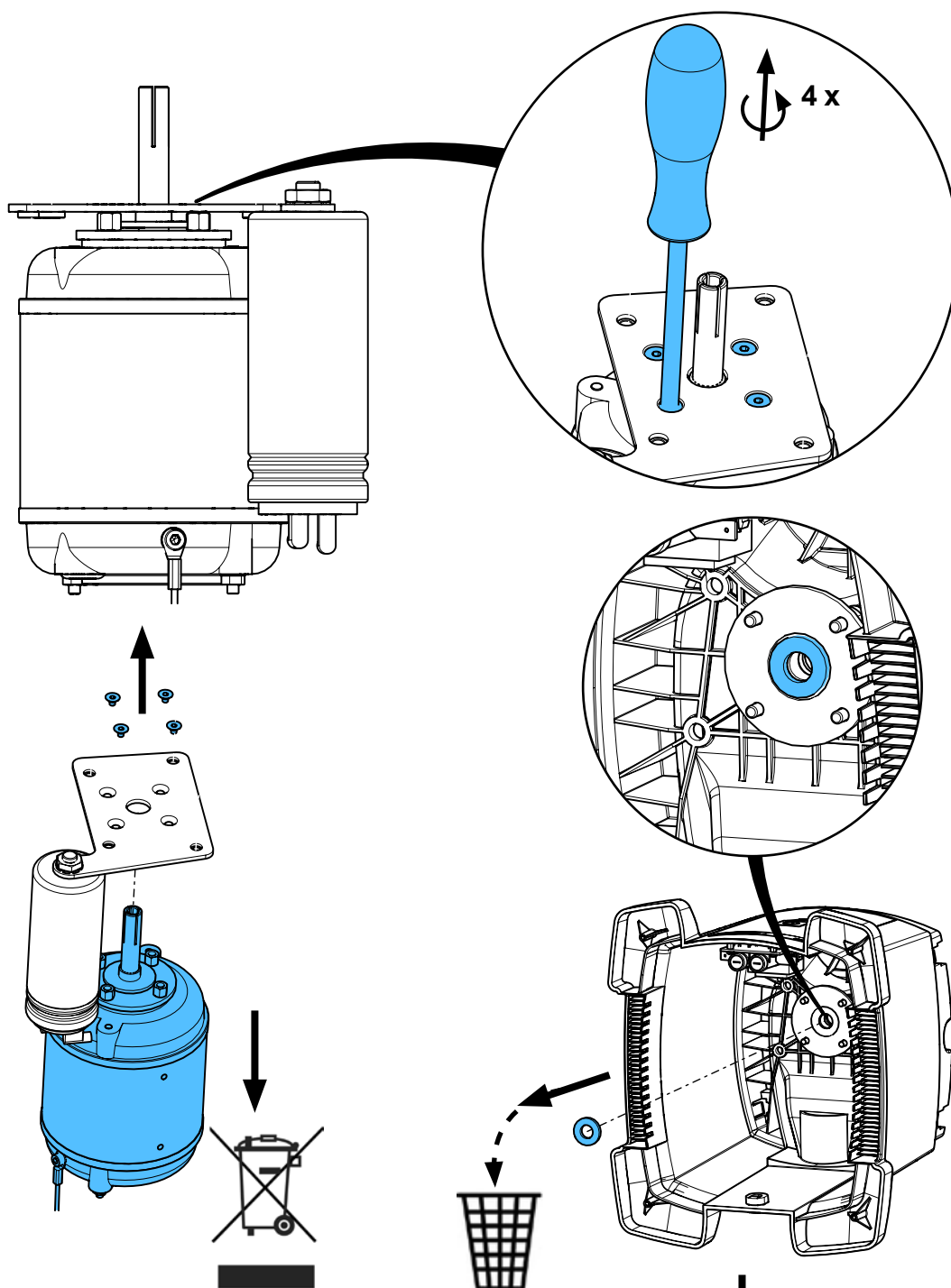


11





12

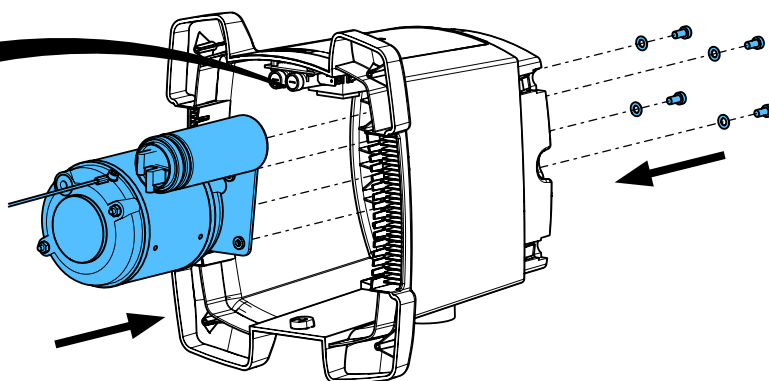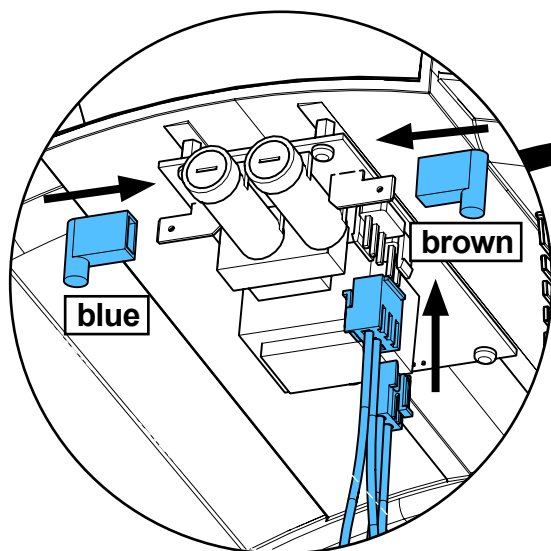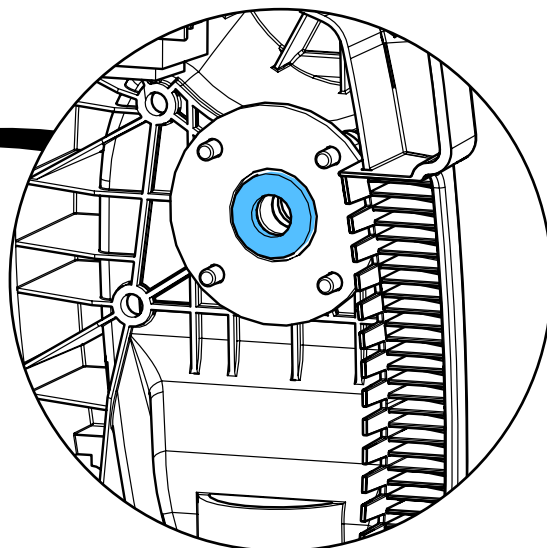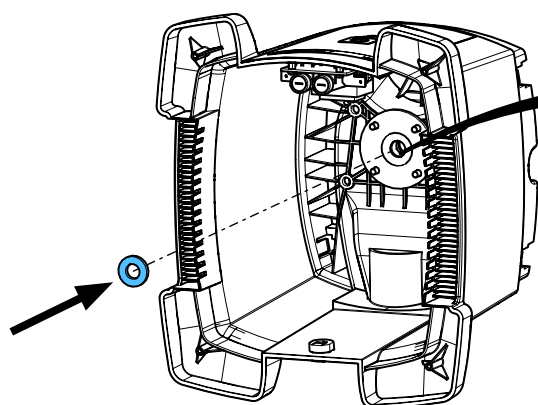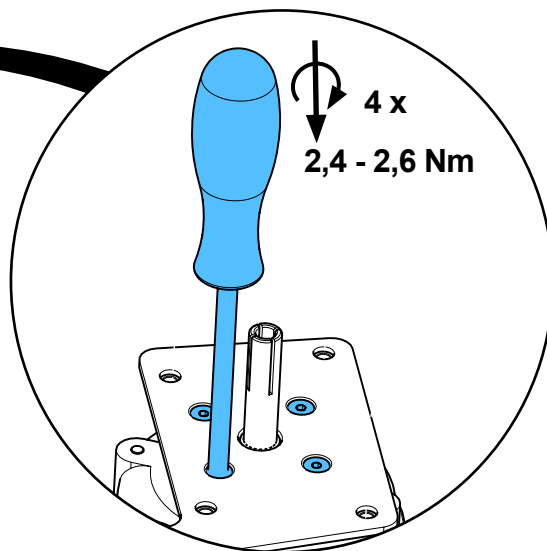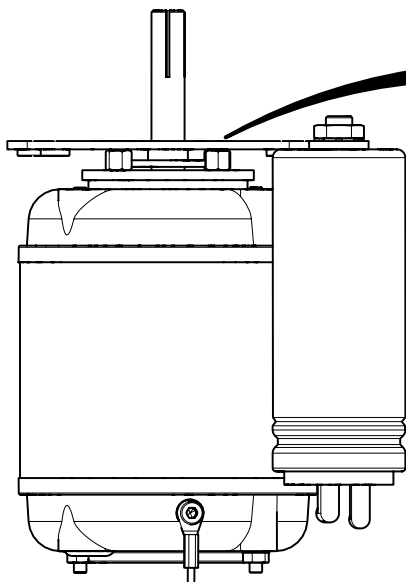


13

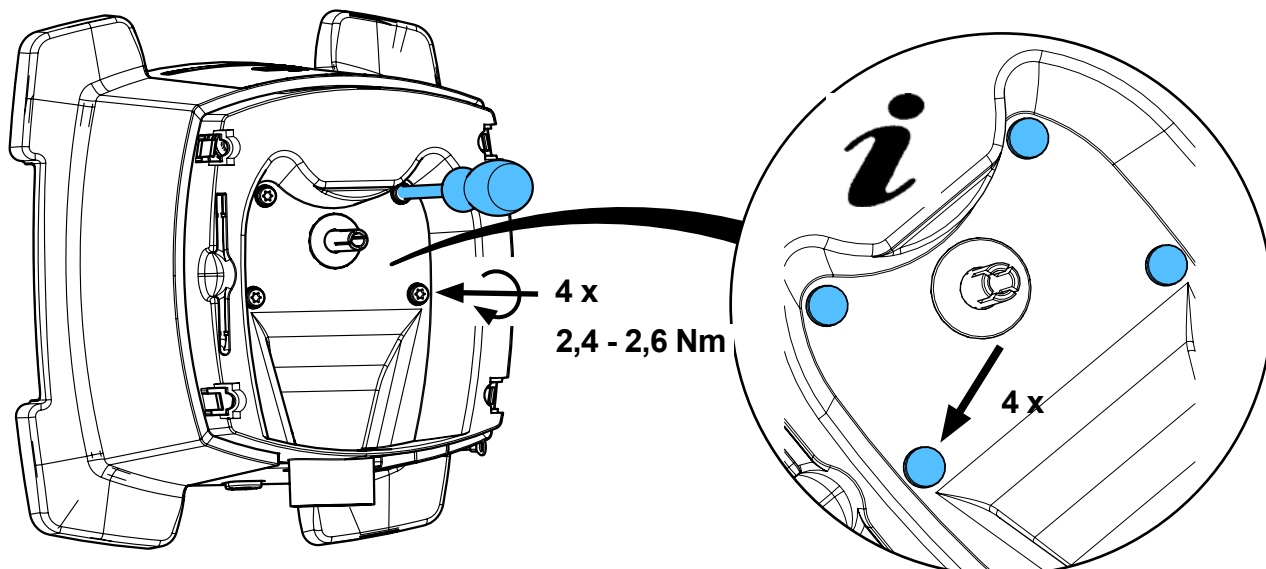



14

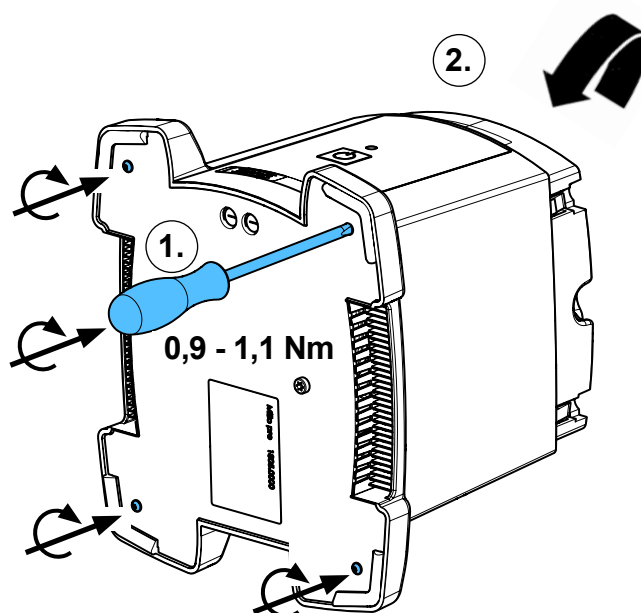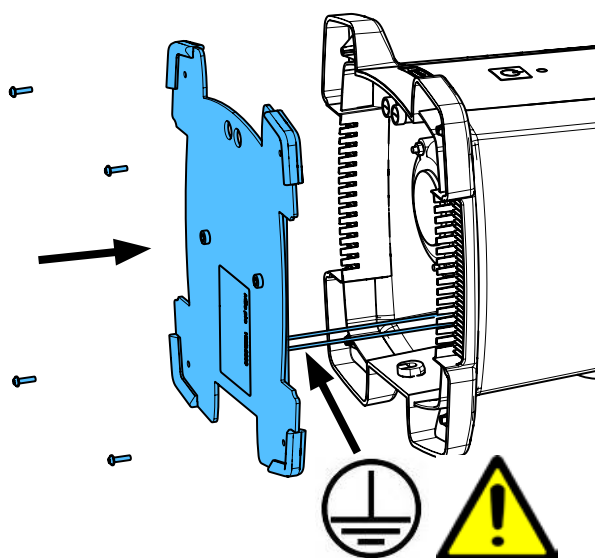



15

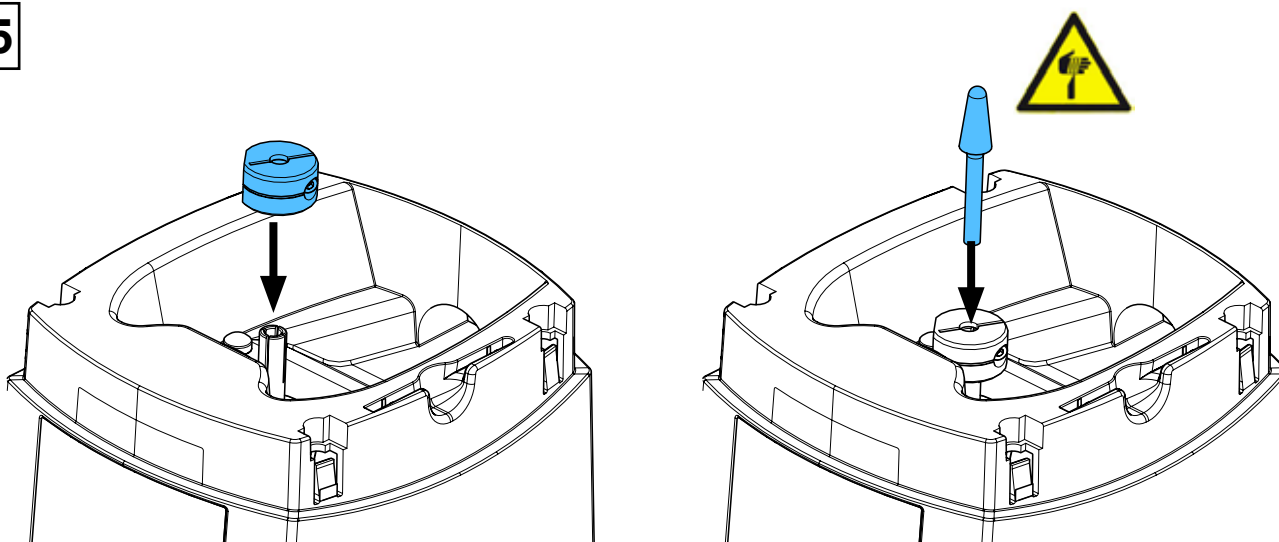




16

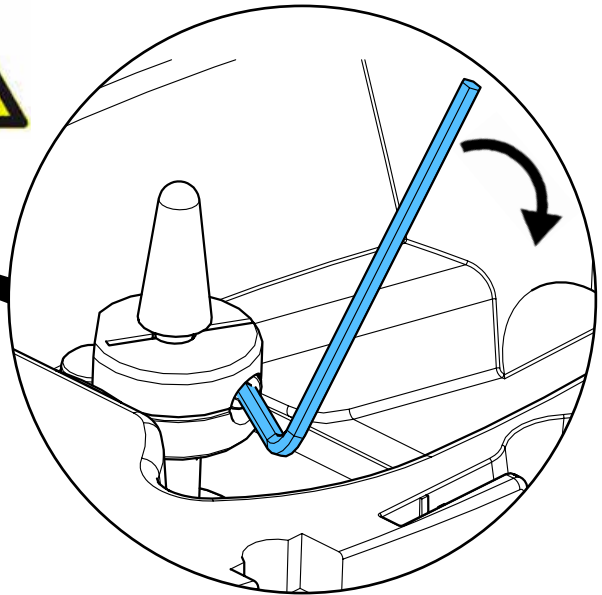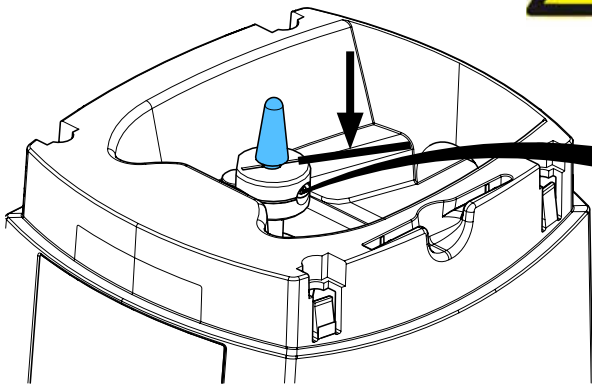

17

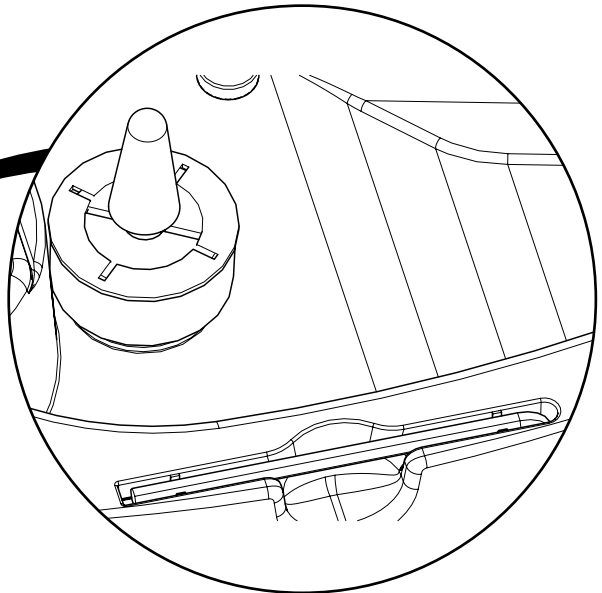
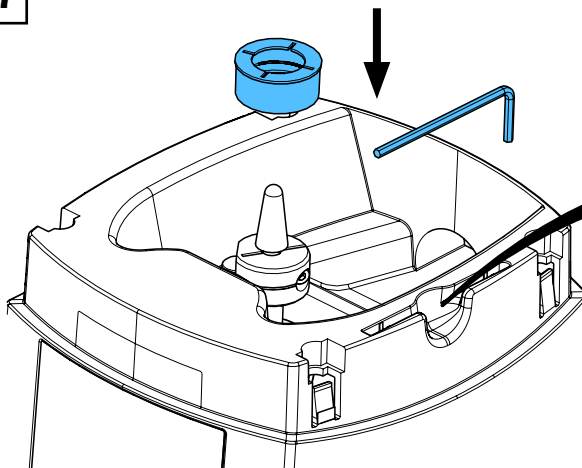


18

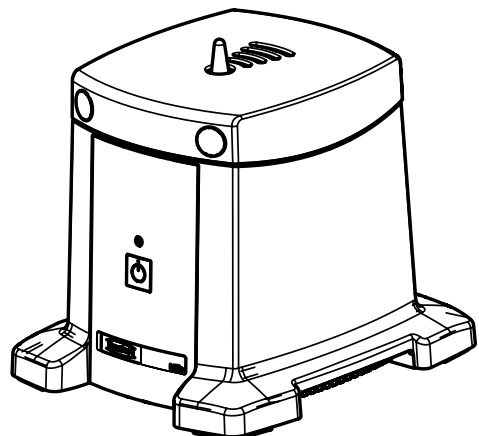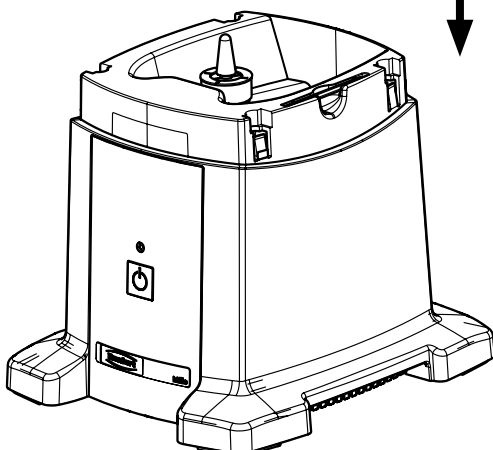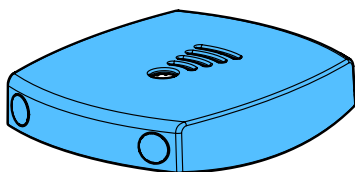


90002 1402 / 90002 1403


Motor • Engine • Moteur • Motore • Motor • Мотор

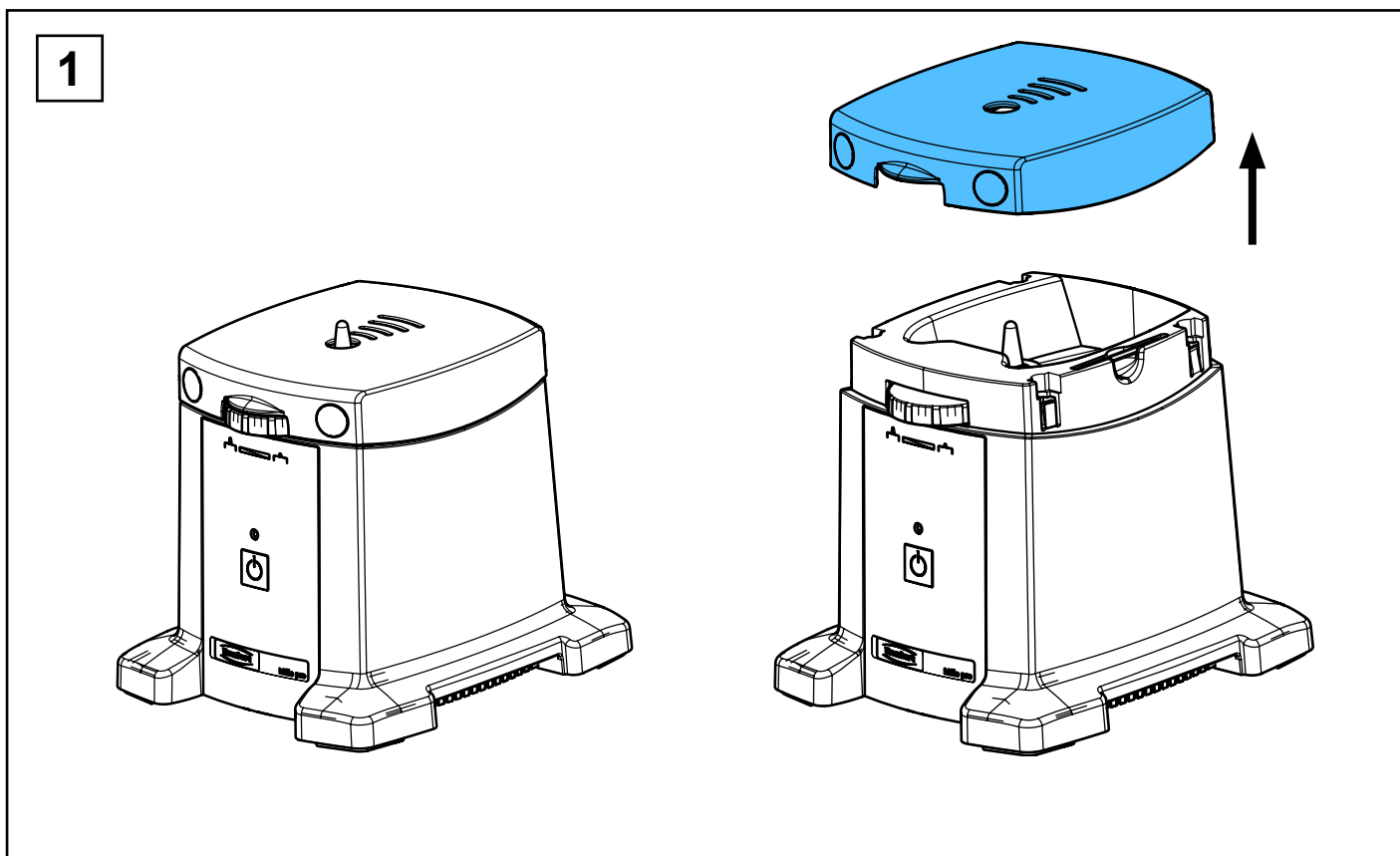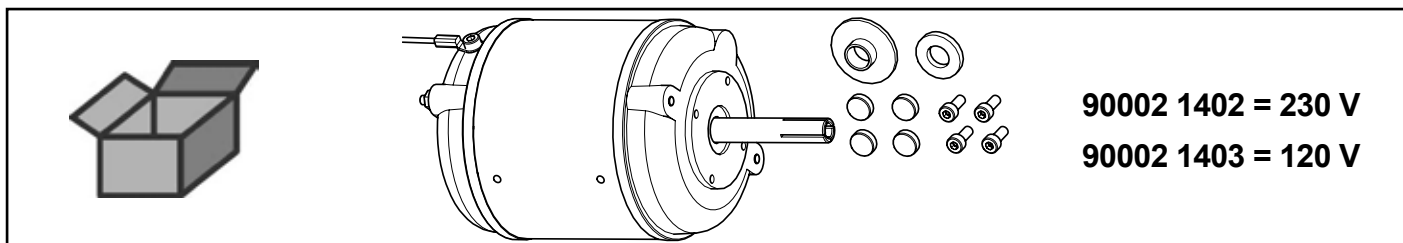


2

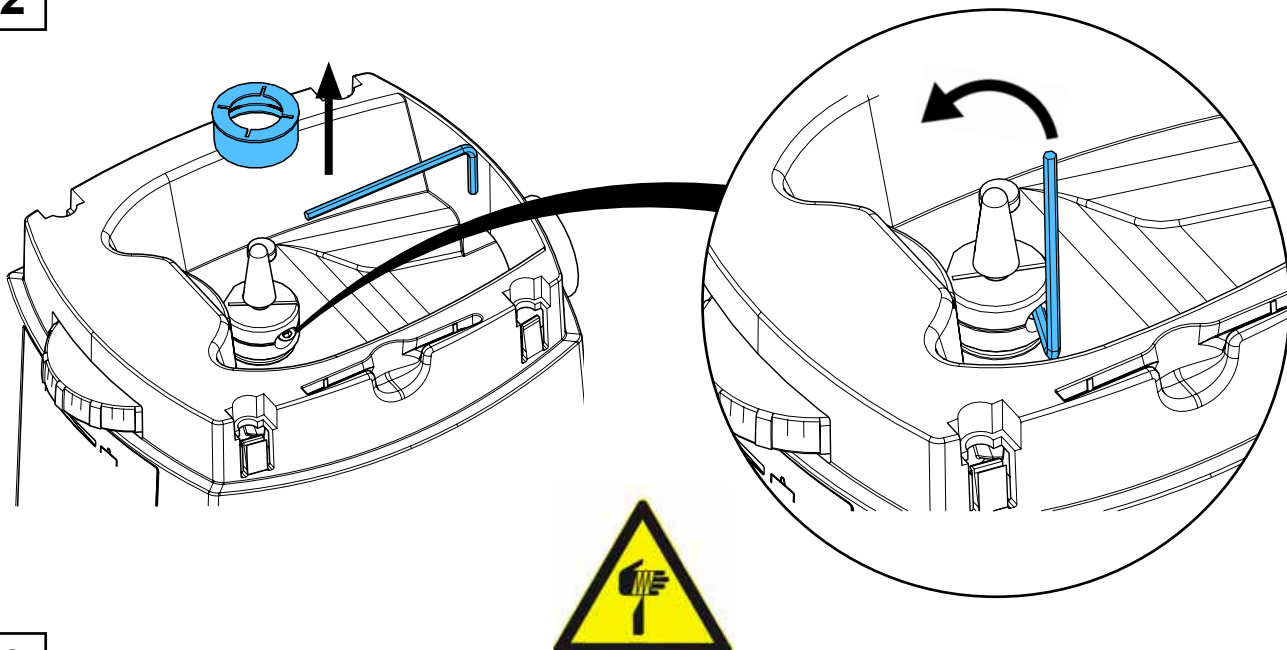



3

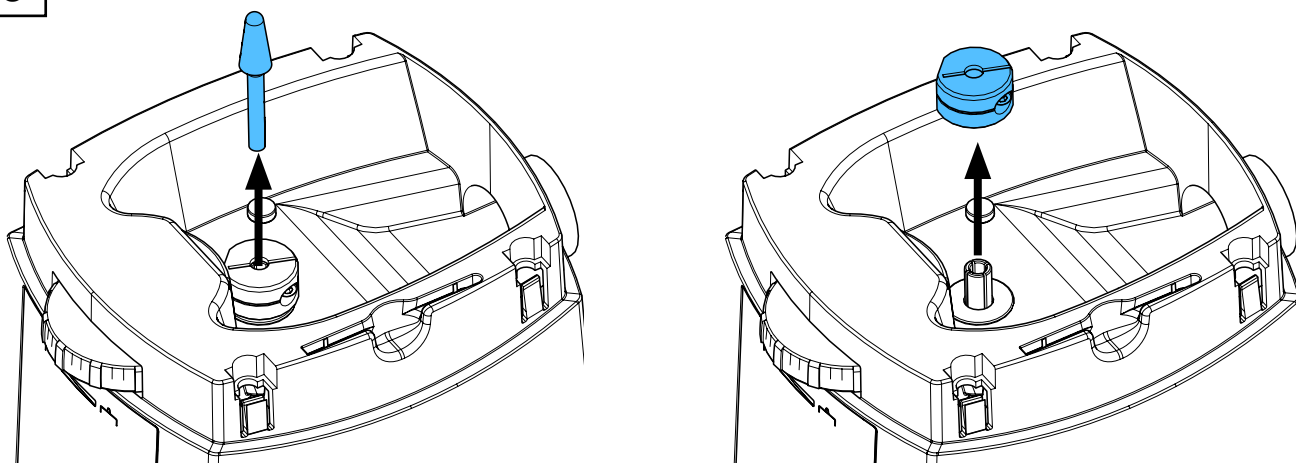



4

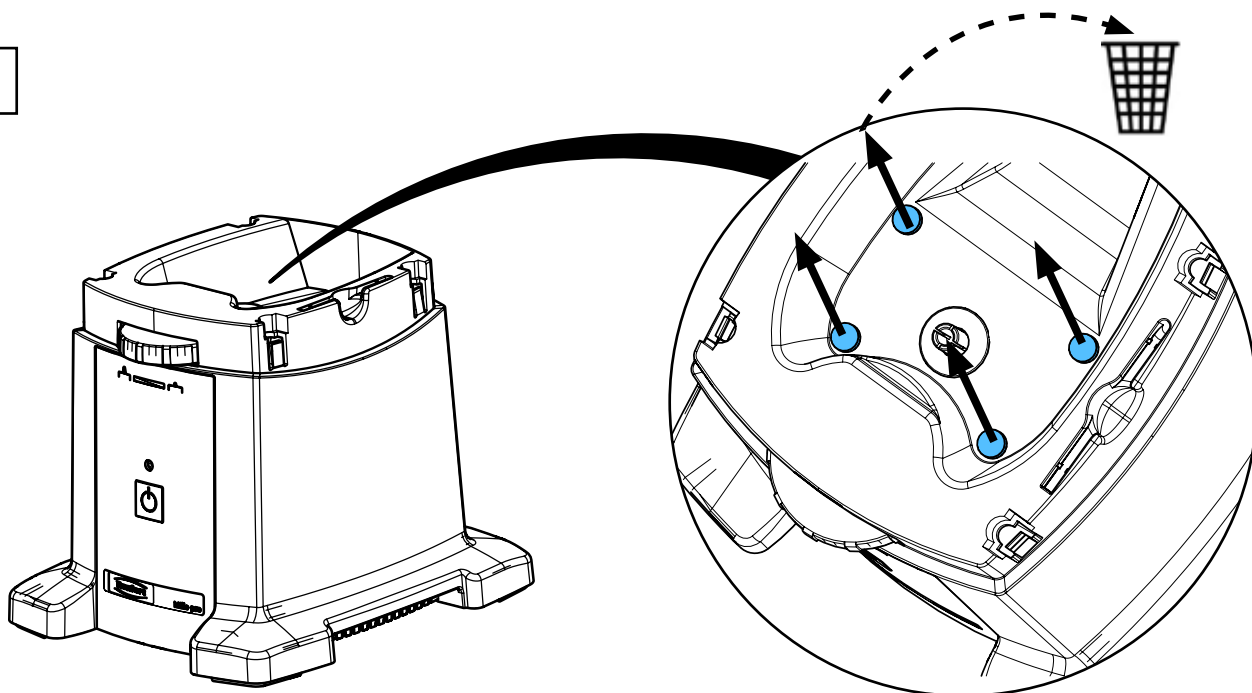



5

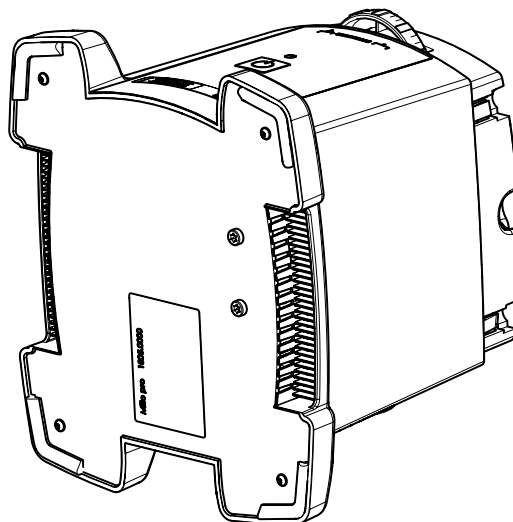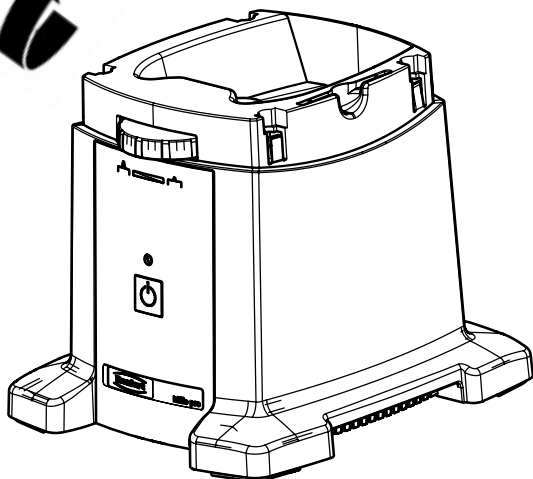


6

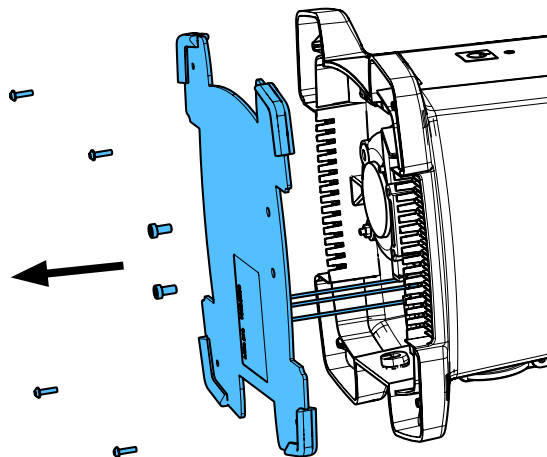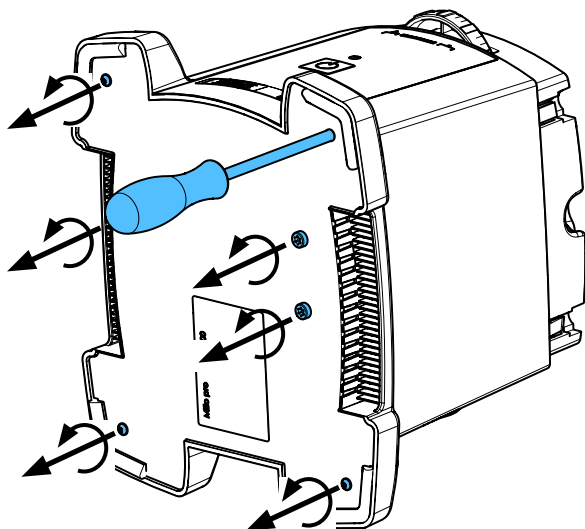


7

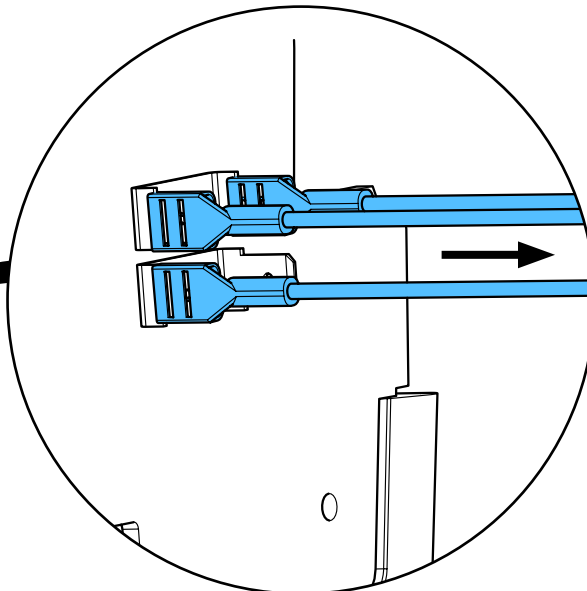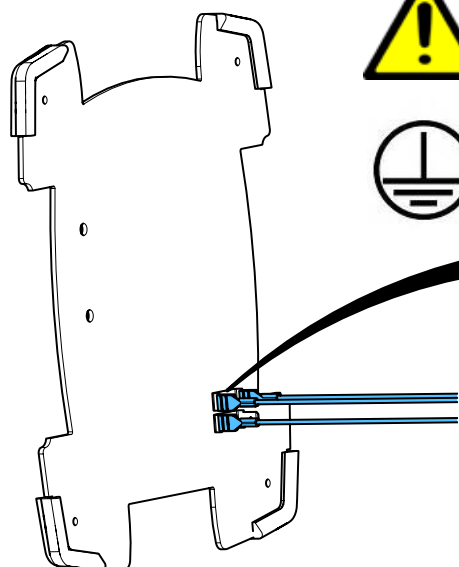


8

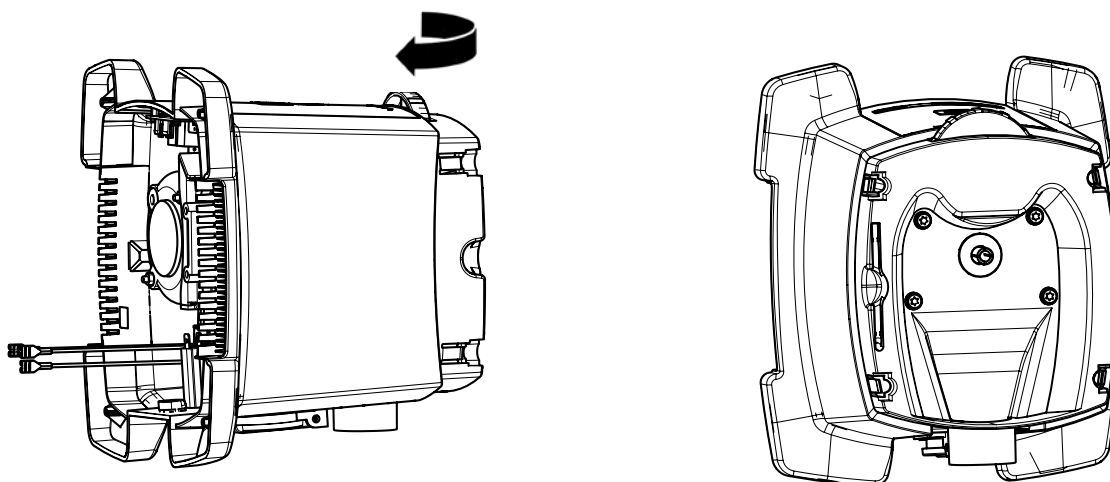


9

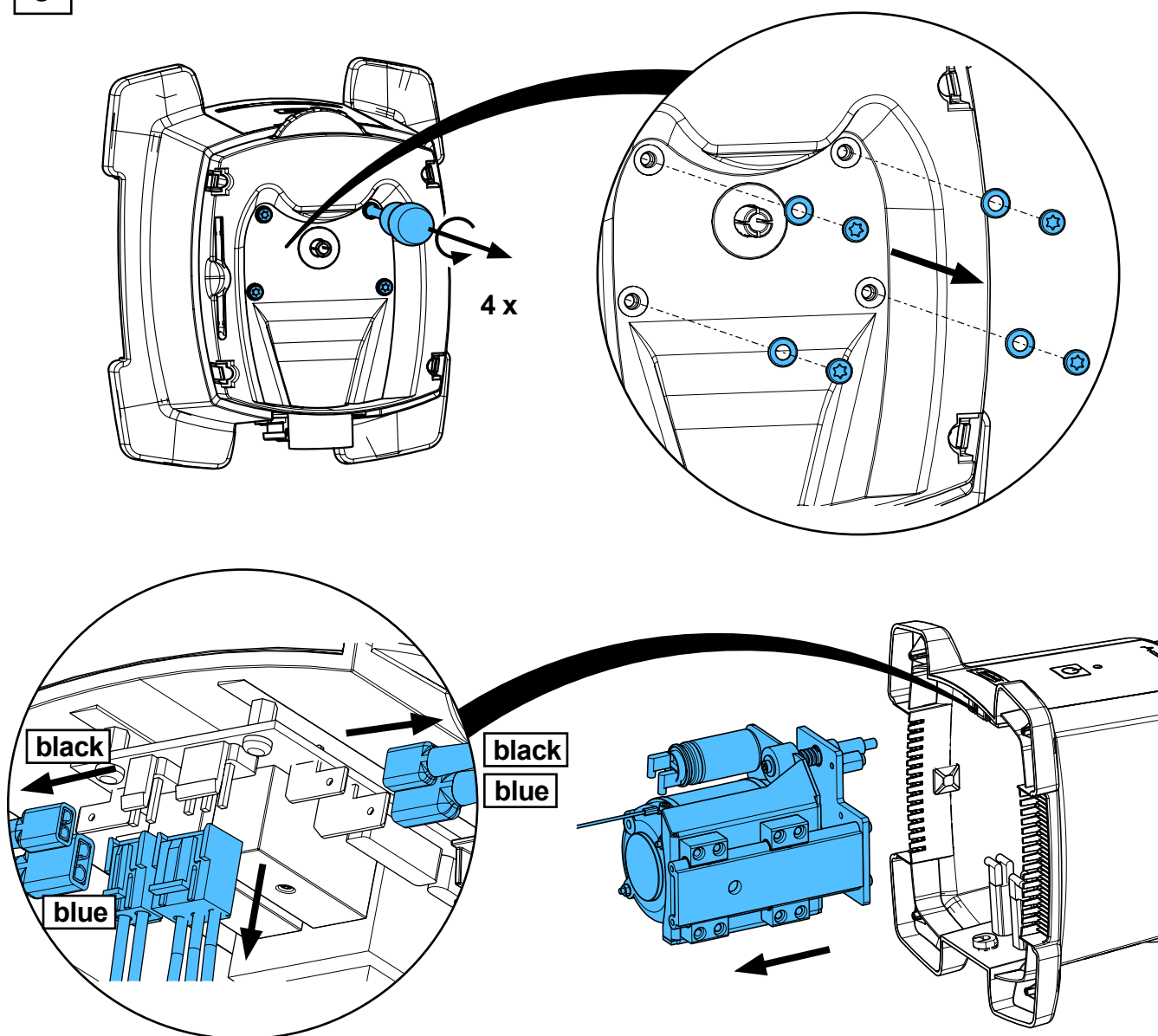


10

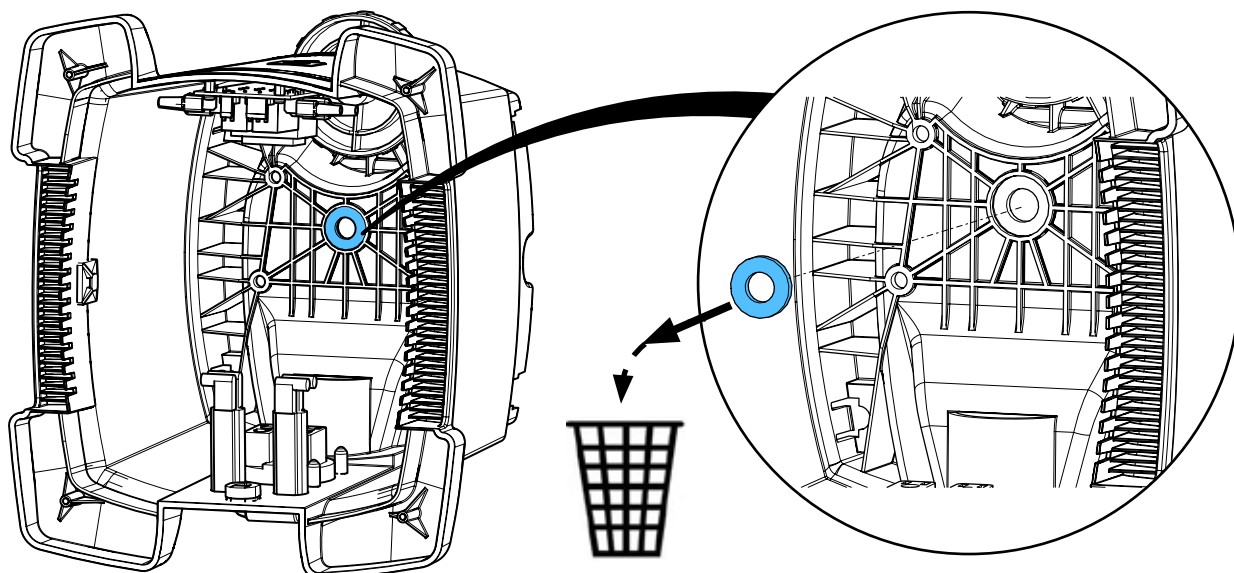


11

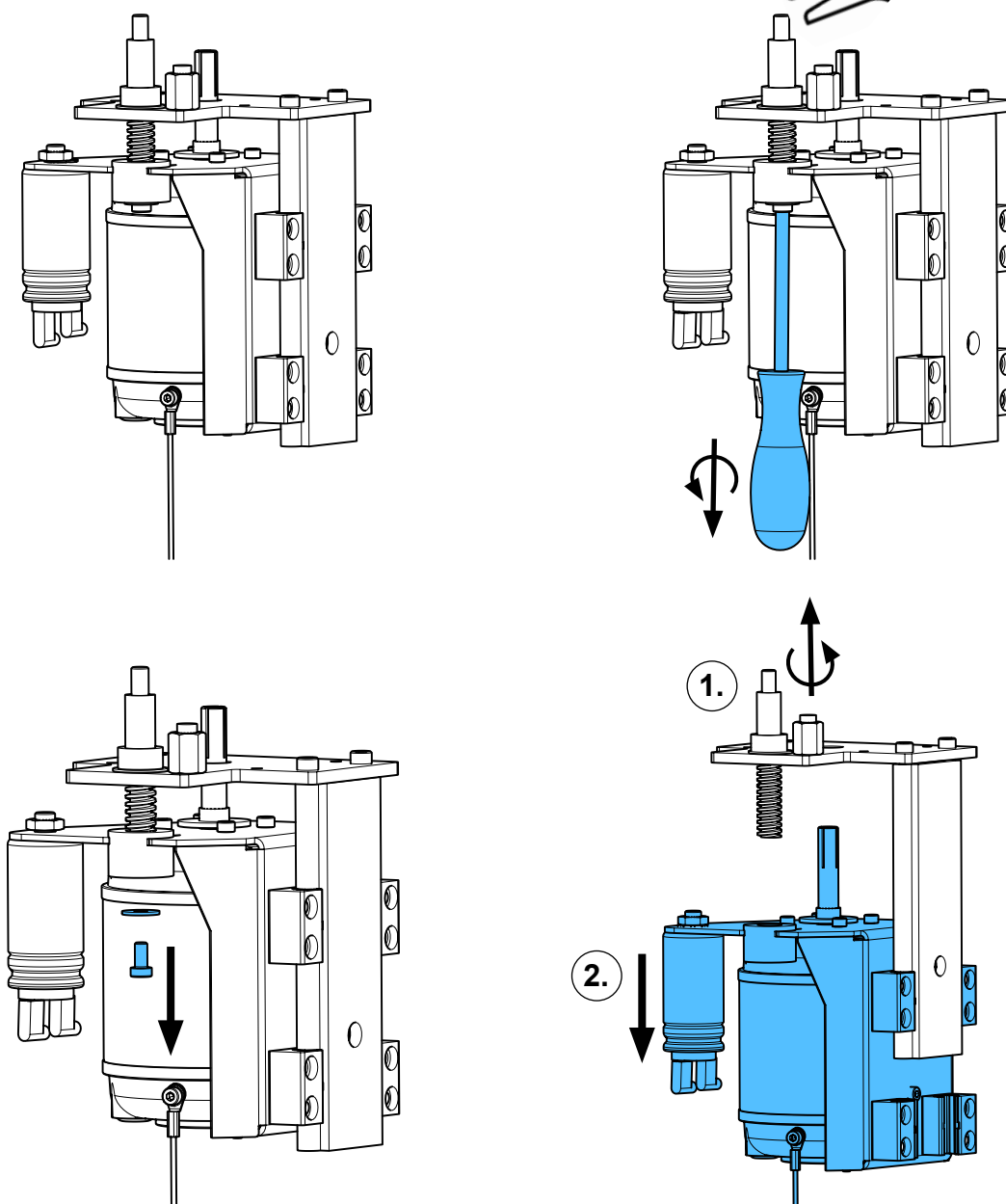


12

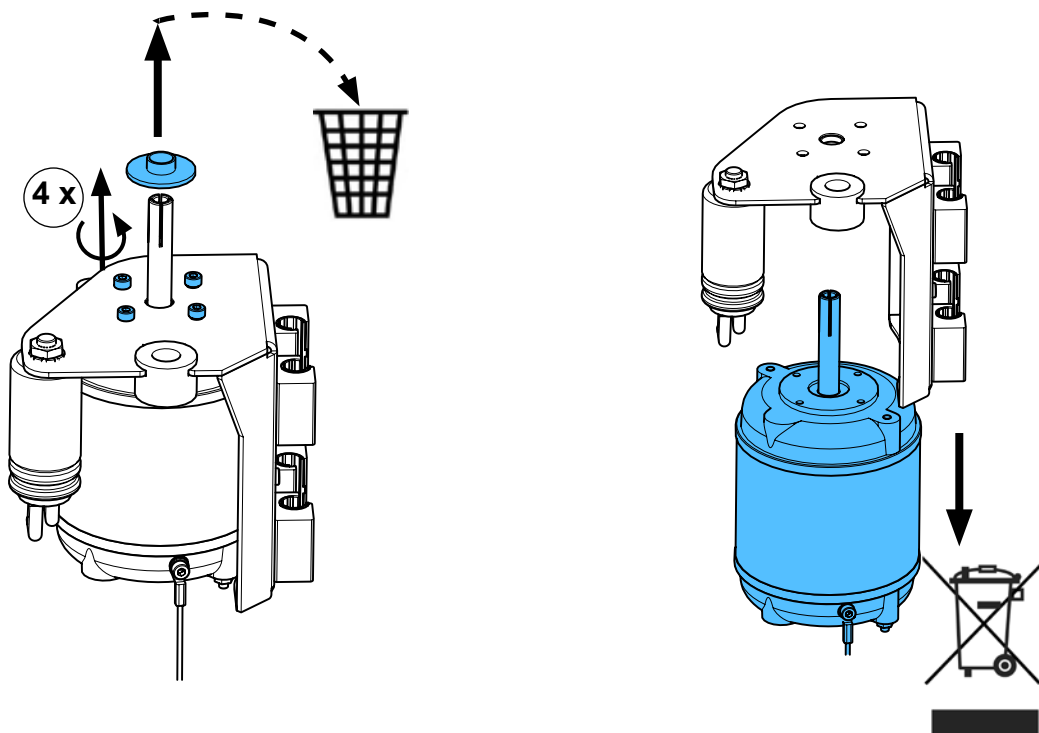






13

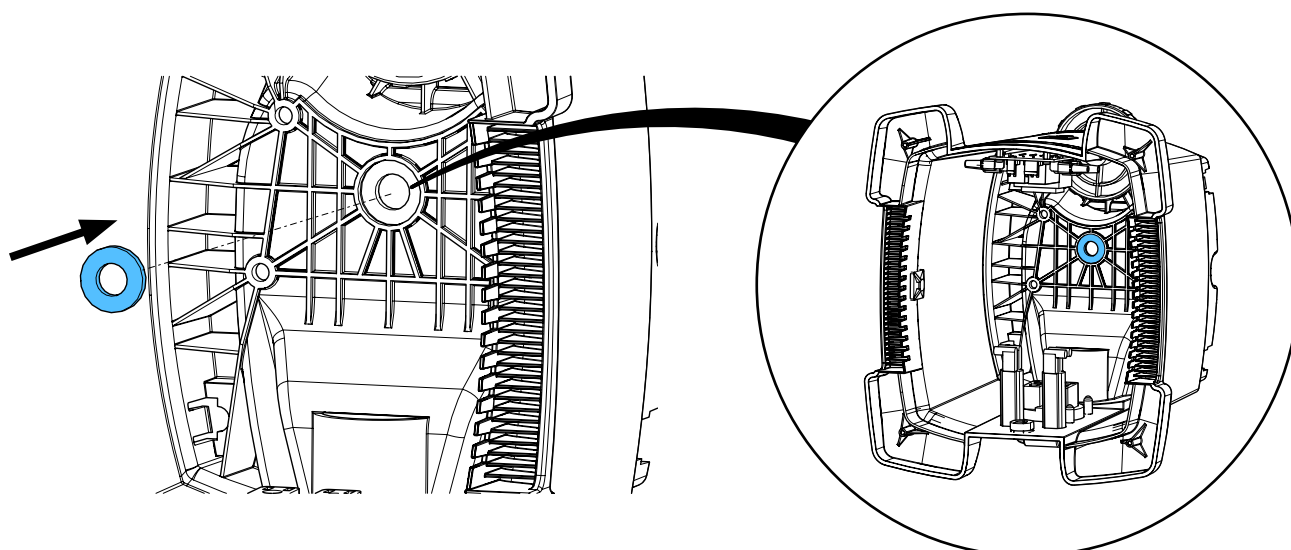


14

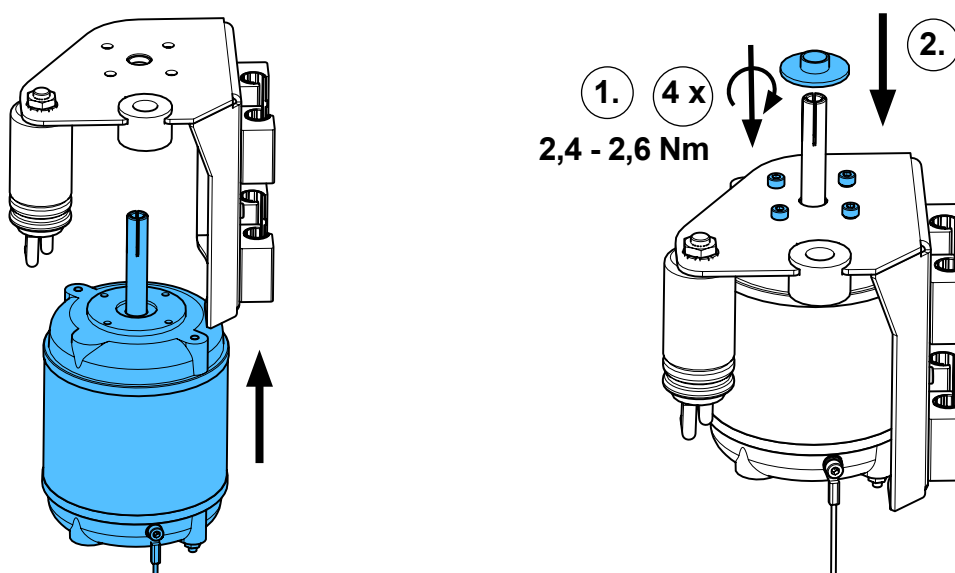


15

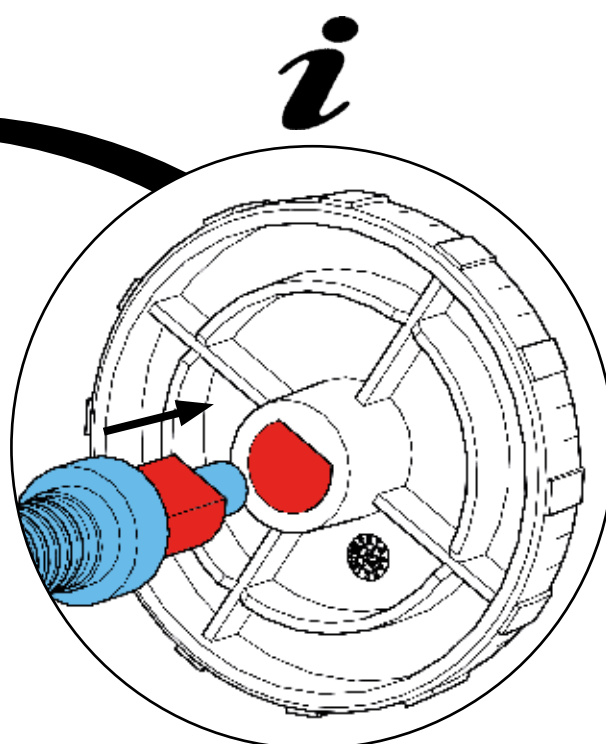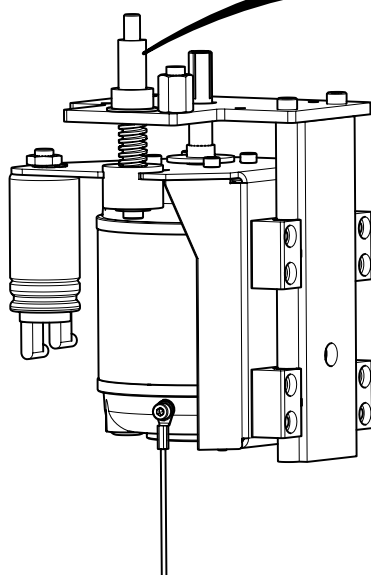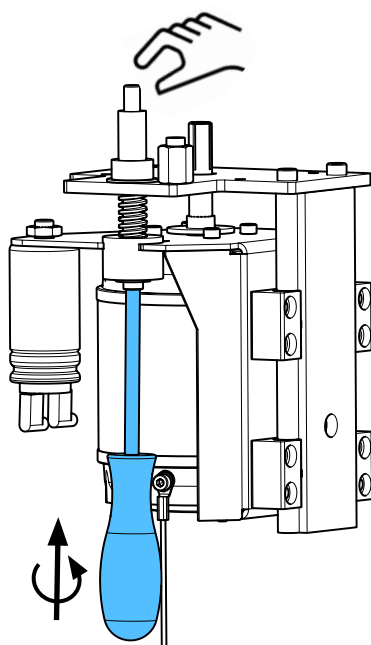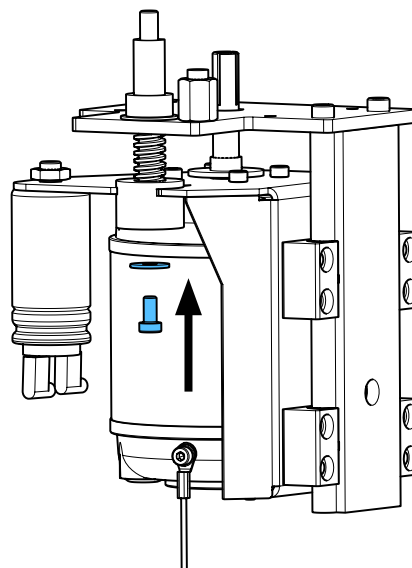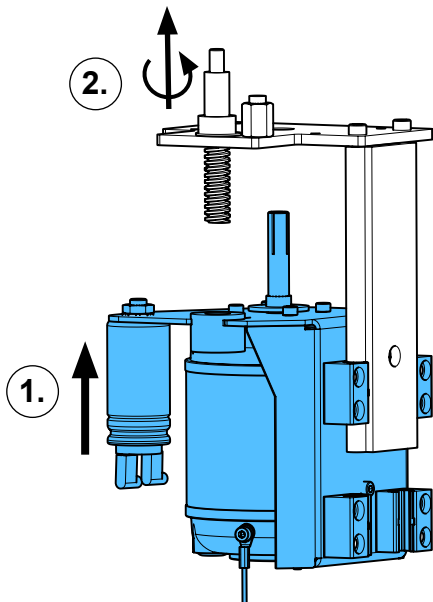


16

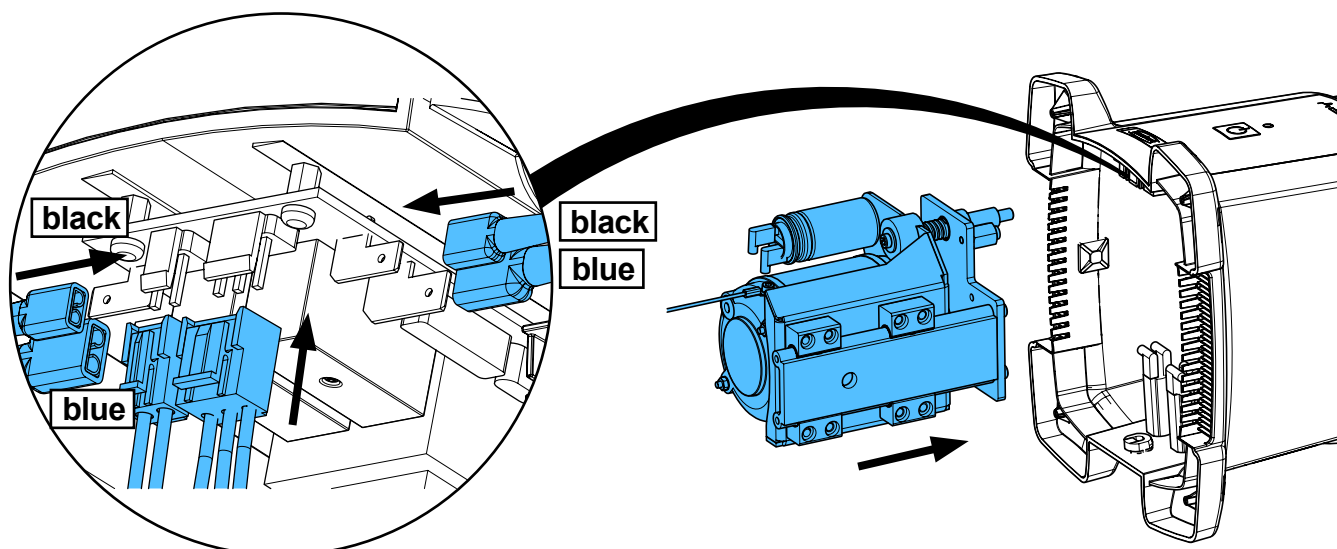


17

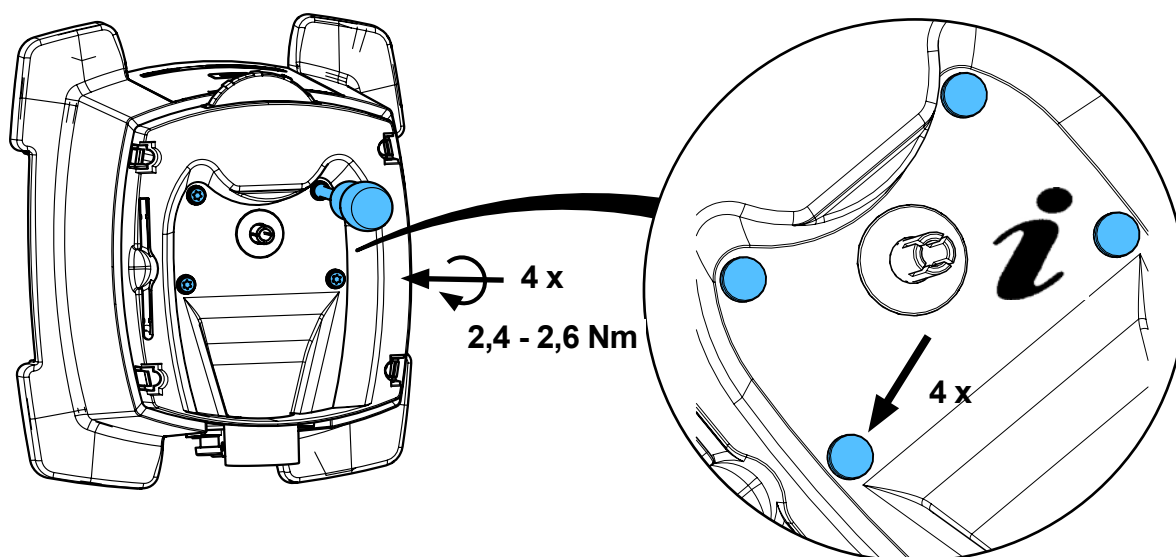


18

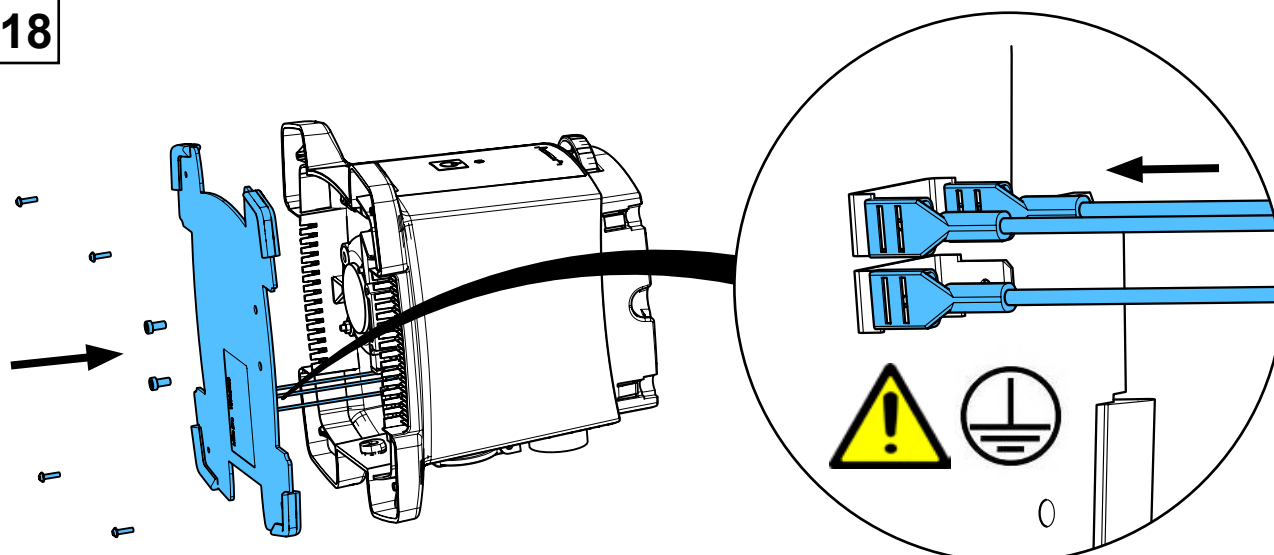


19

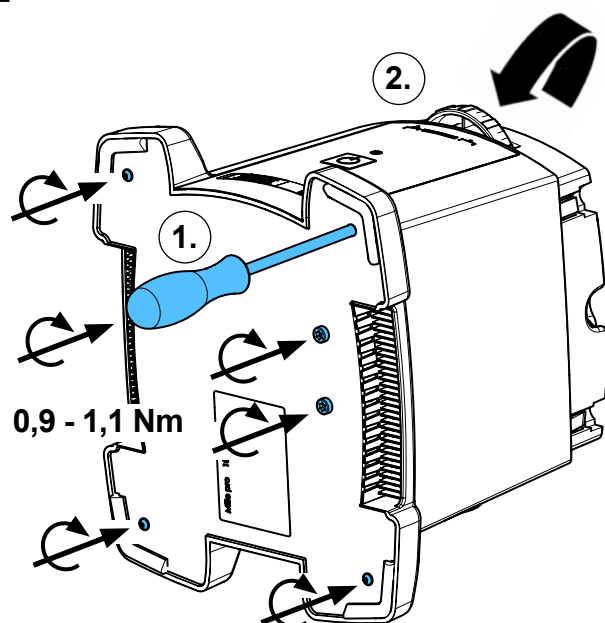


20

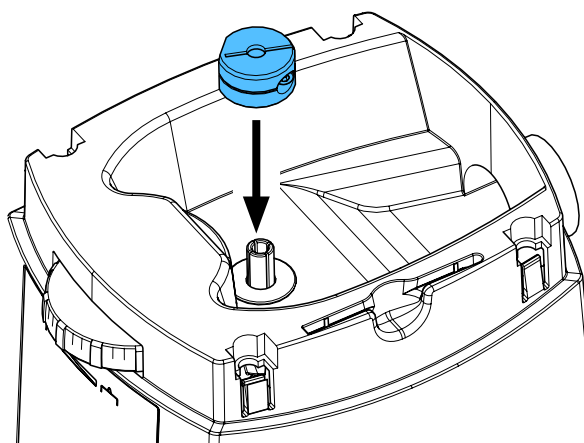


21

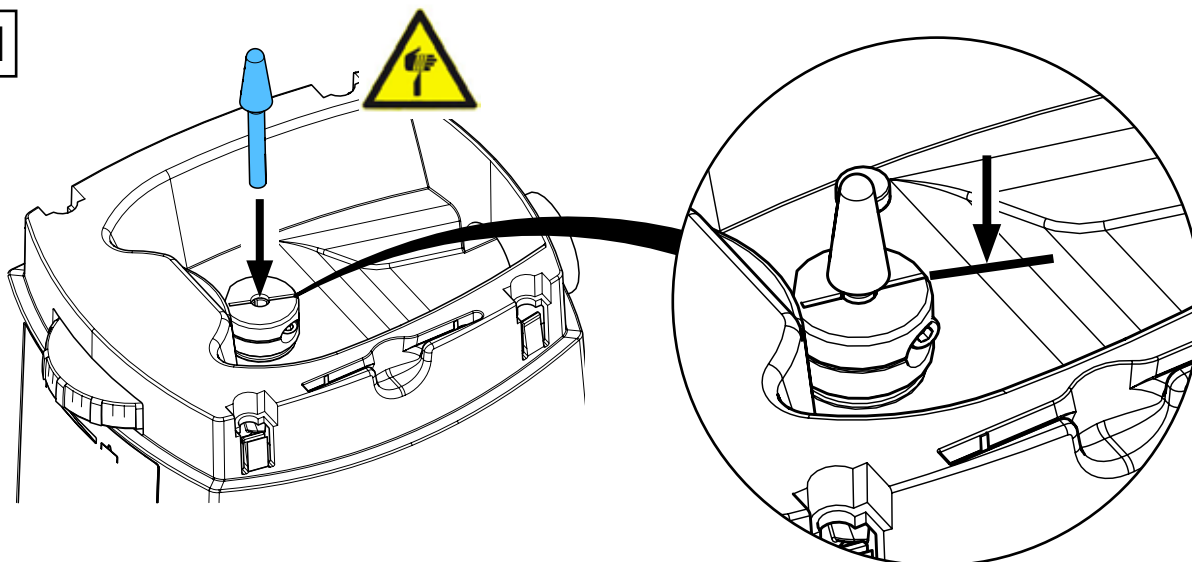


22

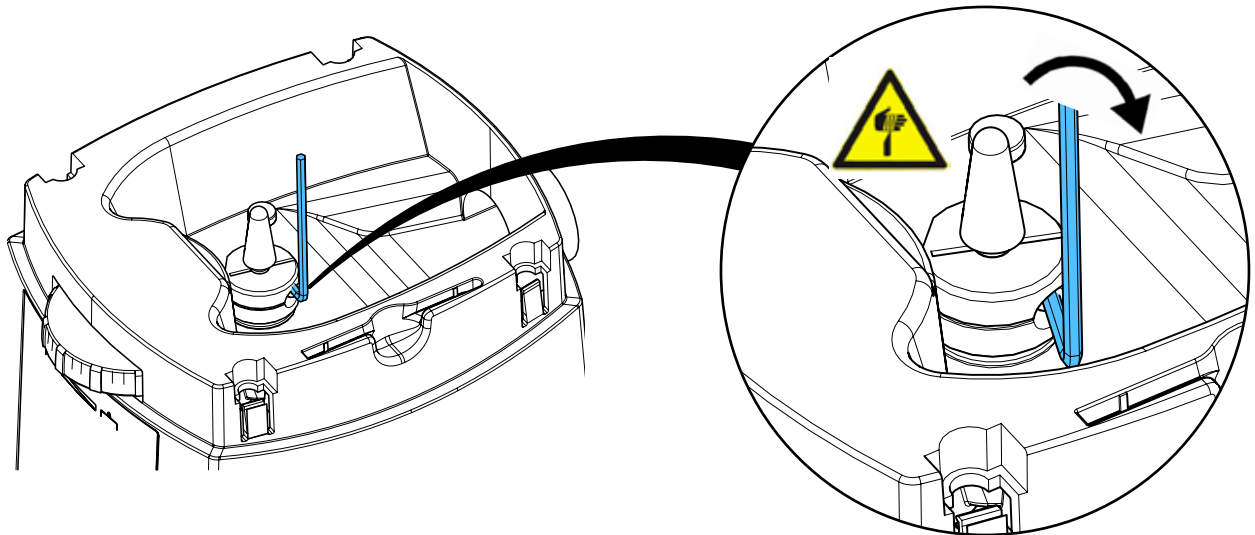

23

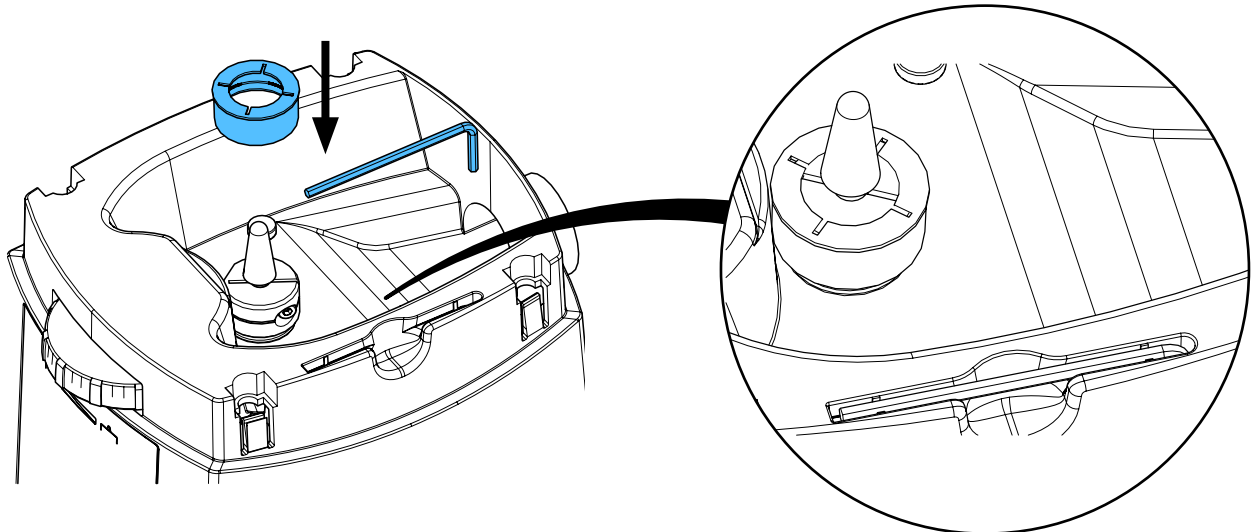


24

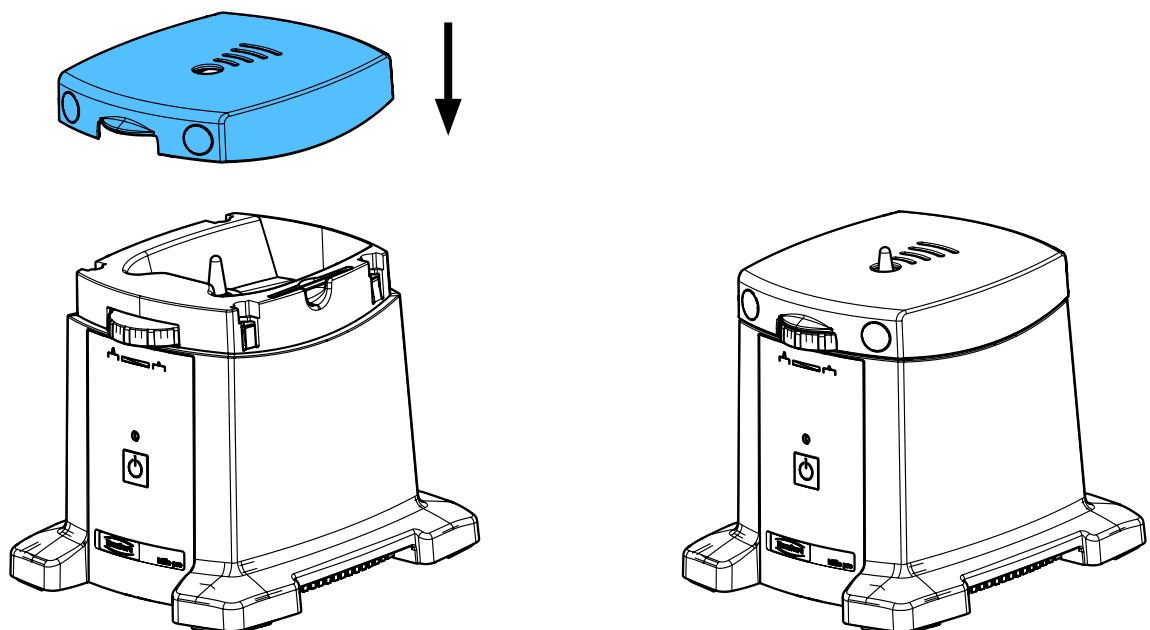


90003 3399 / 90003 3401

Motorkondensator • Engine condenser • Condensateur du moteur
Condensatore del motore • Condensador del motor • Конденсатор мотора

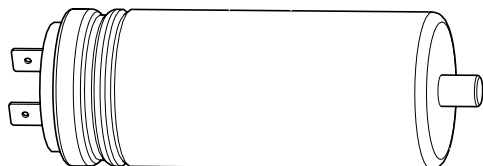


90003 3399 = 230 V

90003 3401 = 120 V



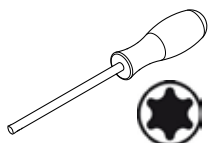


T10
T25

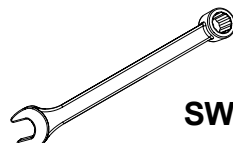


SW 13


1 Millo 1 - 18 / Millo pro 19 - 38

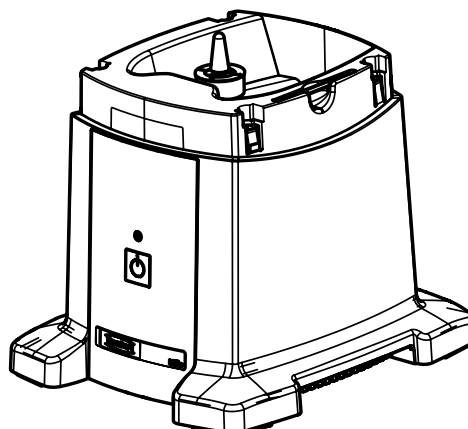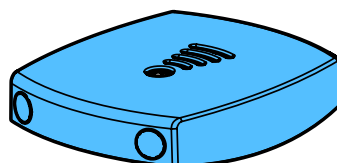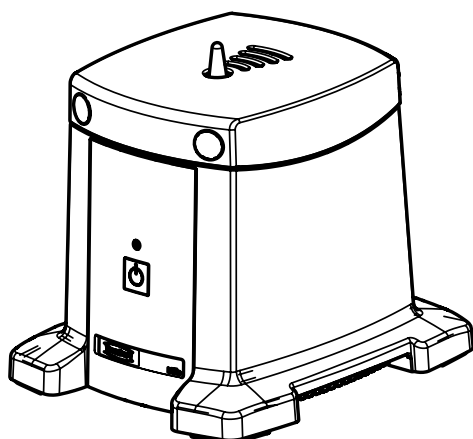


2

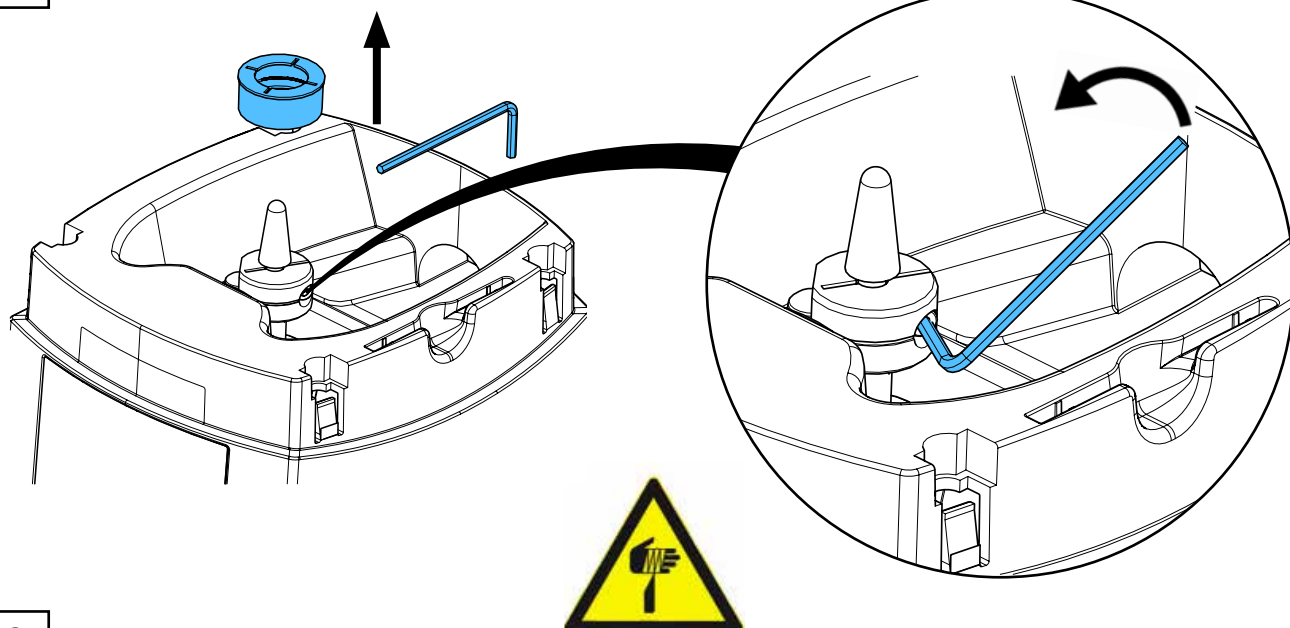


3

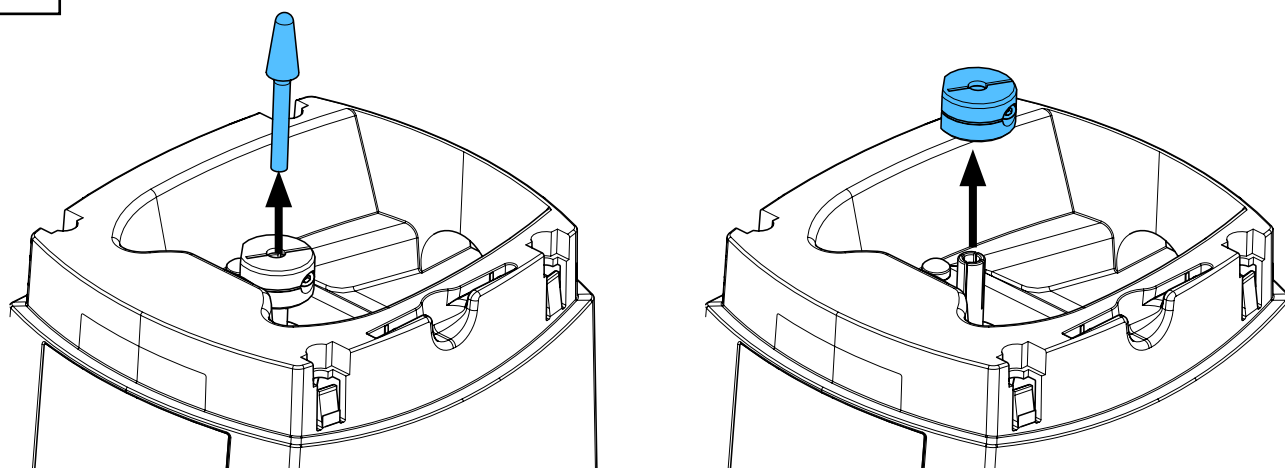


4

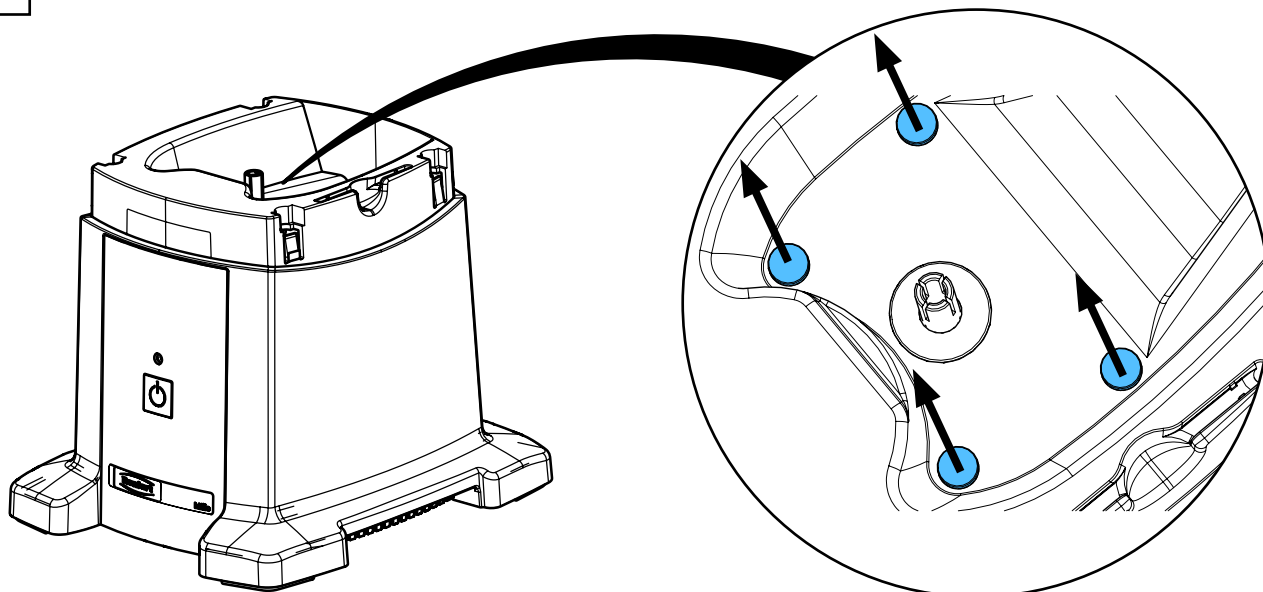


5

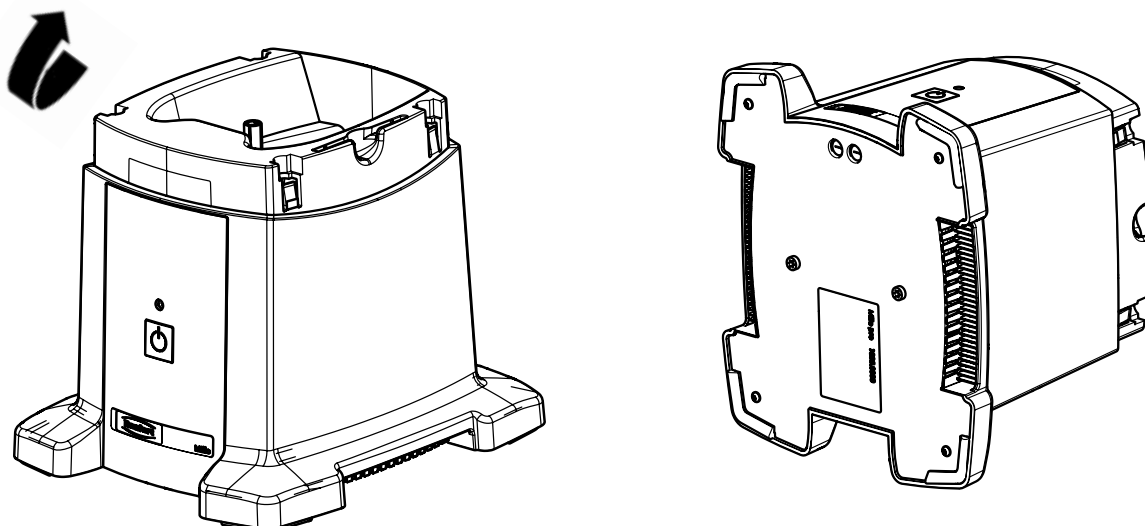


6

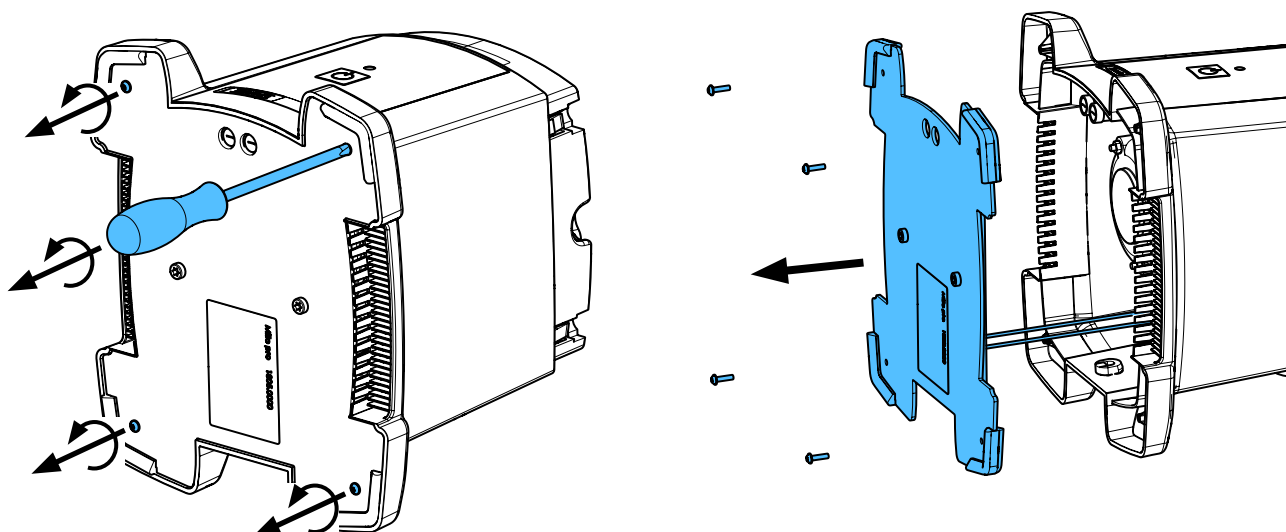


7

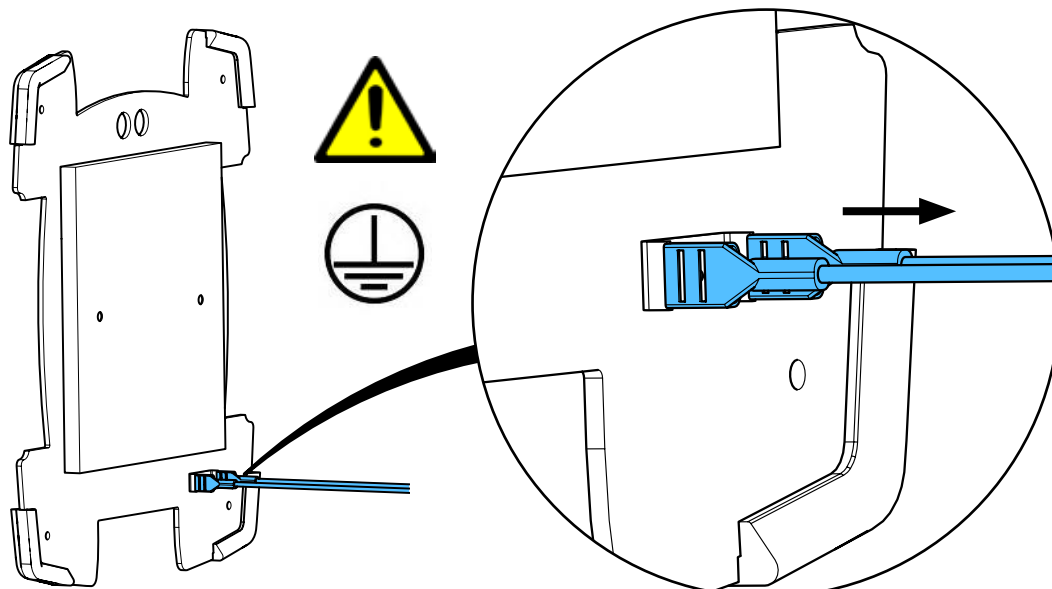



8

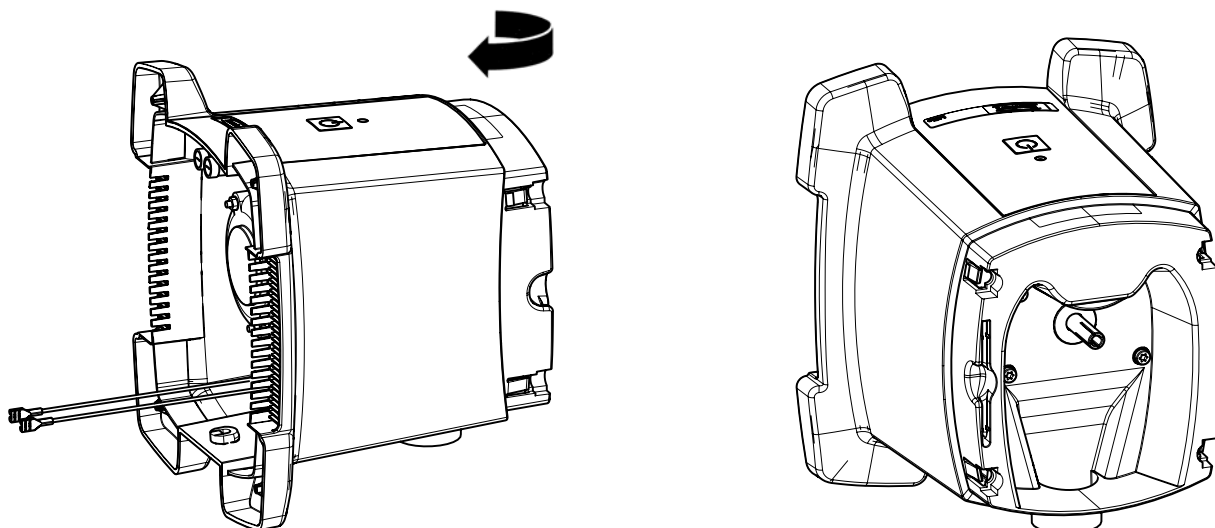


9

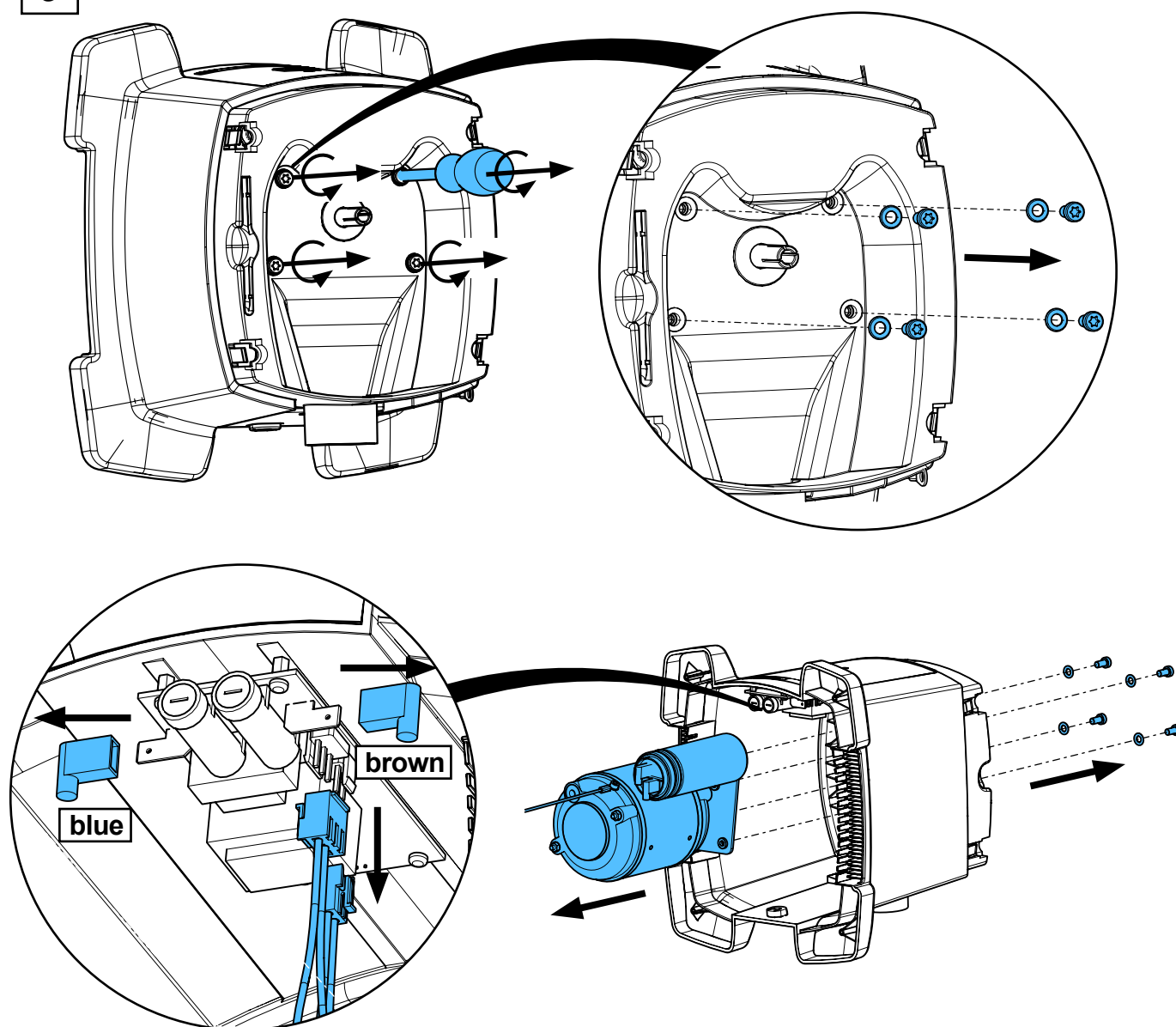


10

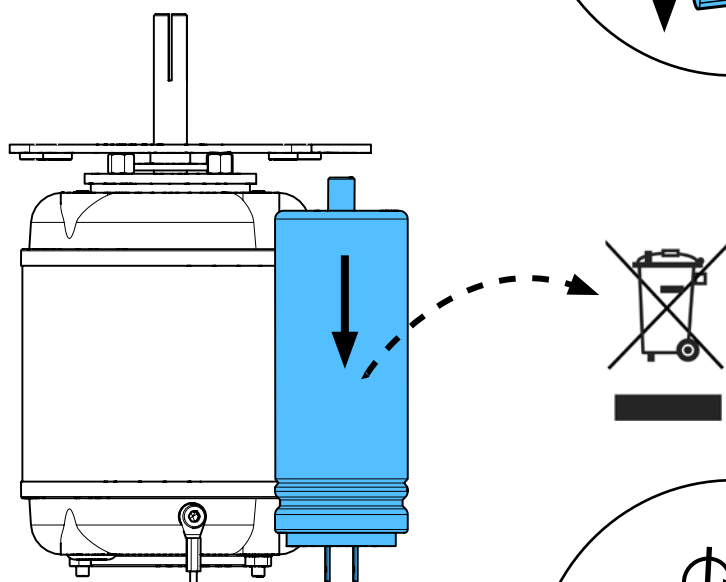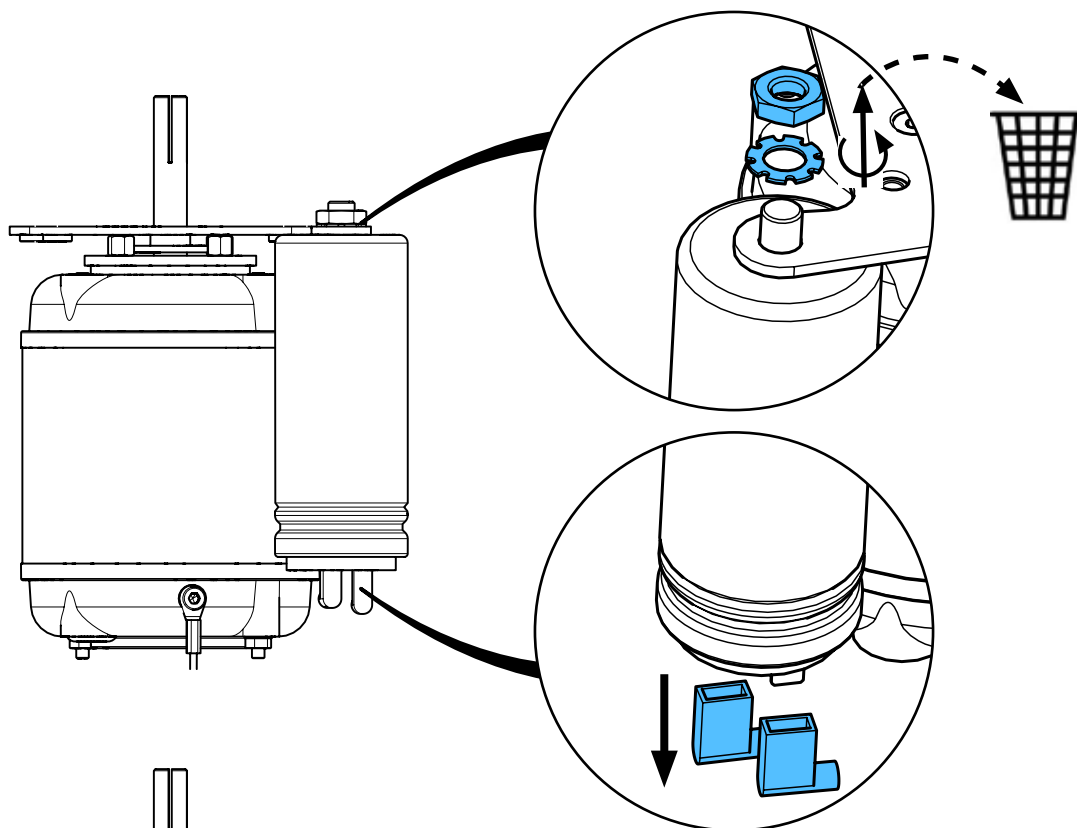


11

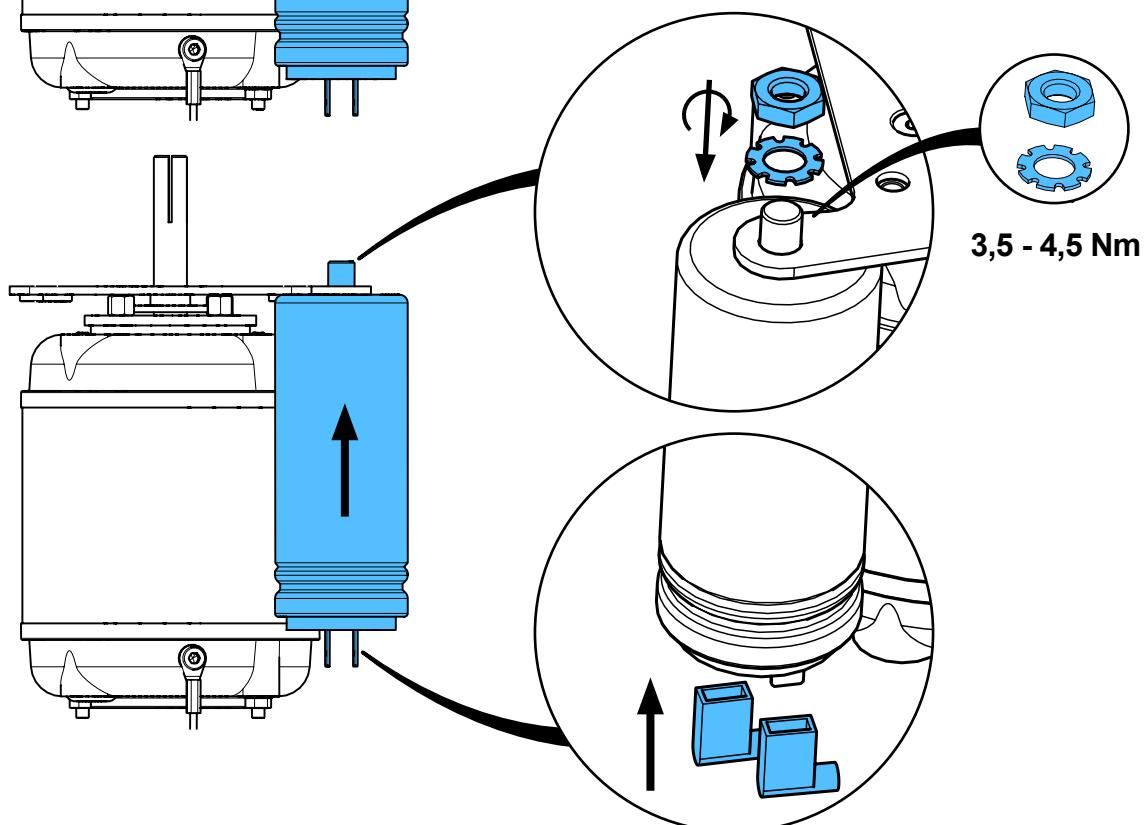


12

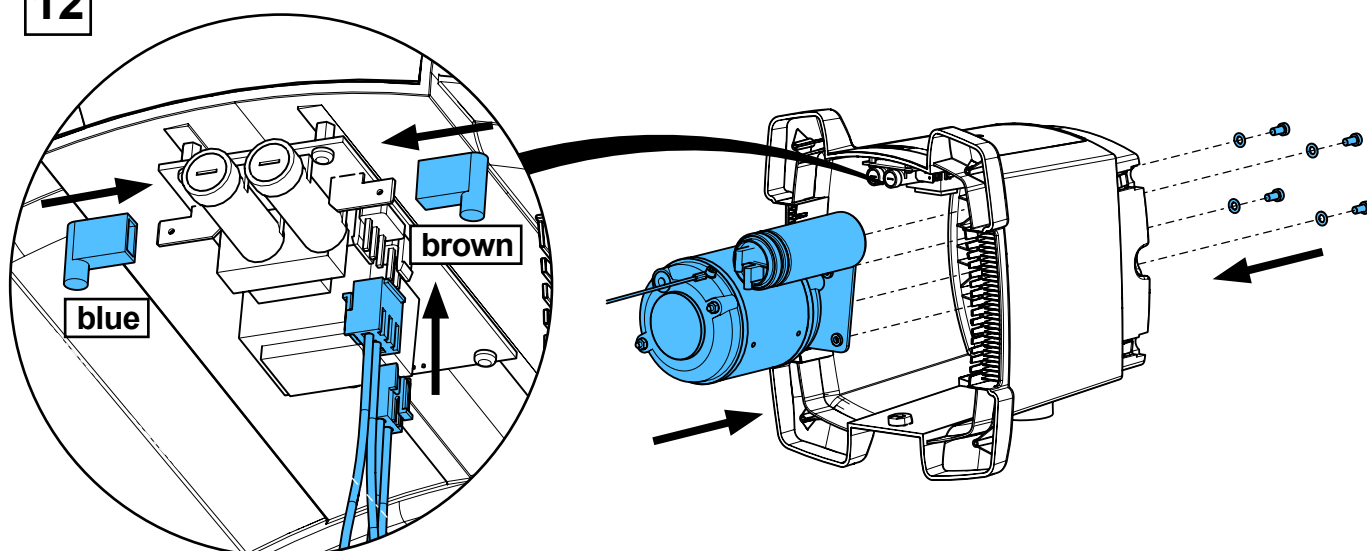



13

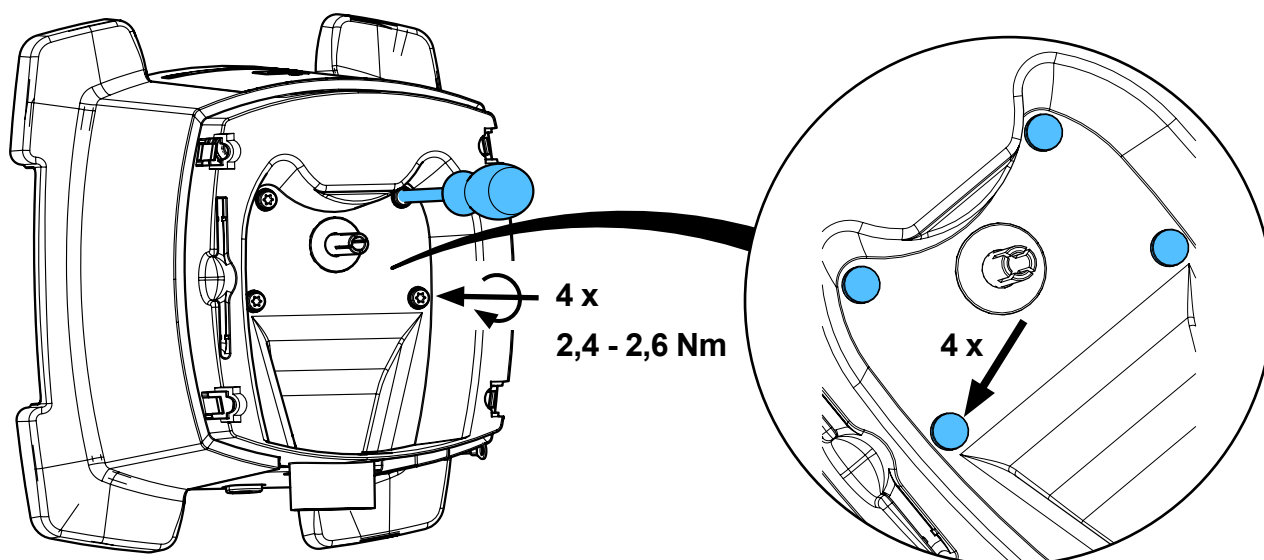



14

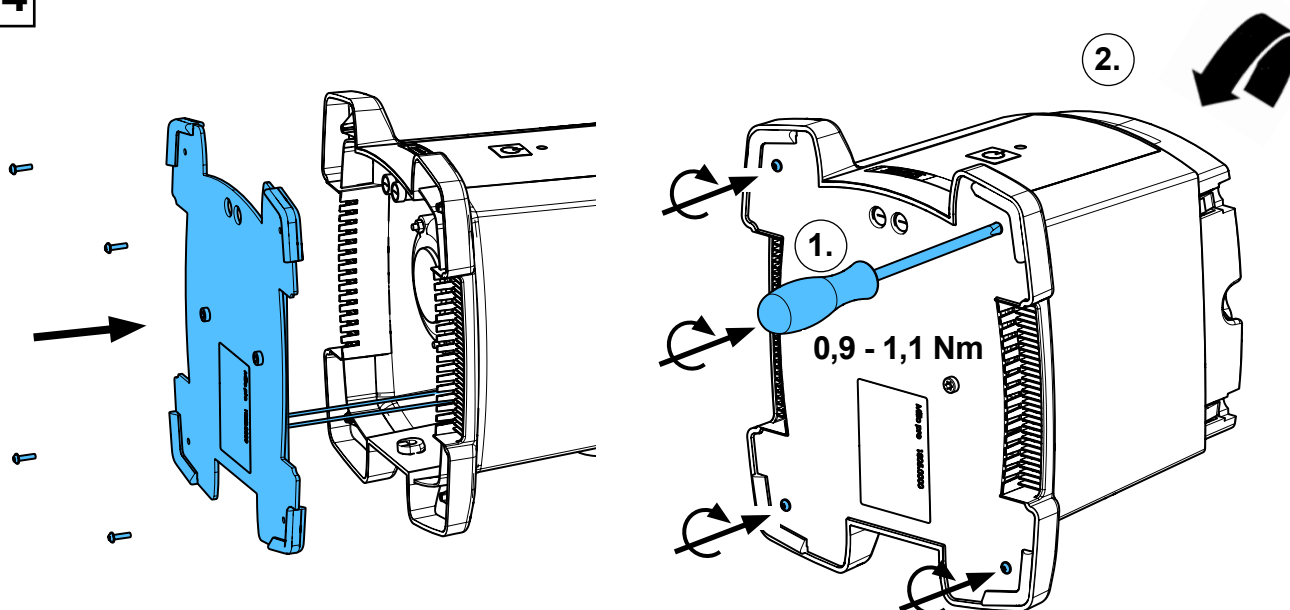



15

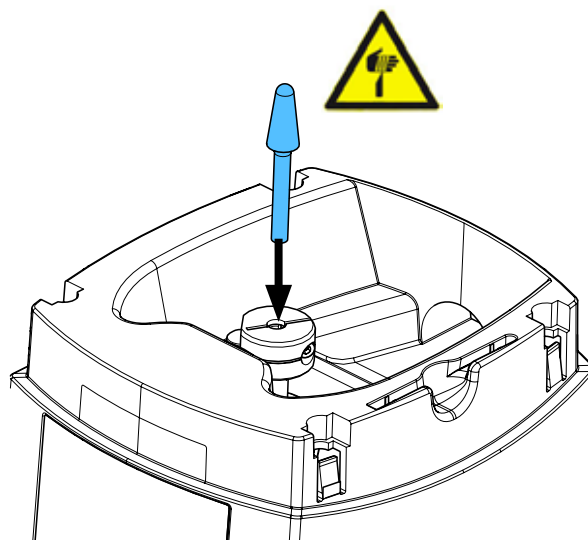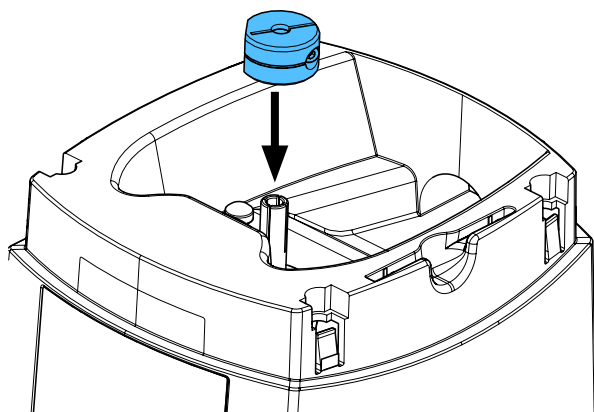


16

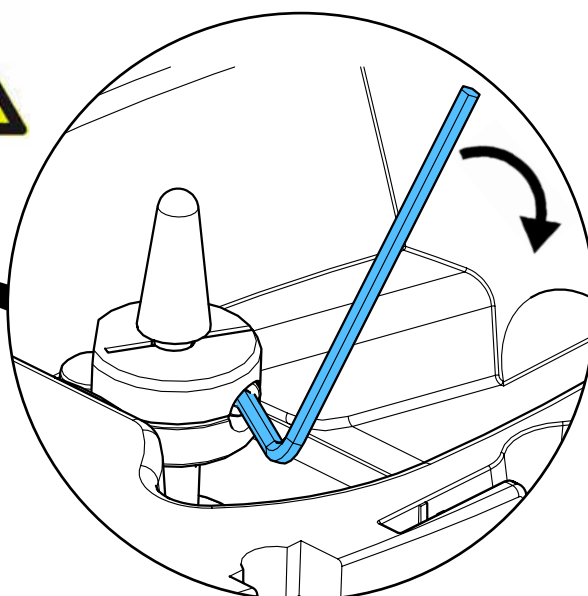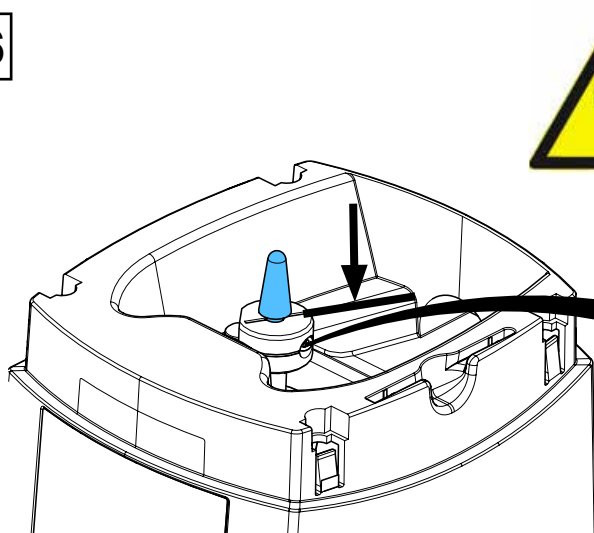



17

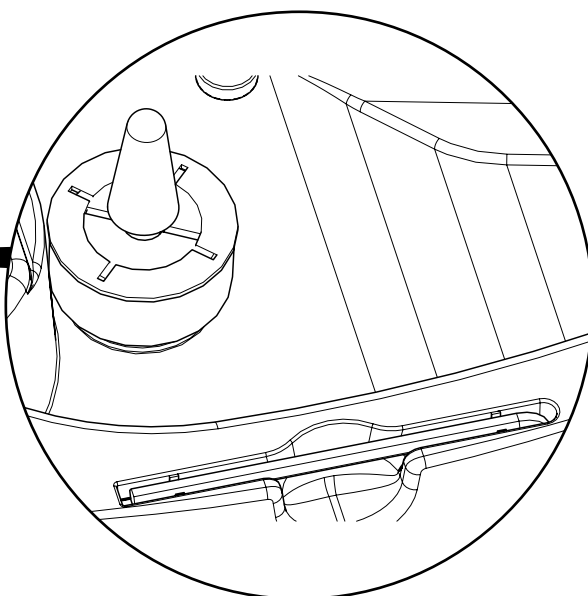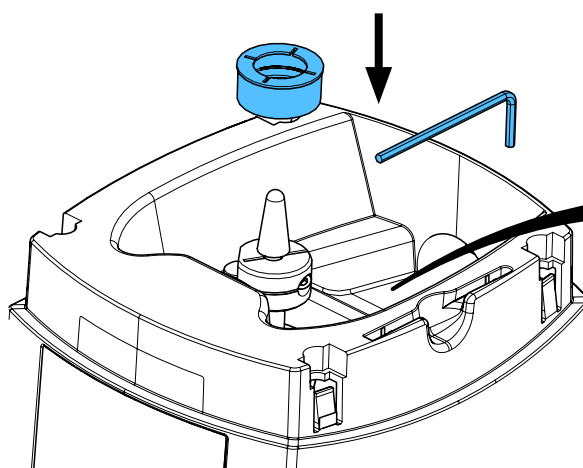



18

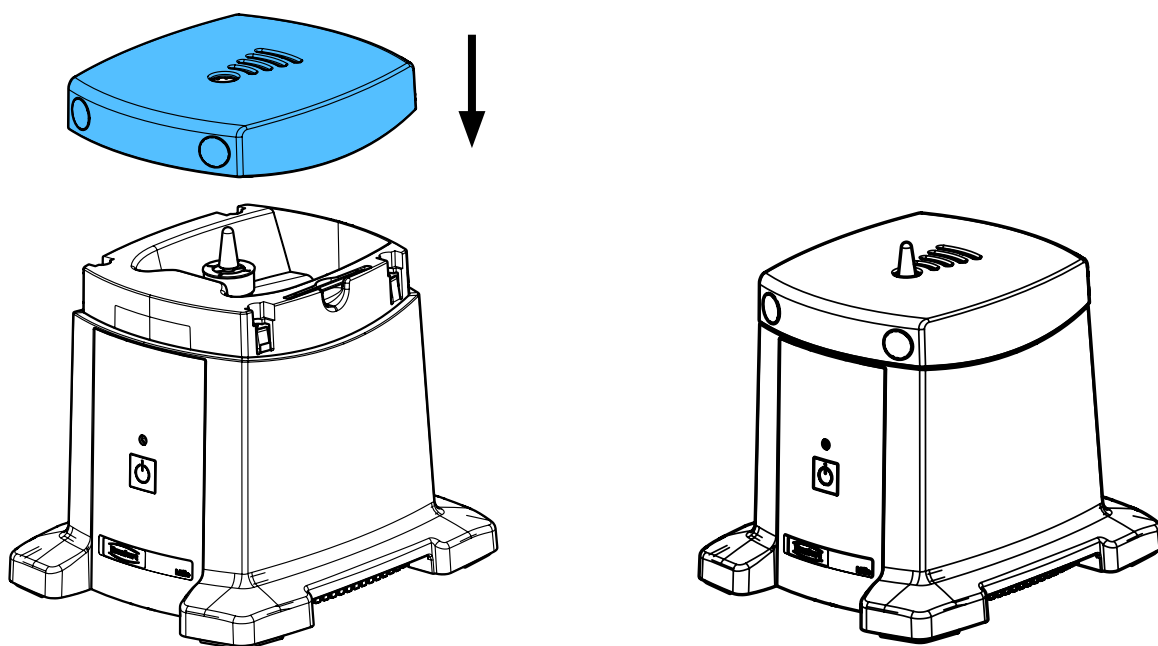



19 Millo pro 19 - 38

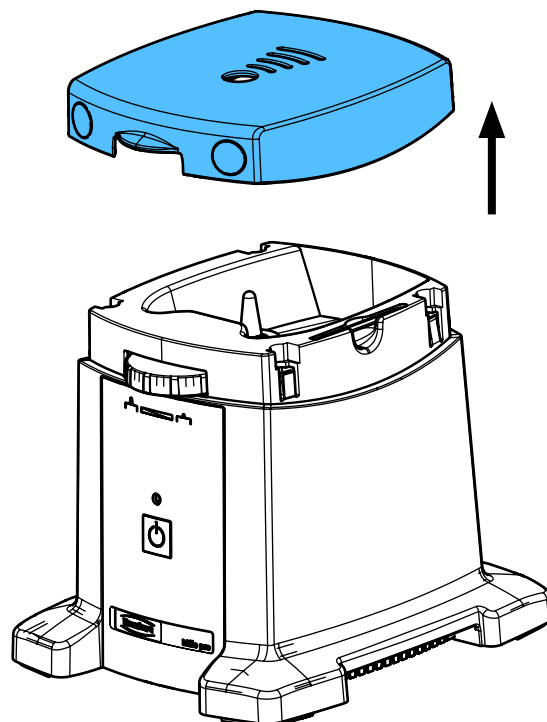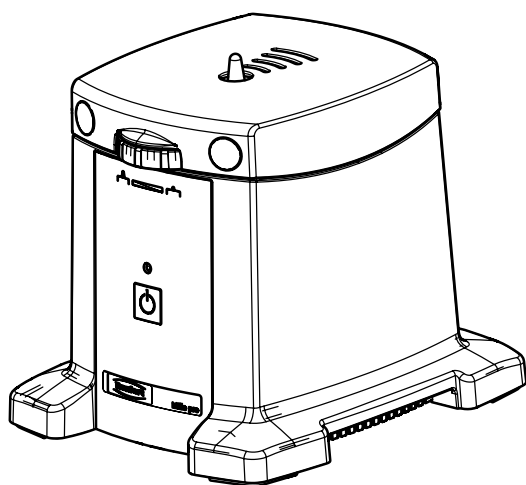


20

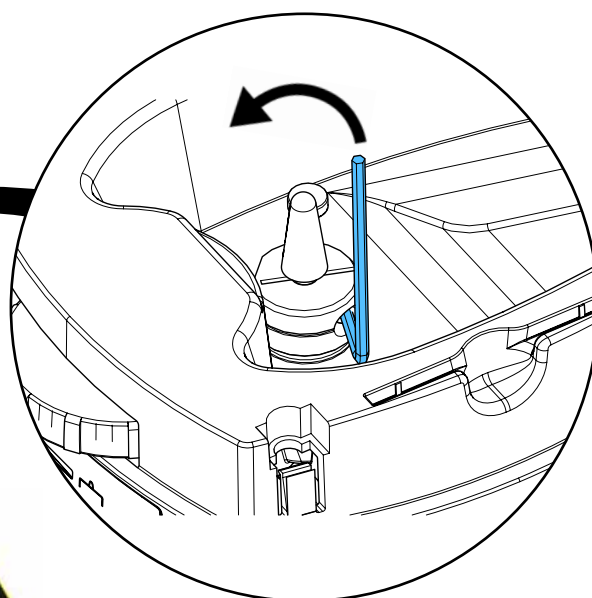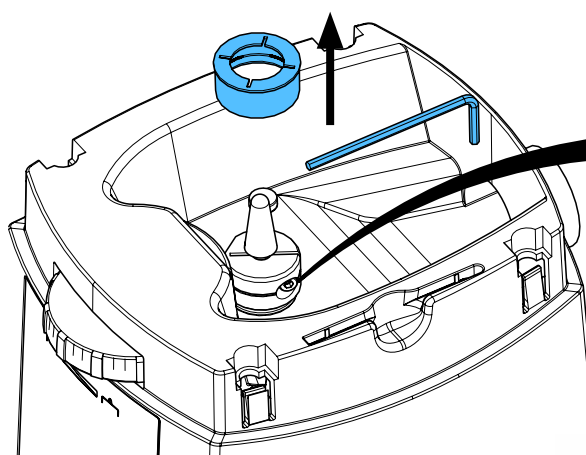


21

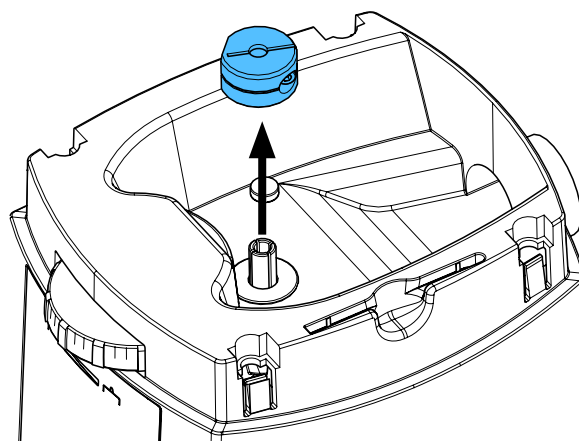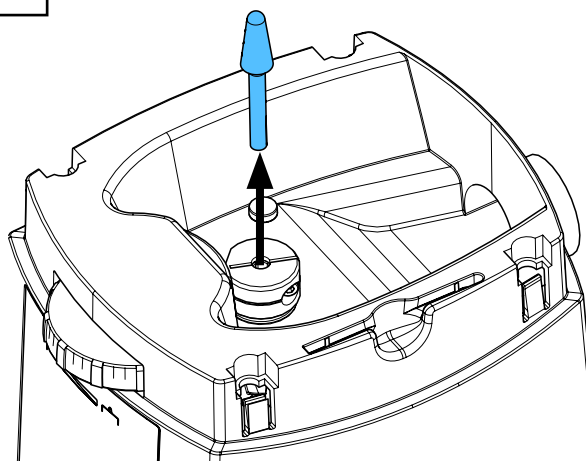


22

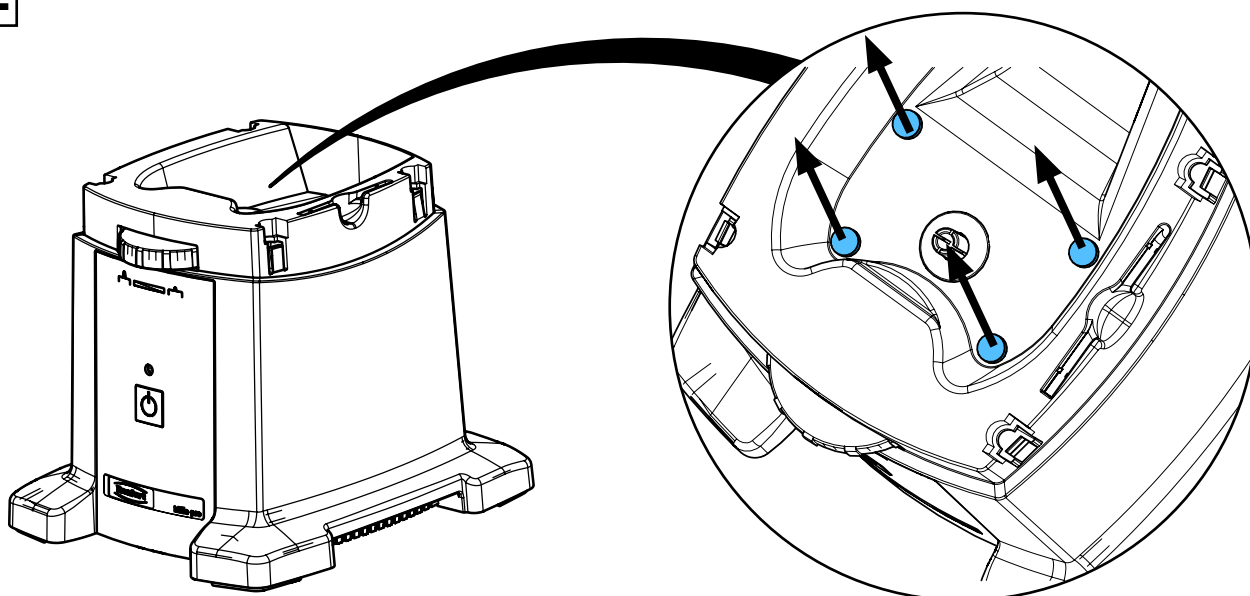


23

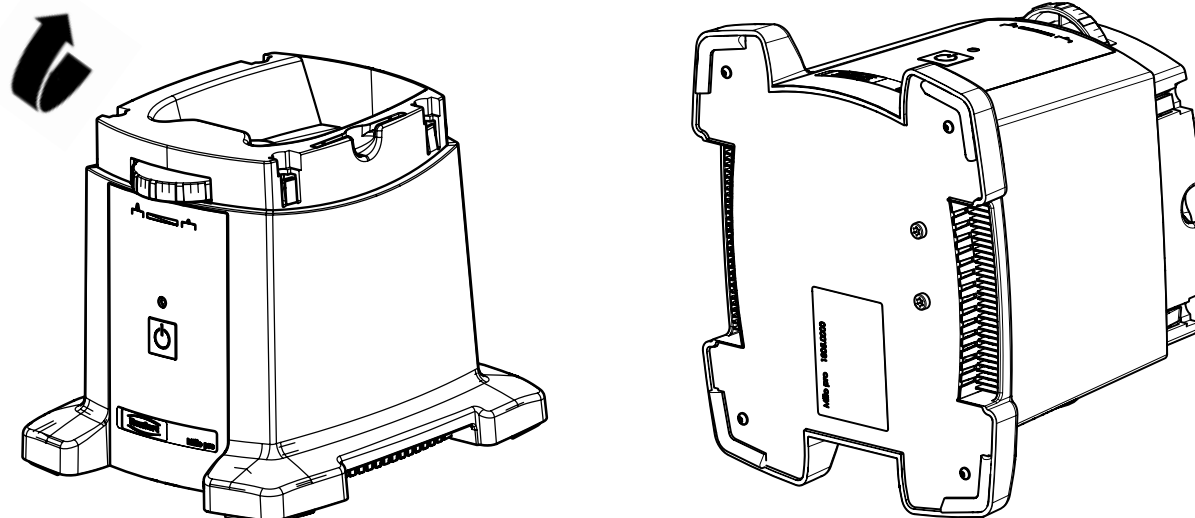


24

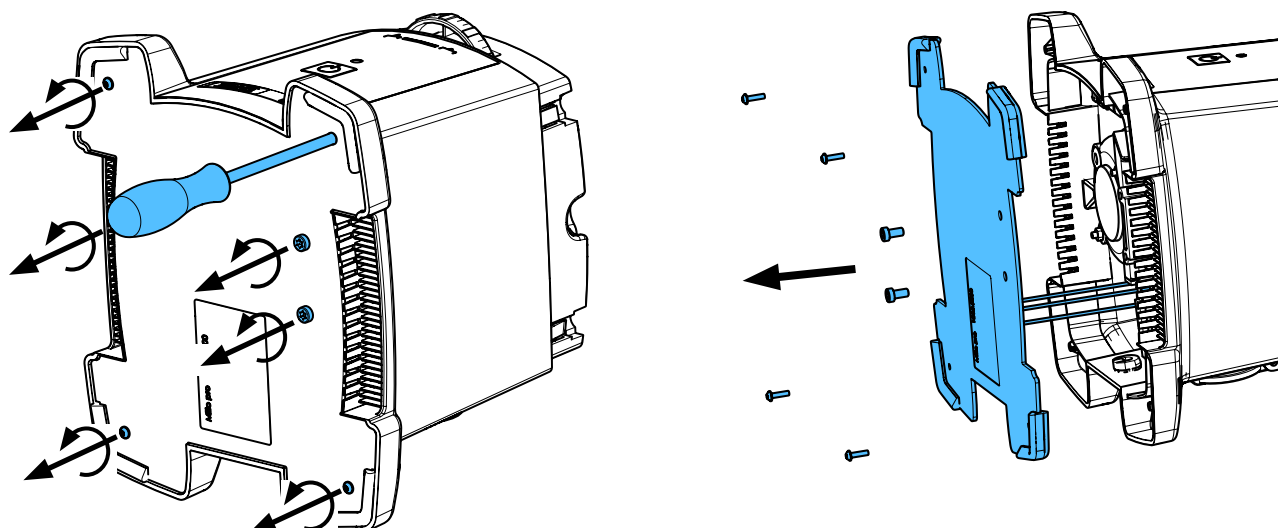


25

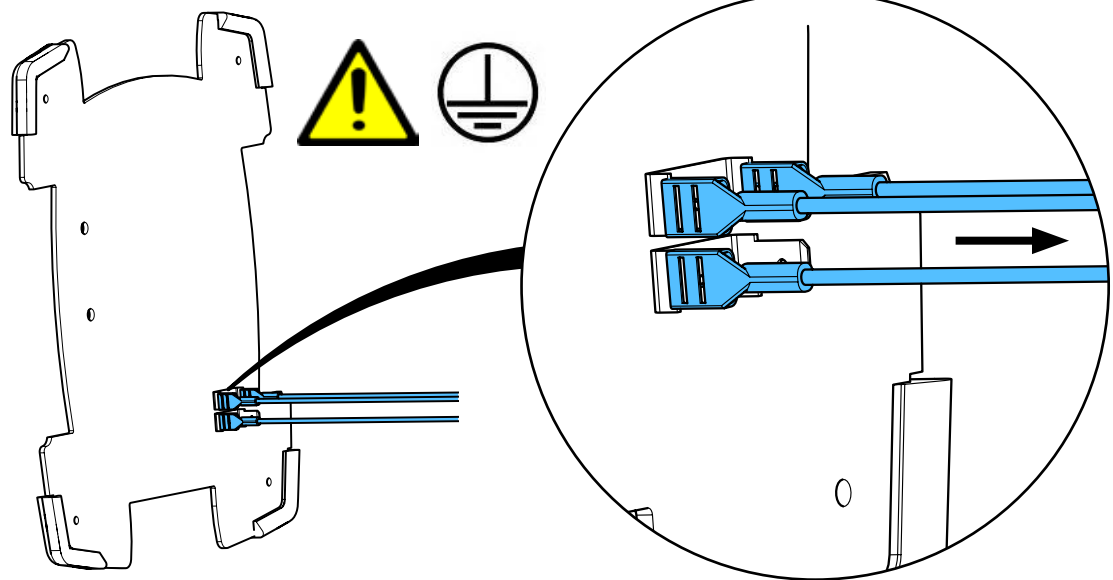


26

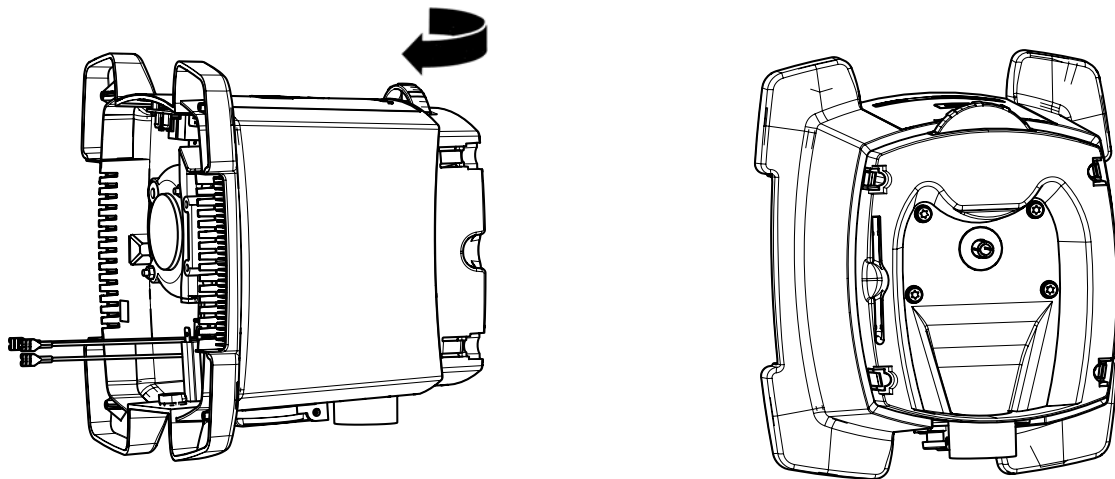


27

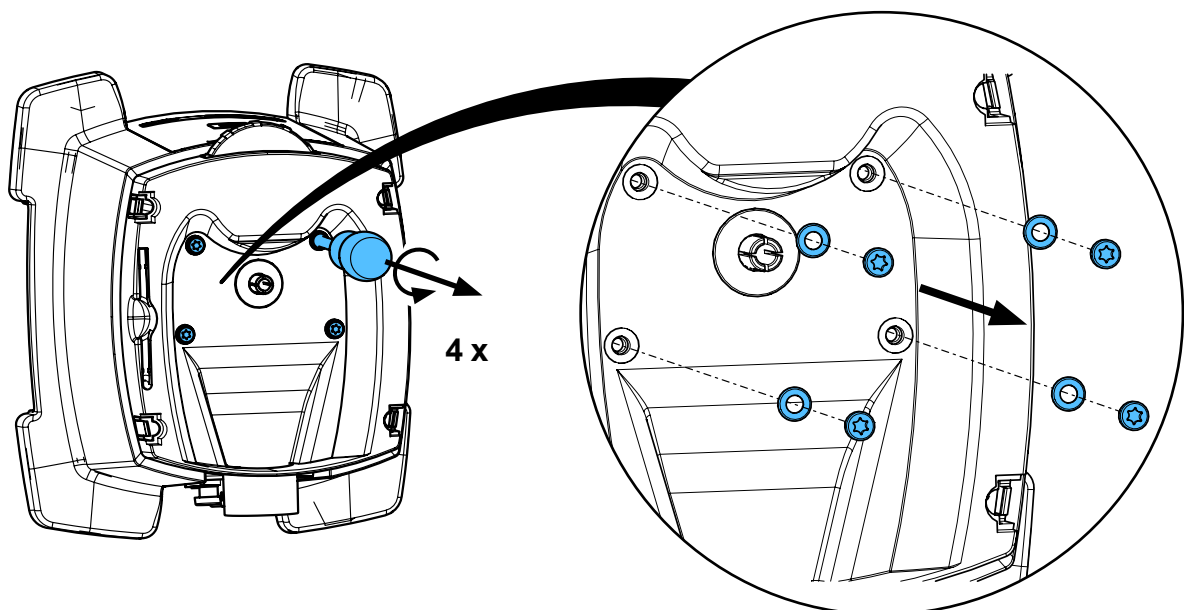


28

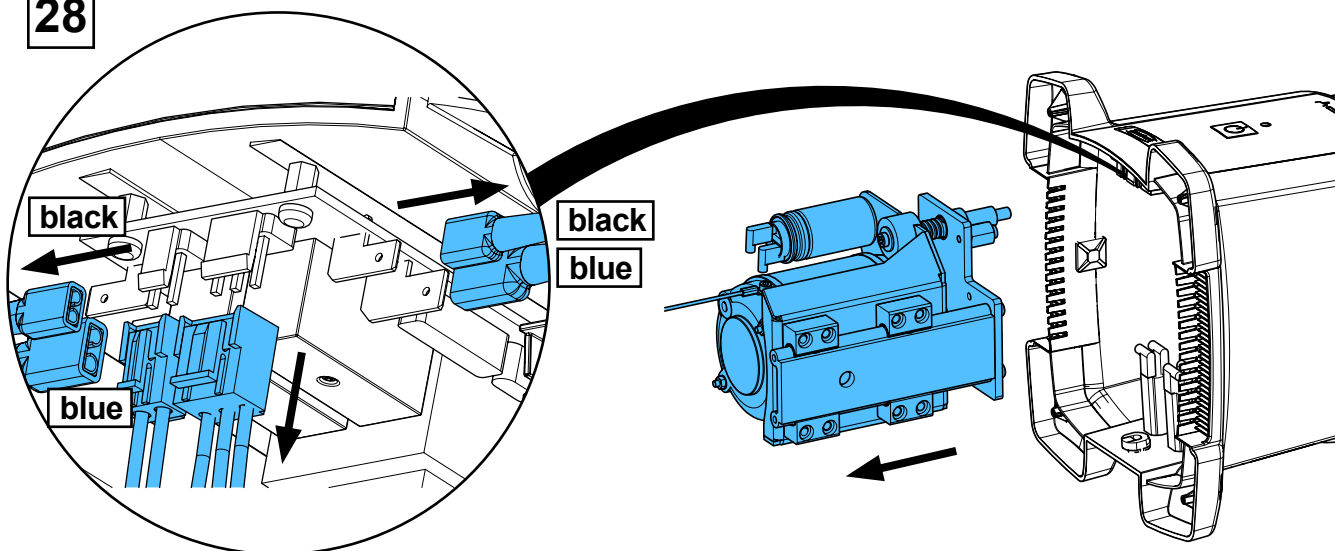


29

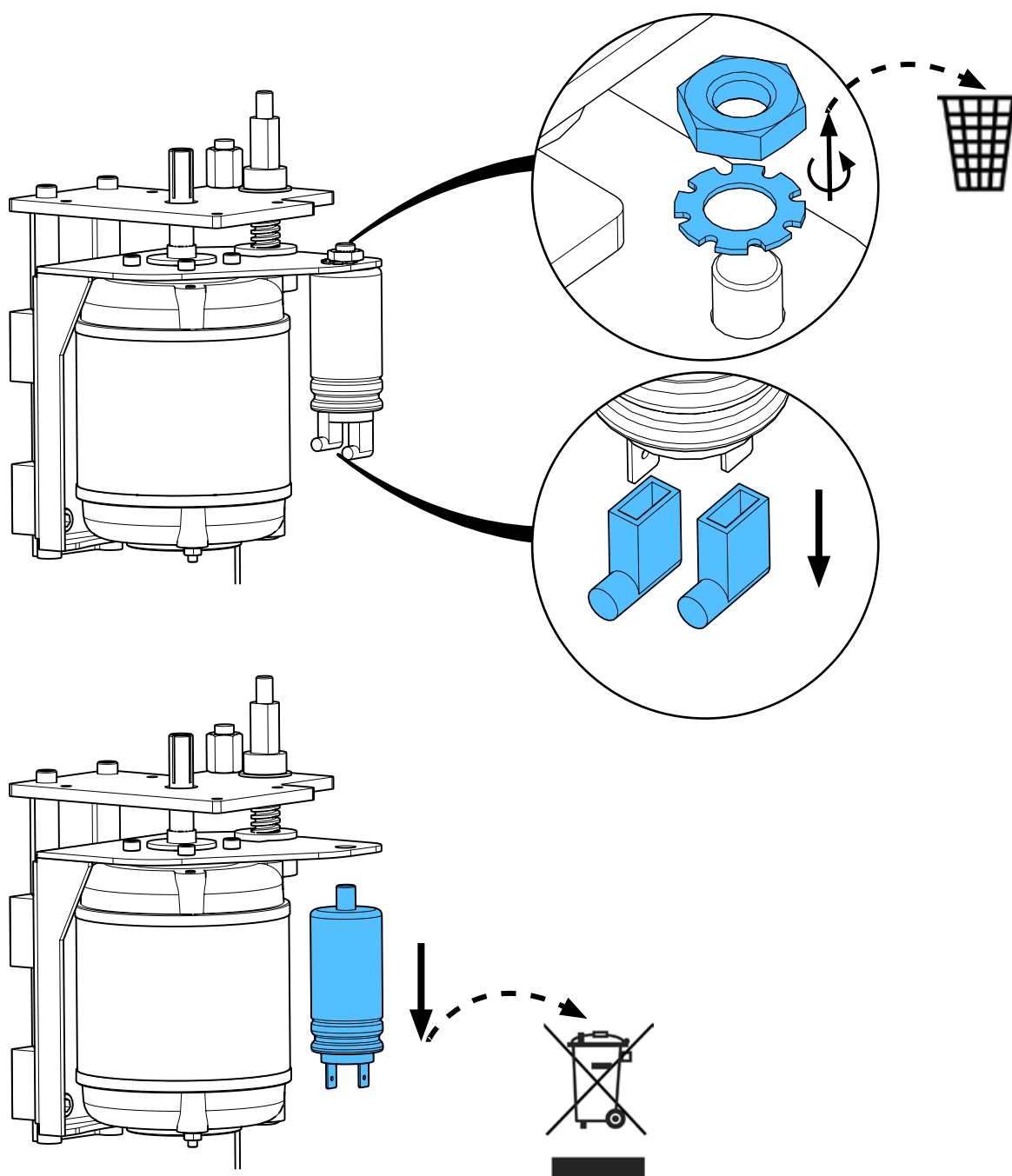


30

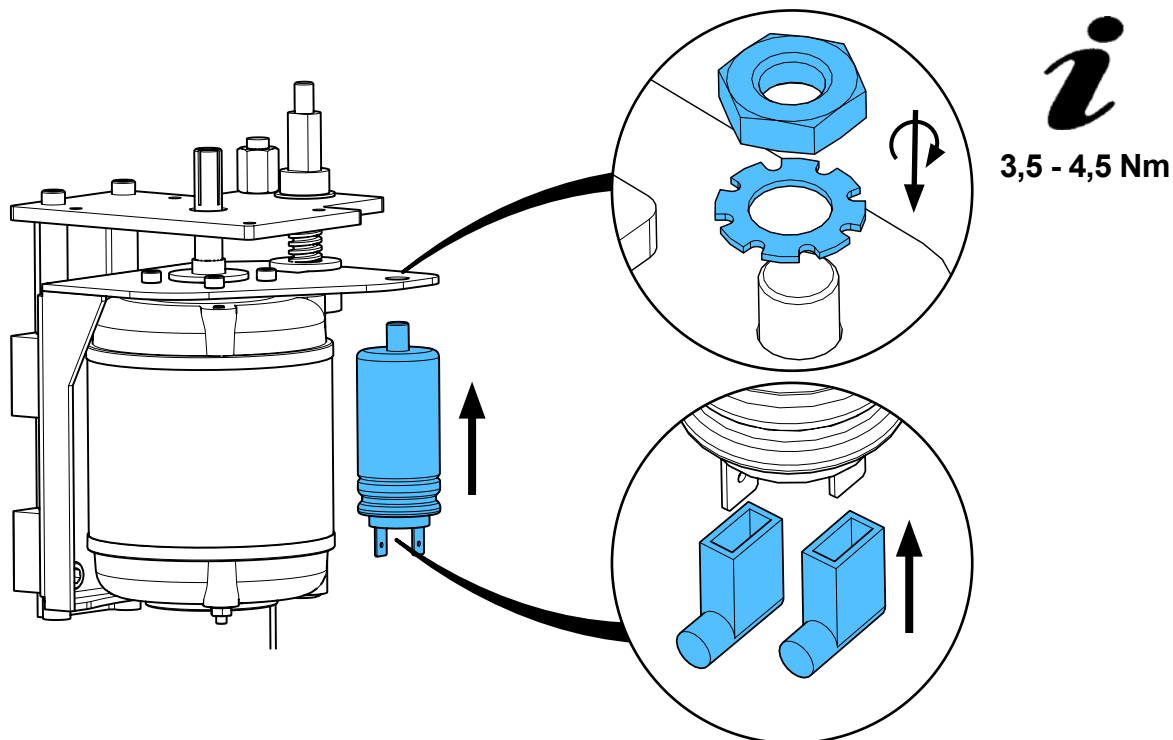


31

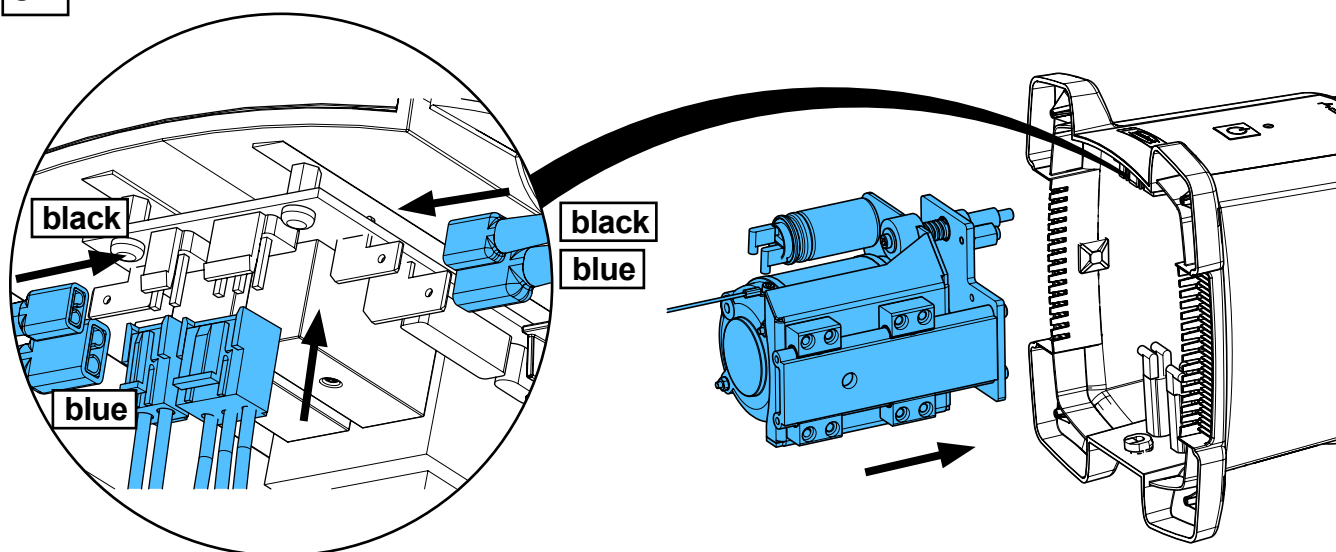


32

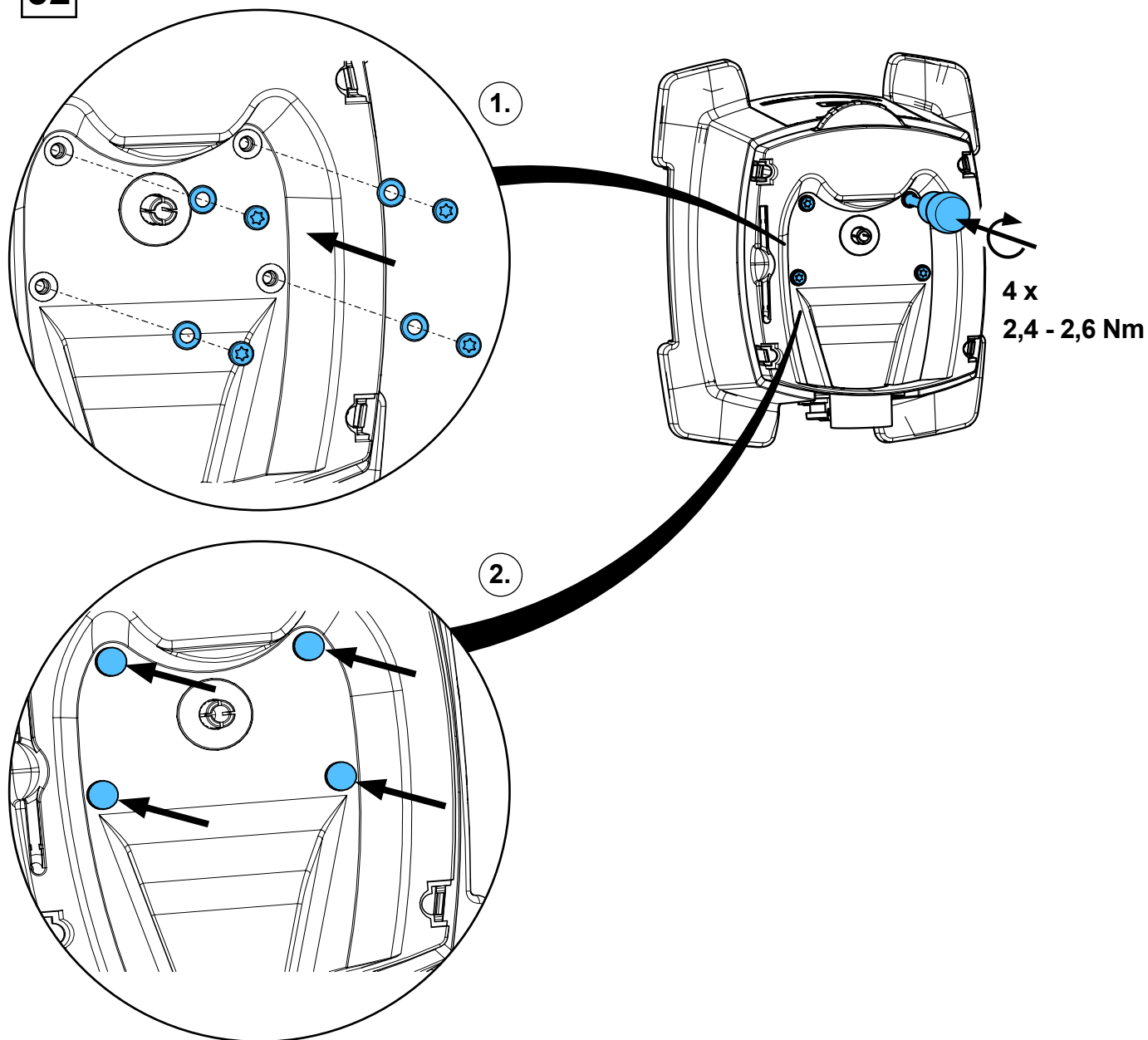



33

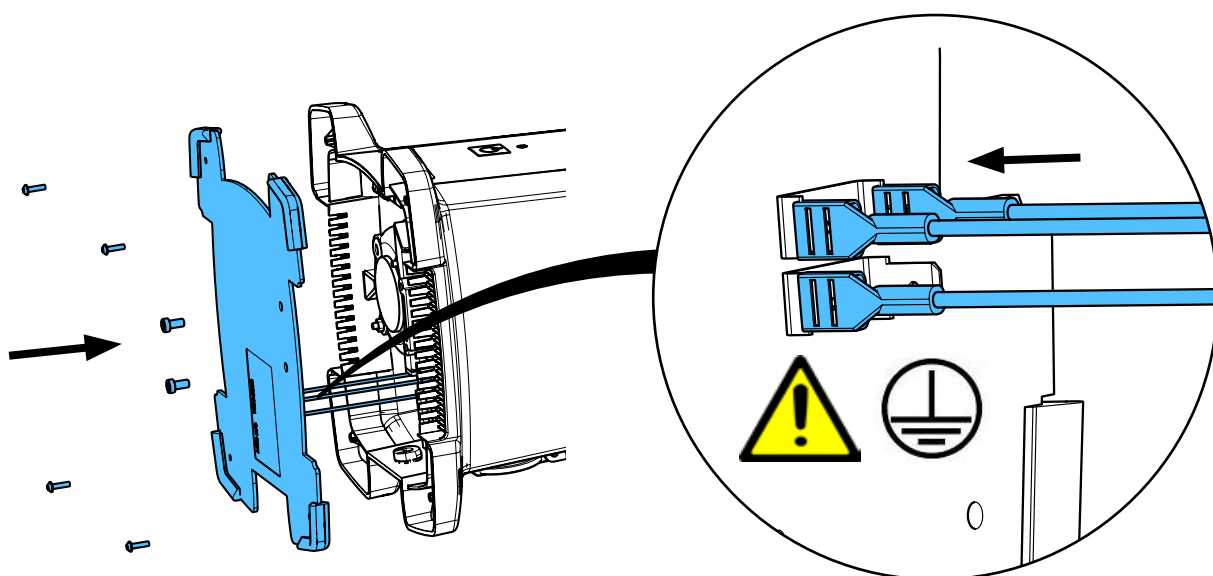


34

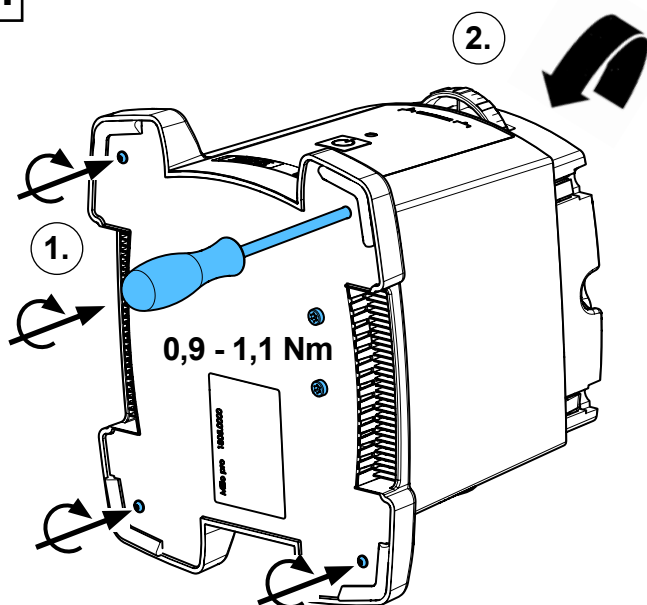


35

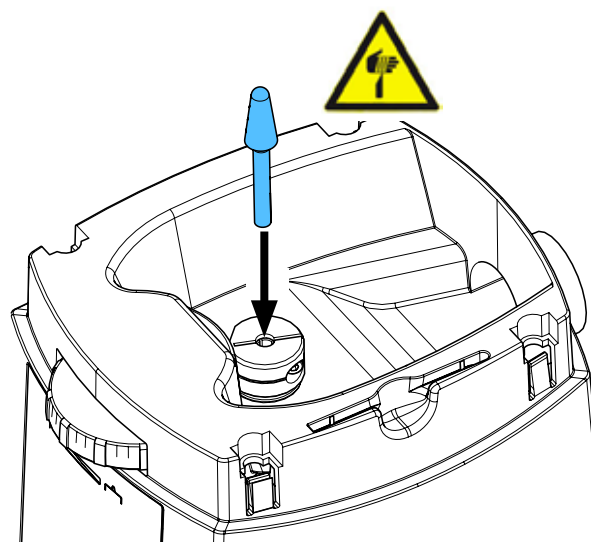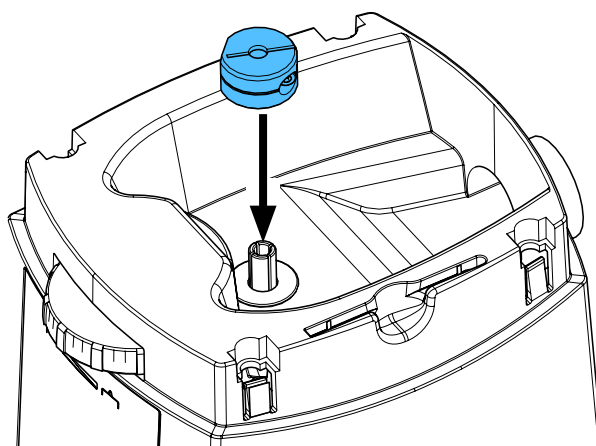


36

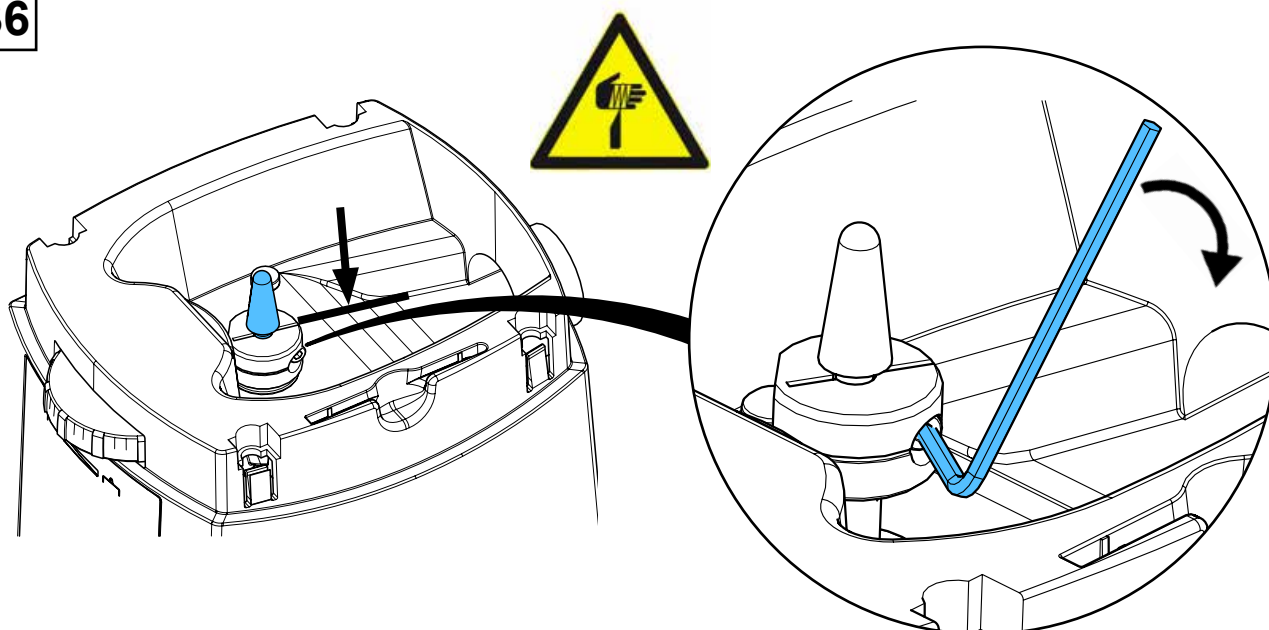

37

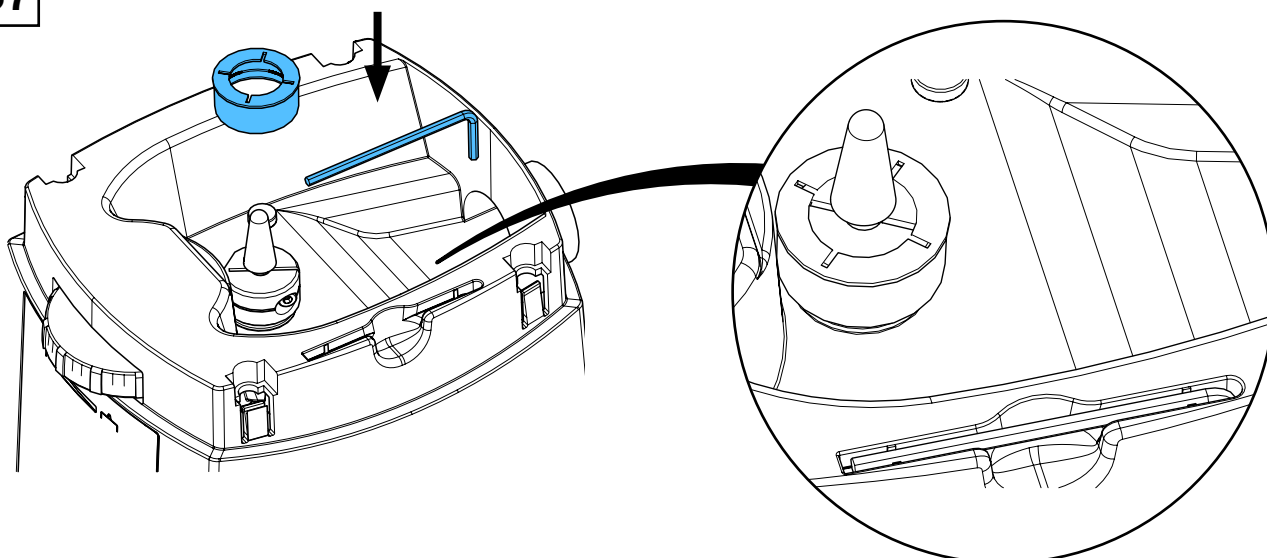

38


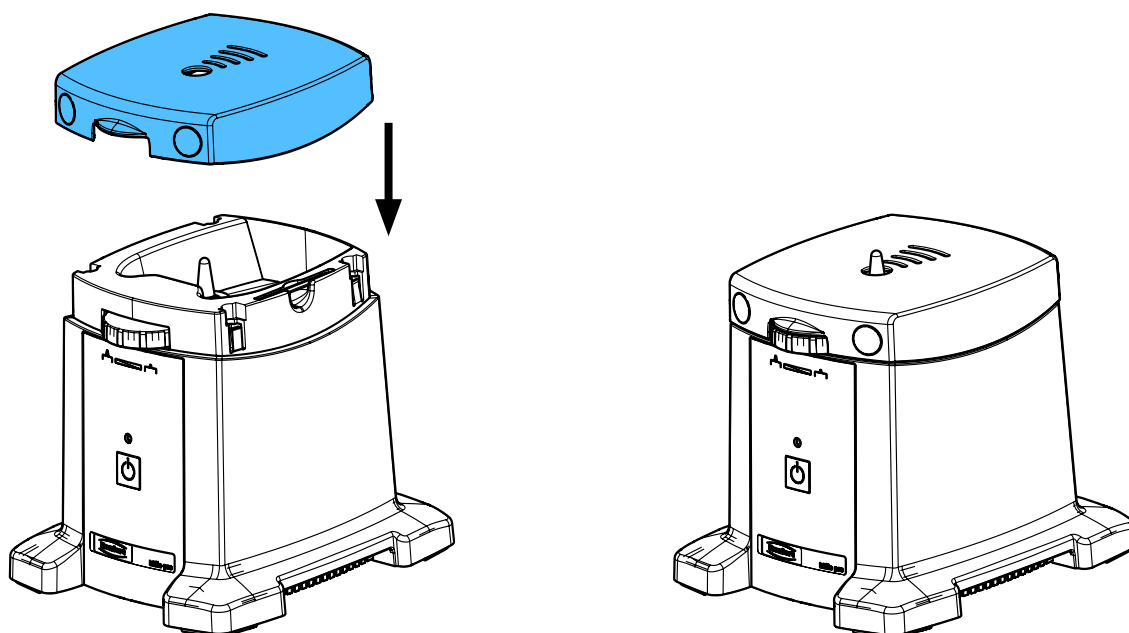


90003 3384

Gerätefüße • Rubber bases • Pieds de l'appareil • Piedini dell'apparecchio
Pies de l'aparato • Ножки прибора





4 x

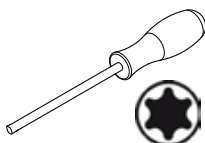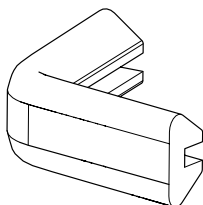



T10

1


Millo / Millo pro

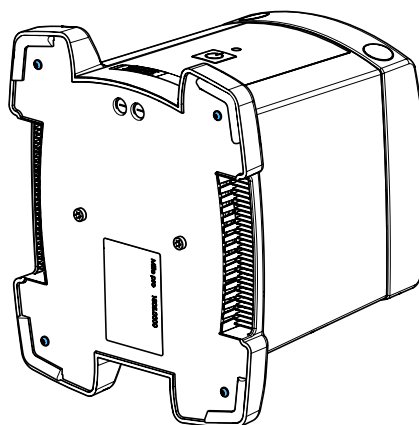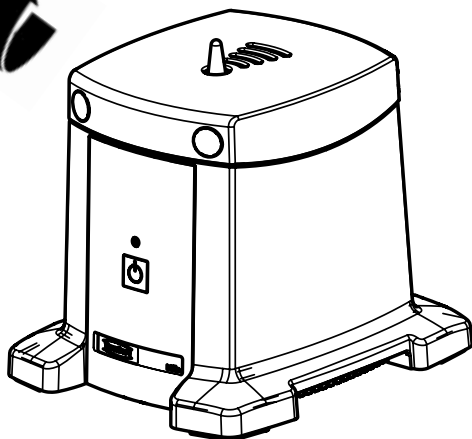


2

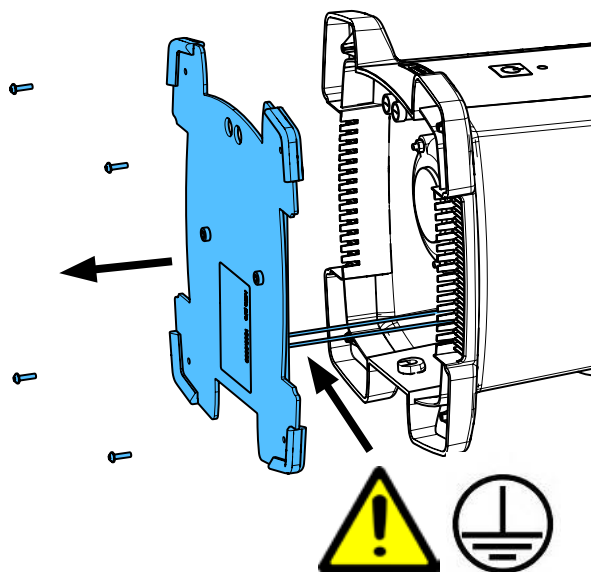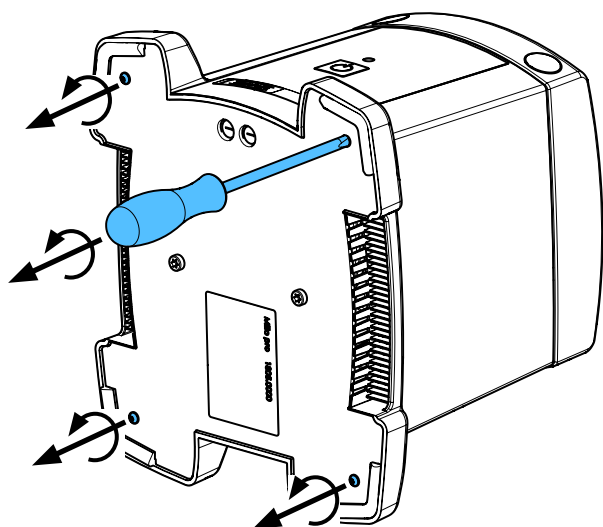


3

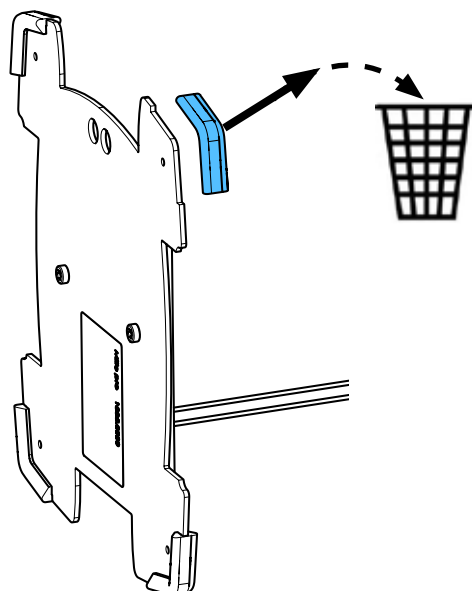



4

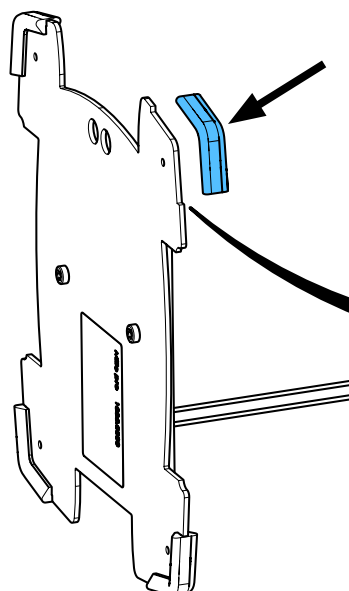


i

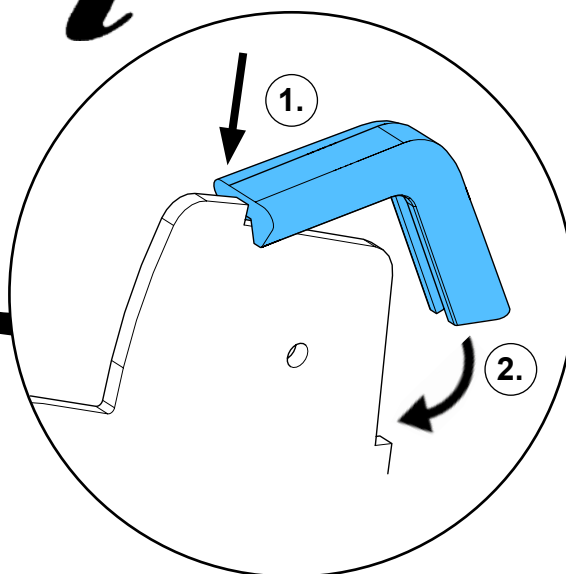


5

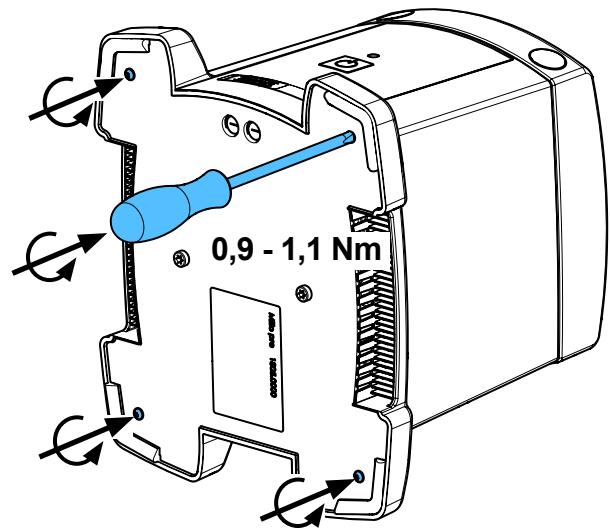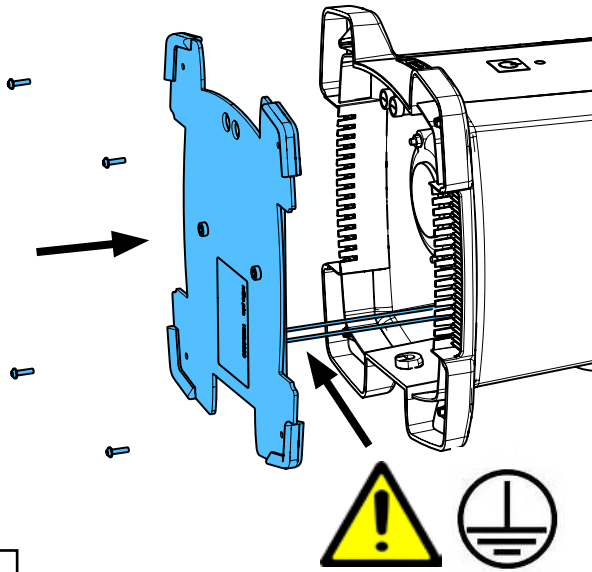

6


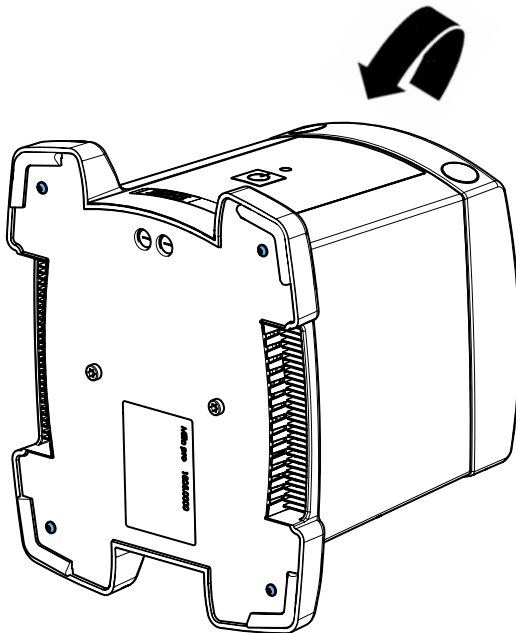


7

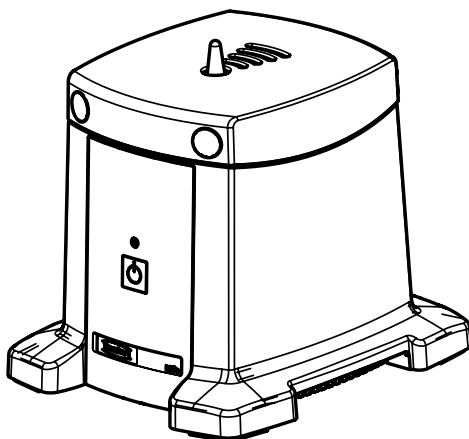


90003 3386


Höhenverstellung kpl • Height adjustment complete • Régleur d'hauteur complet
Regolatore d'altezza completo • Ajusta de altura completo • Установка по высоте

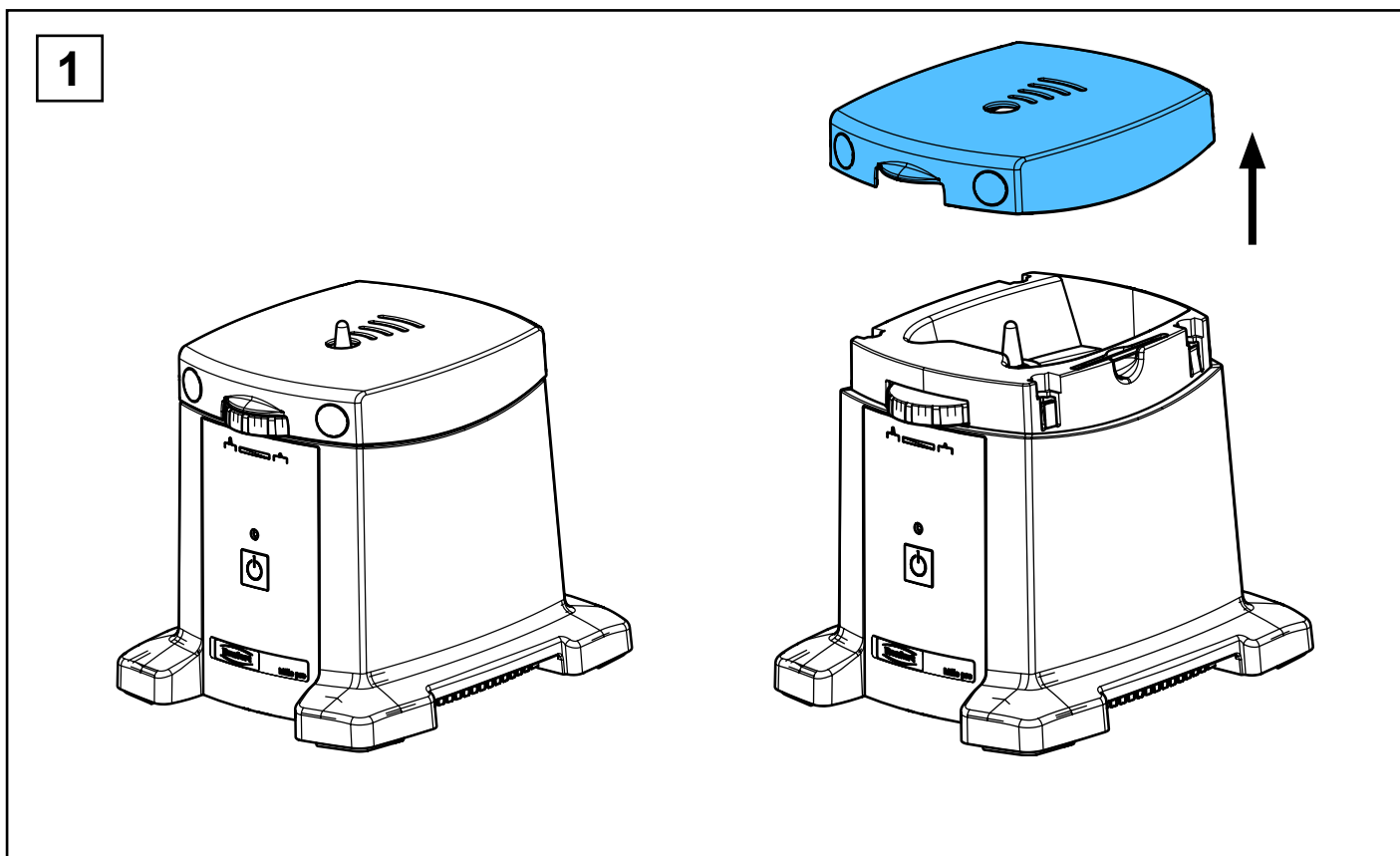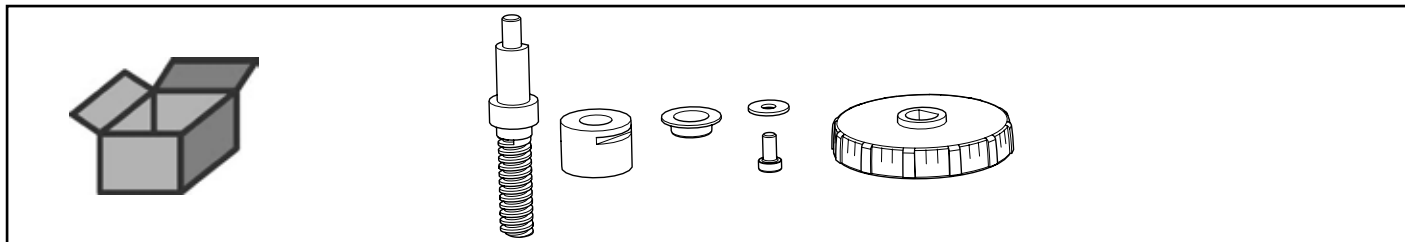


2

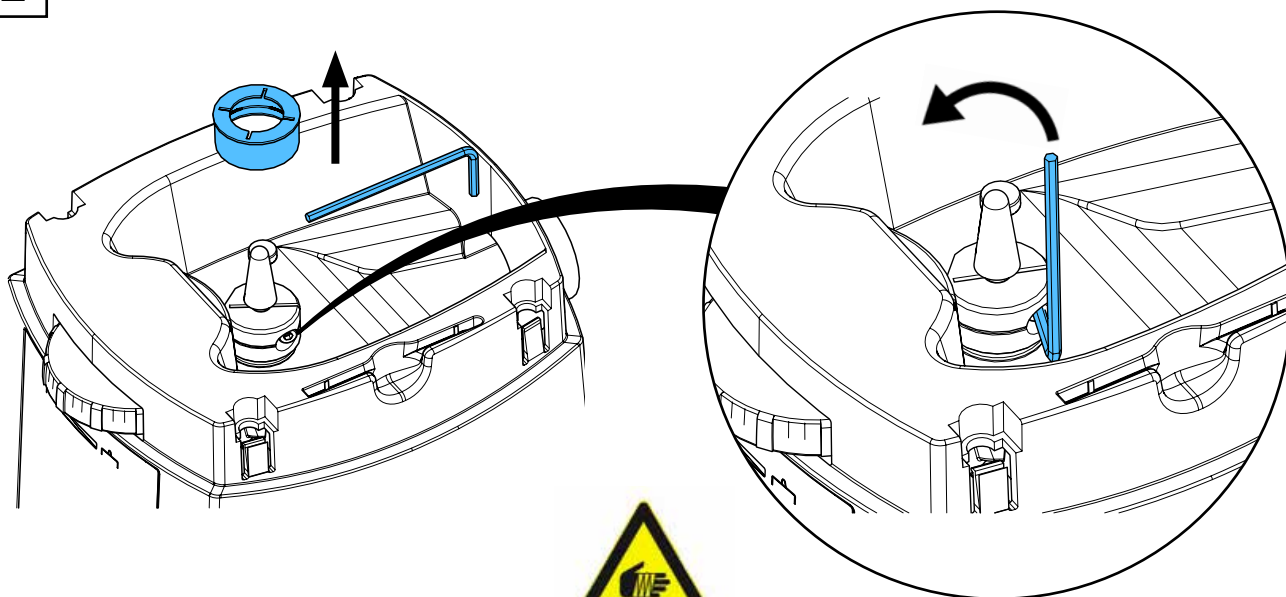



3

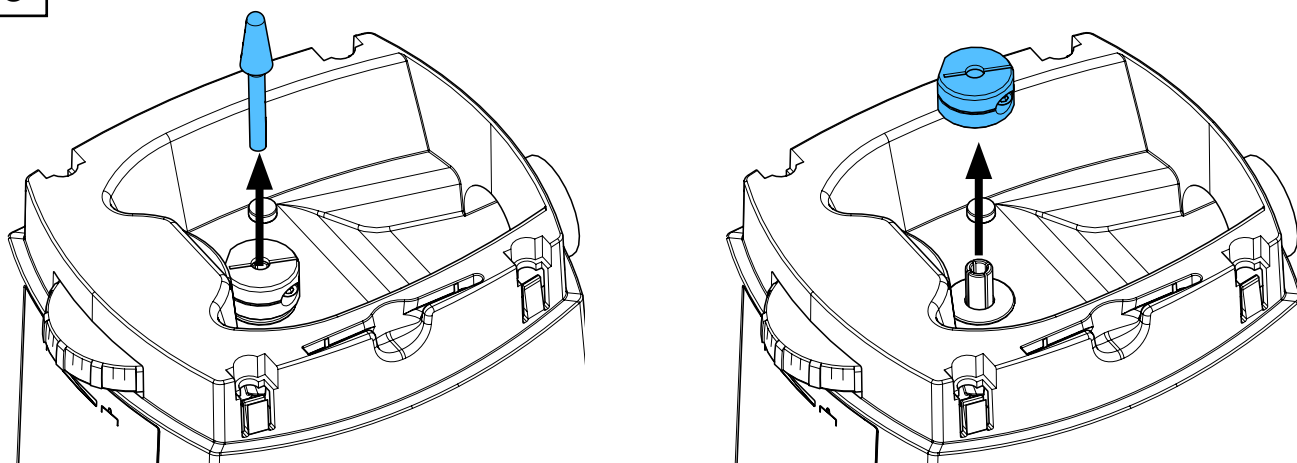



4

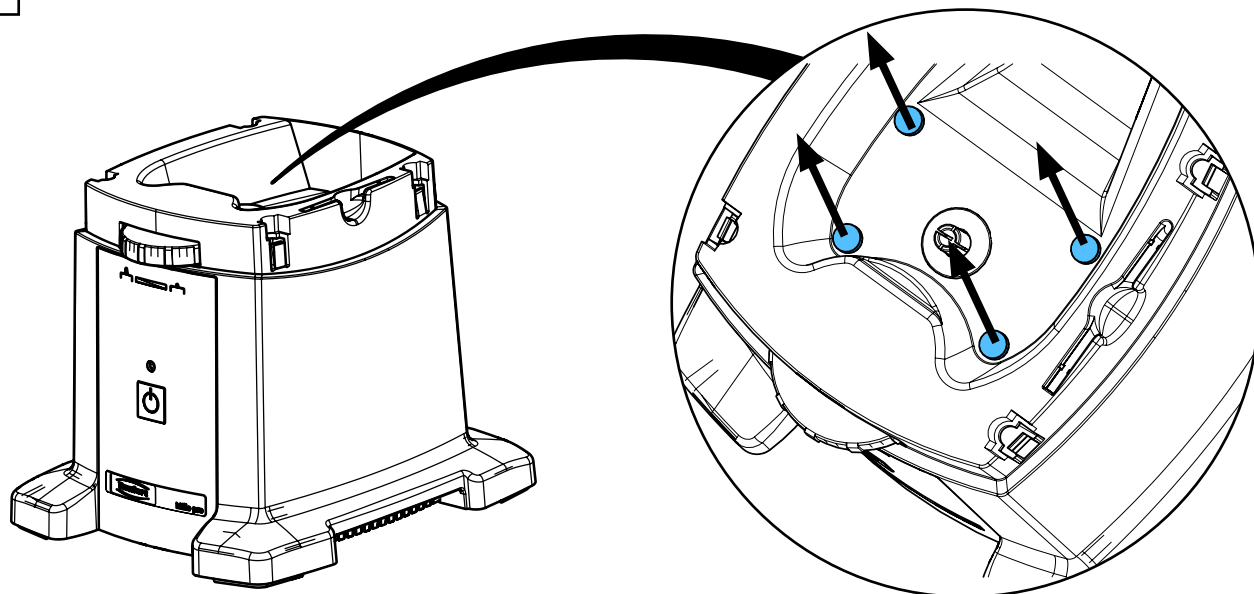



5

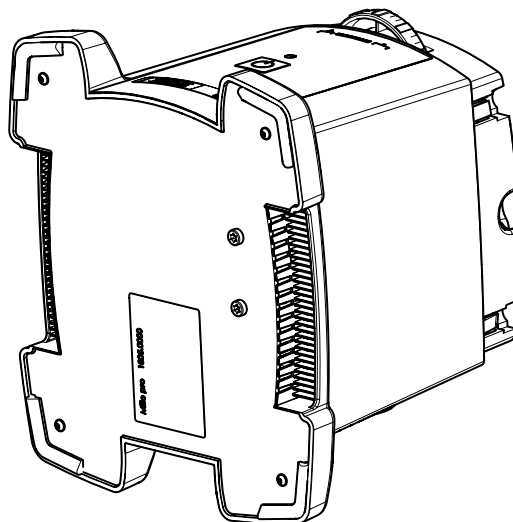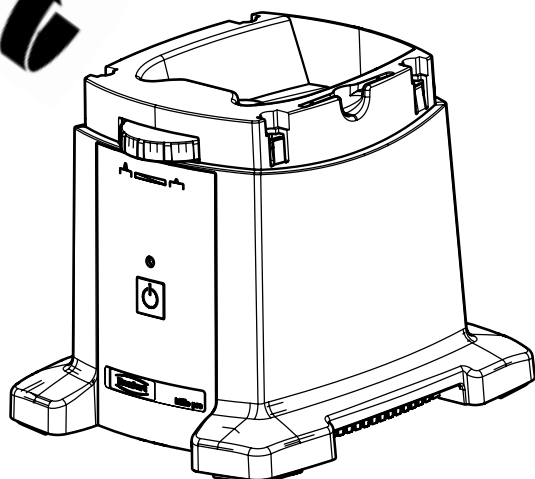


6

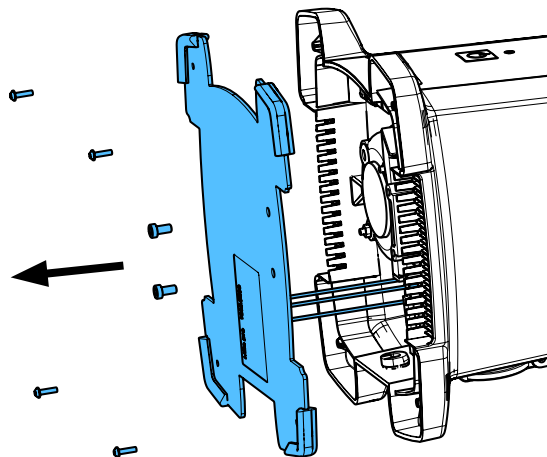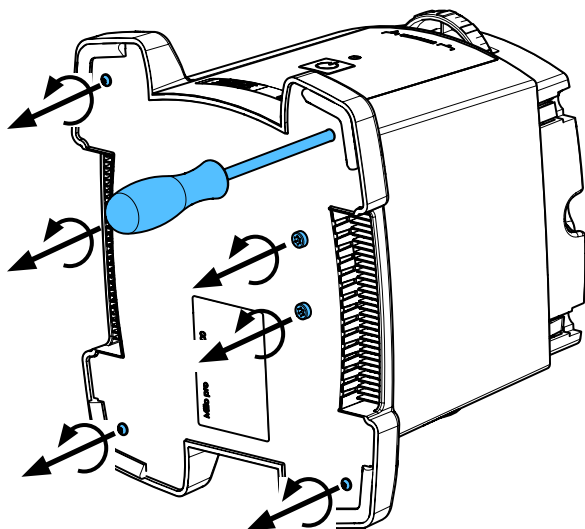


7

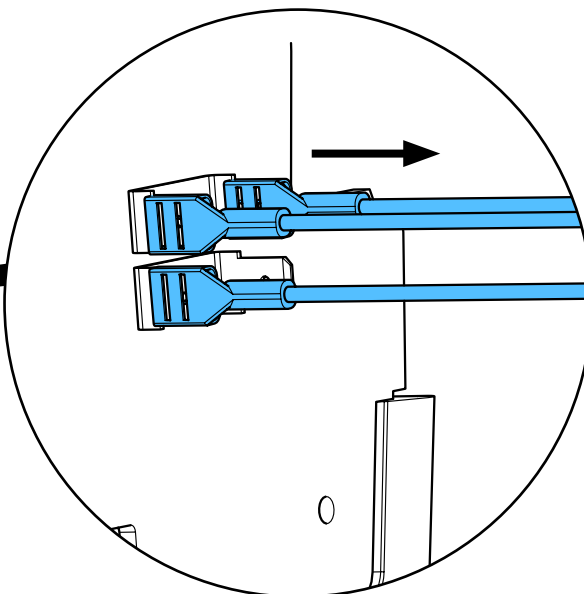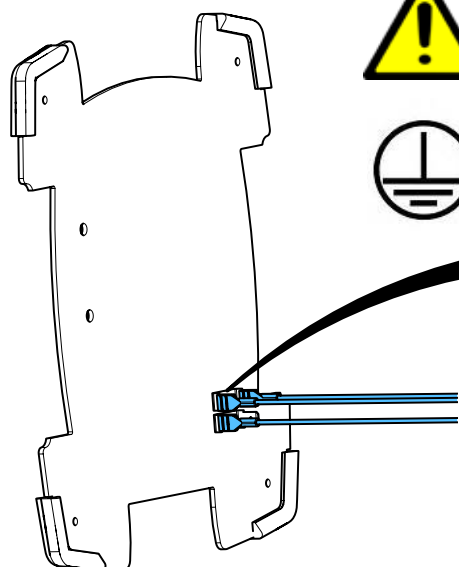


8

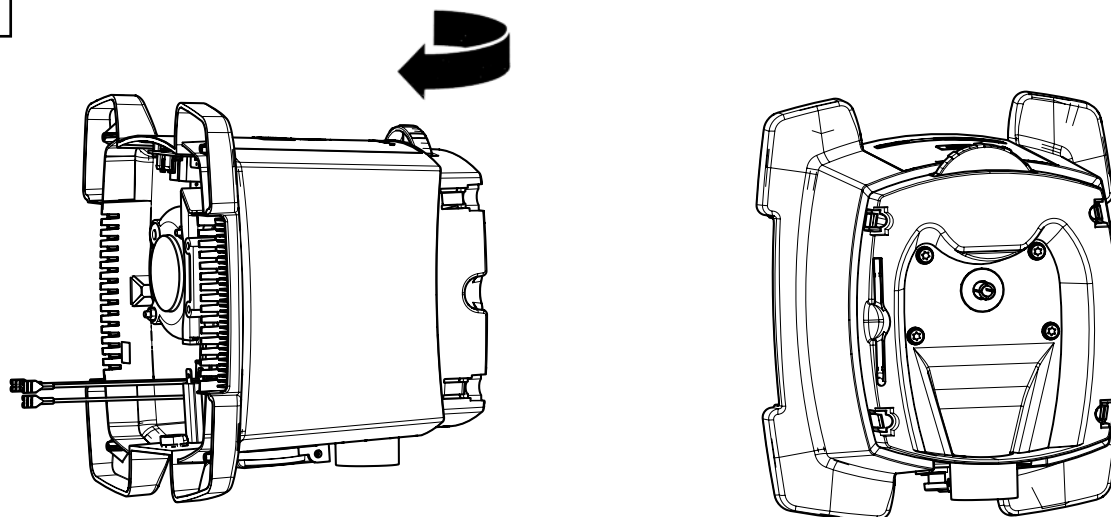





9

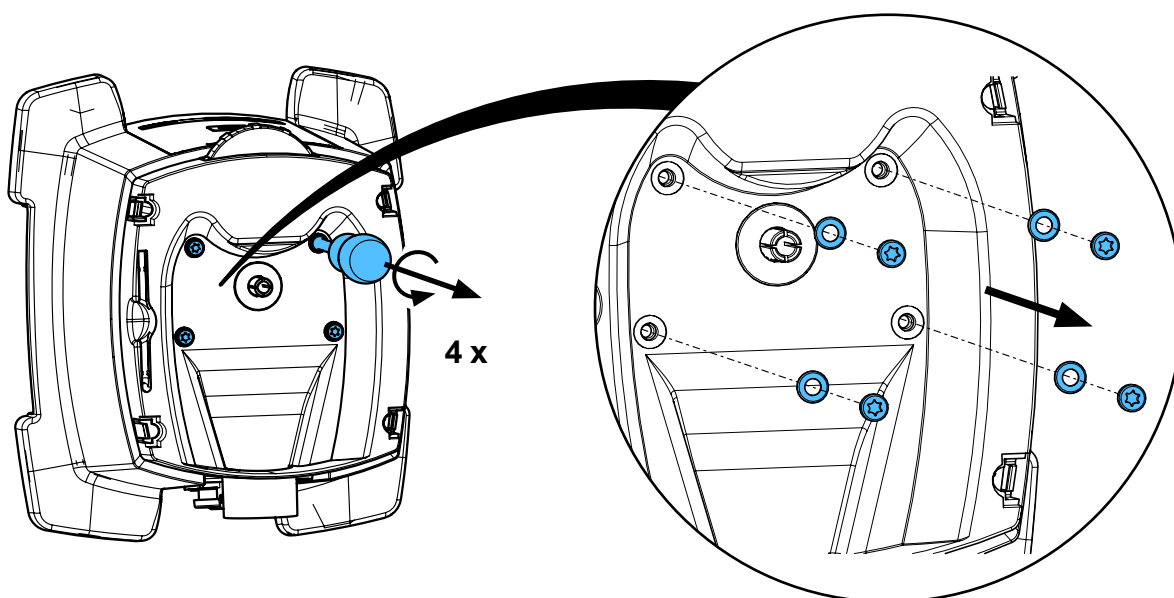


10

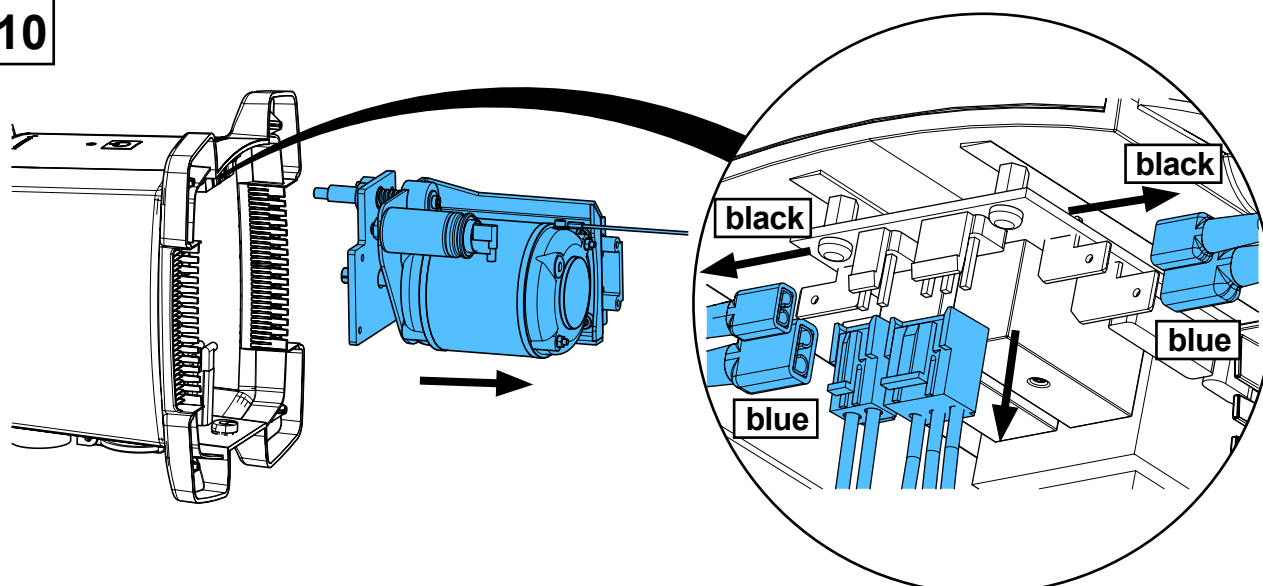


11

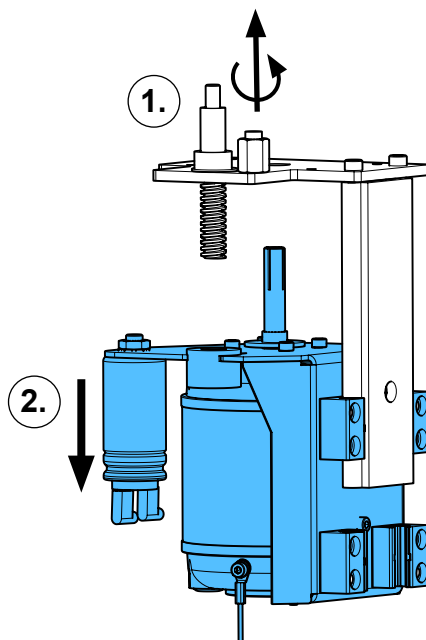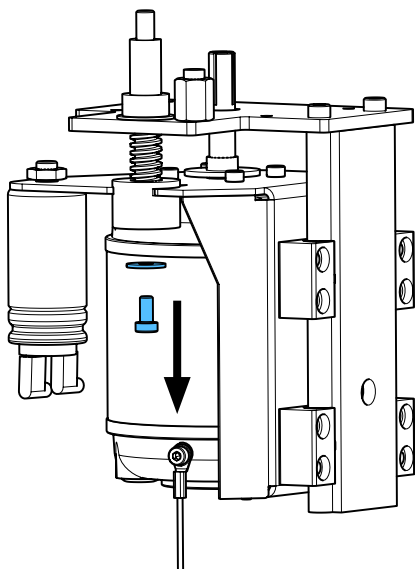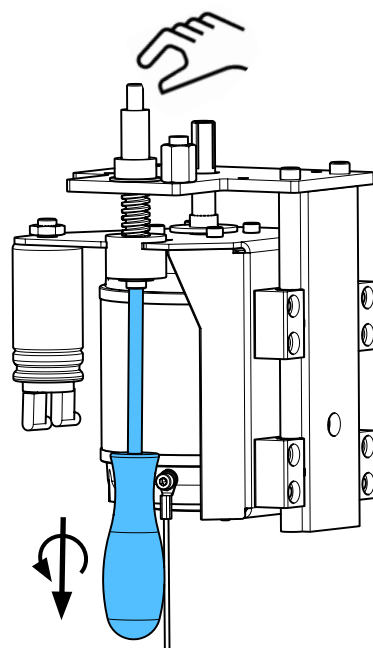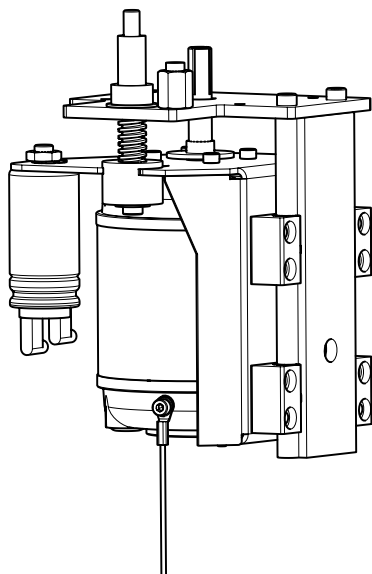


12

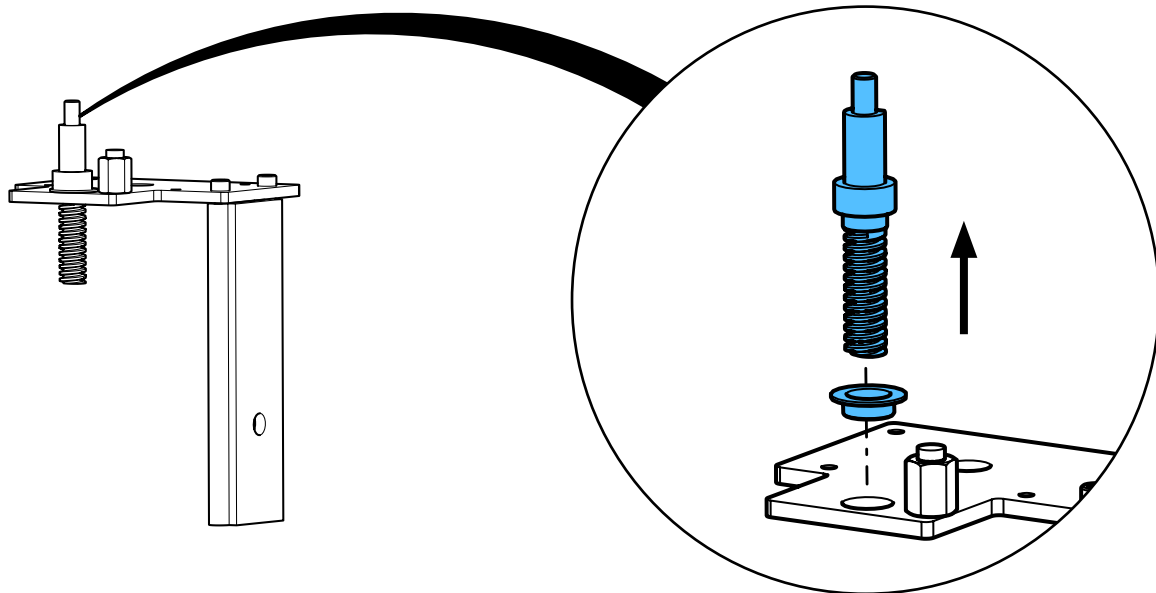



13

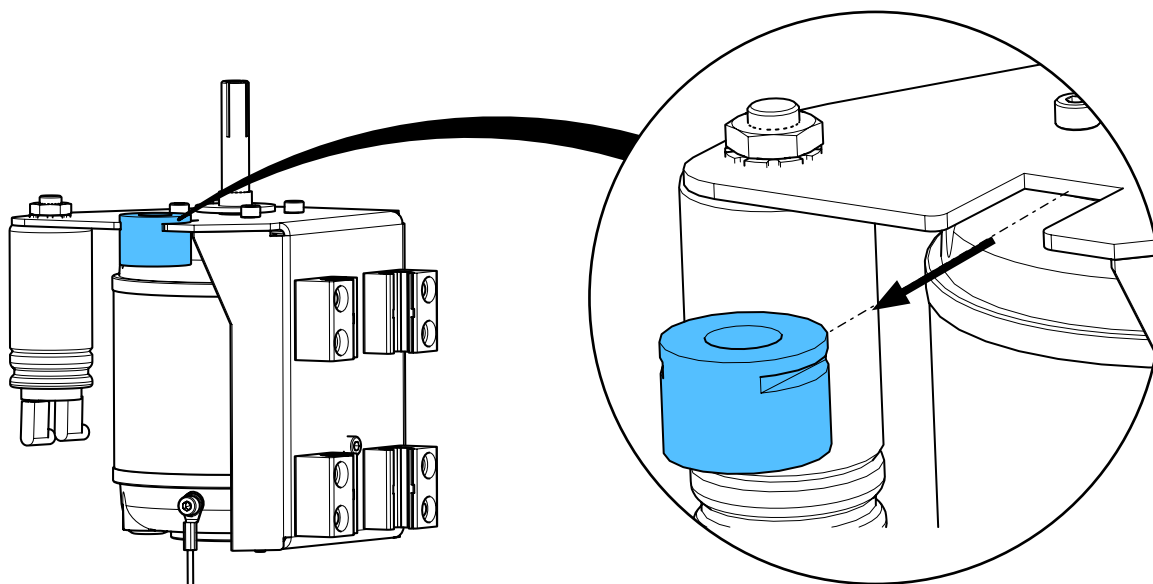





14

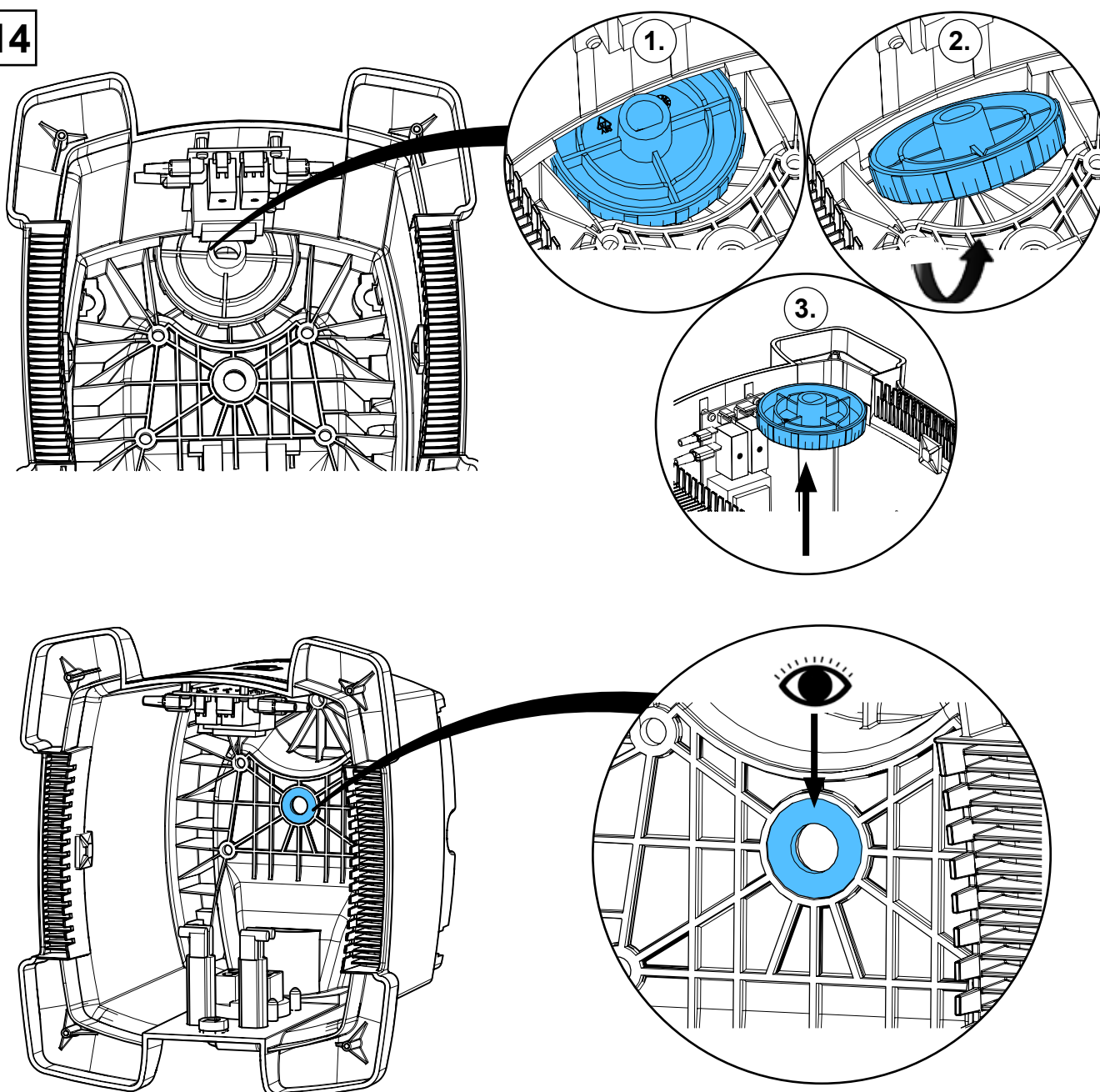



15

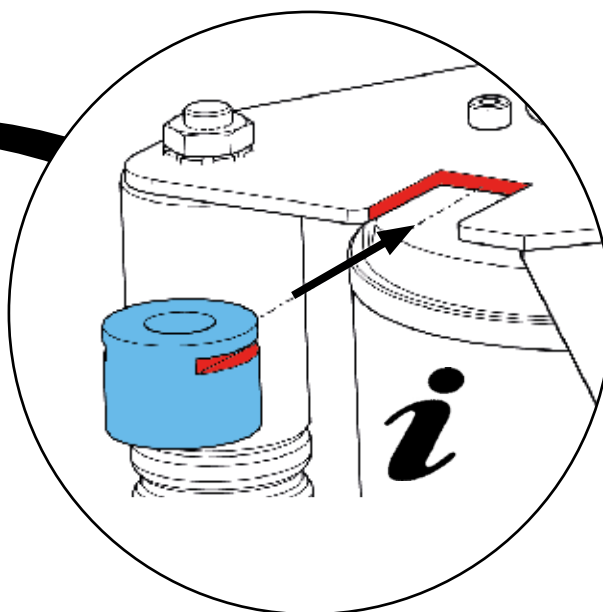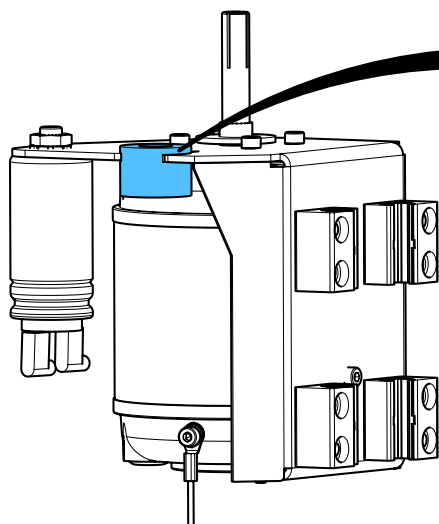


16

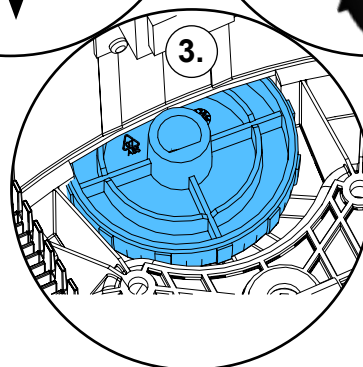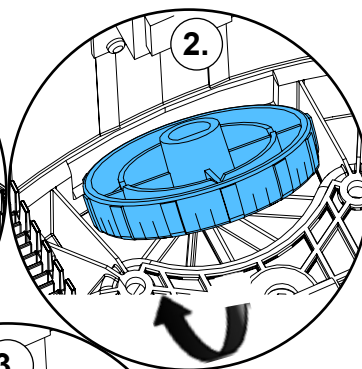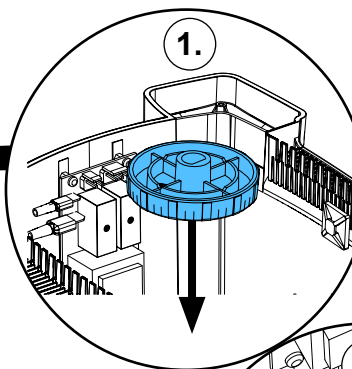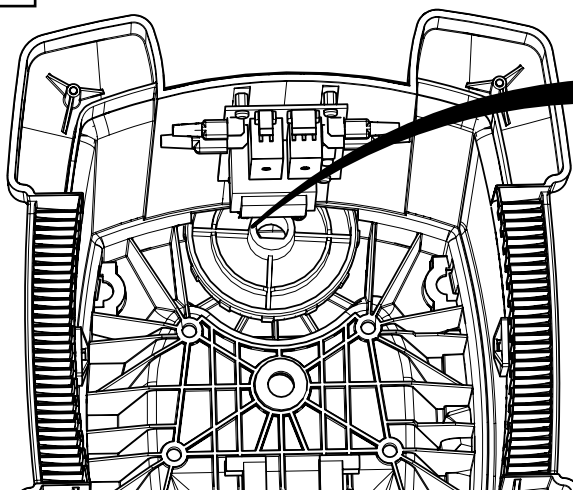




17

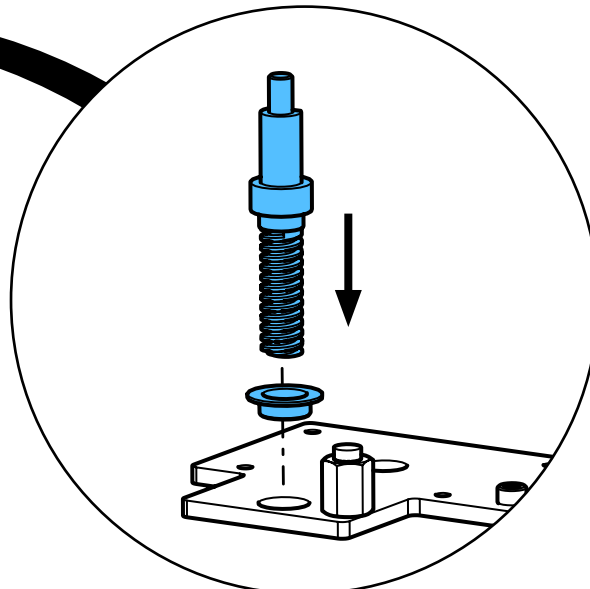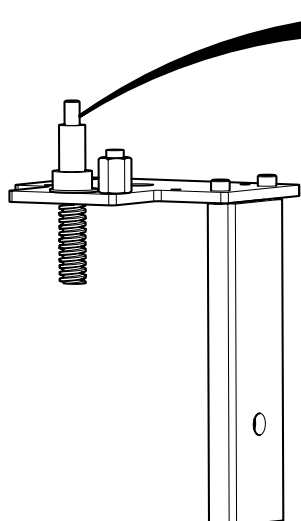


18

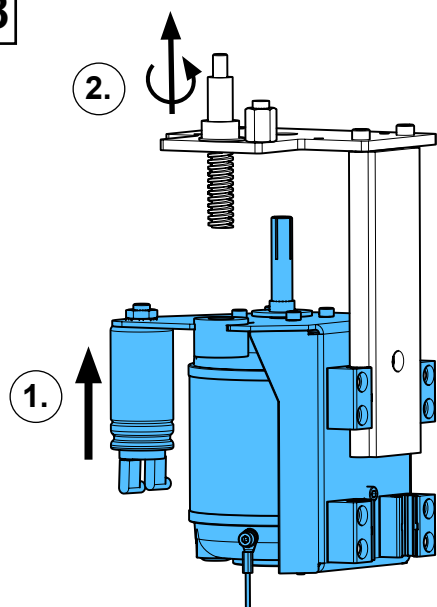


19

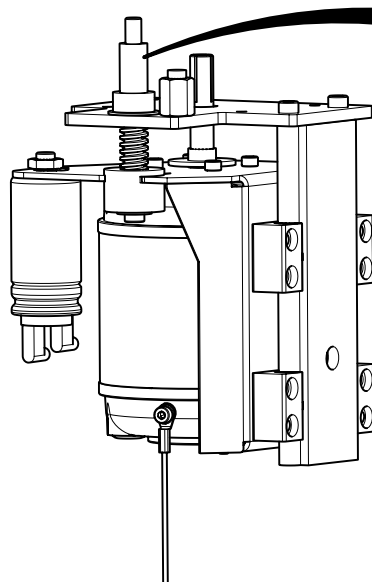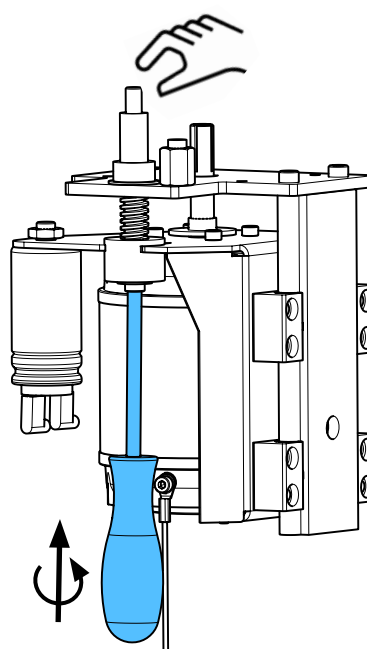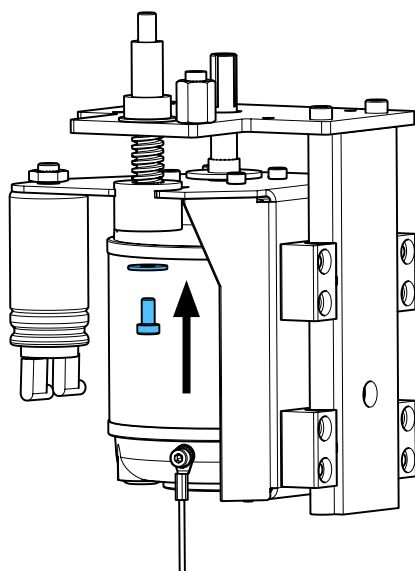


i

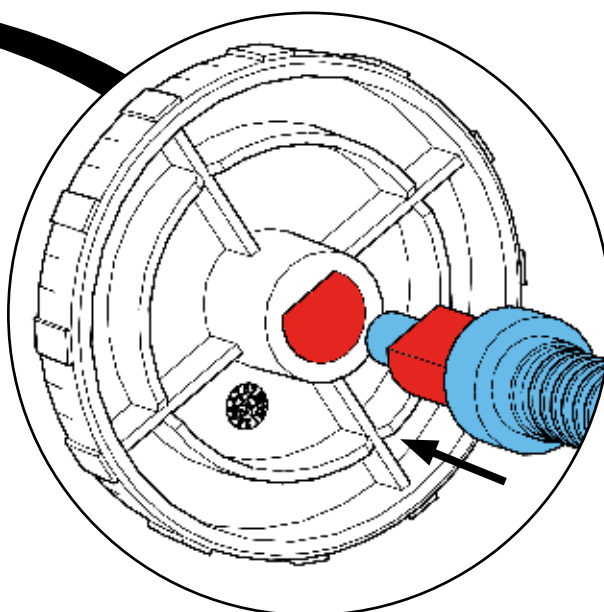


20

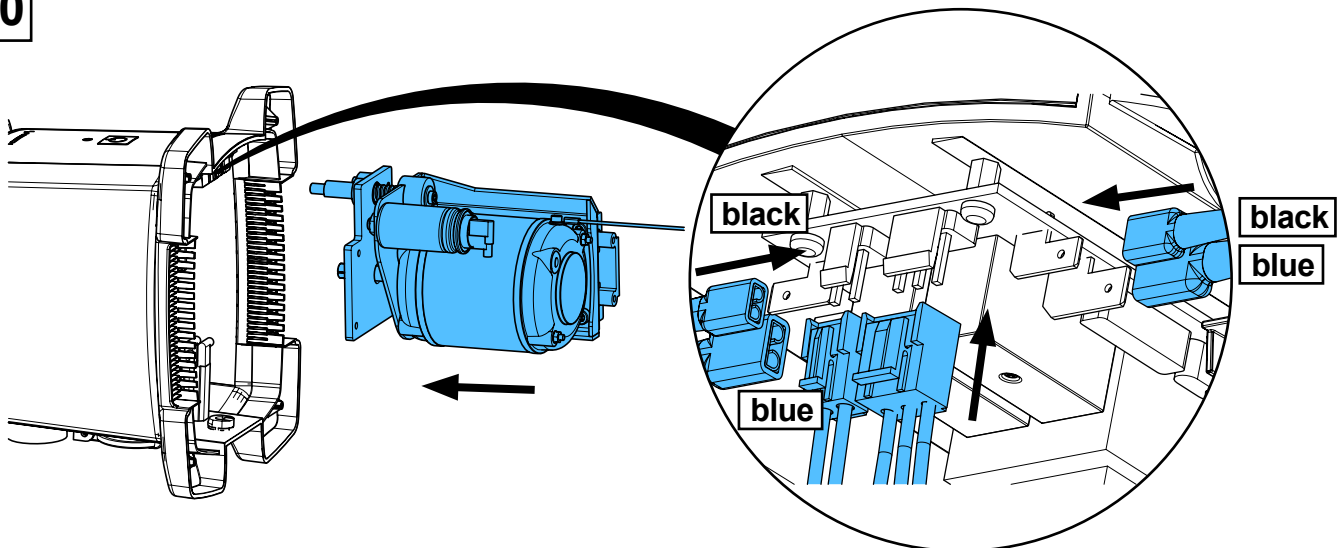


21

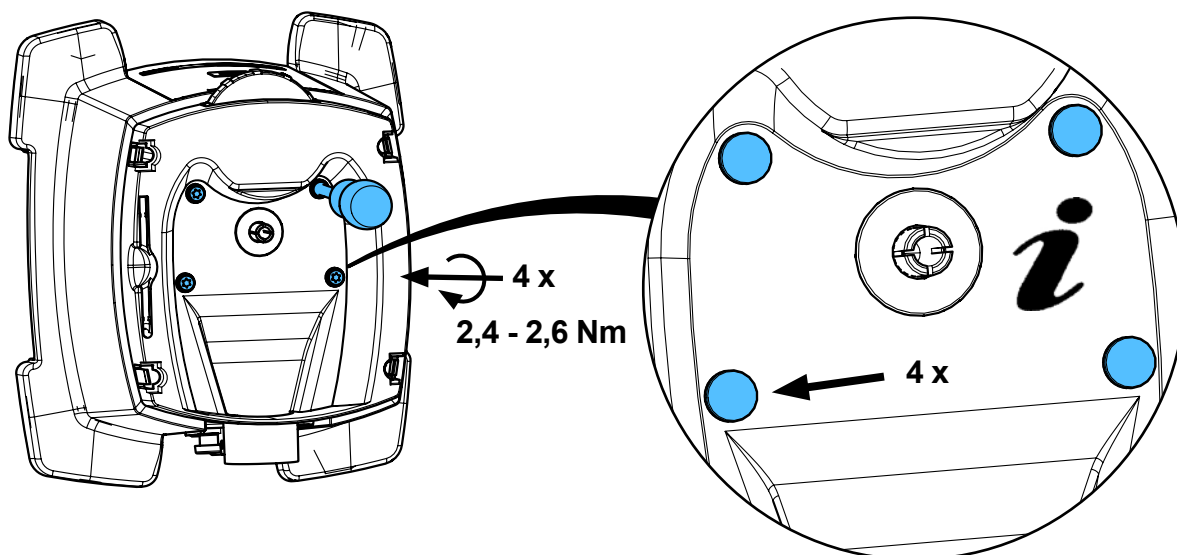


22

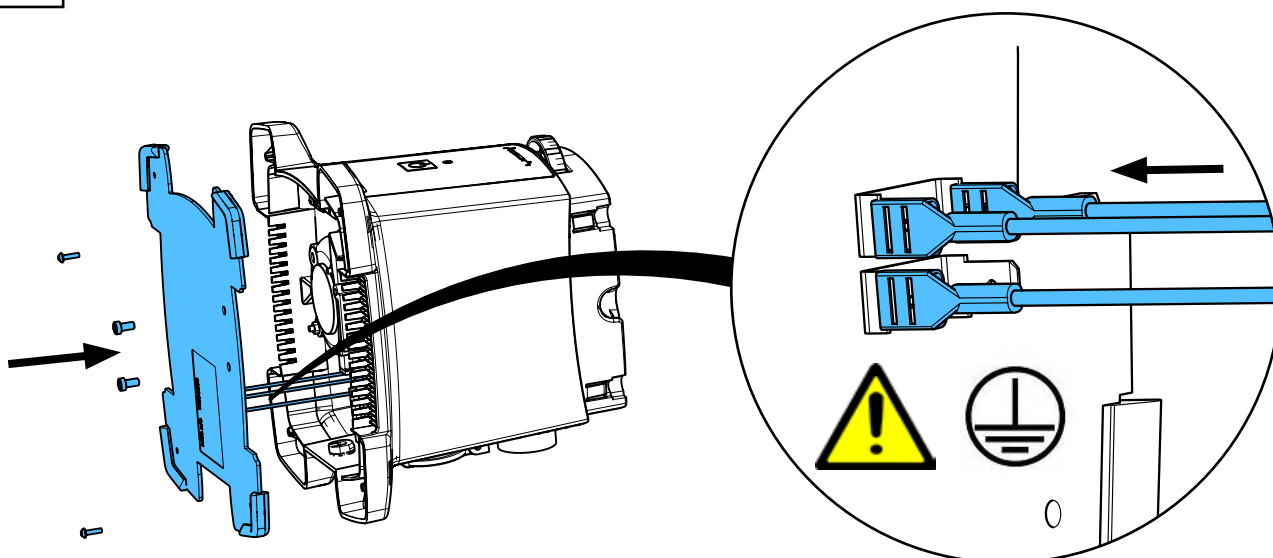


23


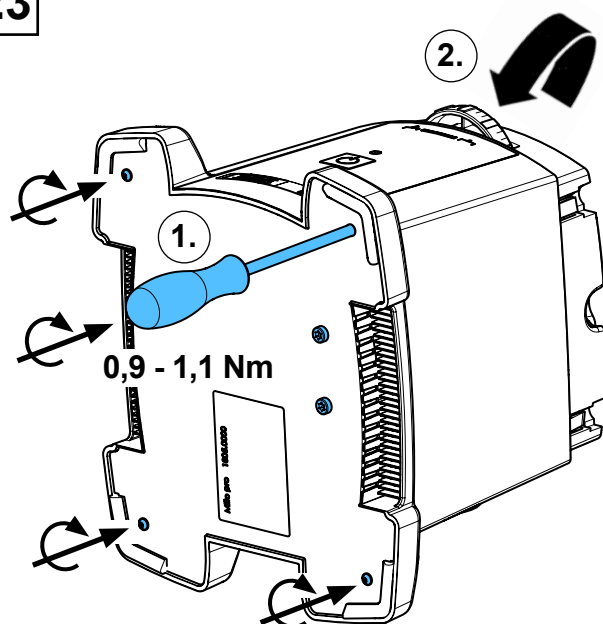


24


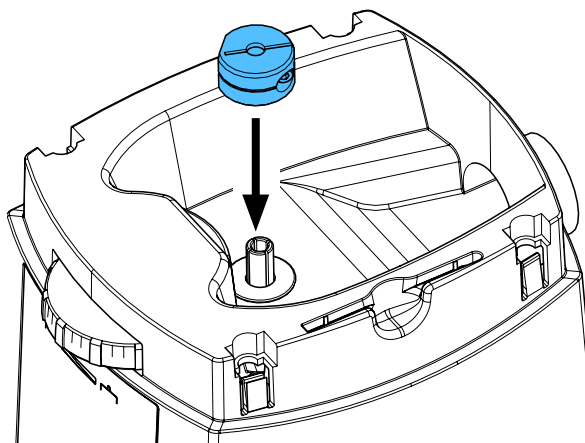


25


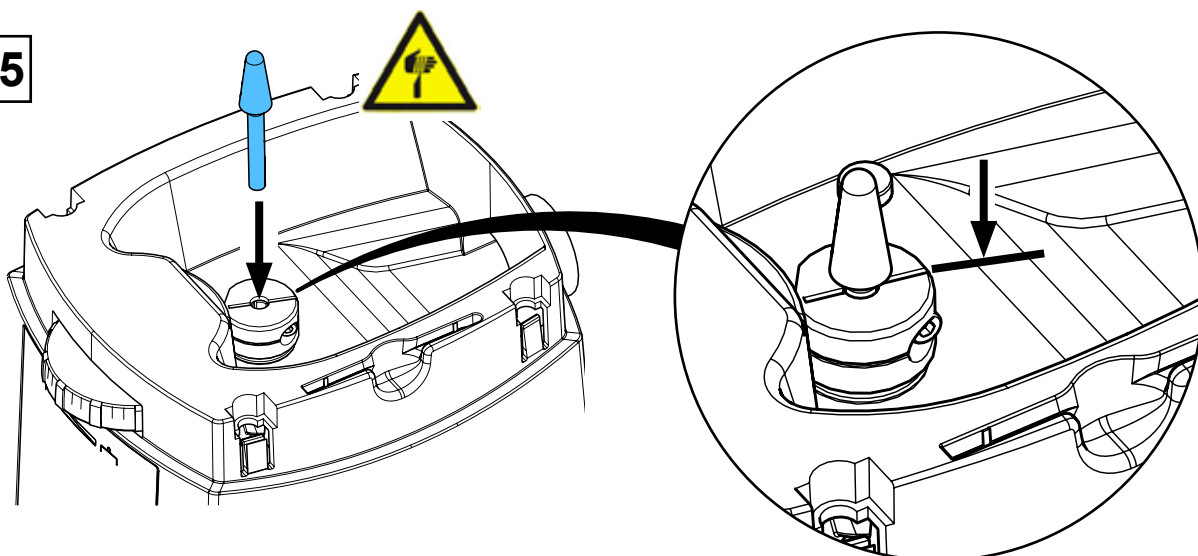


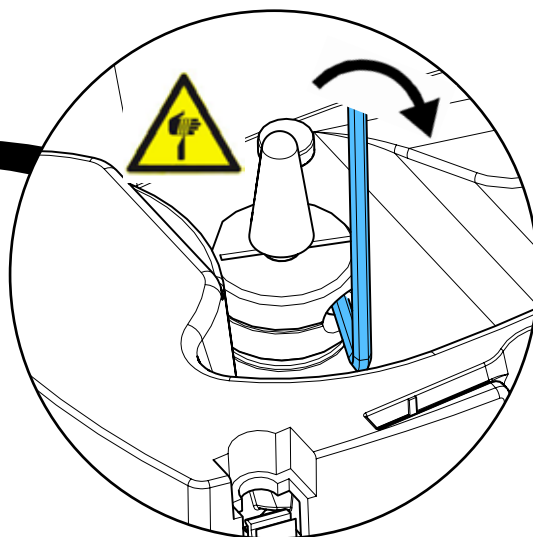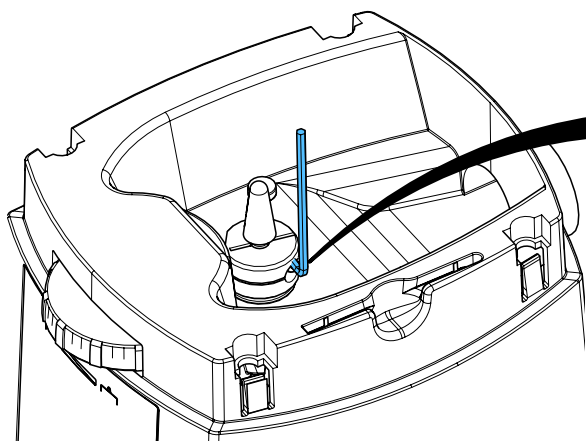

26

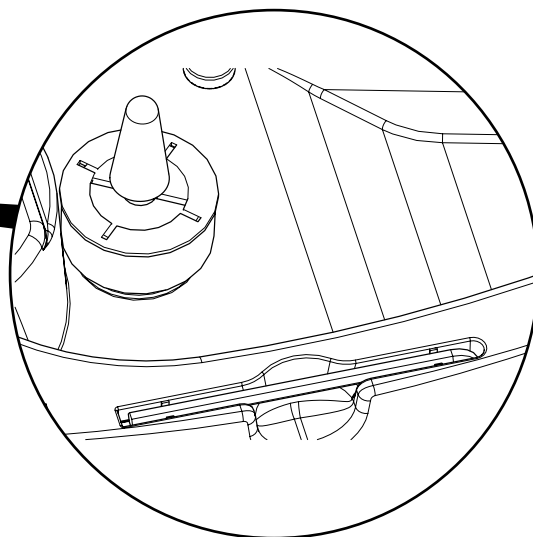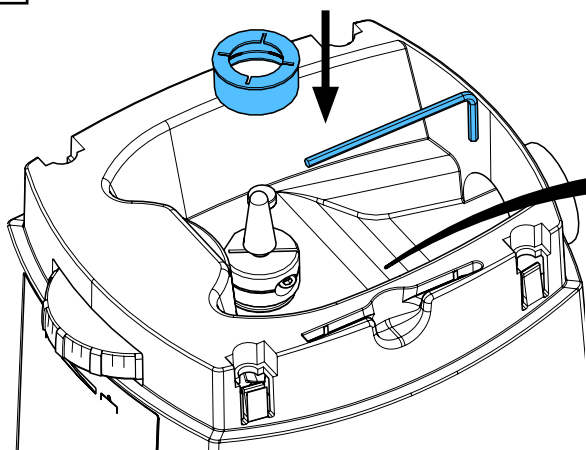

27


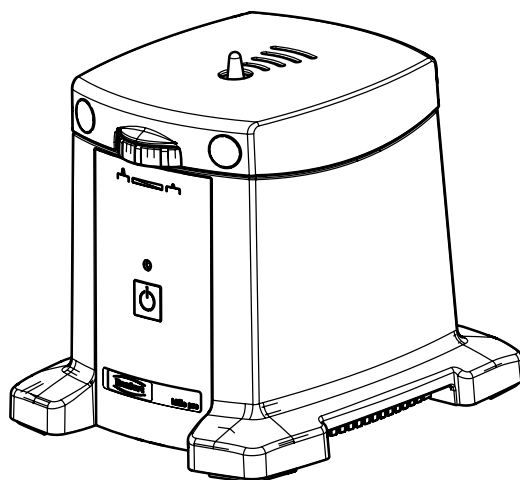
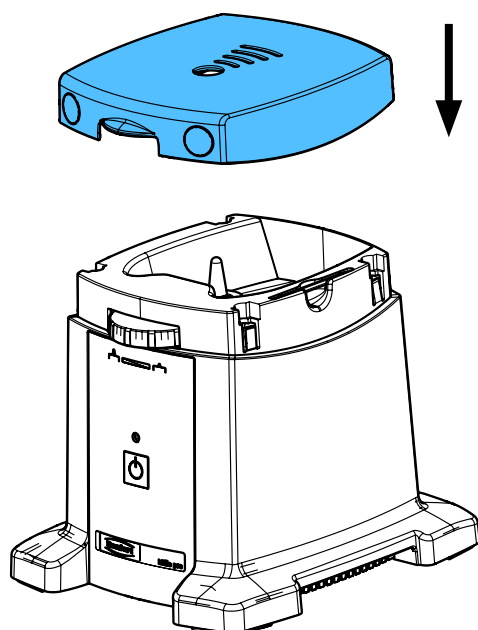


90003 3405



Überstromschutzschalter • Overcurrent circuit breaker

Protecteur contre les surintensités • Protezione da sovratensioni



Protección de sobrecargas • Защитный выключатель при перегрузке тока

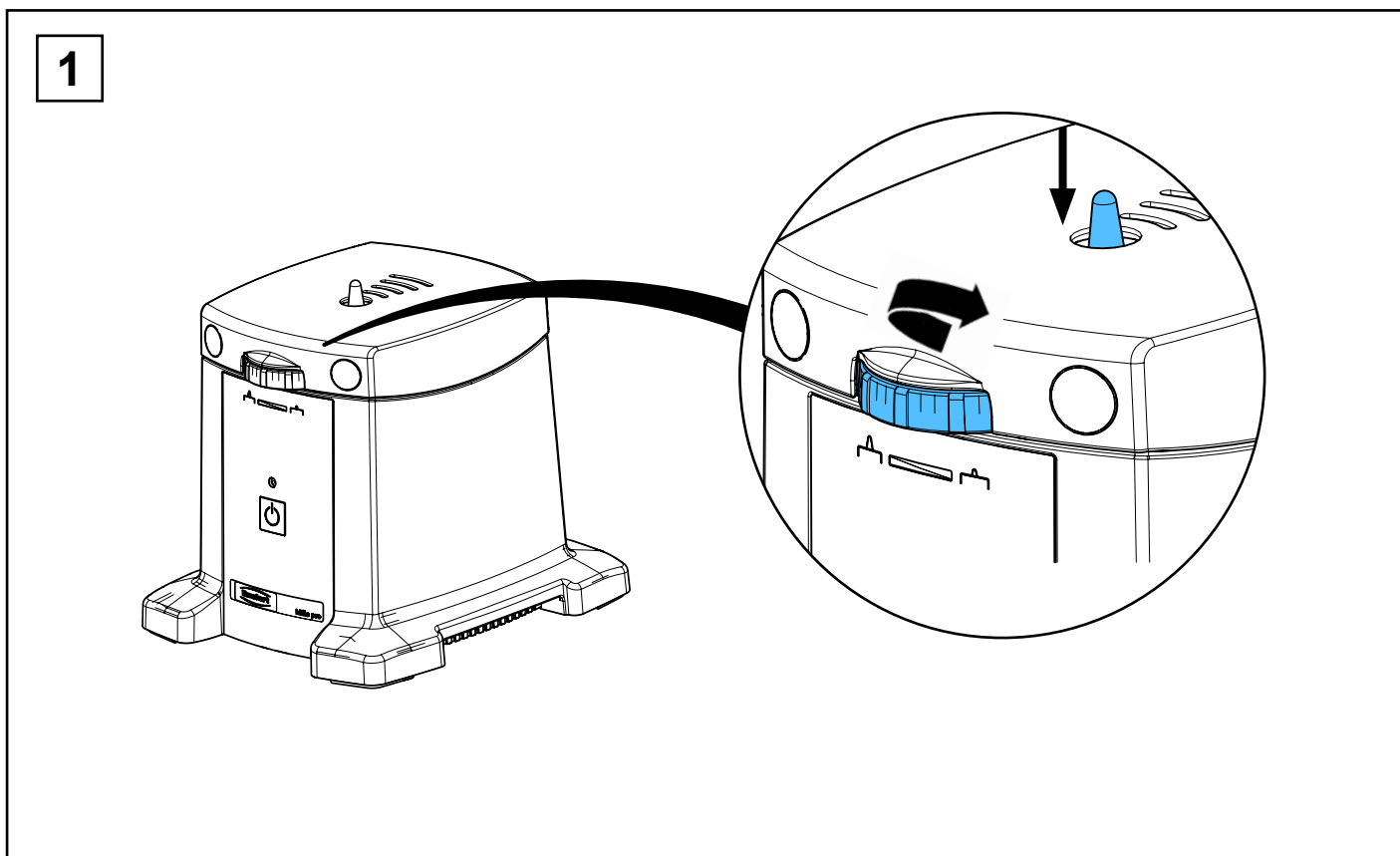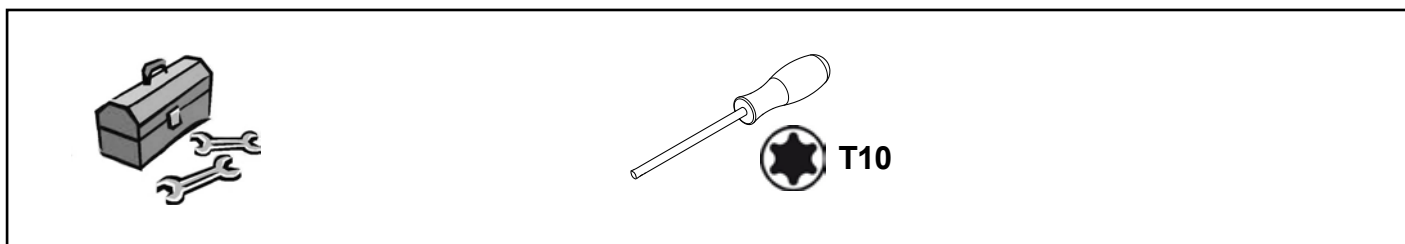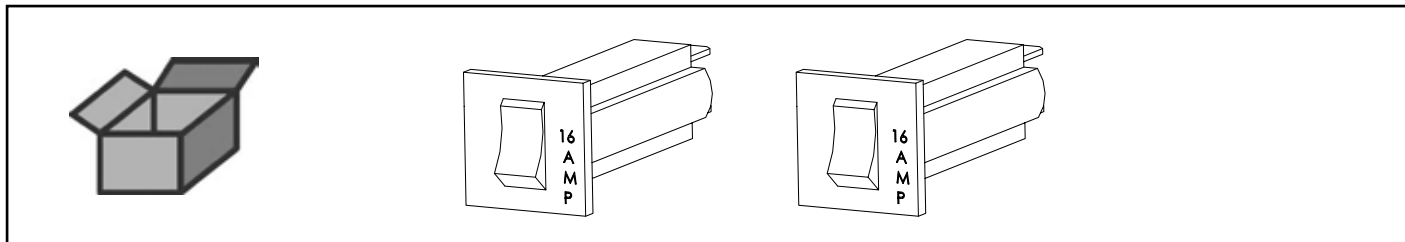



2

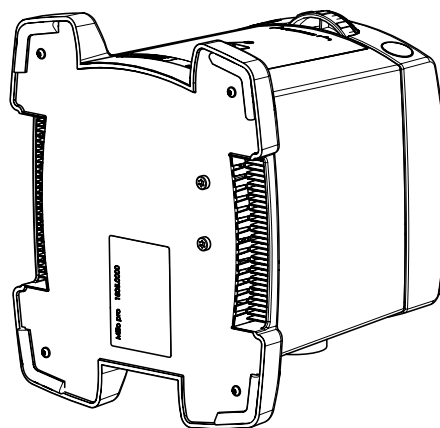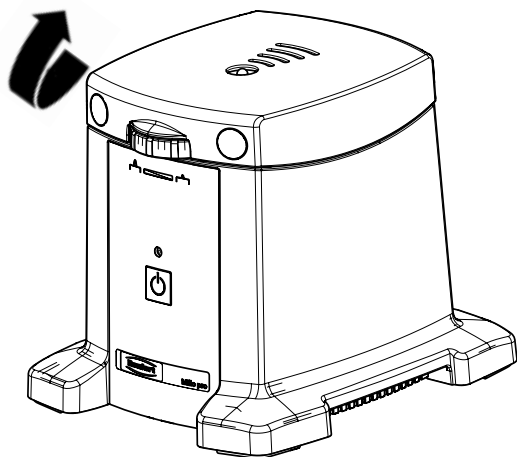



3

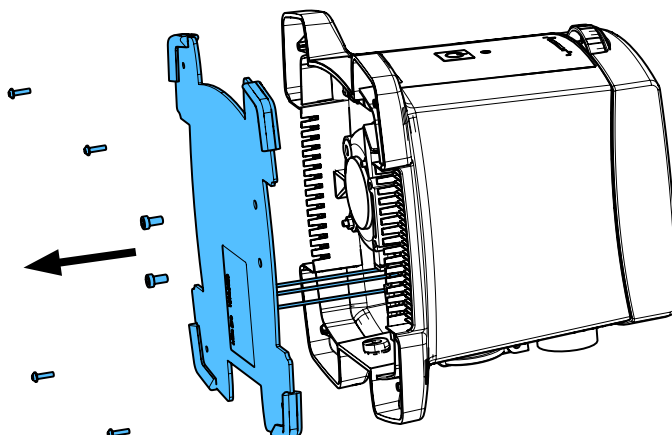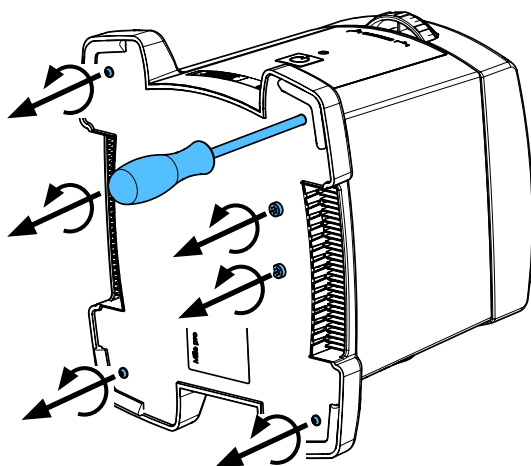





4

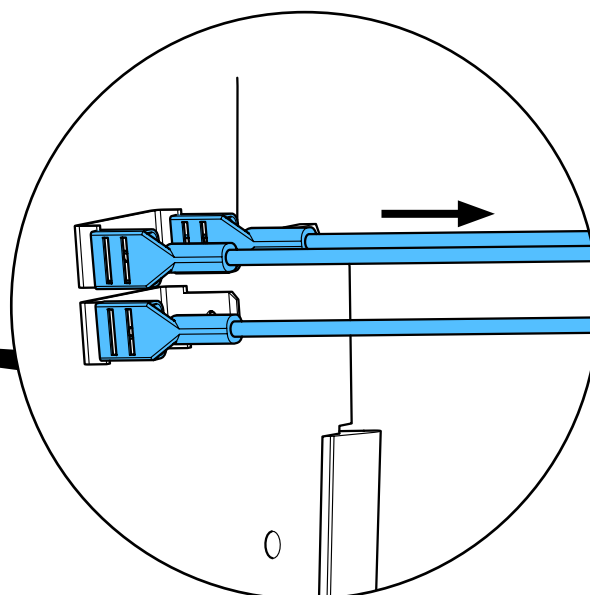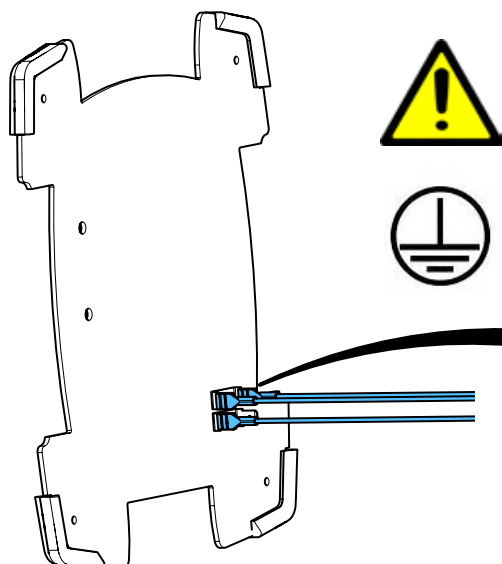



5

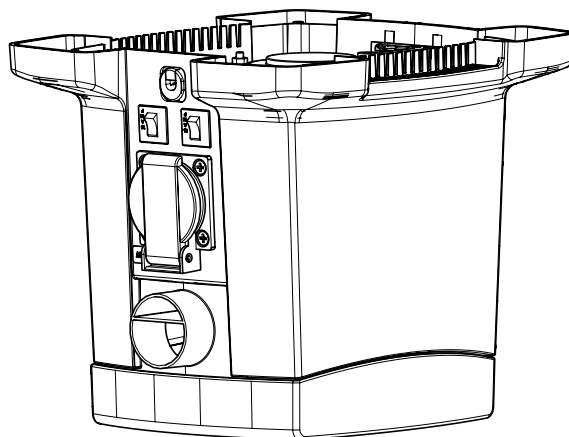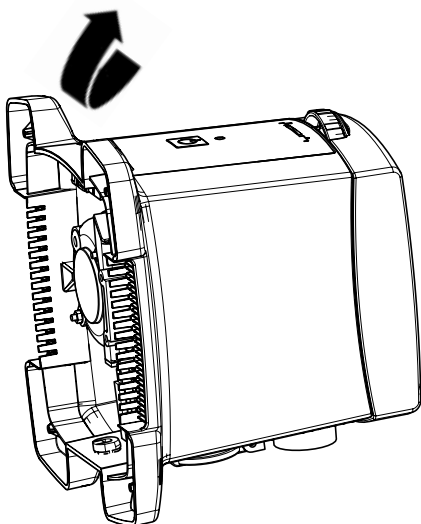


6

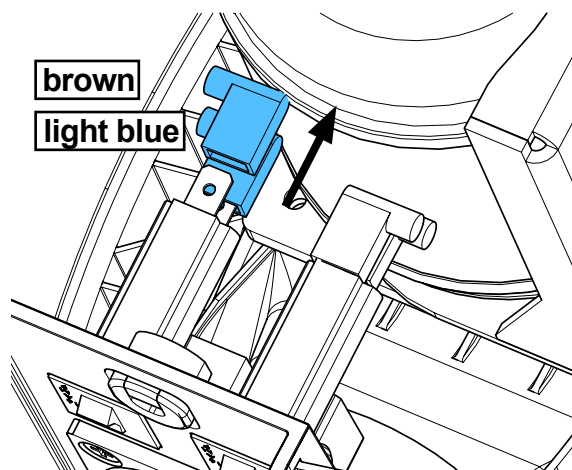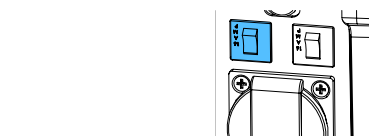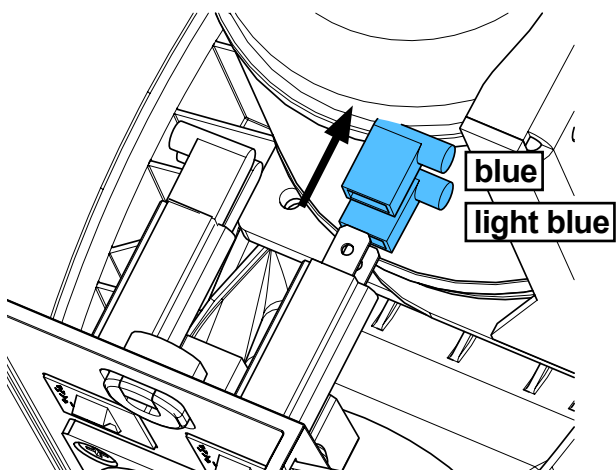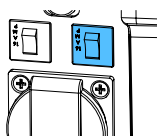


7

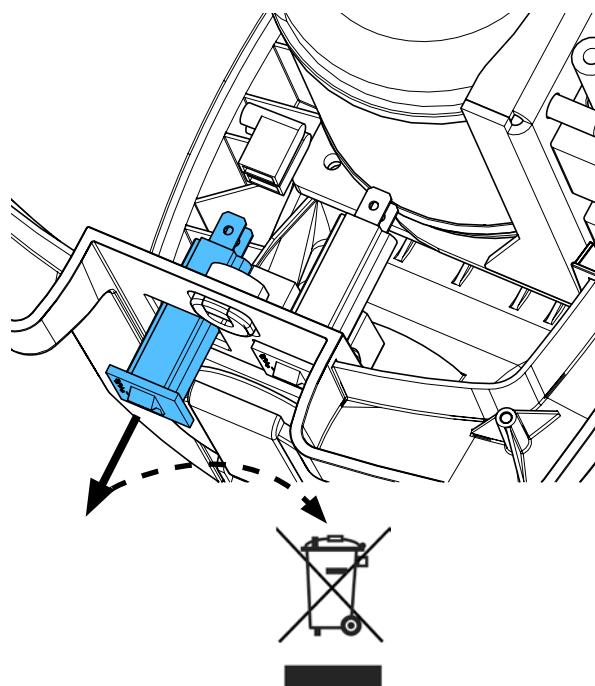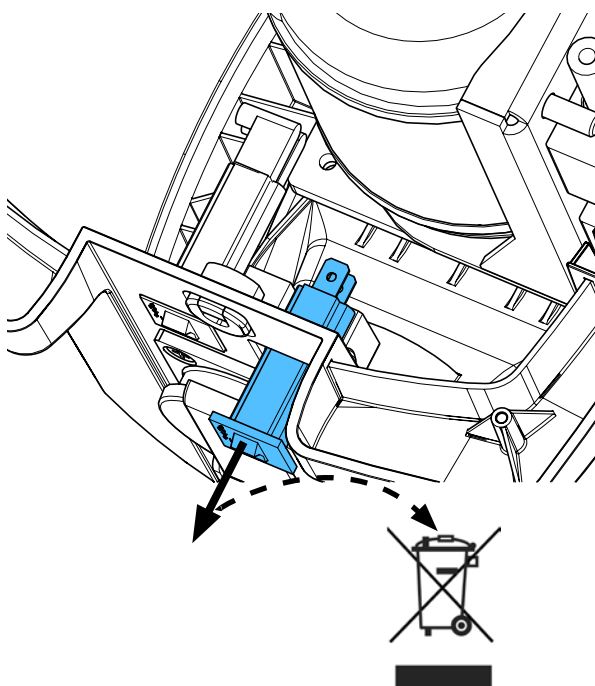


8

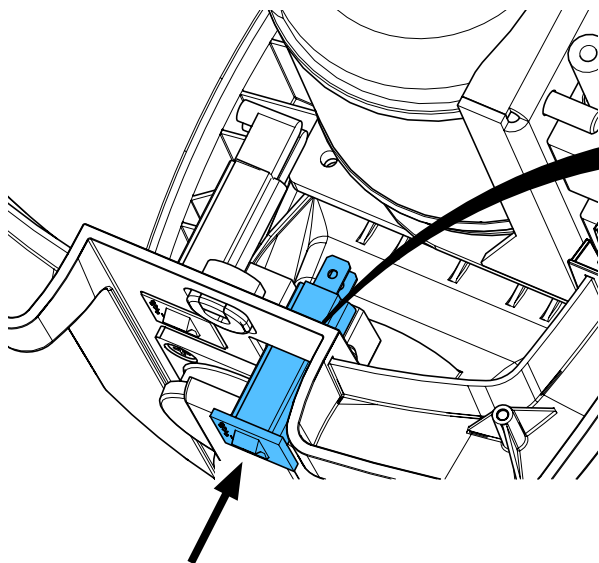




9

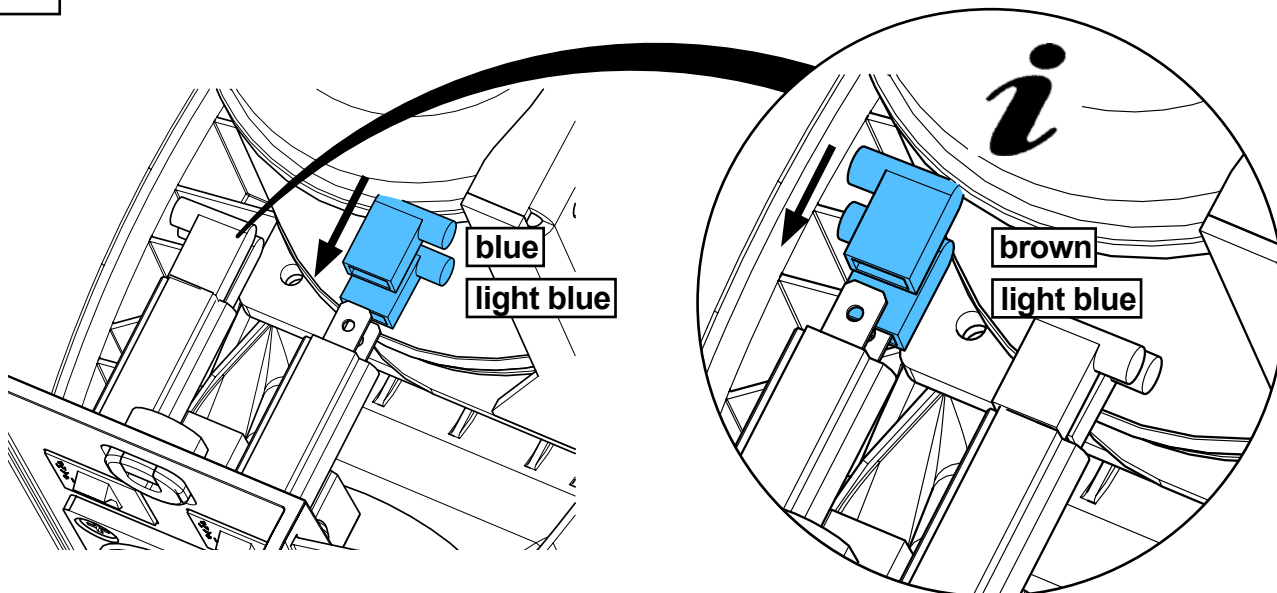



10

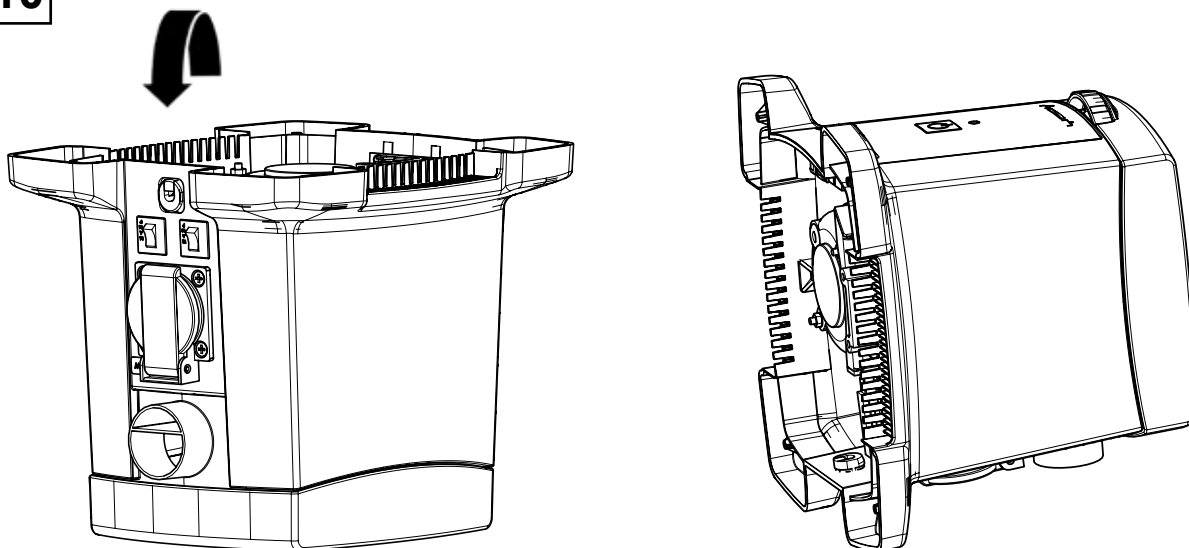

11

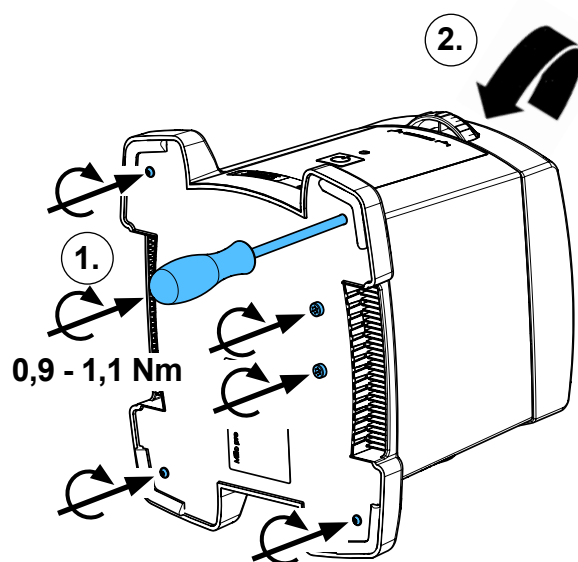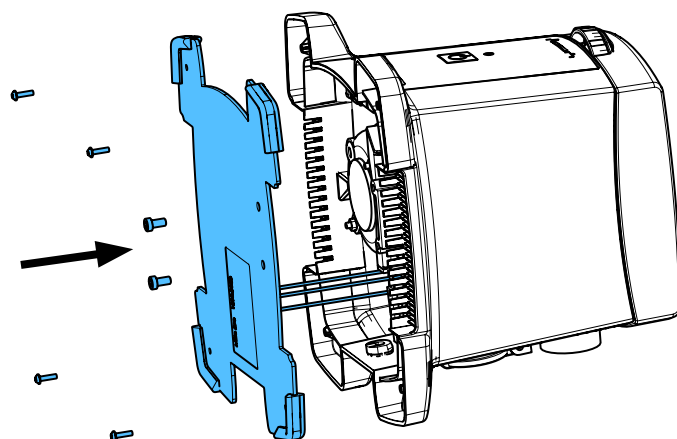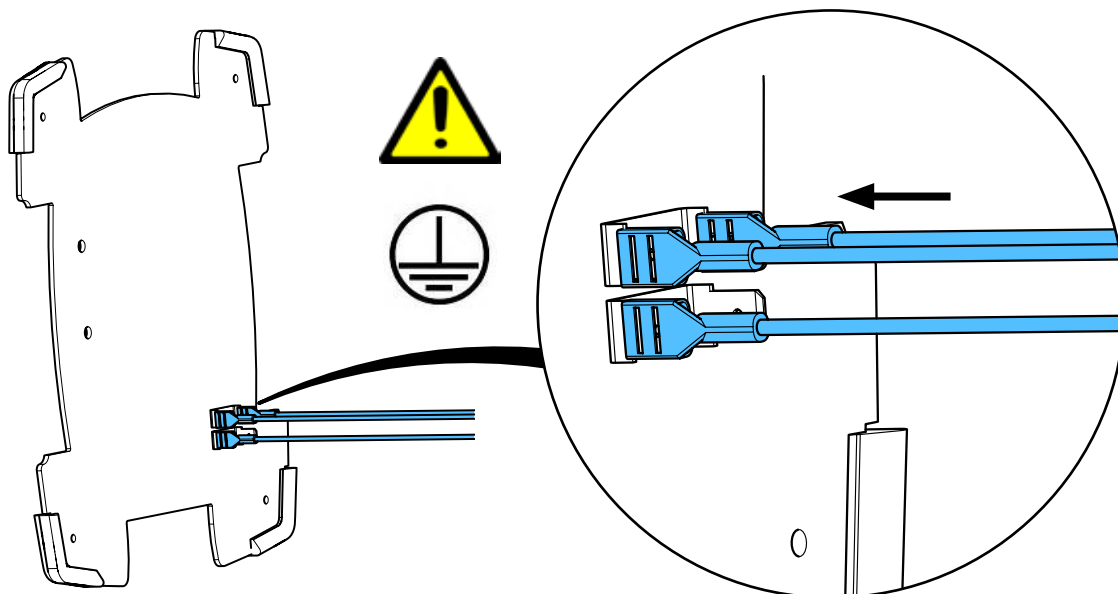

12

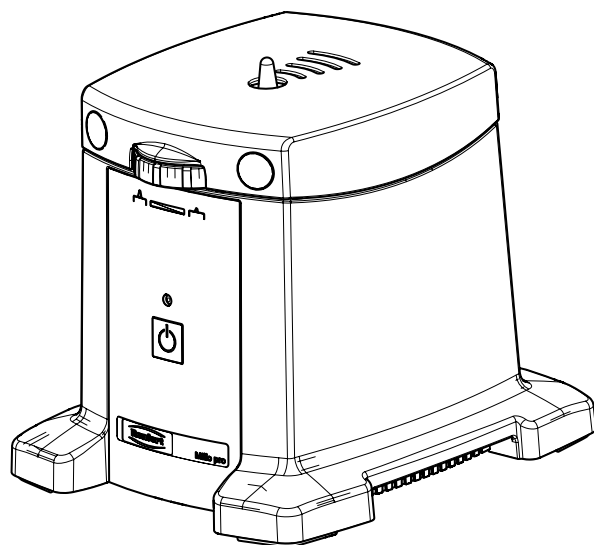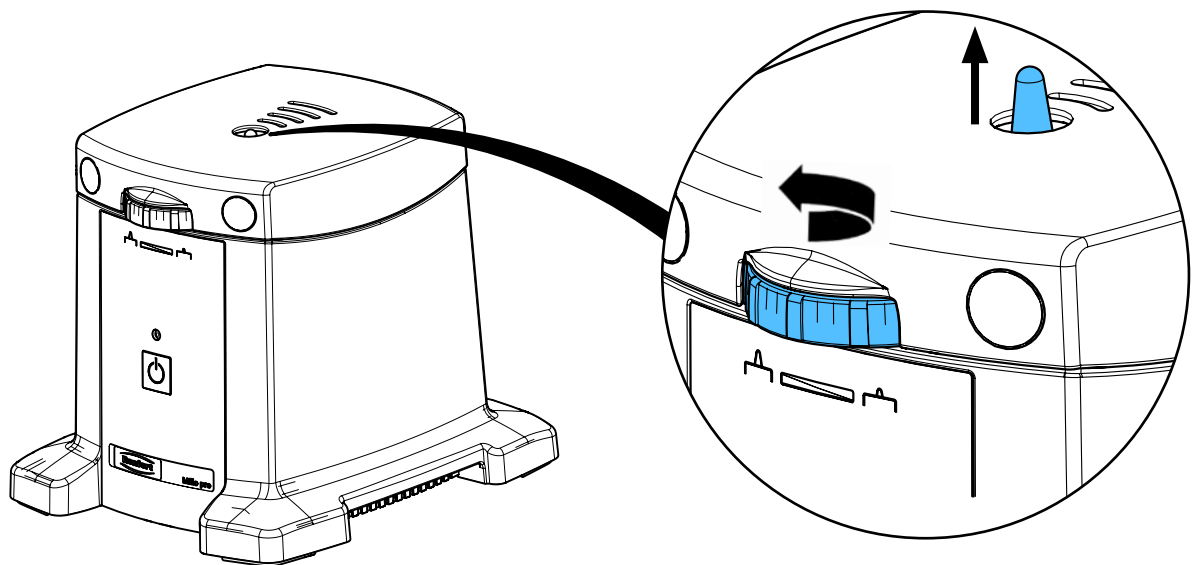

Hochaktuell und ausführlich auf ...
Up to date and in detail at ...
Actualisé et détaillé sous ...
Aggiornato e dettagliato su ...
La máxima actualidad y detalle en ...
Актуально и подробно на ...

www.renfert.com

Renfert GmbH • Untere Gießwiesen 2 • 78247 Hilzingen/Germany
Tel.: +49 7731 82 08-0 • Fax: +49 7731 82 08-70
www.renfert.com • info@renfert.com

Renfert USA • 3718 Illinois Avenue • St. Charles IL 60174/USA
Tel.: +1 6307 62 18 03 • Fax: +1 6307 62 97 87
www.renfert.com • info@renfertusa.com
USA: Free call 800 336 7422