




making work easy

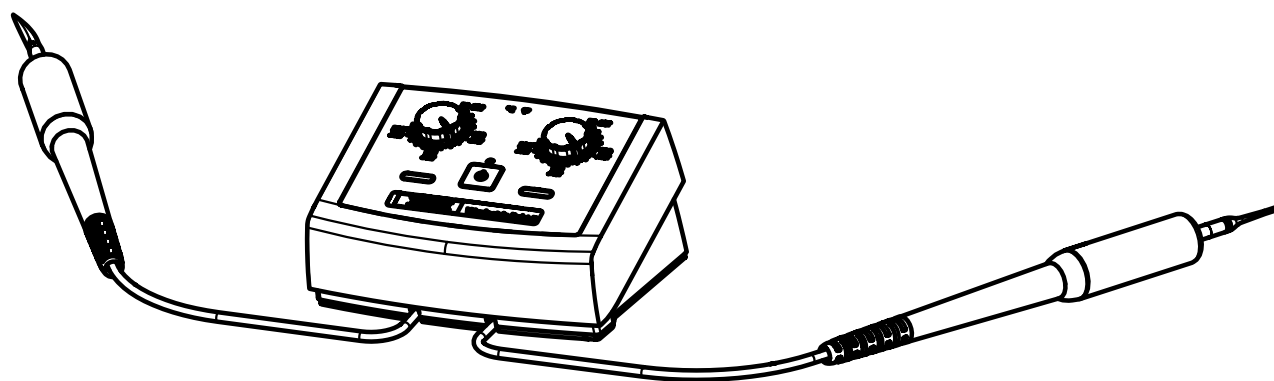

Waxelectric light II

Service Instructions

Made in Germany

21-6921 05122025

Contents

2154 0001 / 2154 0002

Heizgriff • Heating handle • Manche chauffant • Manipolo • Mango • Нагревательная рукоятка

90003 5948

Griffstück Kork • Cork-grip • Manche en liège • Impugnatura di sughero • Empuñadura de corcho • Зажим для пробки

90003 6138 / 90003 3752

Netzteil • Power supply • Poste secteur • Trasformatore • Trasformatore • Блок питания

90003 2962

O-Ring • Retainer ring • Joint thorique • Guarnizioni • Anillo en O • Блок питания

90003 4564

Gehäusefüsse • Rubber bases • Pieds en caoutchouc • Piedini di gomma • Pies de goma • Ножки прибора

90003 5975 / ..5976 / ..8652 / ..8653 / ..8654 / ..8680 / ..2773

Netzkabel • Power cord • Cordon d'alimentation • Cavo di rete • Cable de red Сетевой кабель

90003 5971

Leiterplatte Waxlectric light II • Mains circuit board Waxlectric light II • Circuit imprimé Waxlectric light II • Circuito stampato Waxlectric light II • Circuito impresso Waxlectric light II • Печатная плата Waxlectric light II

90003 5953

Bedienfolie Waxlectric light II • Operating foil Waxlectric light II • Feuille de comande Waxlectric light II • Lamina di comando Waxlectric light II • Lámina de mando Waxlectric light II • Плёнка для панели управления Waxlectric light II





Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу


making work easy

2154 0001 / 2154 0002


Heizgriff • Heating handle • Manche chauffant • Manipolo • Mango •
Нагревательная рукоятка

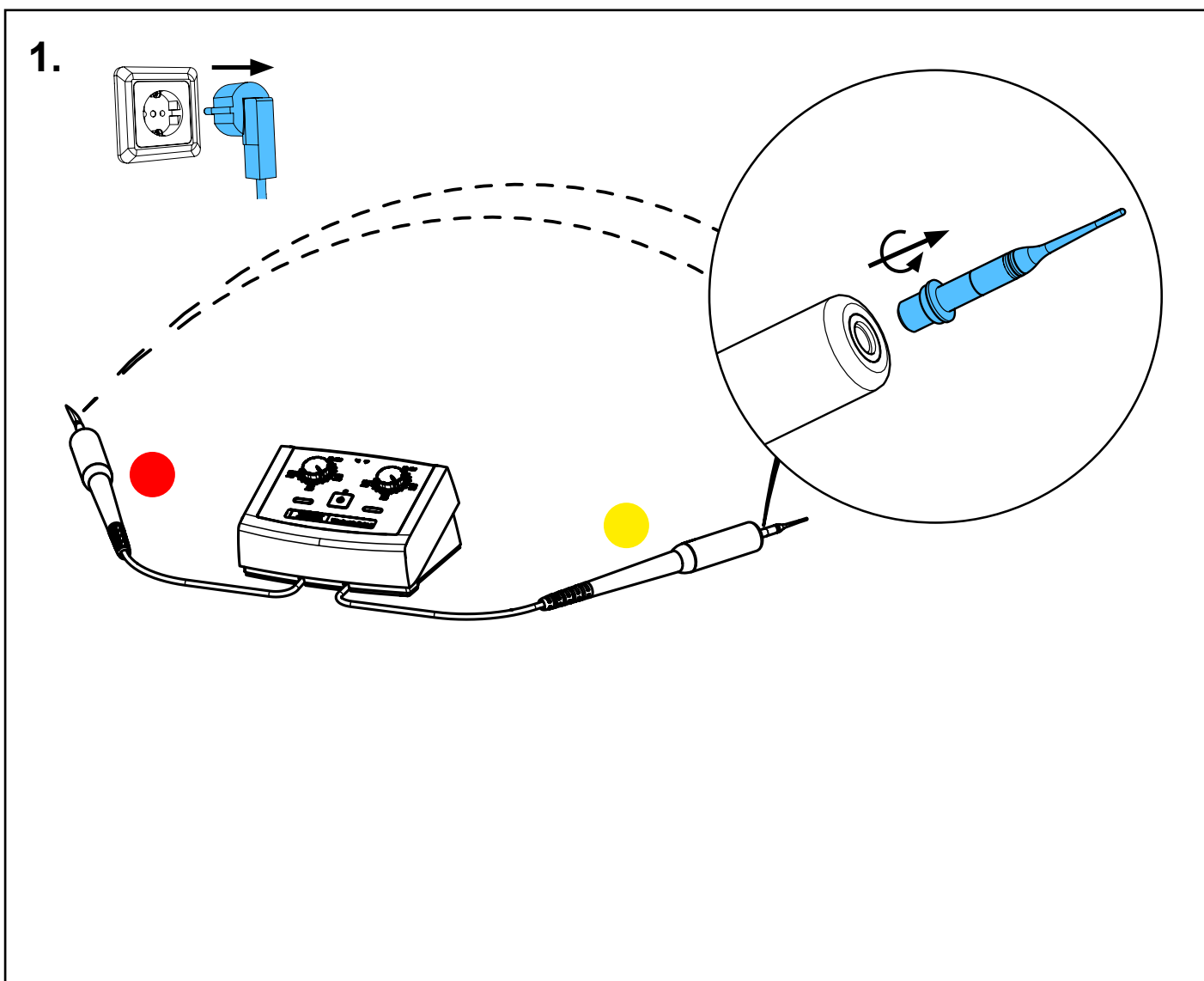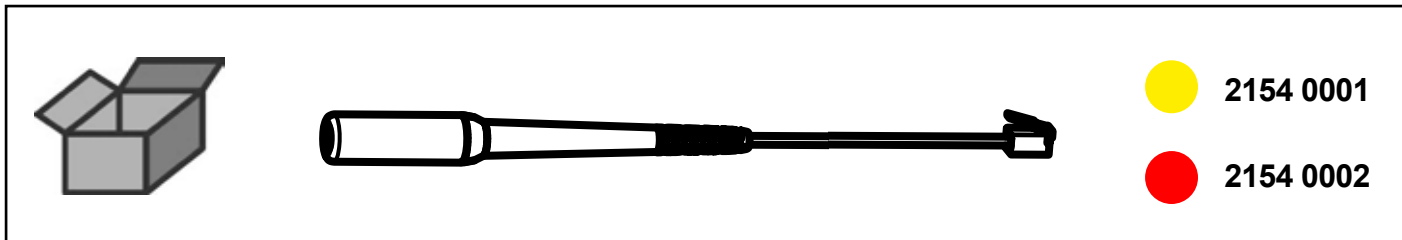


2.

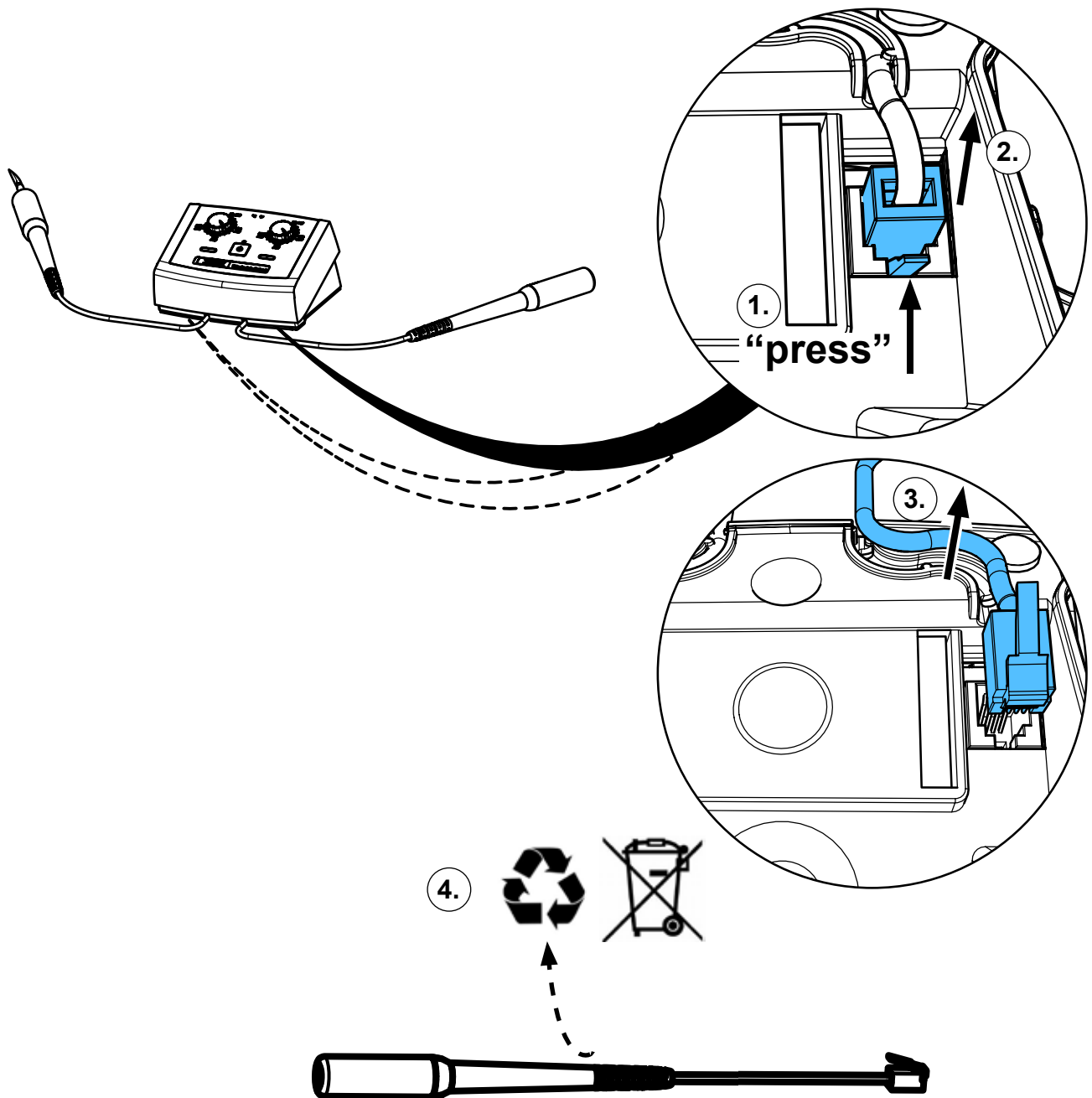


3.

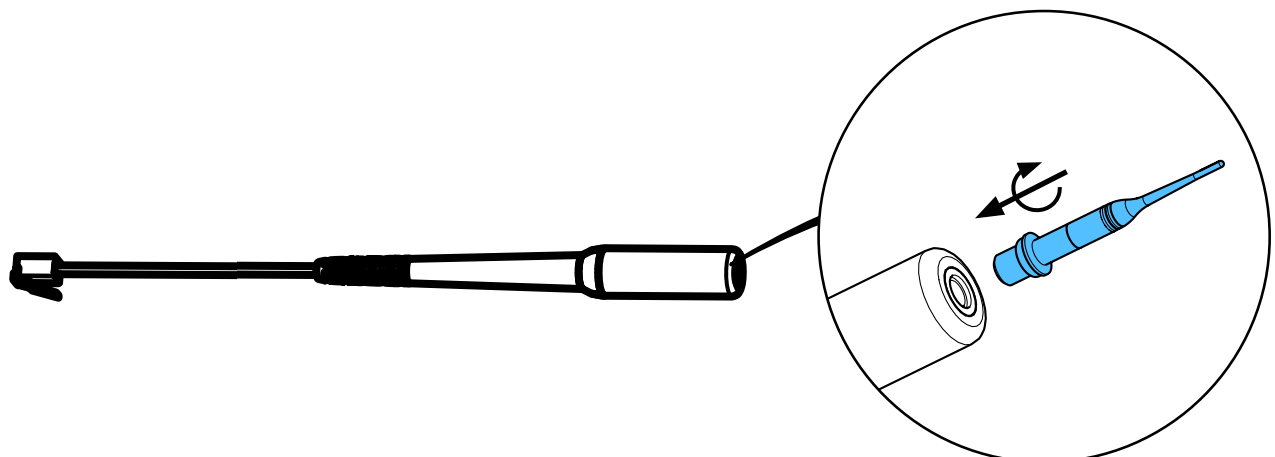

4.

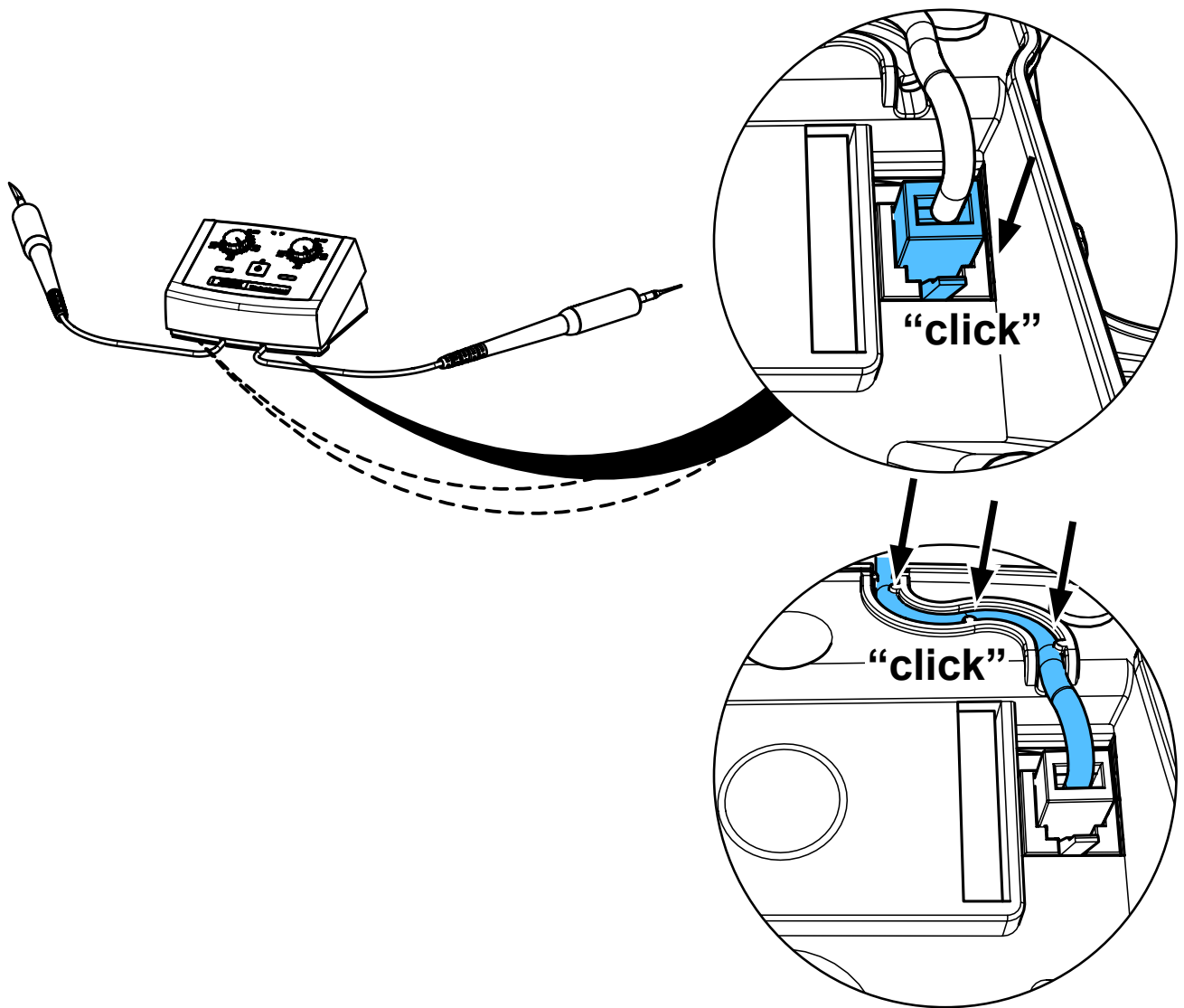

5.

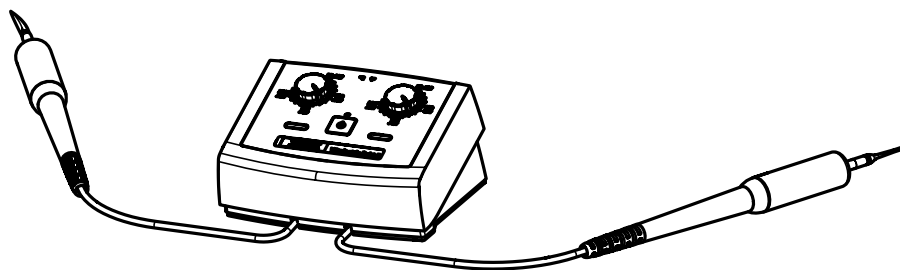




making work easy



Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу

90003 5948


Griffstück Kork • Cork-grip • Manche en liège • Impugnatura di sughero •
Empuñadura de corcho • Зажим для пробки

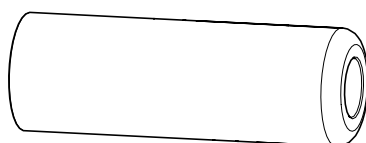


1.

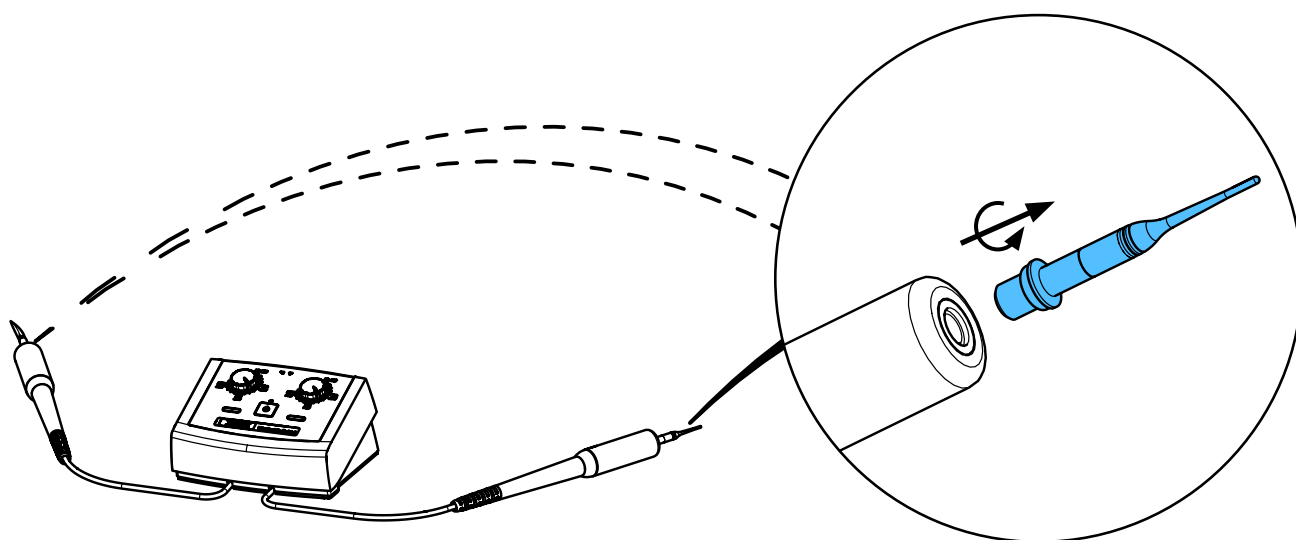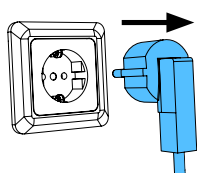


2.

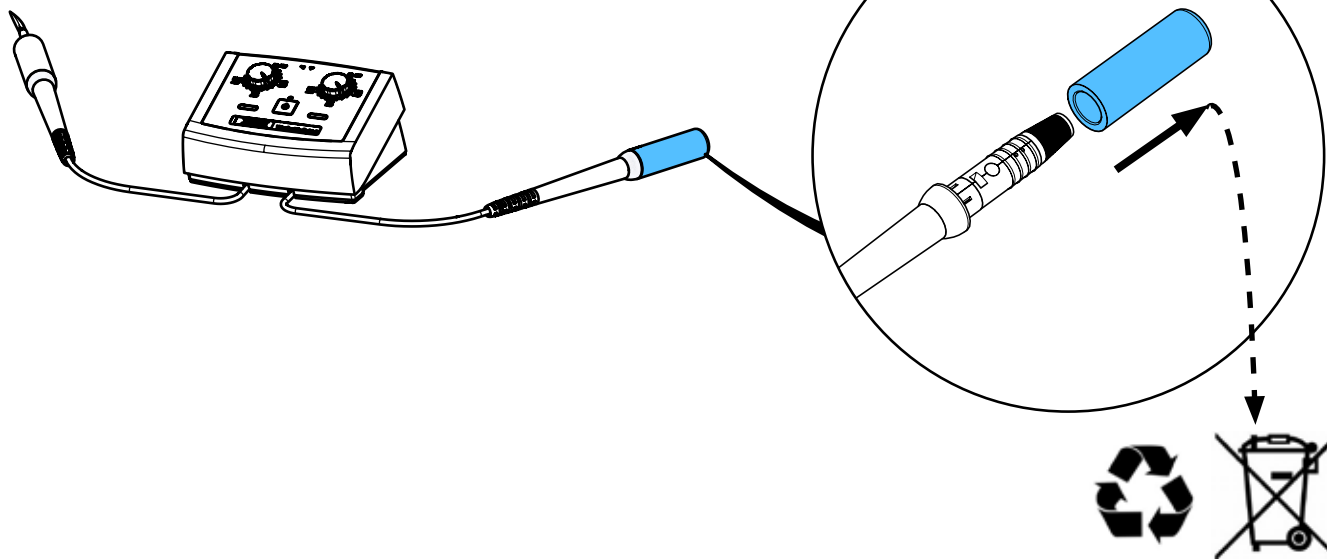

3.

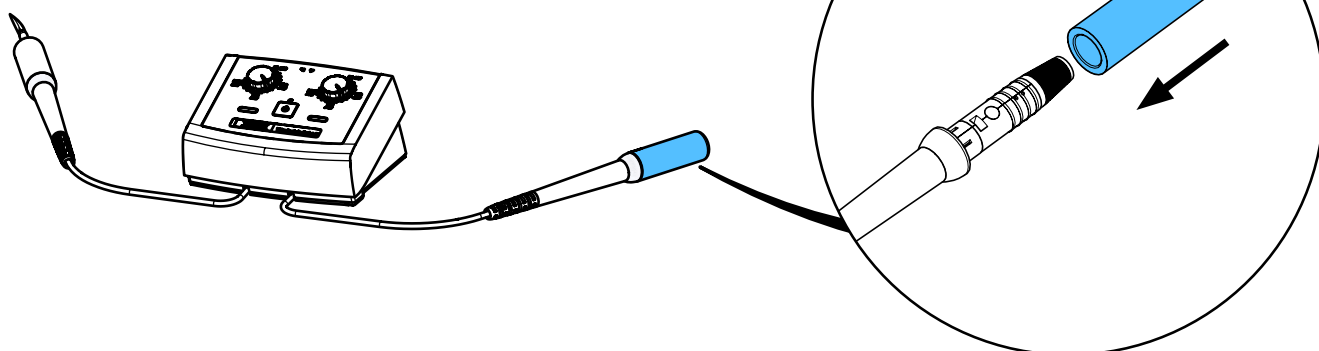

4.

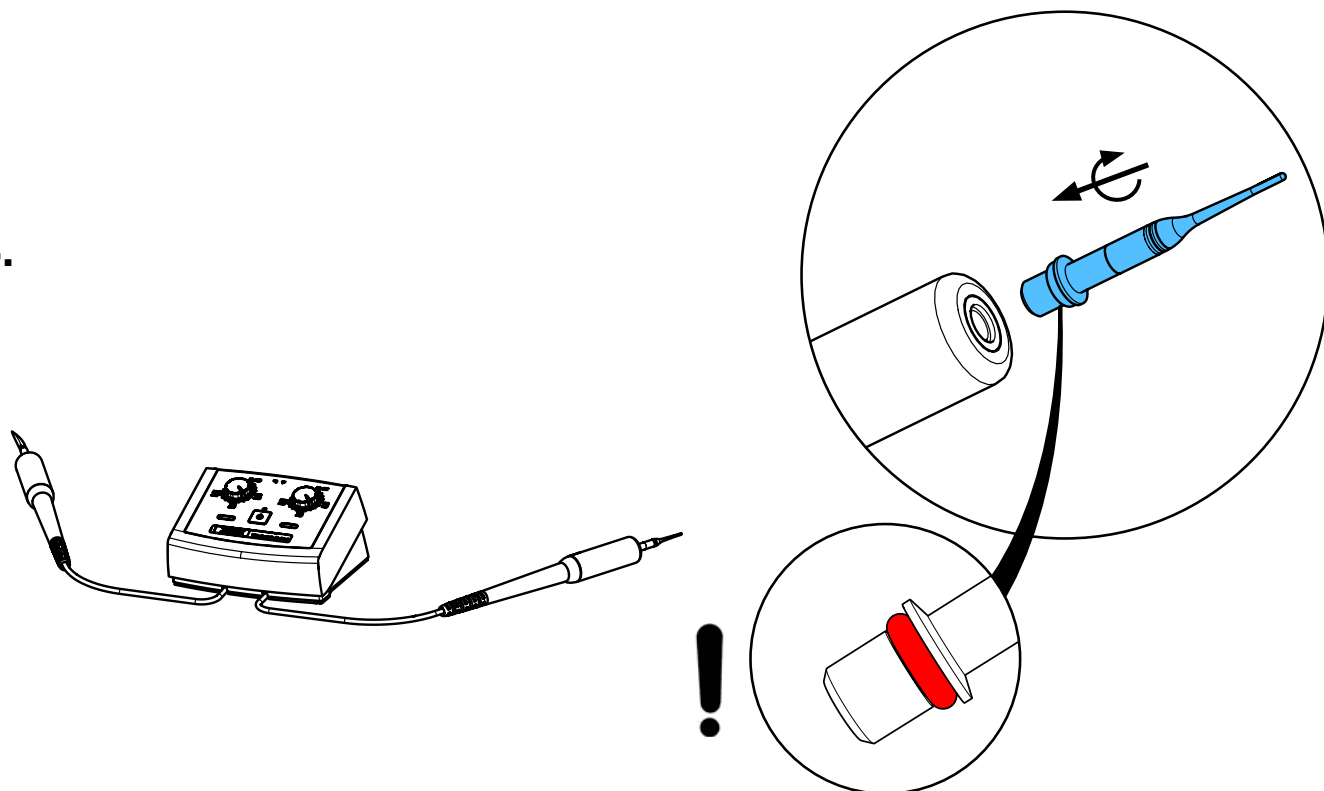





making work easy


Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу

90003 6138 / 90003 3752


Netzteil • Power supply • Poste secteur • Trasformatore • Trasformatore •
Блок питания

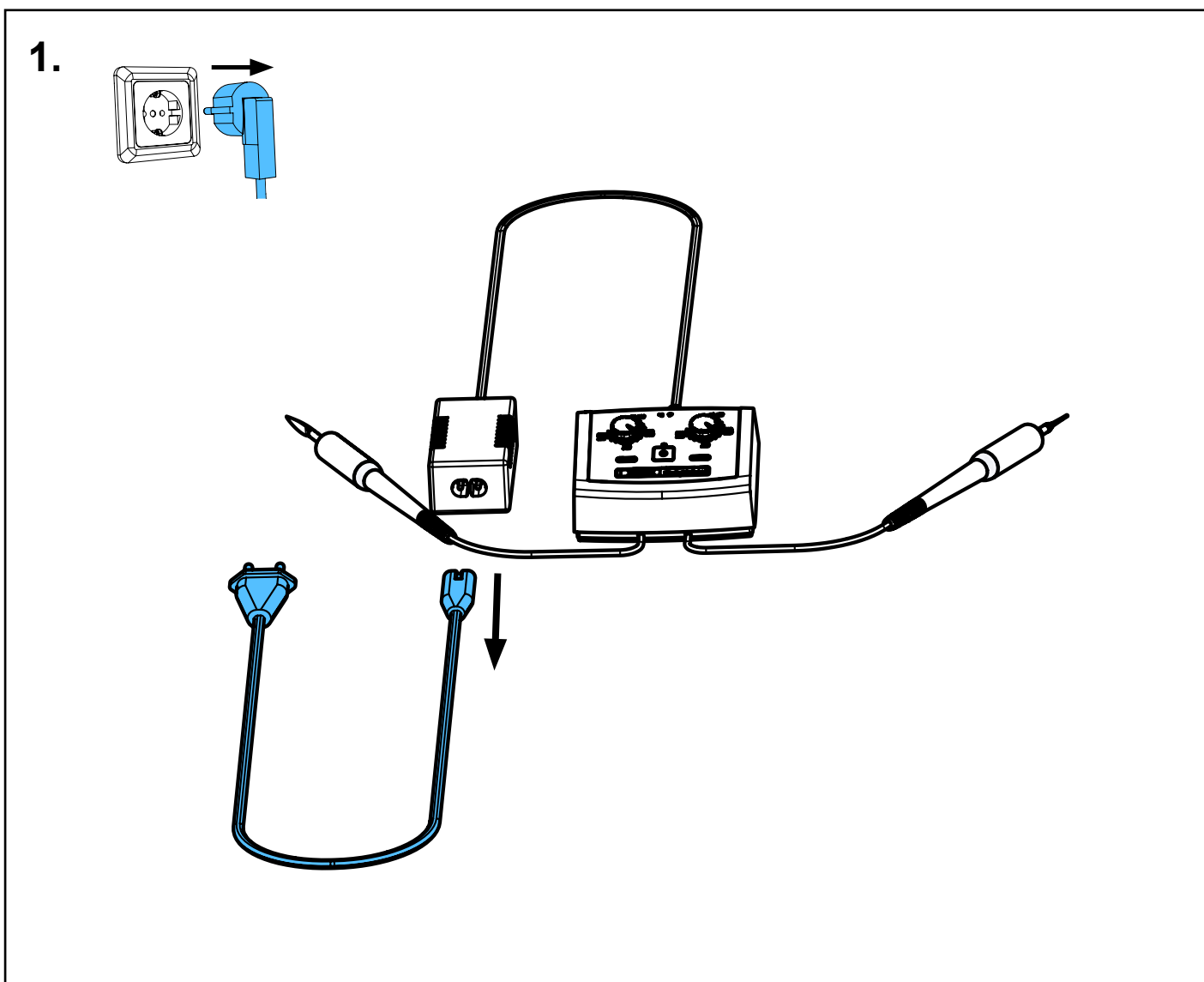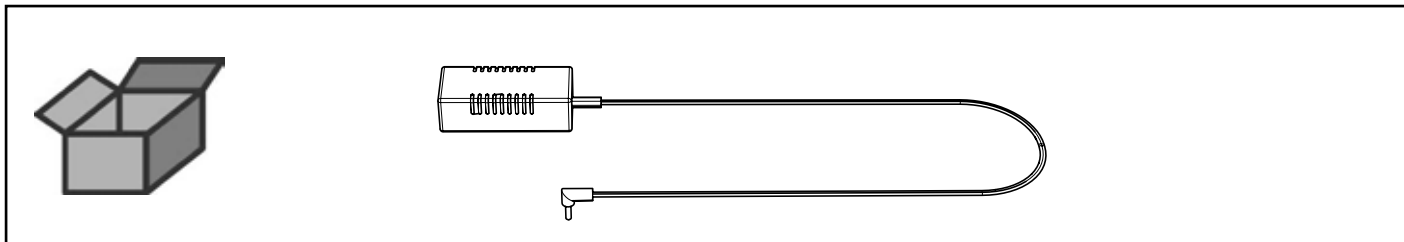


2.

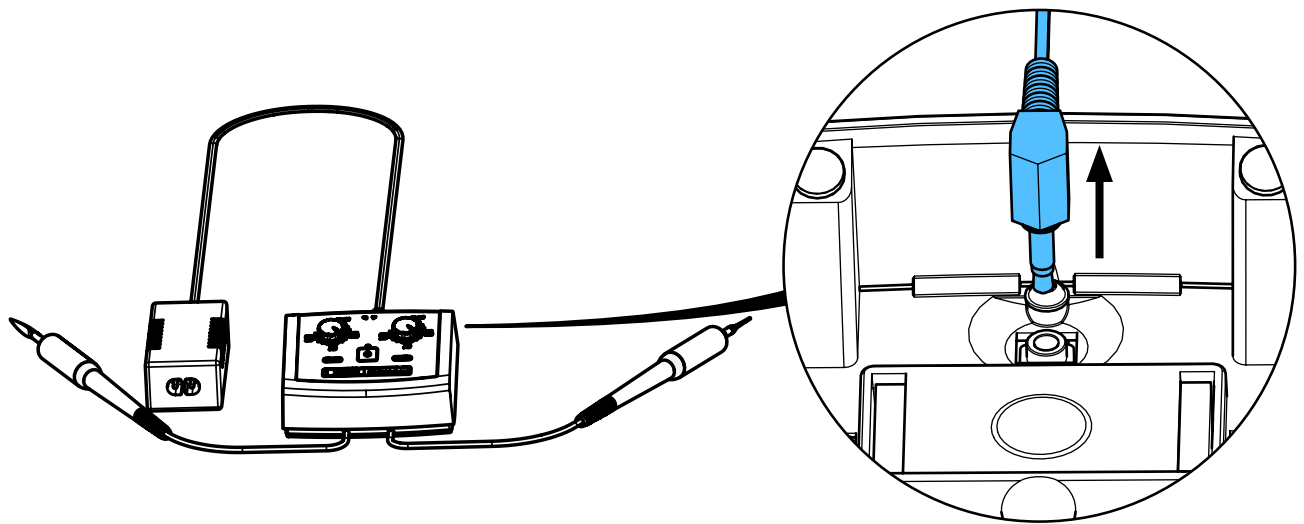


3.

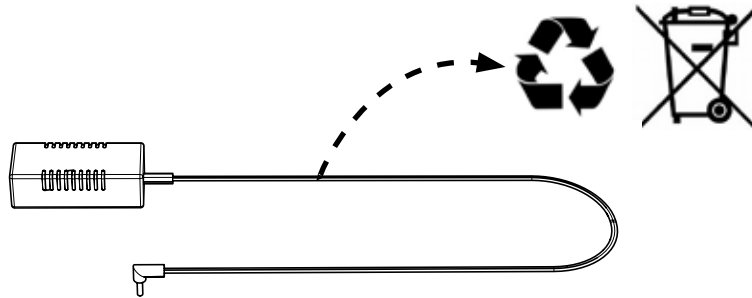

4.

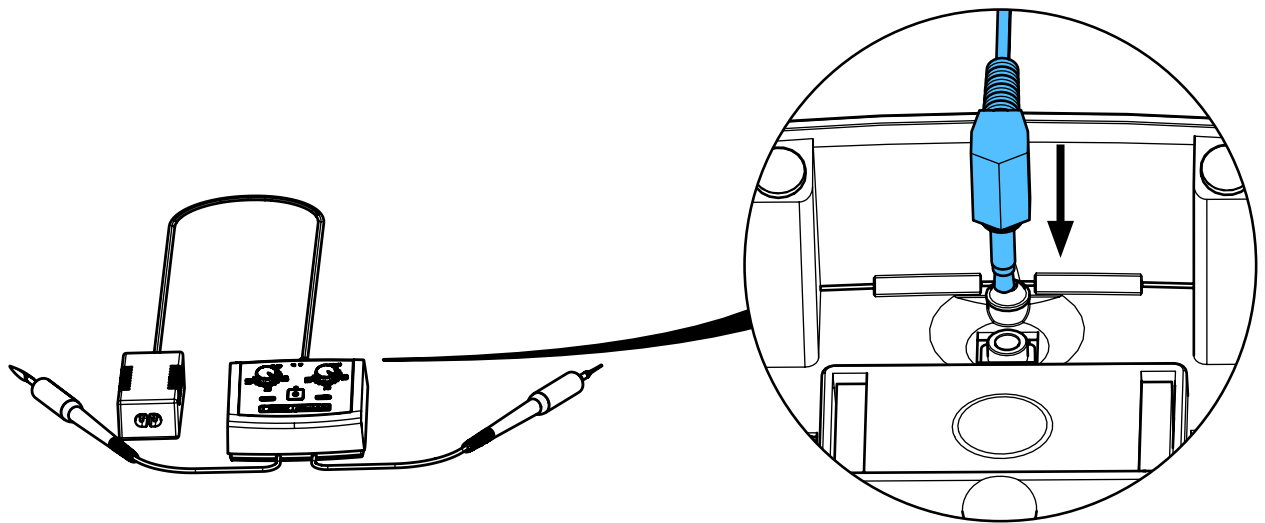

5.

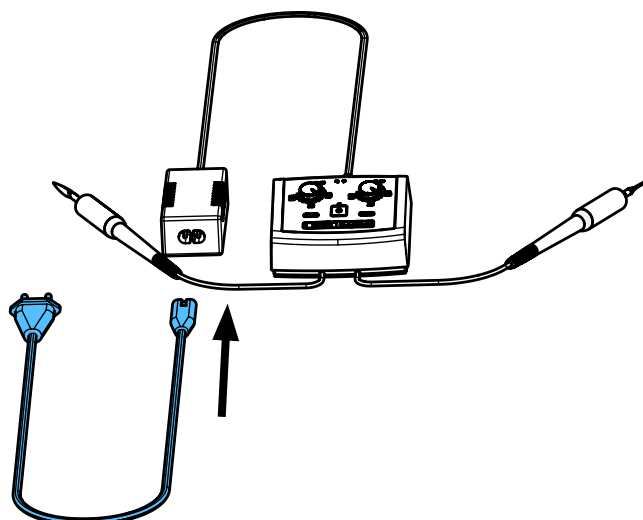



making work easy


Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу

90003 2962



O-Ring • Retainer ring • Joint thorique • Guarnizioni • Anillo en O •
Блок питания




5 x

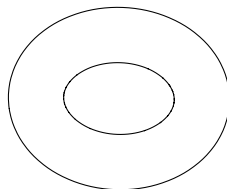


1.

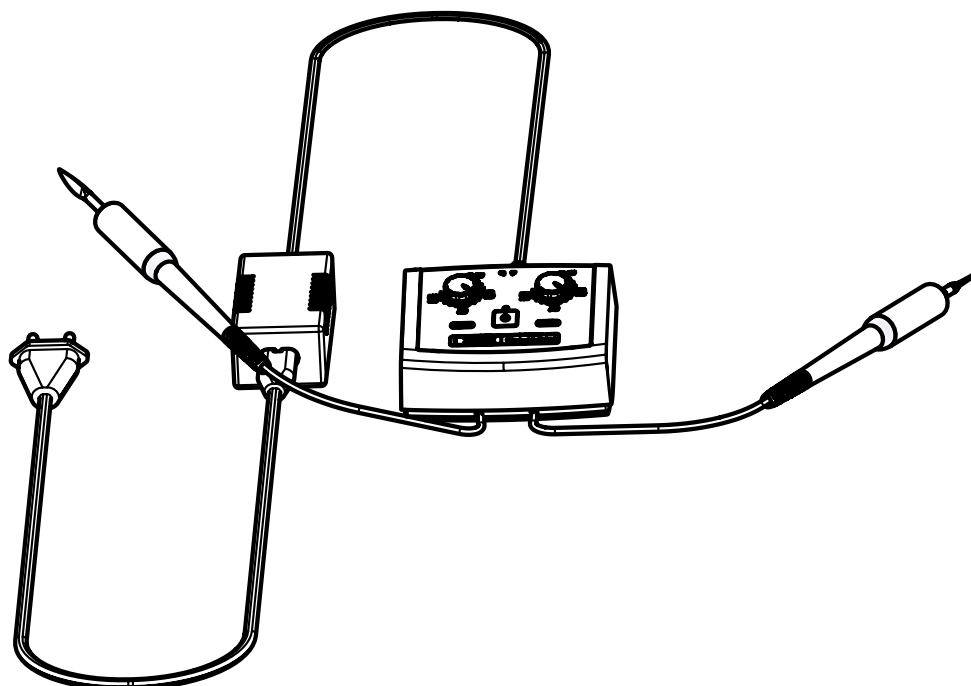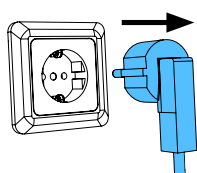


2.

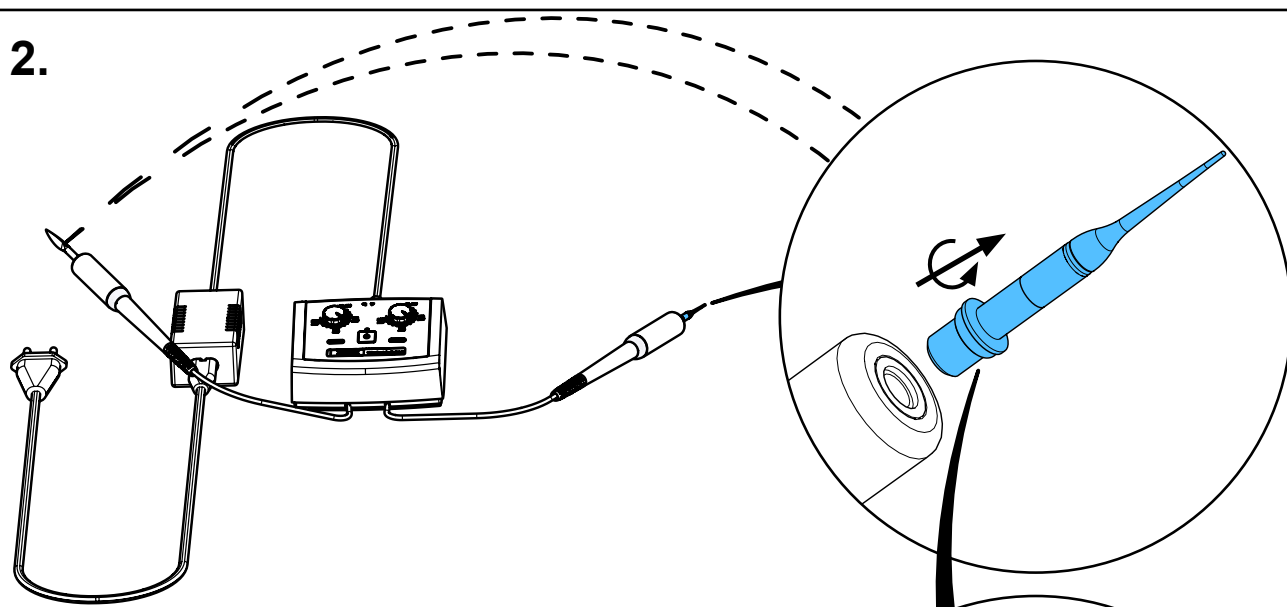


3.

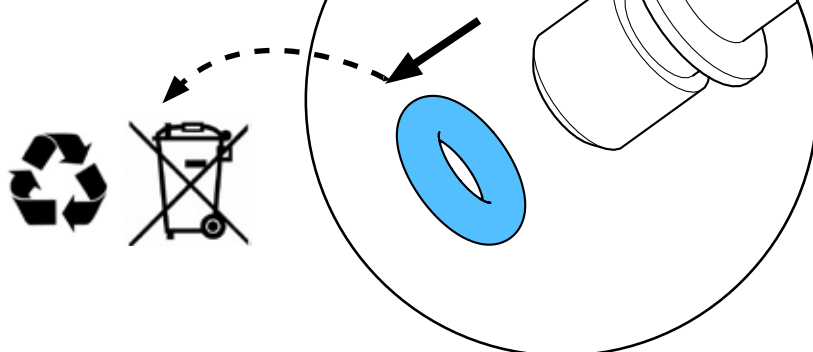

4.

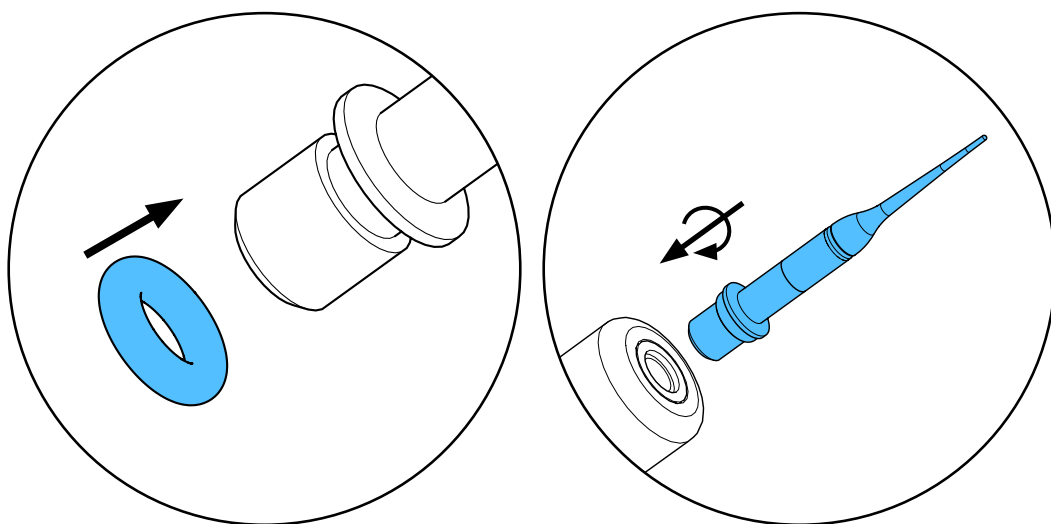

5.

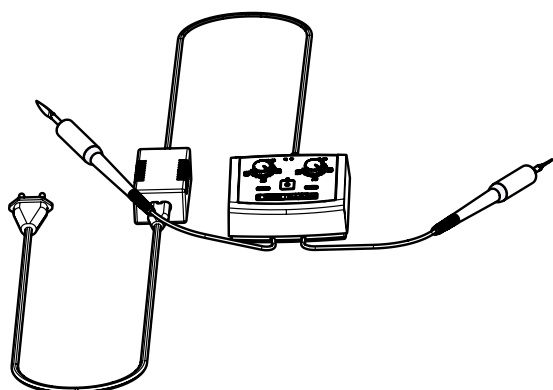




making work easy



Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу

90003 4564


Gehäusefüsse • Rubber bases • Pieds en caoutchouc • Piedini di gomma •
Pies de goma • Ножки прибора

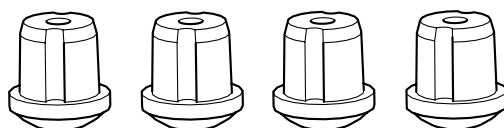


1.

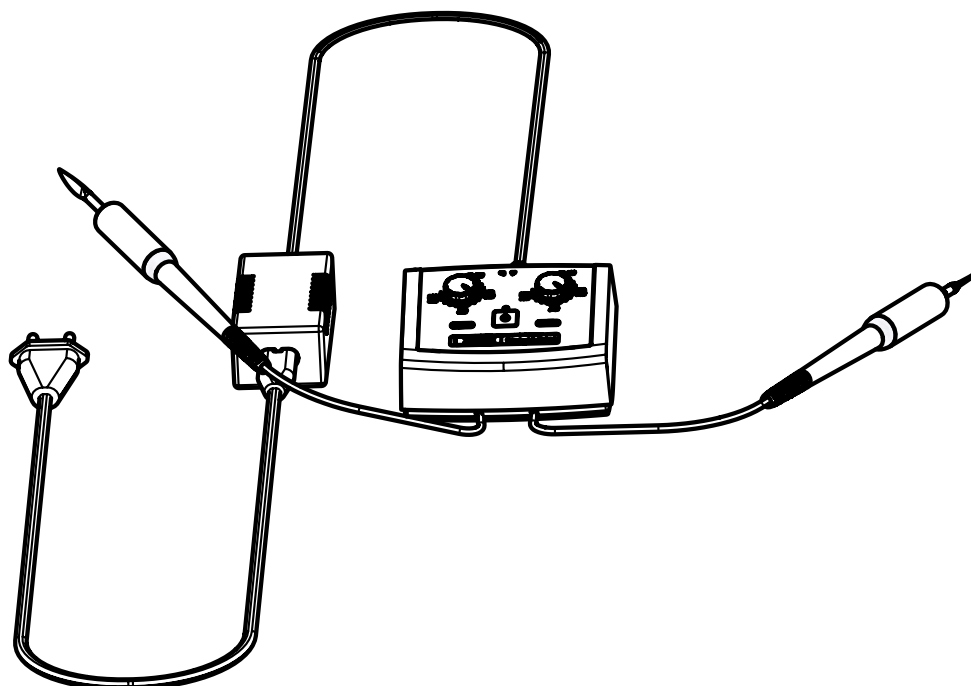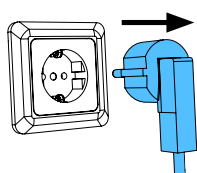


2.

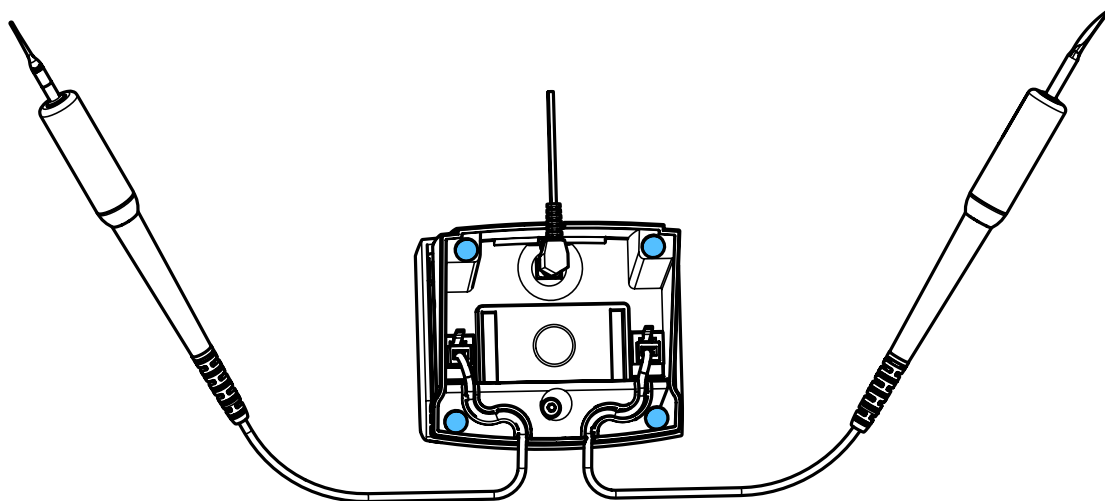


3.

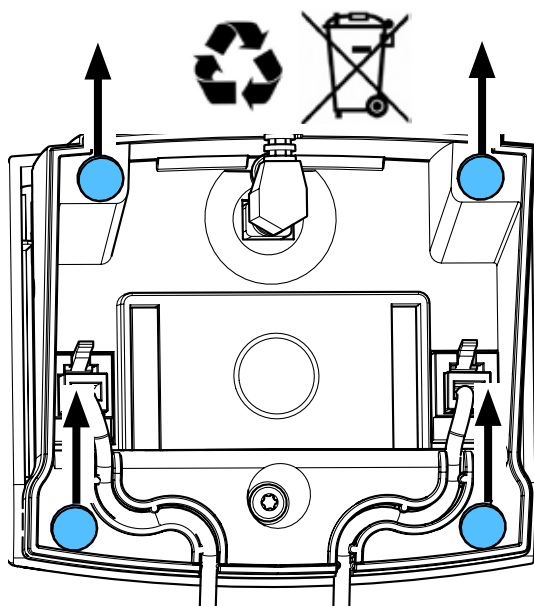

4.

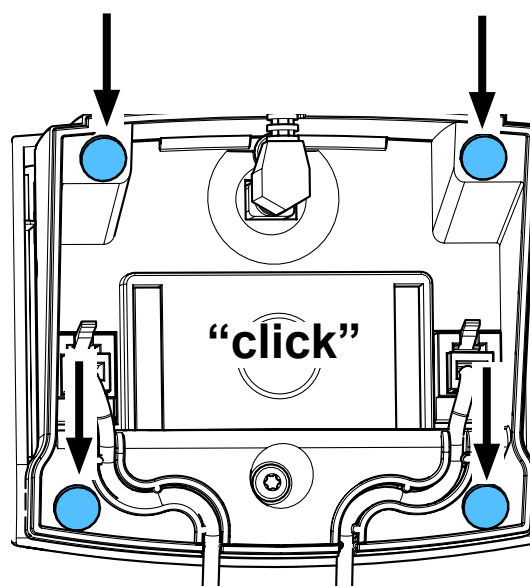

5.

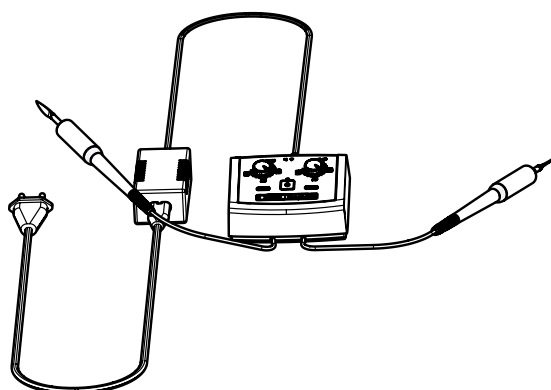





making work easy


Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу

90003 5975/..5976/..8652/..8653/..8654/..8680/..2773


Netzkabel • Power cord • Cordon d'alimentation • Cavo di rete • Cable de red
Сетевой кабель

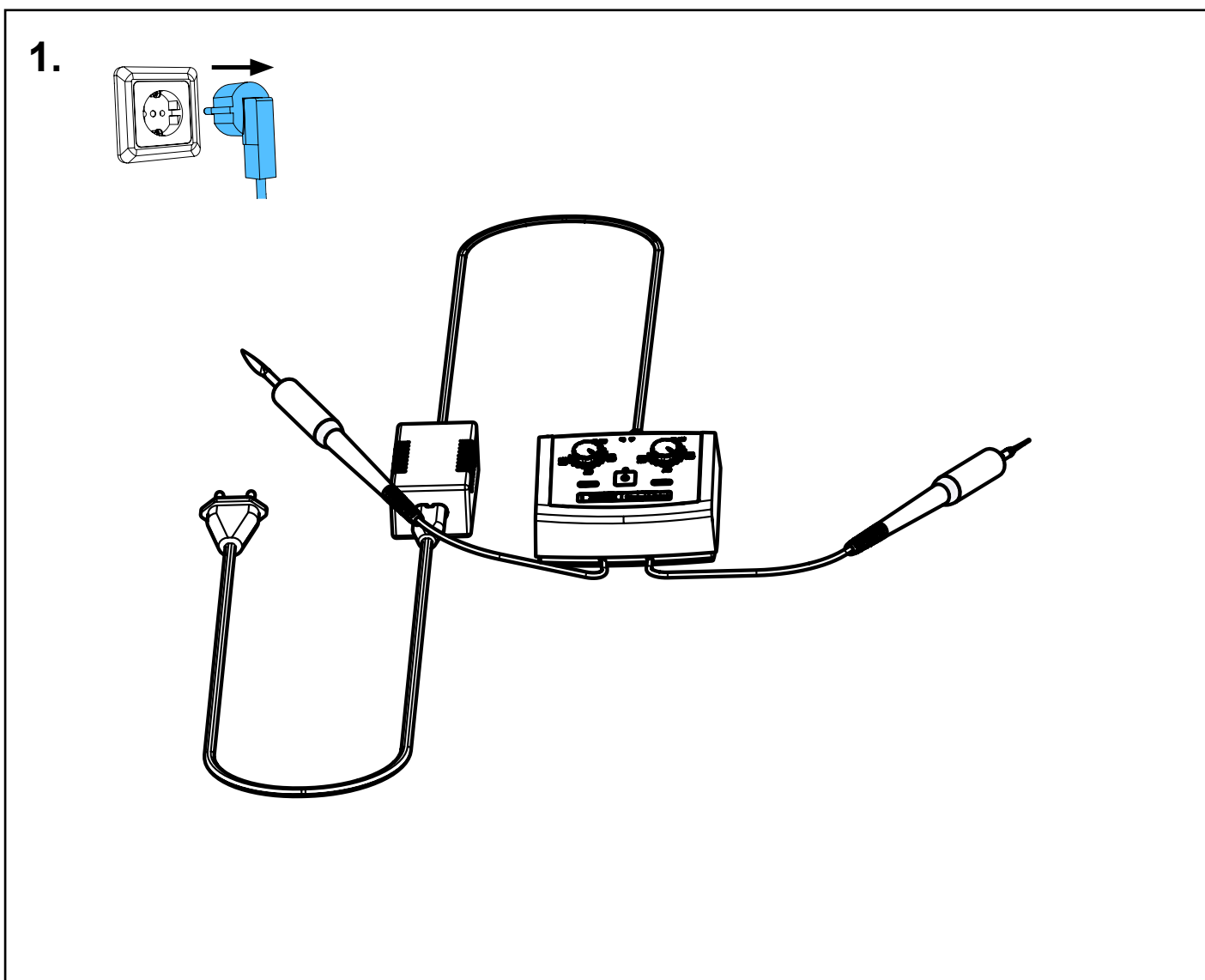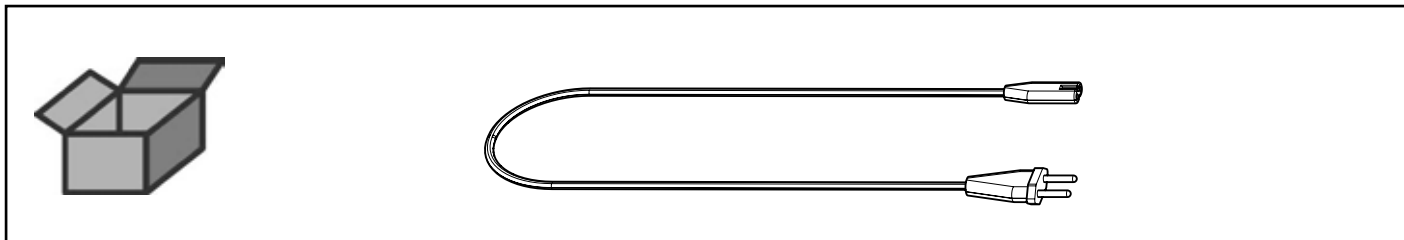


2.

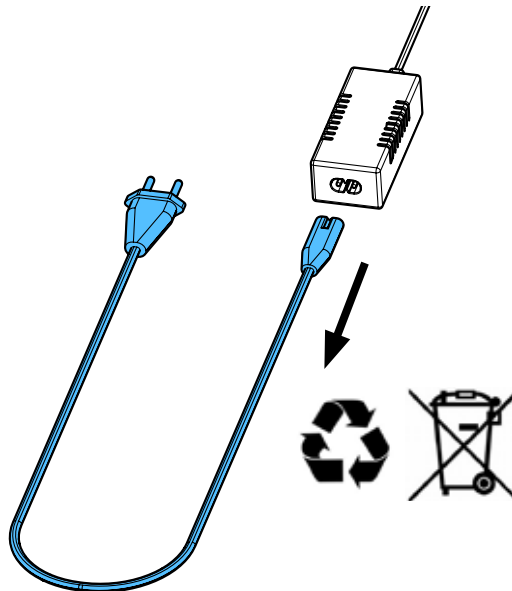

3.

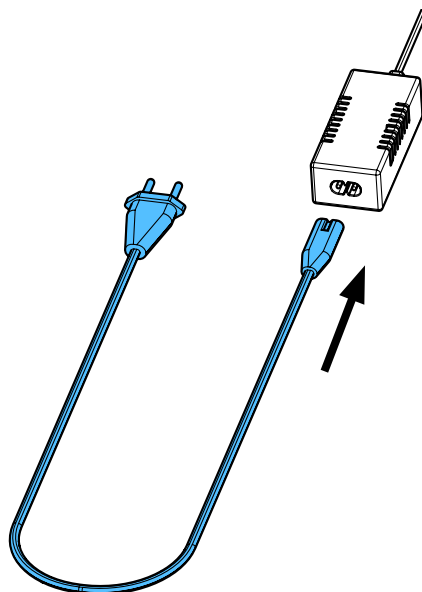

4.

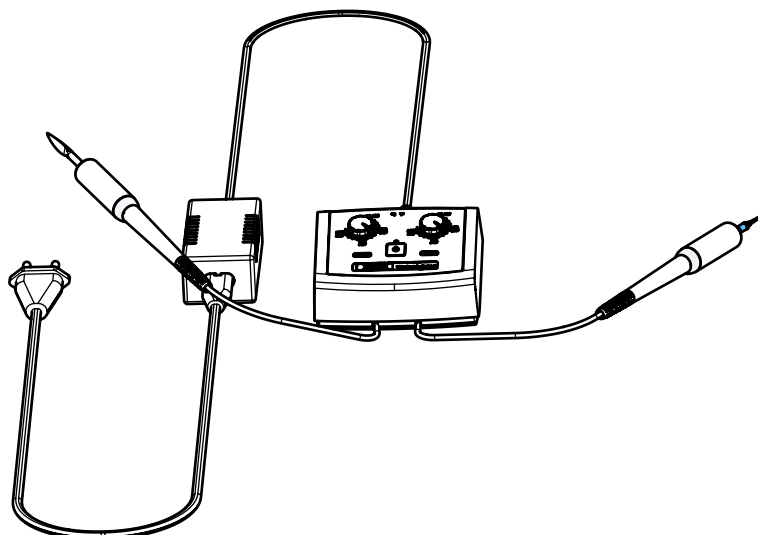






making work easy




Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу

90003 5971




Leiterplatte Waxelectric light II • Mains circuit board Waxelectric light II • Circuit imprimé Waxelectric light II • Circuito stampato Waxelectric light II • Circuito impreso Waxelectric light II • Печатная плата Waxelectric light II

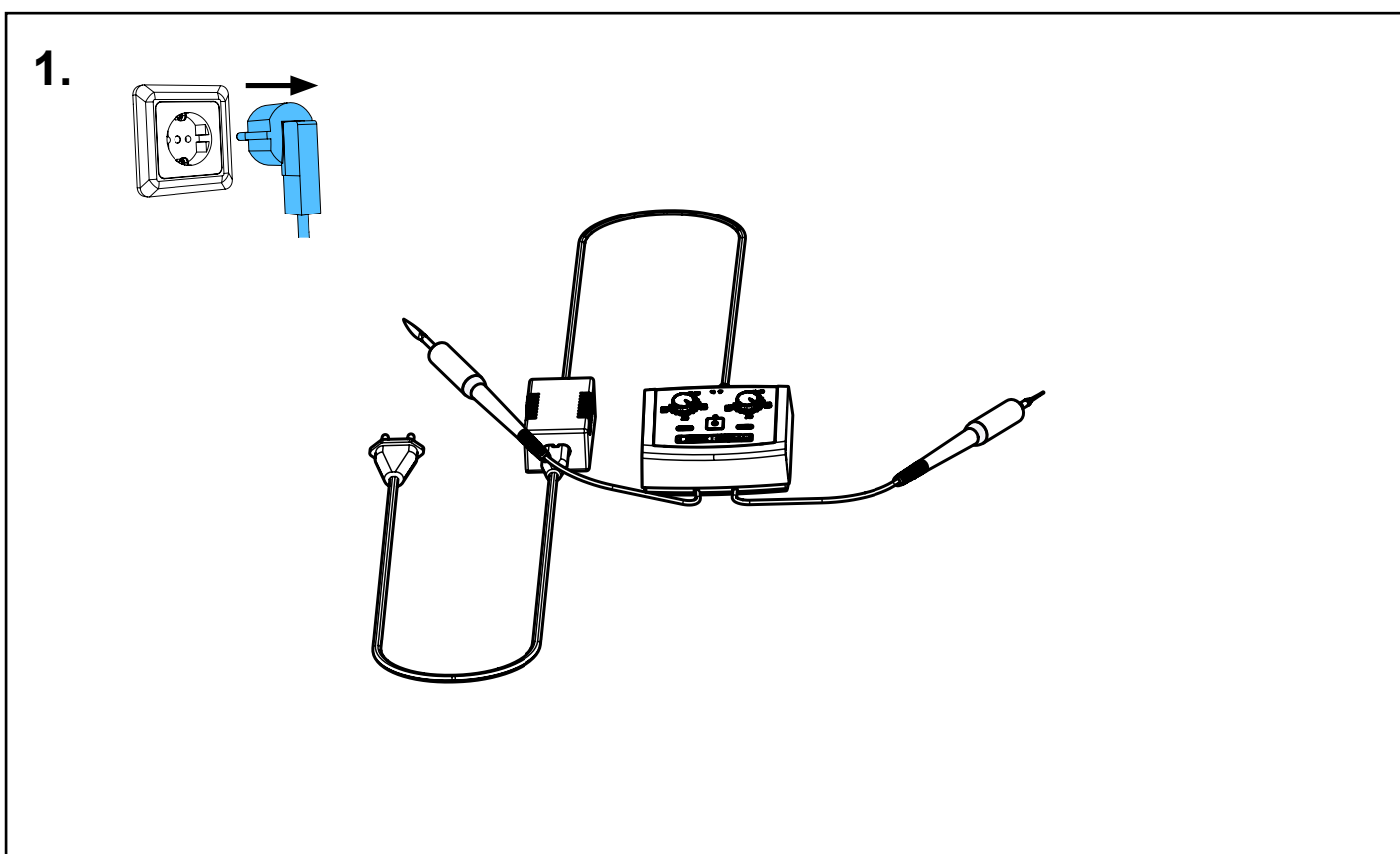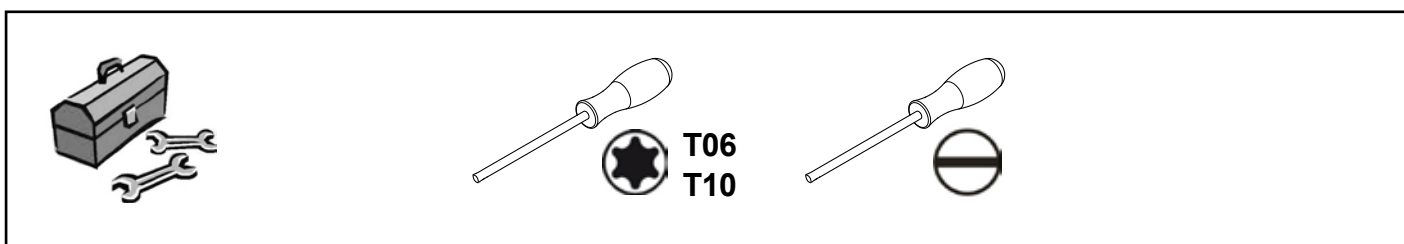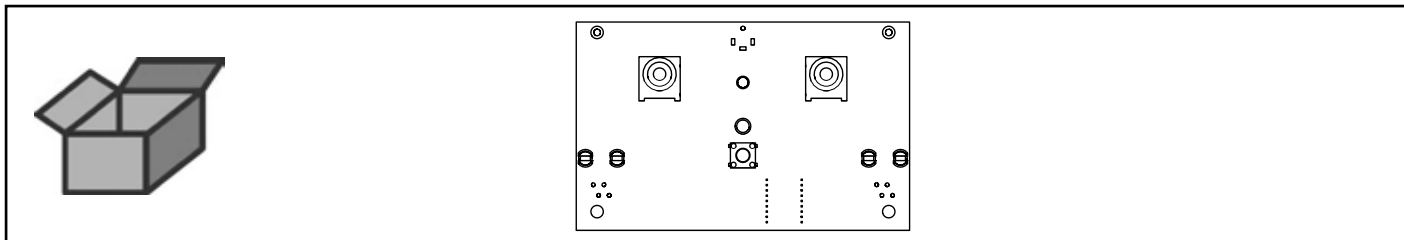


2.

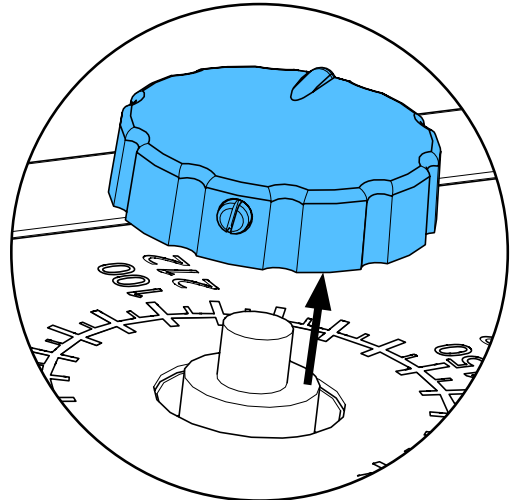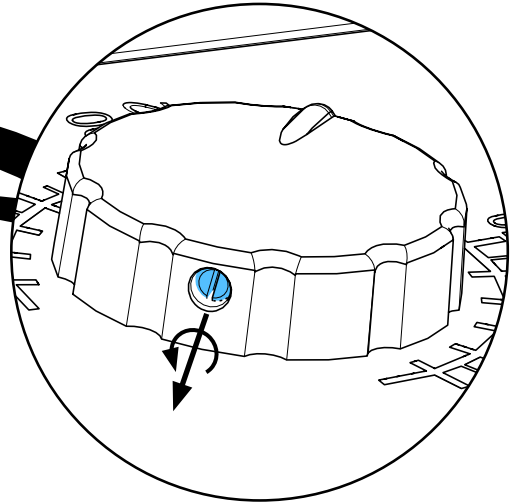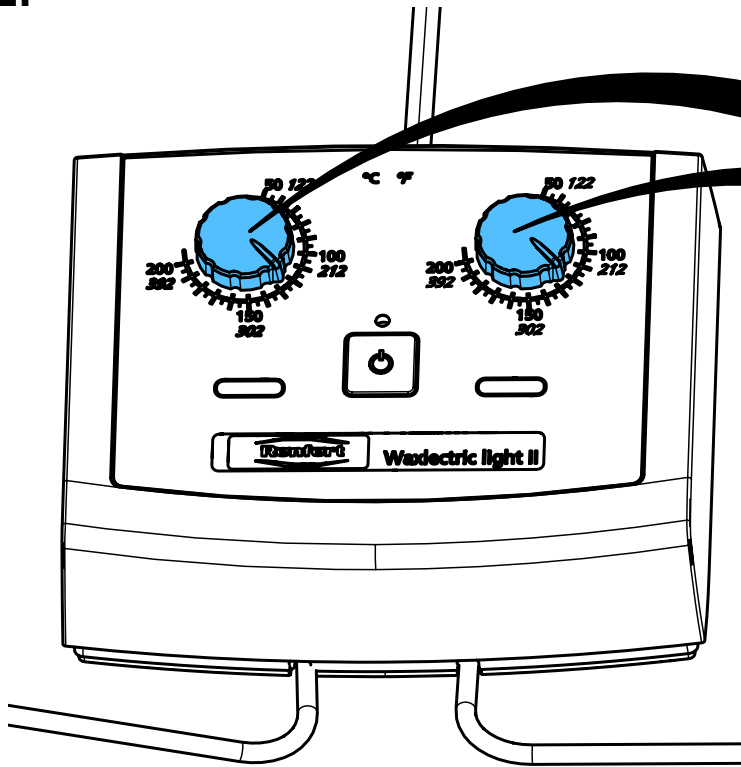


3.

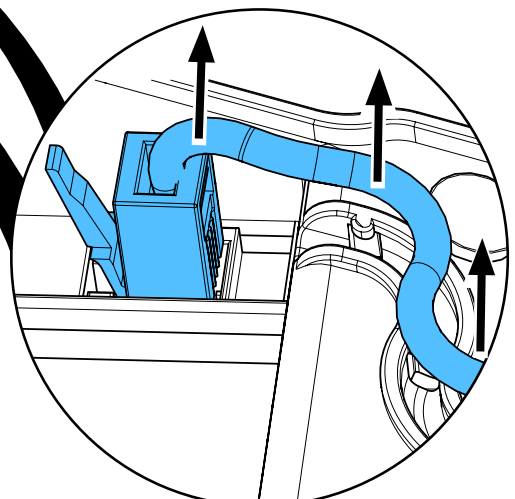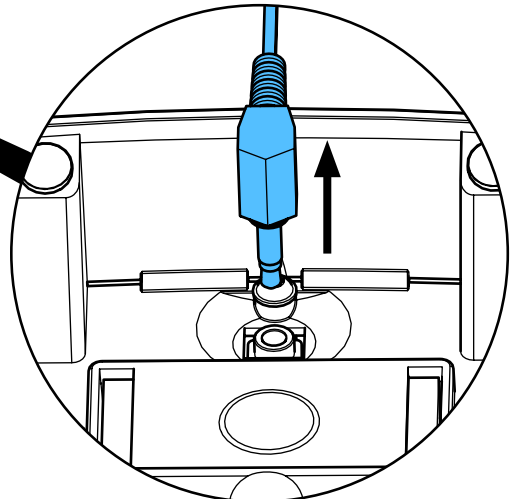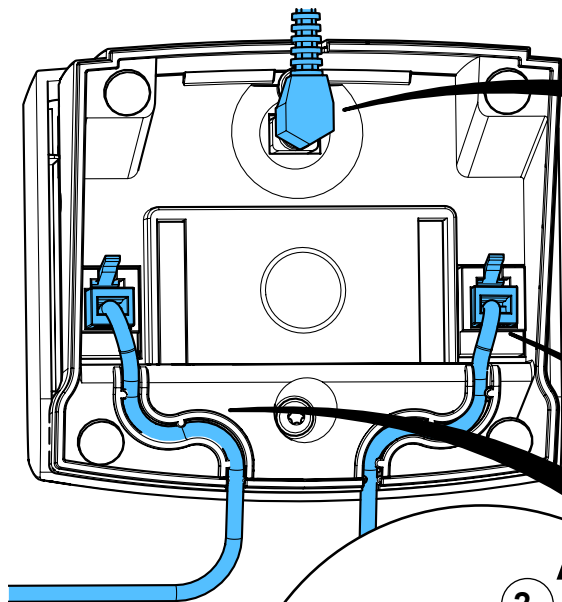


4.

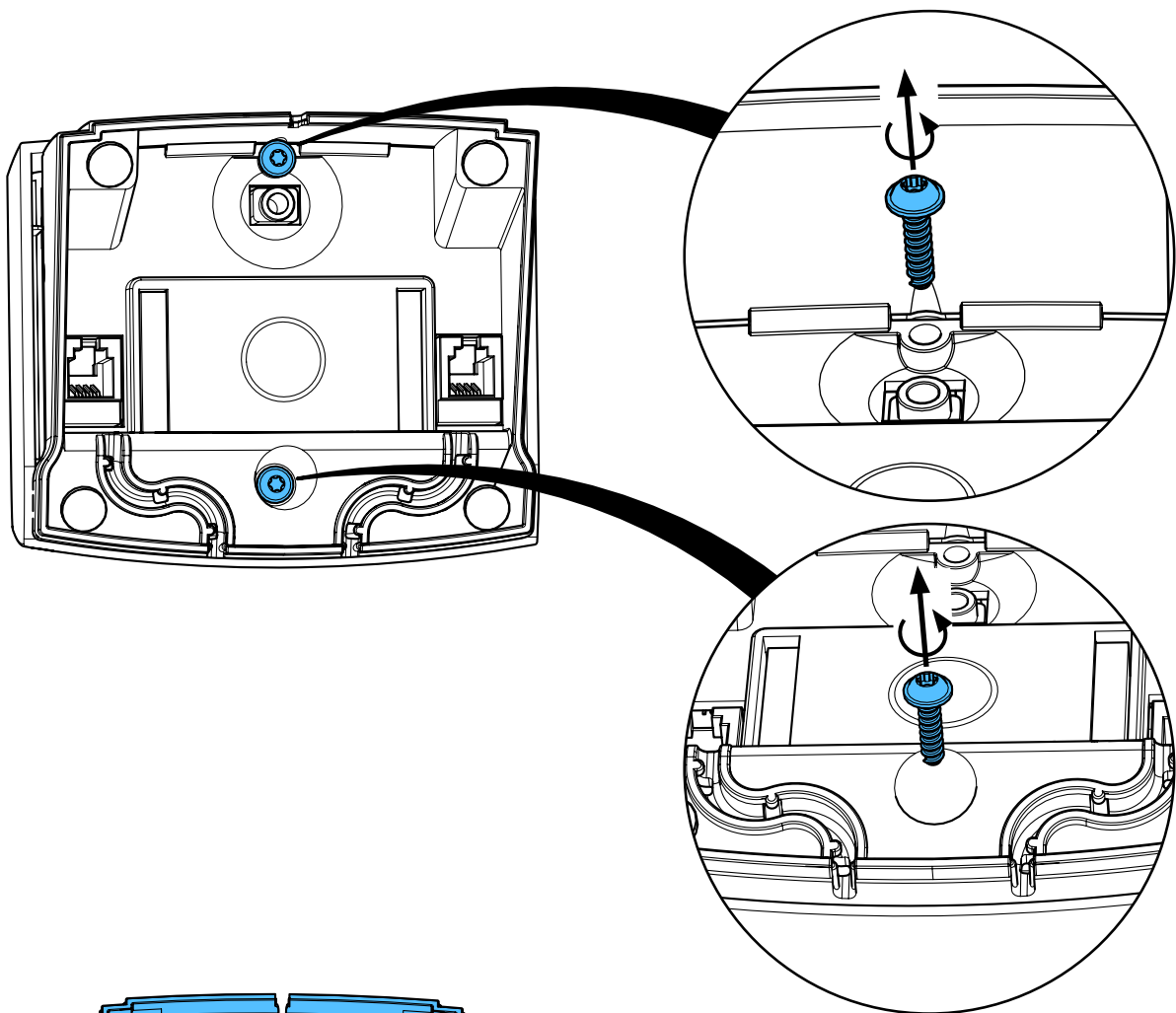


5.

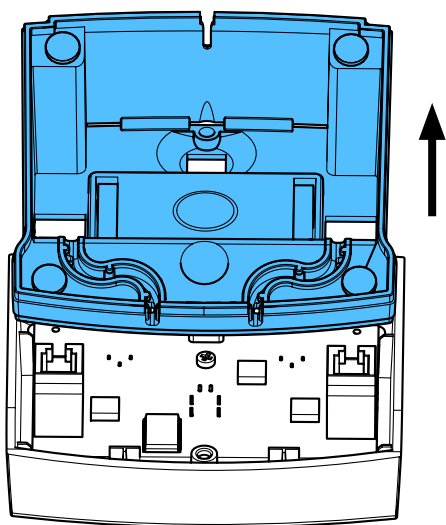


6.

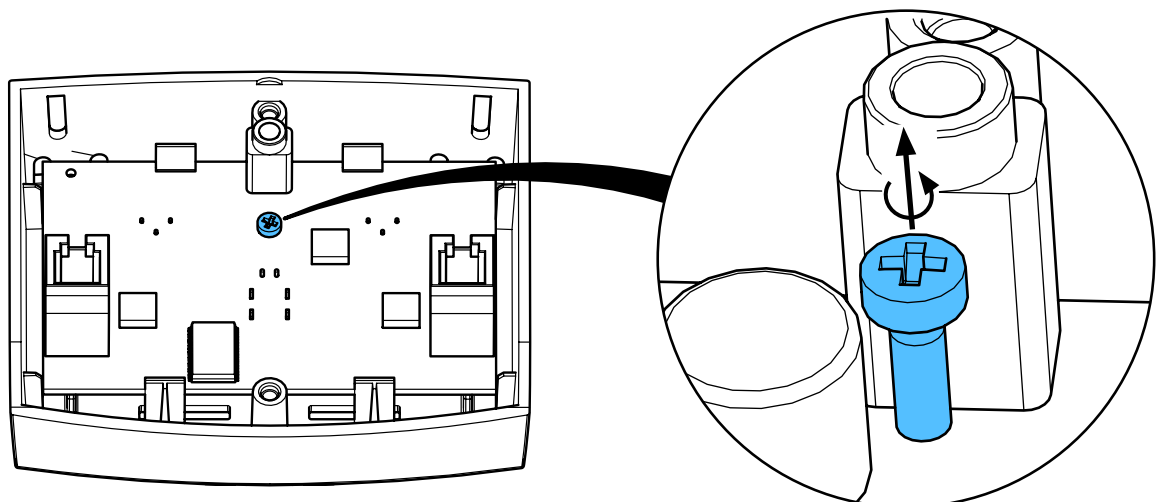


7.

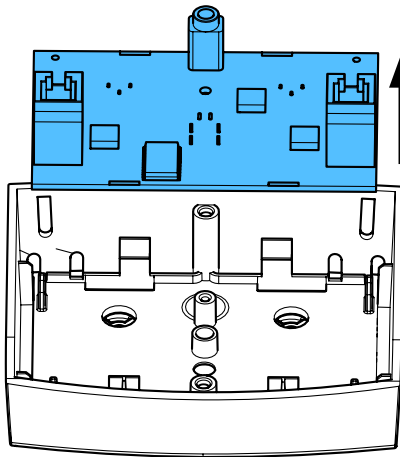


i

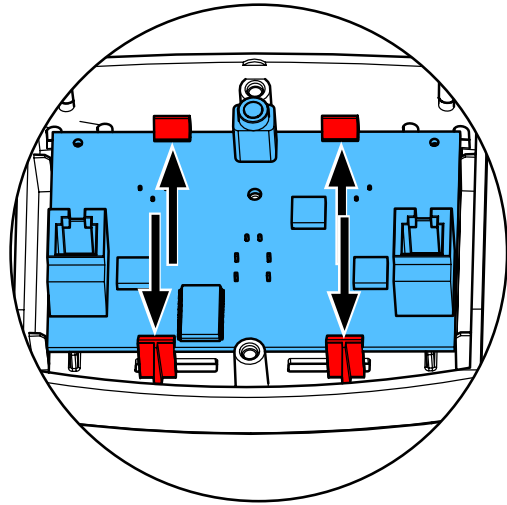



8.

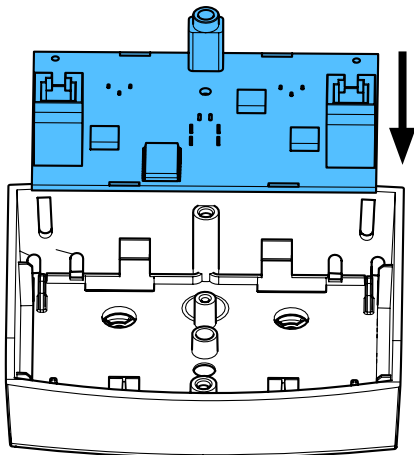


i

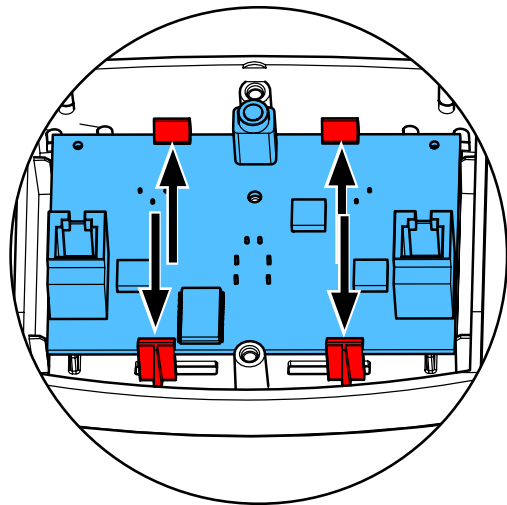


9.

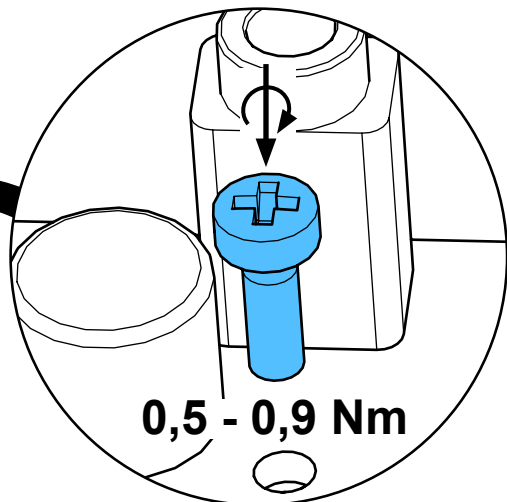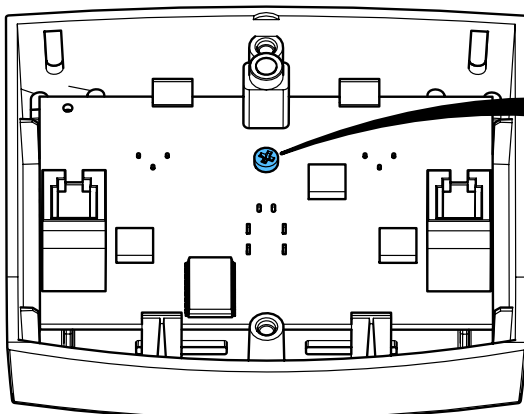


10.

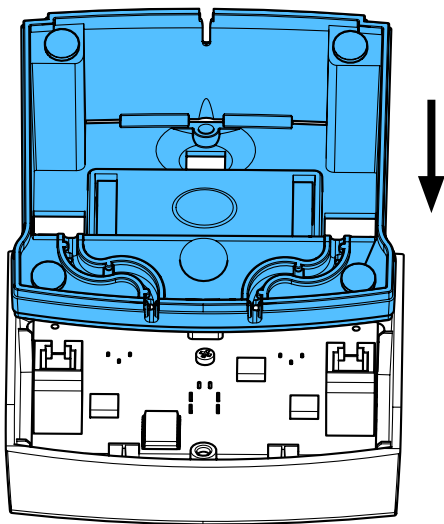


11.

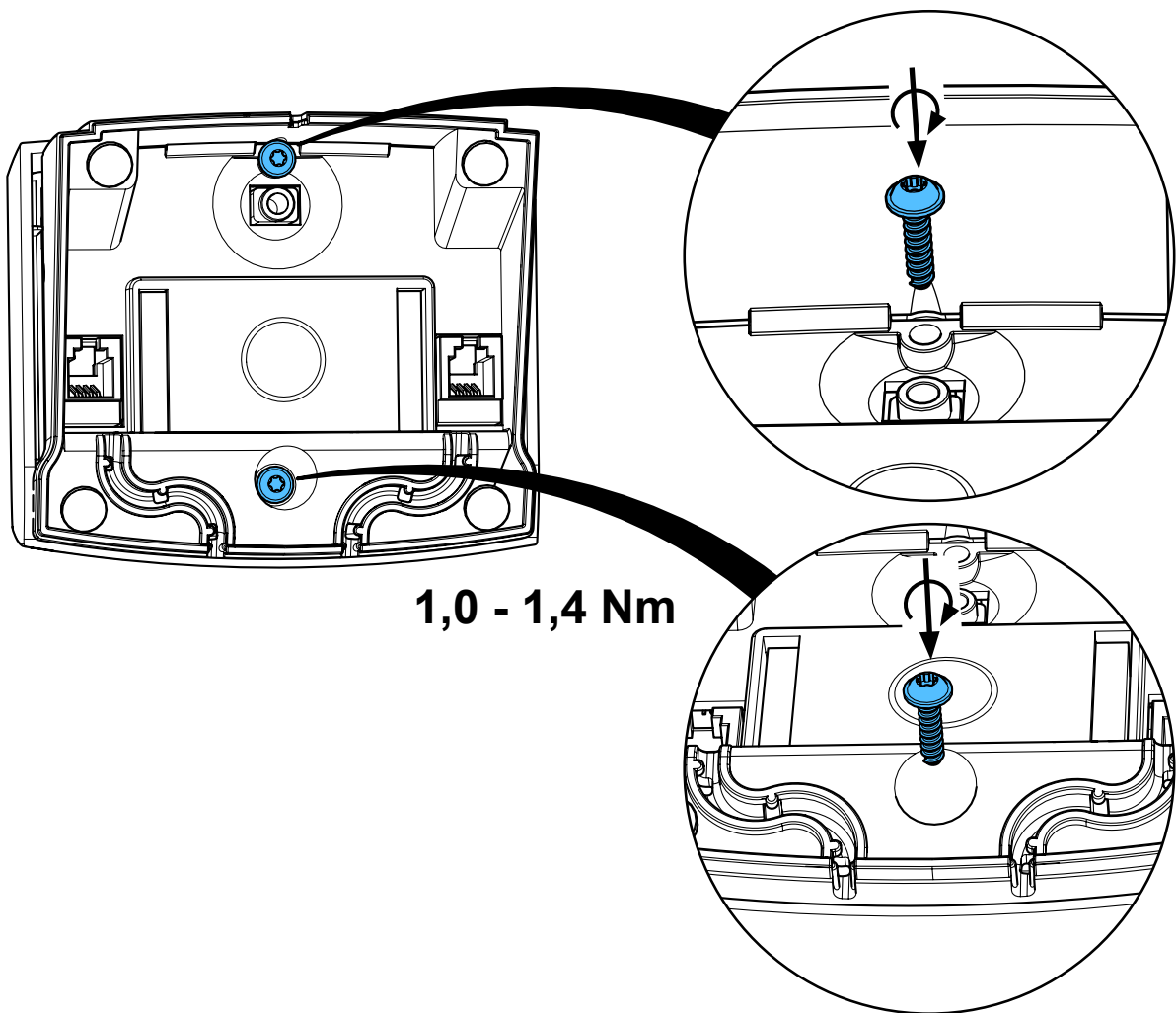

12.




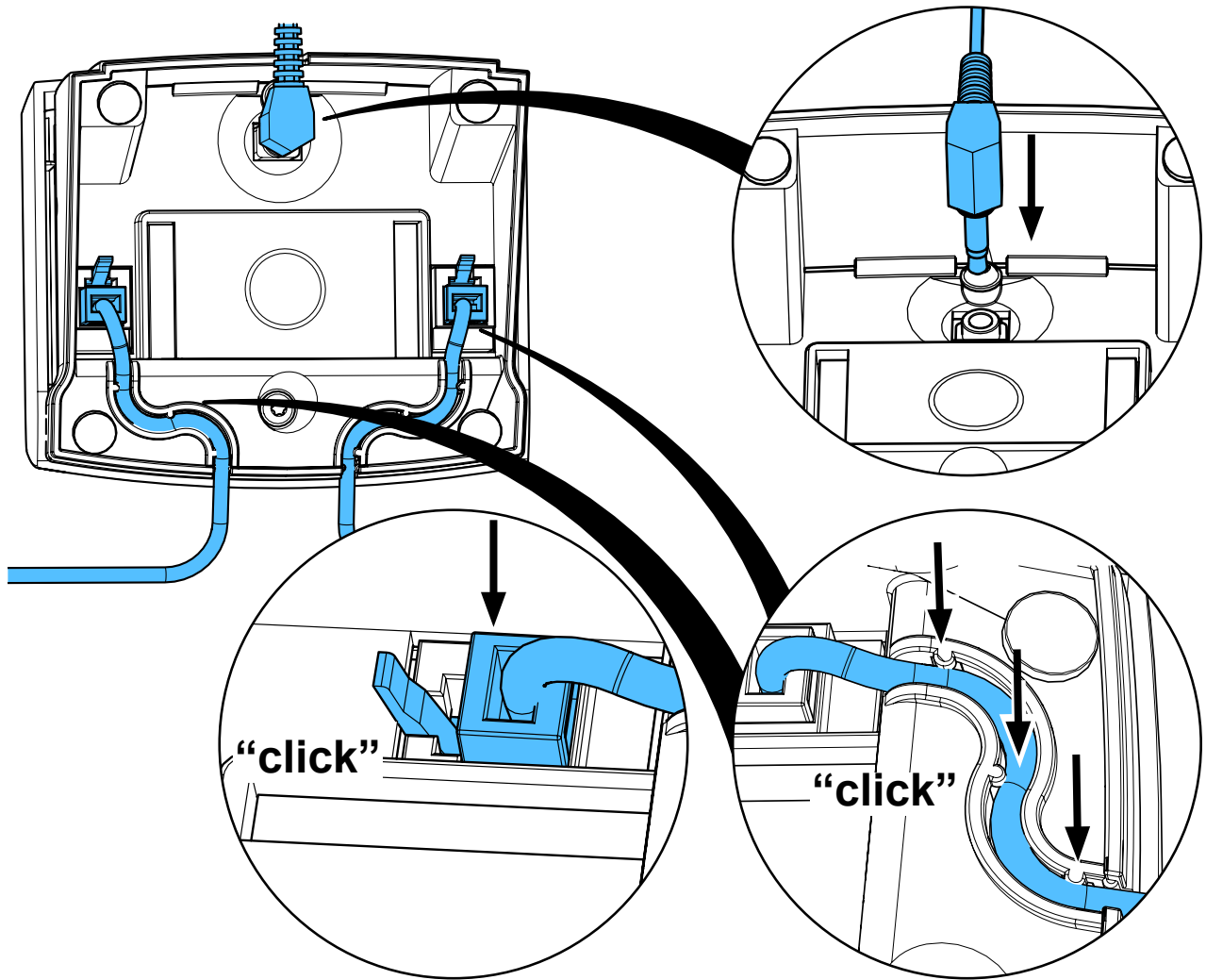


13.

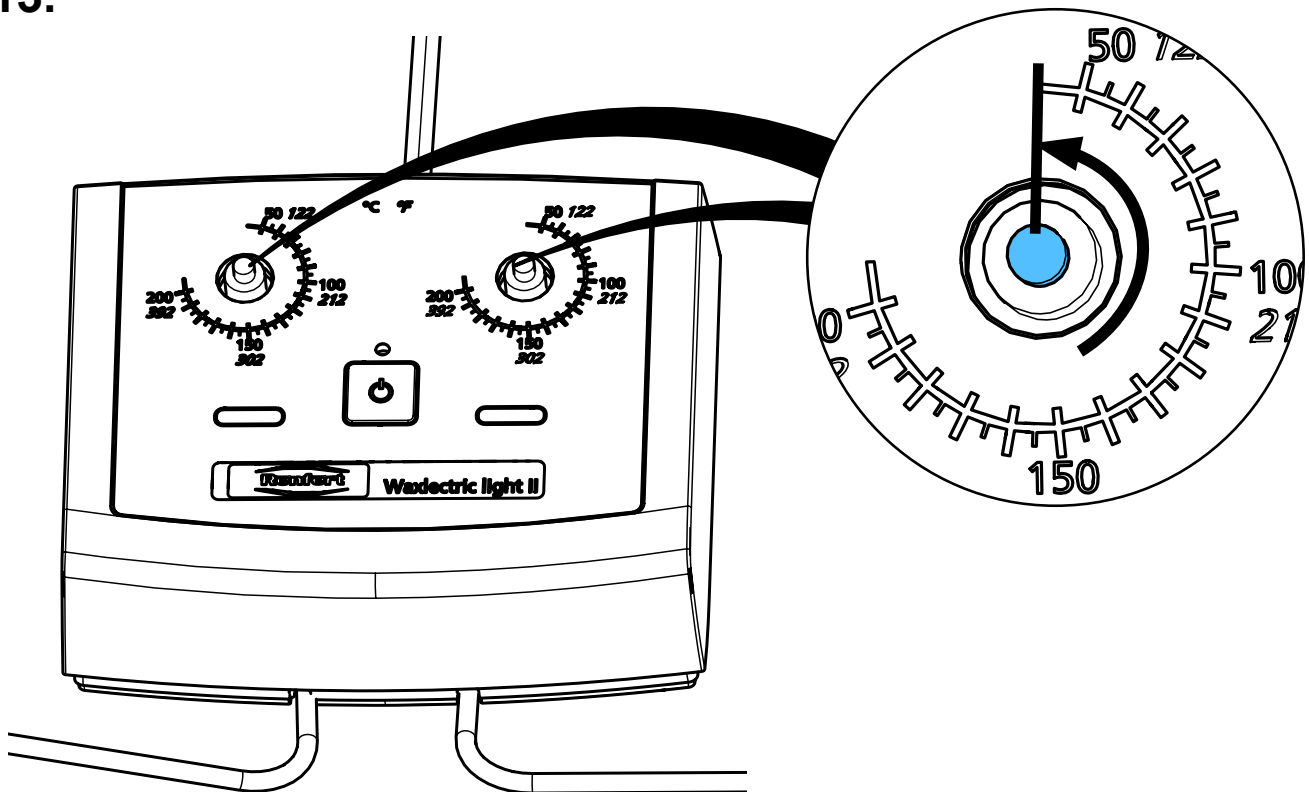


14.

i

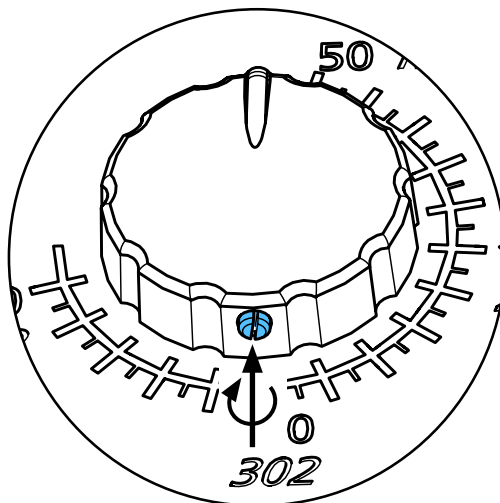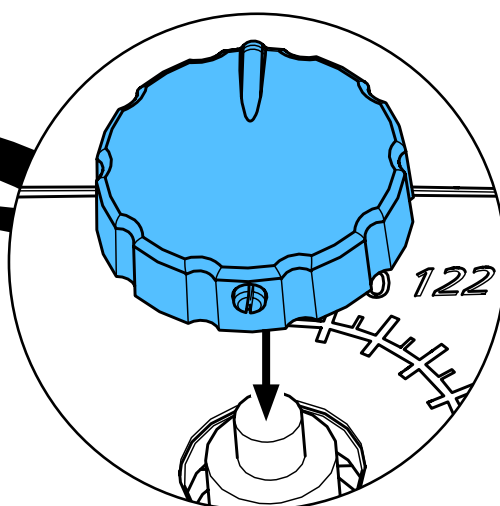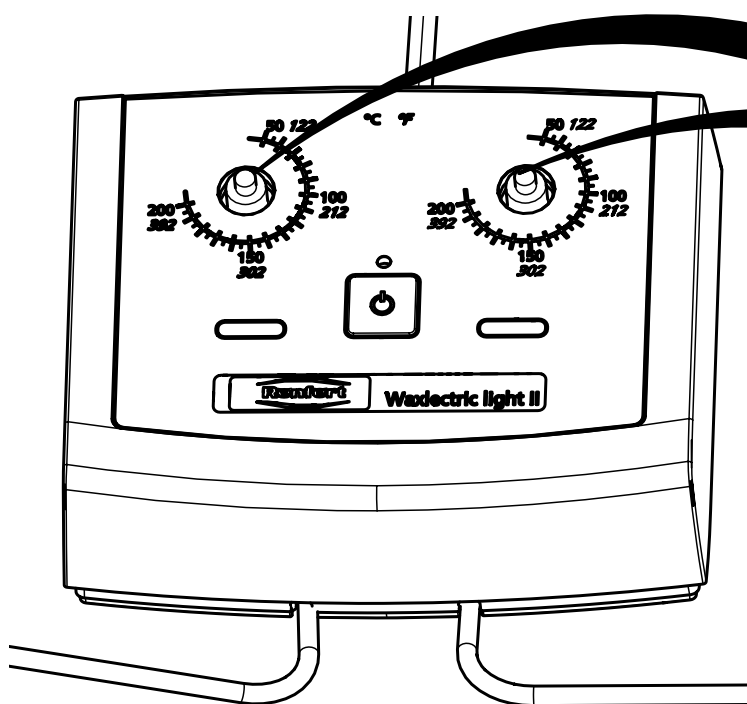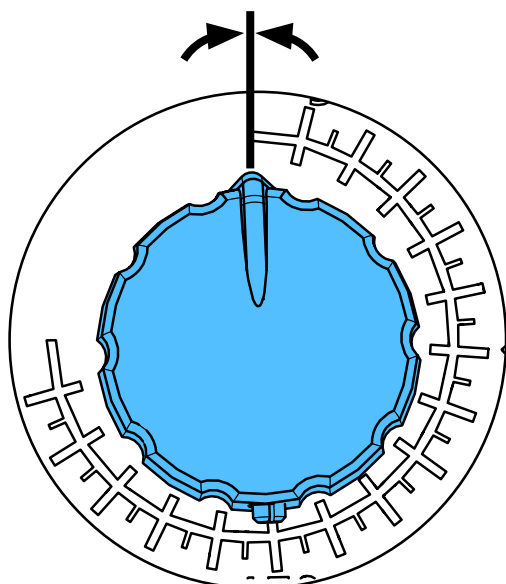

15.

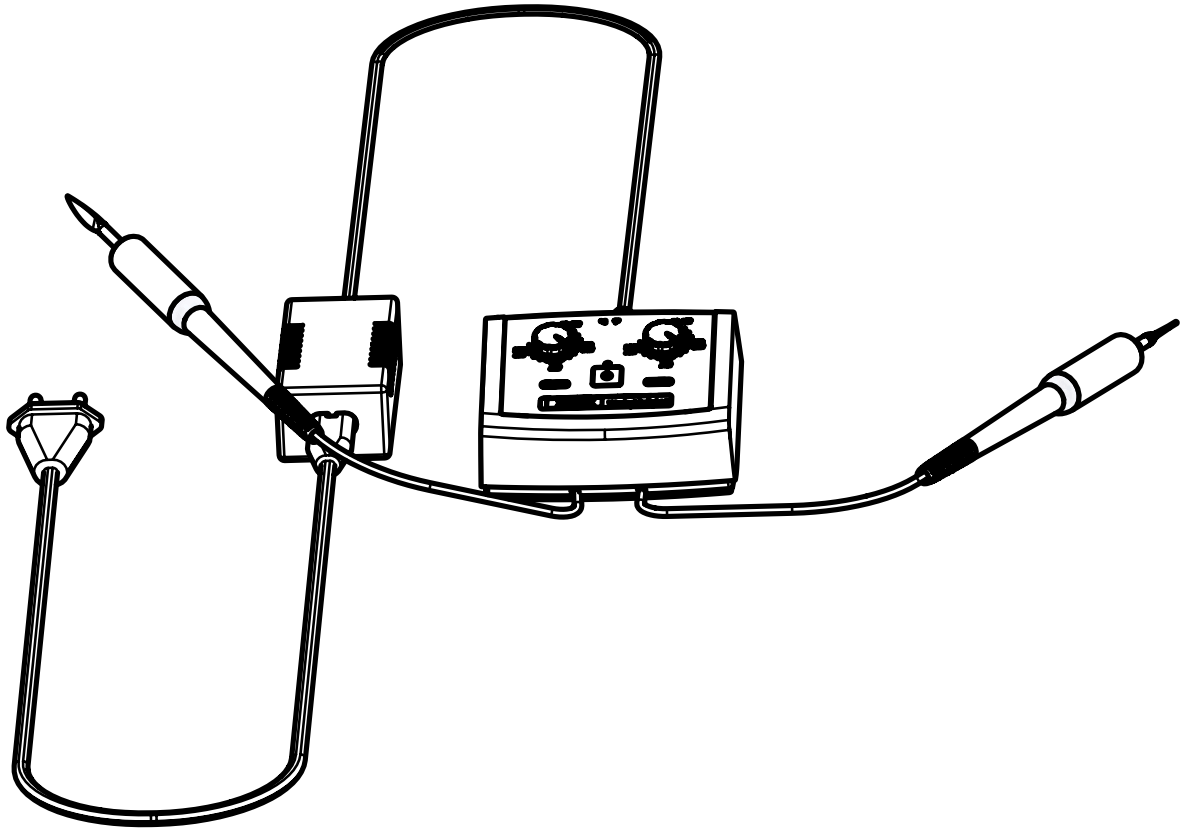






making work easy


Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу

90003 5953


Bedienfolie Waxelectric light II • Operating foil Waxelectric light II • Feuille de
comande Waxelectric light II • Lamina di comando Waxelectric light II • Lámina de
mando Waxelectric light II • Плёнка для панели управления Waxelectric light II

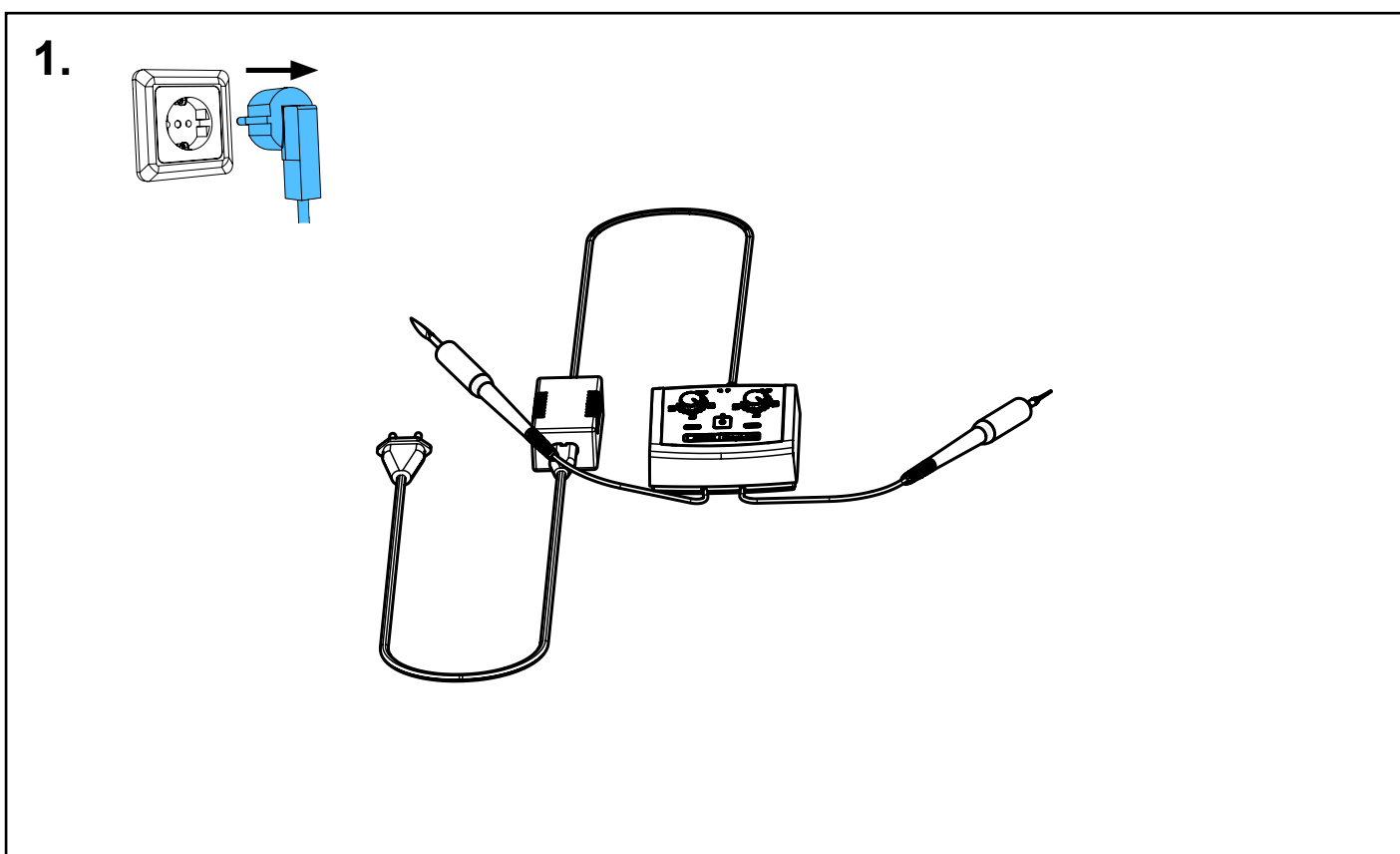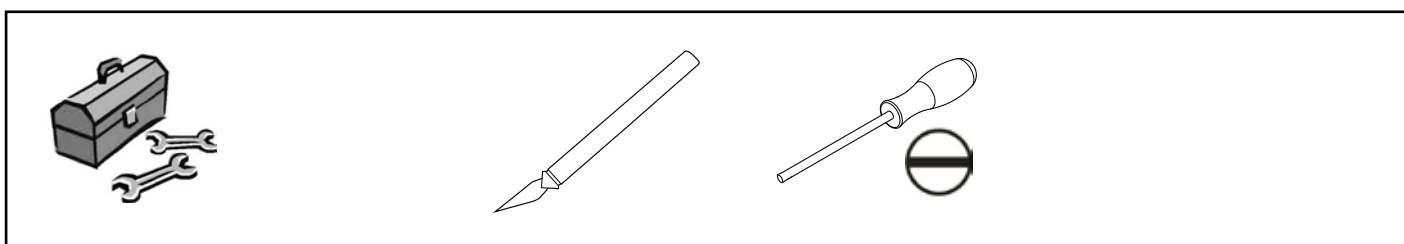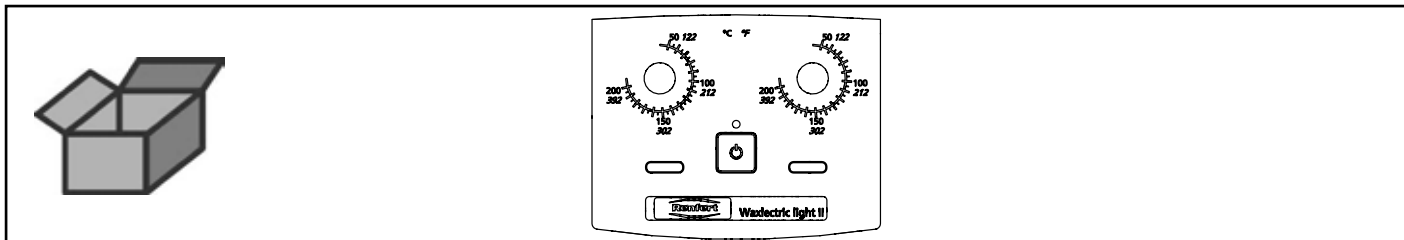


2.

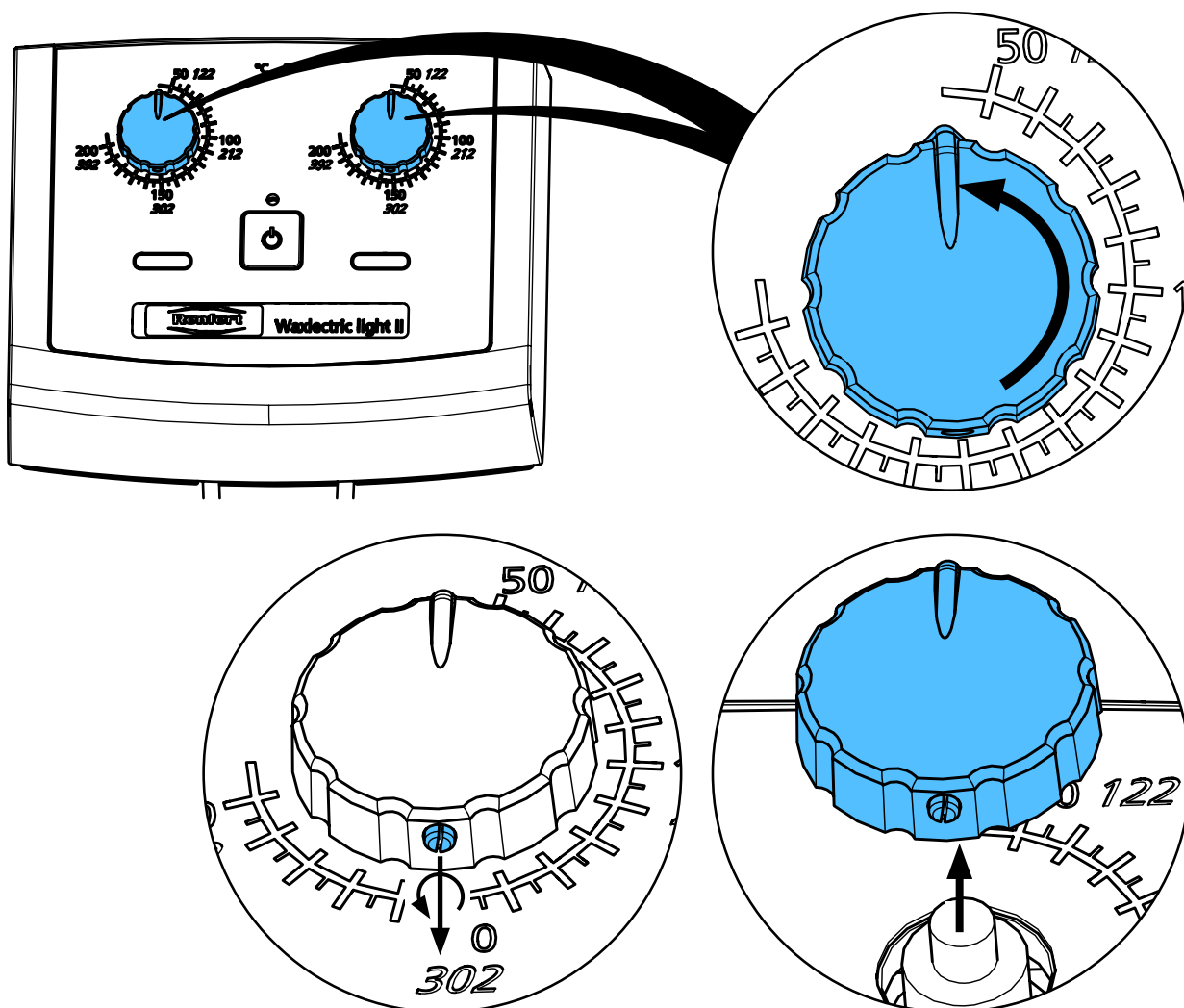


3.

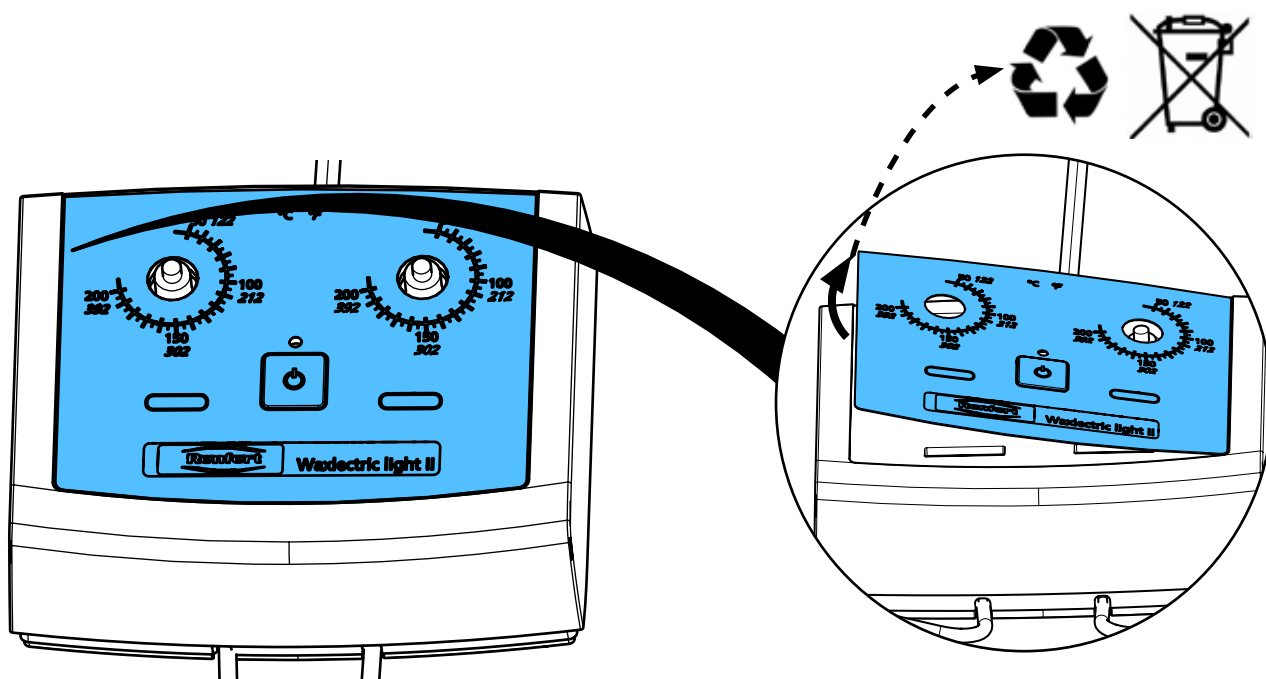


4.

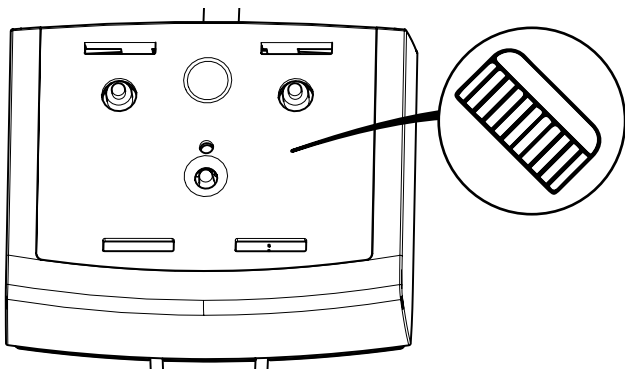


5.

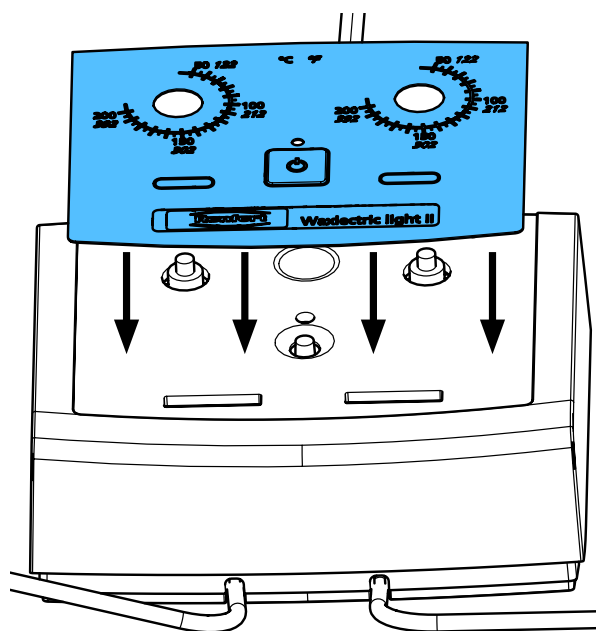

6.

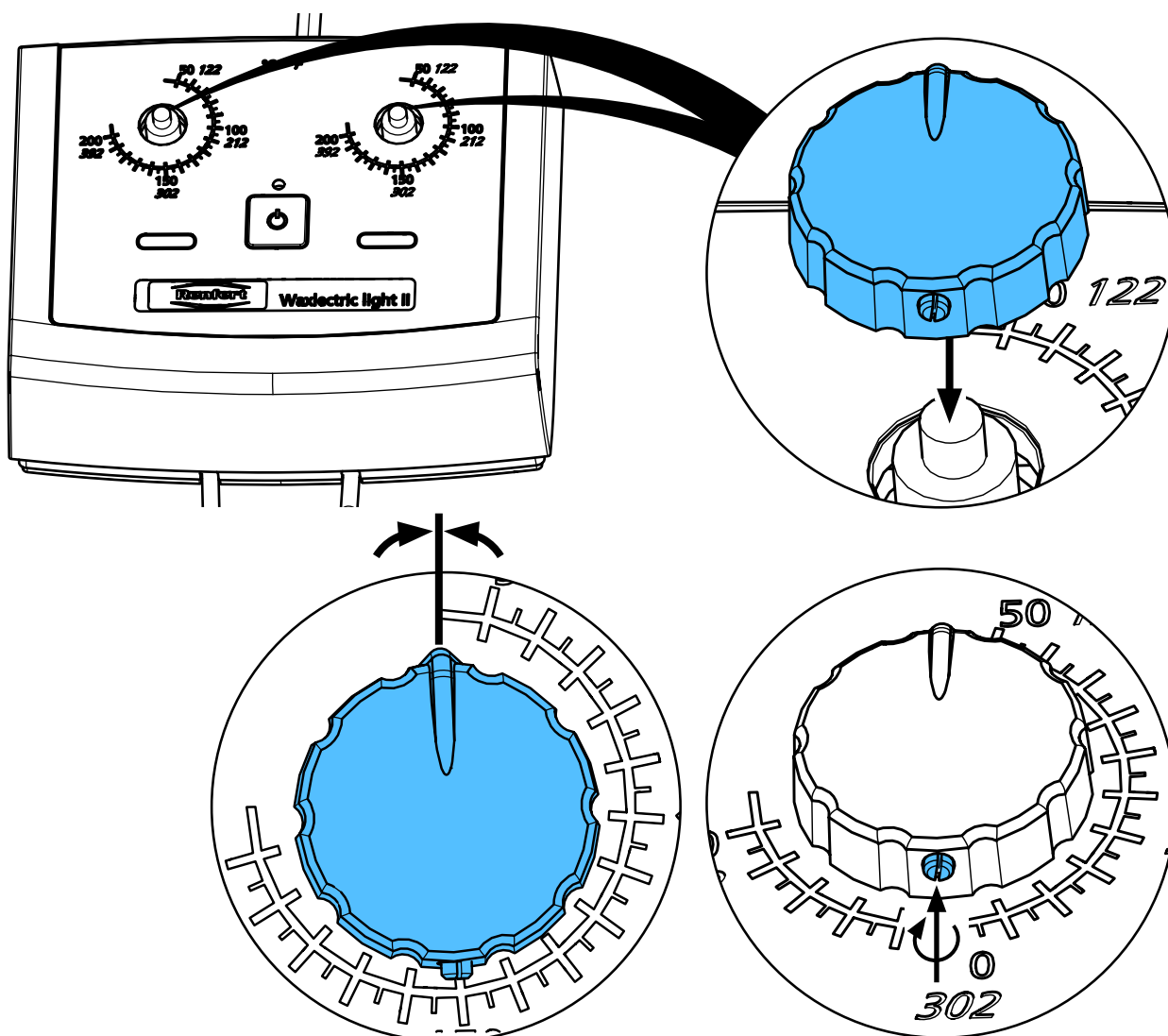

7.

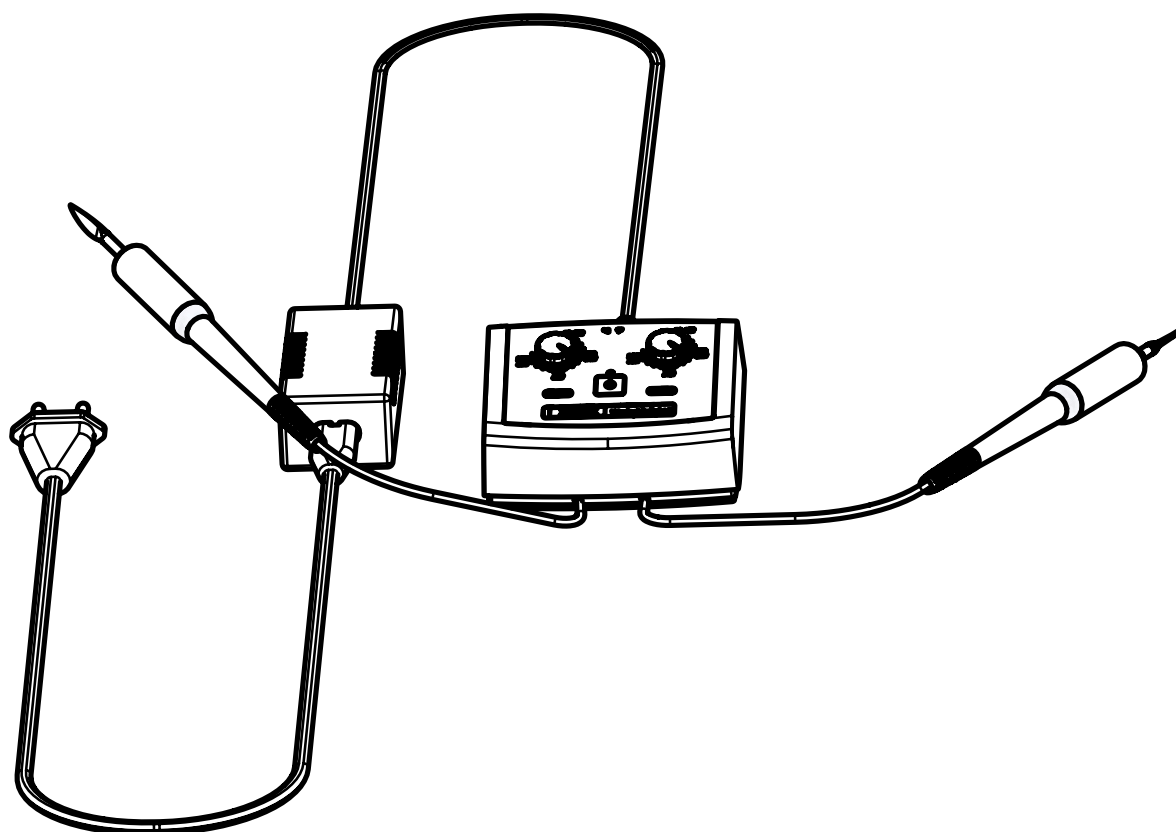

**Hochaktuell und ausführlich auf ...
Up to date and in detail at ...
Actualisé et détaillé sous ...
Aggiornato e dettagliato su ...
La máxima actualidad y detalle en ...
Актуально и подробно на ...**

www.renfert.com

Renfert GmbH • Untere Gießwiesen 2 • 78247 Hilzingen/Germany
Tel.: +49 7731 82 08-0 • Fax: +49 7731 82 08-70
www.renfert.com • info@renfert.com

Renfert USA • 3718 Illinois Avenue • St. Charles IL 60174/USA
Tel.: +1 6307 62 18 03 • Fax: +1 6307 62 97 87
www.renfert.com • info@renfertusa.com
USA: Free call 800 336 7422